

**Renfert**

making work easy



# BASIC master

Service Instructions

Made in Germany

21-6912 16082019

# Contents

## 90002 1428

Lampe kpl. • Lamp complete • Lampe complète • Lampada completa  
Lámpara completa • Светильник полный

## 90003 6659

Lampenabdeckung • Cover for lamp • Protection pour lampe • Protezione per la lampada  
Cobertura de lámpara • Крышка для светильника

## 90003 6121 /...6122 /...5938 /...2769 /...6070 /...2767 /...2768

Netzkabel • Power cord • Cordon d'alimentation • Cavo di rete • Cable de red  
Сетевой кабель

## 2958 0050 / 2958 0250

Zusatztank • Additional tank • Silo supplémentaire • Serbatoio supplementare  
Depósito adicional • Дополнительные резервуары

## 90003 5801

Ersatztankdeckel • Lid f.replacement tank • Couvercle p. silo rec. • Coperchio p.dep. ric.  
Тара p.depósito de rec. • Крышка запасного бачка

## 90002 1431

Filterpatrone • Filter cartridge • Cartouche de filtres • Cartuccia dei filtri con raccordo  
Cartucho de filtración • Патрон фильтра

## 90003 3035

O-Ring DIN 3771 82x4 NBR 55 • O-ring DIN 3771 82x4 NBR 55 • Joint thorique DIN 3771 82x4  
NBR 55 • Guarnizione DIN 3771 82x4 NBR 55 • Anillo en O DIN 3771 82x4 NBR 55  
Кольцо круглого сечения DIN 3771, 82x4 NBR 55

## 92947 0002

Kleinteileset Basic • Hardware Basic • Petit materiel d.montage Basic • Kit di mi-nuteria Basic  
Juego de piezas pequeñas Basic • Комплект мелких деталей

## 90003 6330

Bodengitter • Base grid • Grille de fond • Griglia di fondo • Rejilla de fondo • Решетка для дна

## 90003 4751

Handlochmanschetten • Hand sleeves • Manchettes • Manicotto • Puños • Манжета серая

## 90003 6339

Handstulpenring • Ring for hand sleeves • Anneau pour gants • Ghiera per mani-cotto  
Anillo para puños • Кольцо манжеты

## 90002 1435

Glasscheibe beschichtet • Coated glass screen • Vitre revêtu • Vetro rivestito  
Cristal con revestimiento • Запасное стекло с покрытием

## 90003 6311

Glasscheibe unbeschichtet • Uncoated glass screen • Vitre non revêtu  
Vetro sen-za rivestimento • Cristal sin revestimiento • Запасное стекло без покрытия

## 92947 0001

Kantenschutzprofil • Edge protection • Bordure de protection • Profilo protettivo  
Tira de protección • Краевой защитный профиль

## 90002 0921

Fußschalter pneumatisch • Pneumatic foot switch • Commande à pieds pneuma-tique  
Comando a pedale pneumatico • Pedal interruptor neumático • Педаль пневматическая

## 90003 6598

EingangsfILTER • Inlet filter • Filtre d'entrée • Filtro d'entrata • Filtro de entrada входной фильтр

**90003 4104**

Belüftungsklappe • Ventilating flap • Clapet de ventilation • Copertura per foro di aerazione  
Charapaleta de ventilación • Вентиляционный клапан

**92900 0050 / 92900 0500**

IT Strahlschlauch inkl. Handstück • IT blasting hose incl. handpiece • IT tuyau de sablage avec  
pièce á main • IT tubo incl. manipolo • IT tubo incl. manipolo • Струйный шланг IT

**90003 6353**

Dosiernippel 25-70 µm steckbar • Plug-in dosing nipple 25-70 µm • Raccord de dosage 25-70 µm  
enfichable • Nipplo di dosaggio 25-70 µm innestabile • Boquilla de dosificación 25-70 µm en-  
chufable • дозирующего сопла 25 мкм - 70 мкм

**90003 6354**

Dosiernippel 70-250µm steckbar • Plug-in dosing nipple 70-250µm • Raccord de dosage 70-250  
µm enfichable • Nipplo di dosaggio 70-250 µm innestabile • Boquilla de dosificación 70-250 µm  
enchufable • дозирующего сопла 70 мкм - 250 мкм

**92947 0003**

Seitenteile Basic • Side panels Basic • Pièces latérales Basic • Parti laterali Basic  
Piezas laterales Basic • боков́а я часть Basic

**90003 8600**

Durchstrahlschutz • Blast-through protection • Protection contre perforation • Pro-tezione contro  
perforazione • Protección contra perforación • Защита от сквозного прохождения струи песка

**90015 0864**

PU-Schlauch 6 x 3,9 mm • PU-hose 6 x 3.9 mm • Tuyau flexible en polyuréthane 6 x 3,9 mm  
Tubo flessibile in poliuretano 6 x 3,9 mm • Tubo flexible en poliuretano 6 x 3,9 mm  
PU-шланга 6 x 3,9 мм

**90003 5800**

Ersatztankmantel Basic • Replacement tank case Basic • Boítier du silo Basic  
Involucro deposito Basic • Envoltura depósito Basic • Корпус запасного бачка Basic

**90003 6386**

Mischkammer • Mixing chamber • Chambre de malaxage • Camera di miscela  
Camara de mezcla • Смесительная камера

**2 1119**

Manometer mit Verschraubung • Manometer with screw fitting • Manomètre avec raccord  
Manometro con filettatura • Manómetro con atornilladura • Манометр с резьбовым  
соединением

**90002 1551**

Tankwahlschalter 2-fach • 2-fold tank selector switch • Commutateur de silo 2-voies • Selettore  
di serbatoio 2-vie • Conmutador de depósito doble • Переключатель на 2 бачка

**90003 6413**

Druckregler • Pressure regulator • Régulateur de pression • Regolatore di pressione  
Regulador de presión • Регулятор давление








Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу


making work easy

90002 1428


Lampe kpl. • Lamp complete • Lampe complète • Lampada completa  
Lámpara completa • Светильник полный

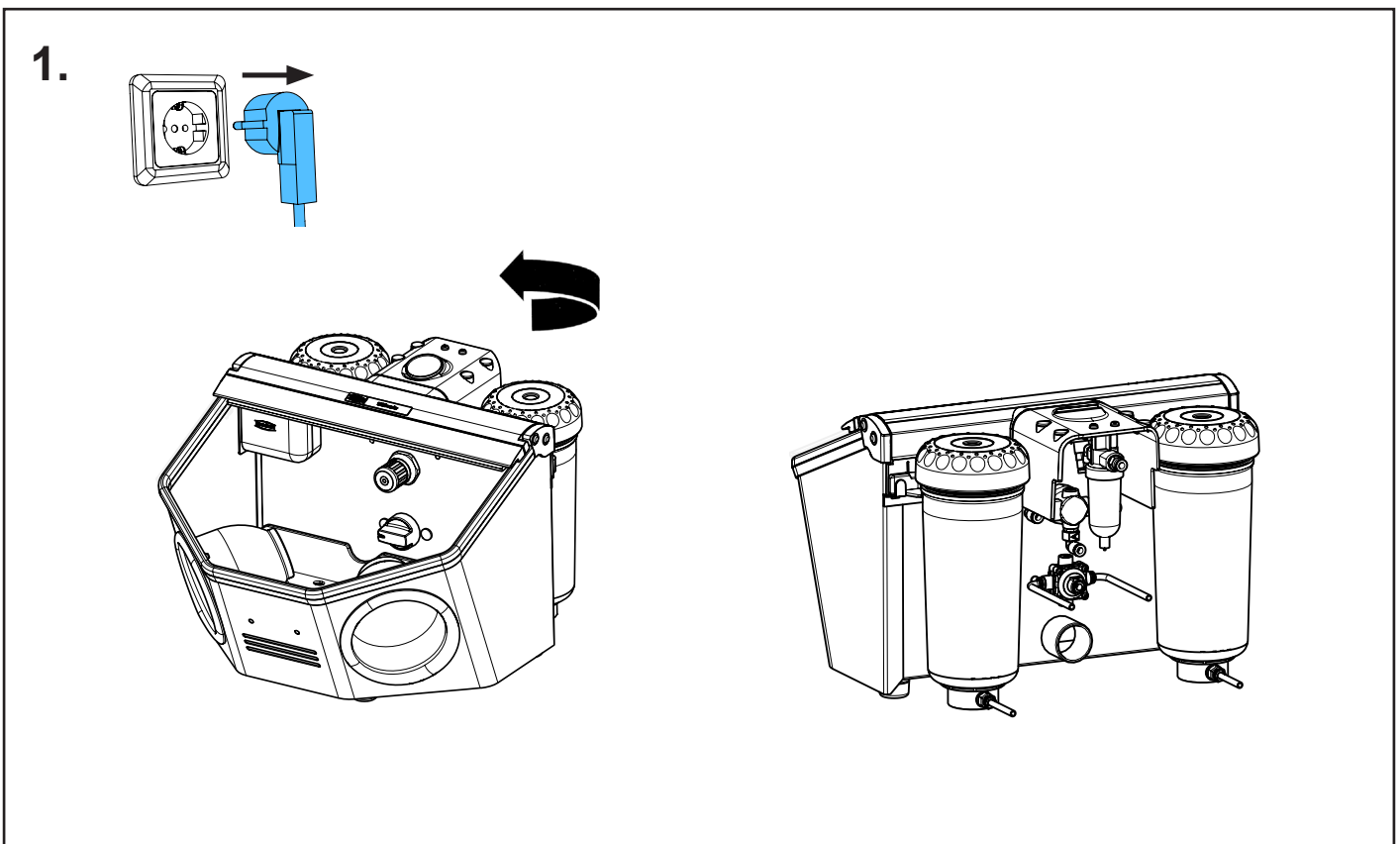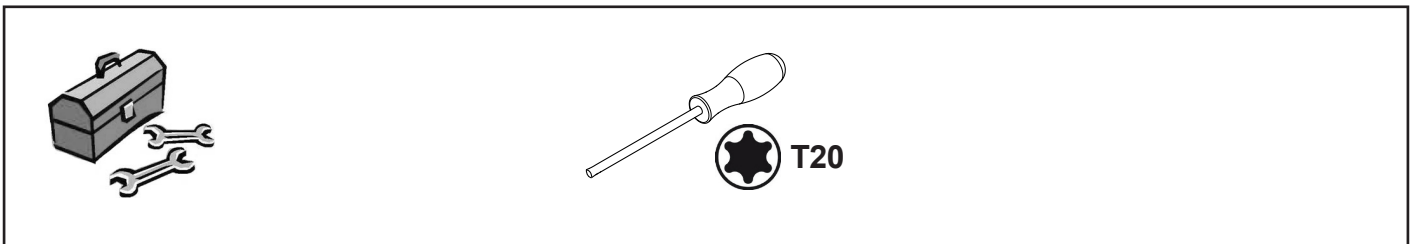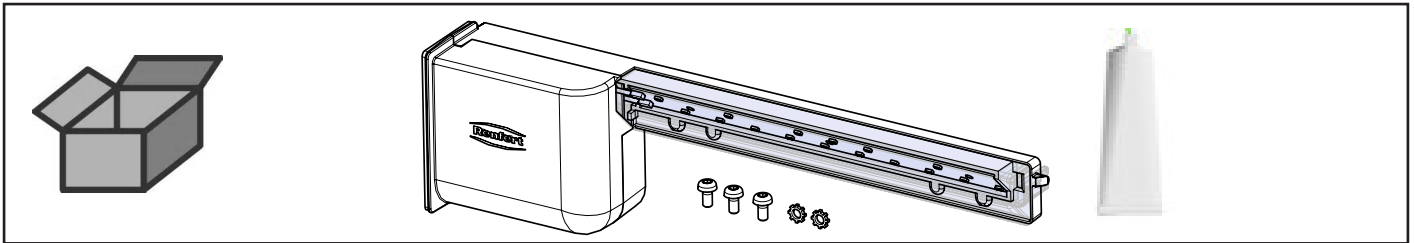


2.

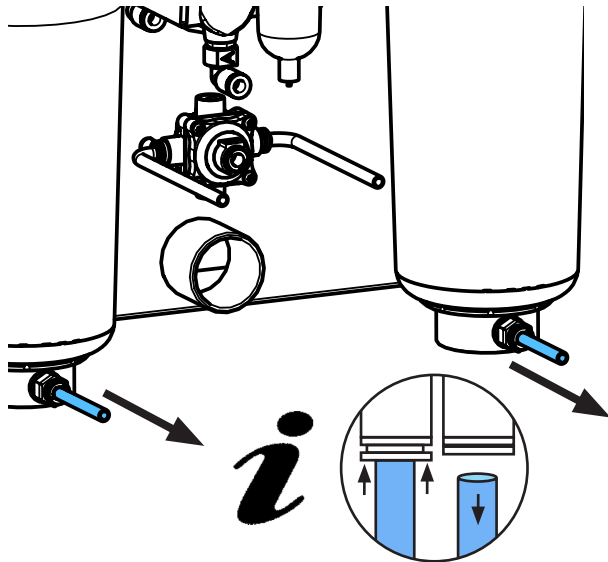


3.

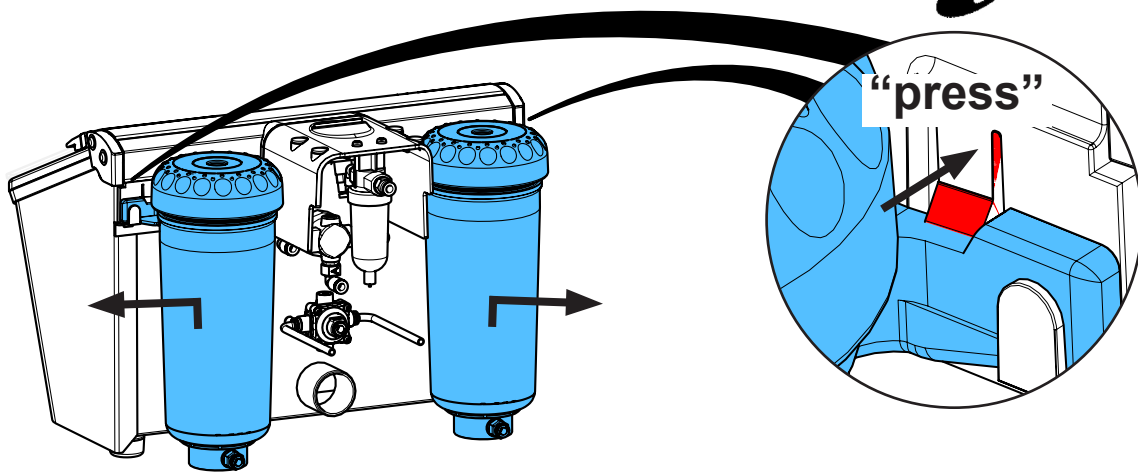


4.

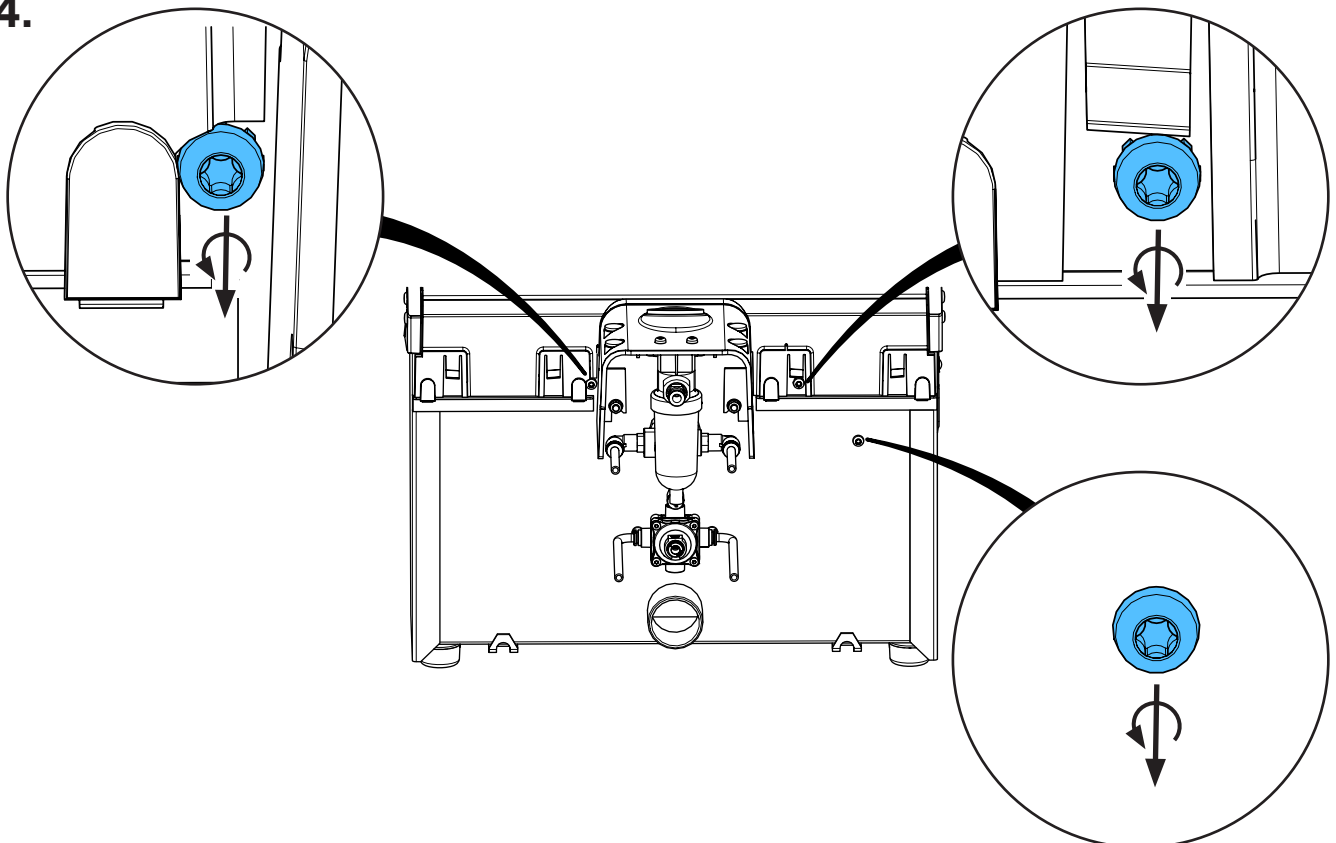


5.

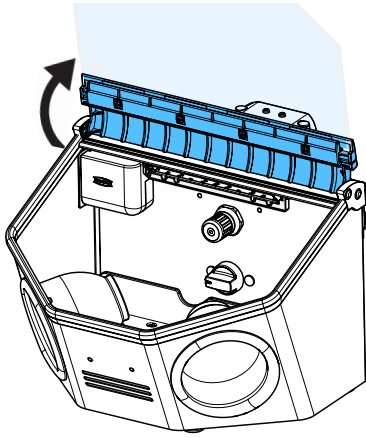


6.

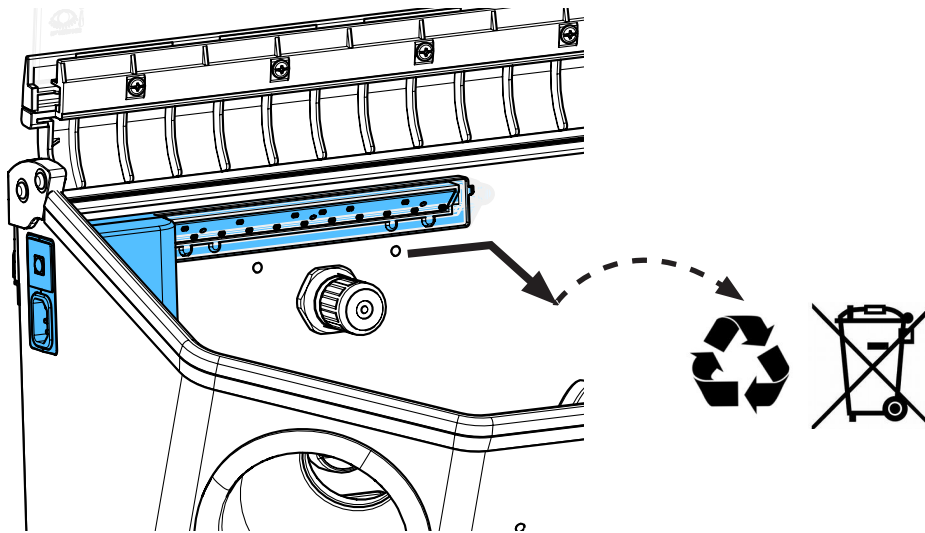


7.

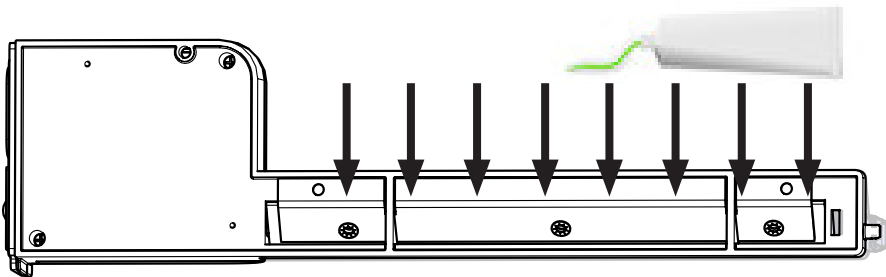

8.

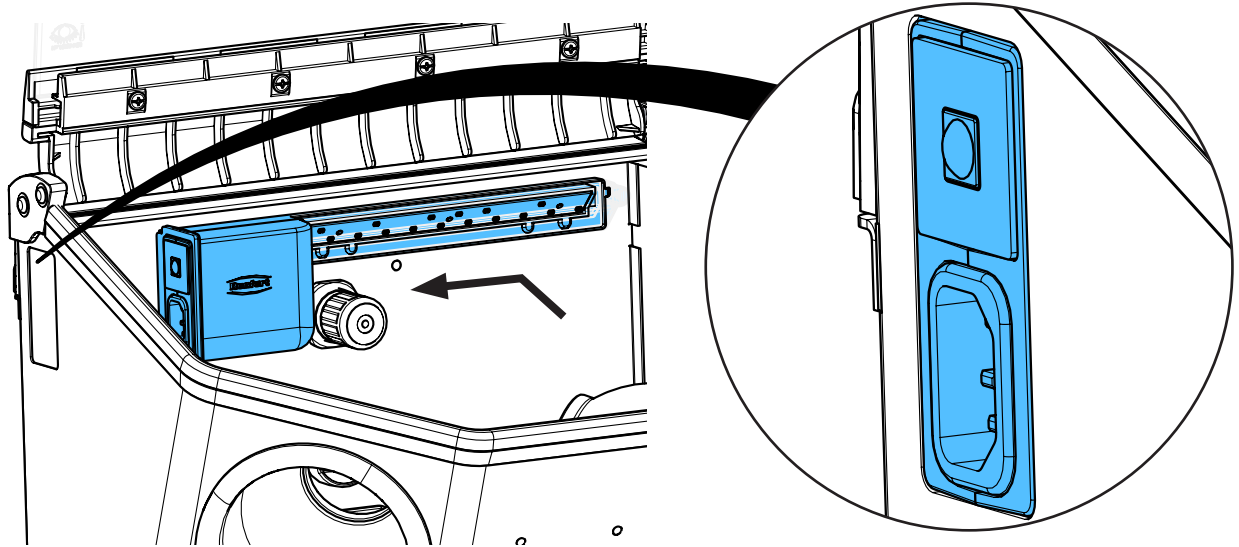


9.

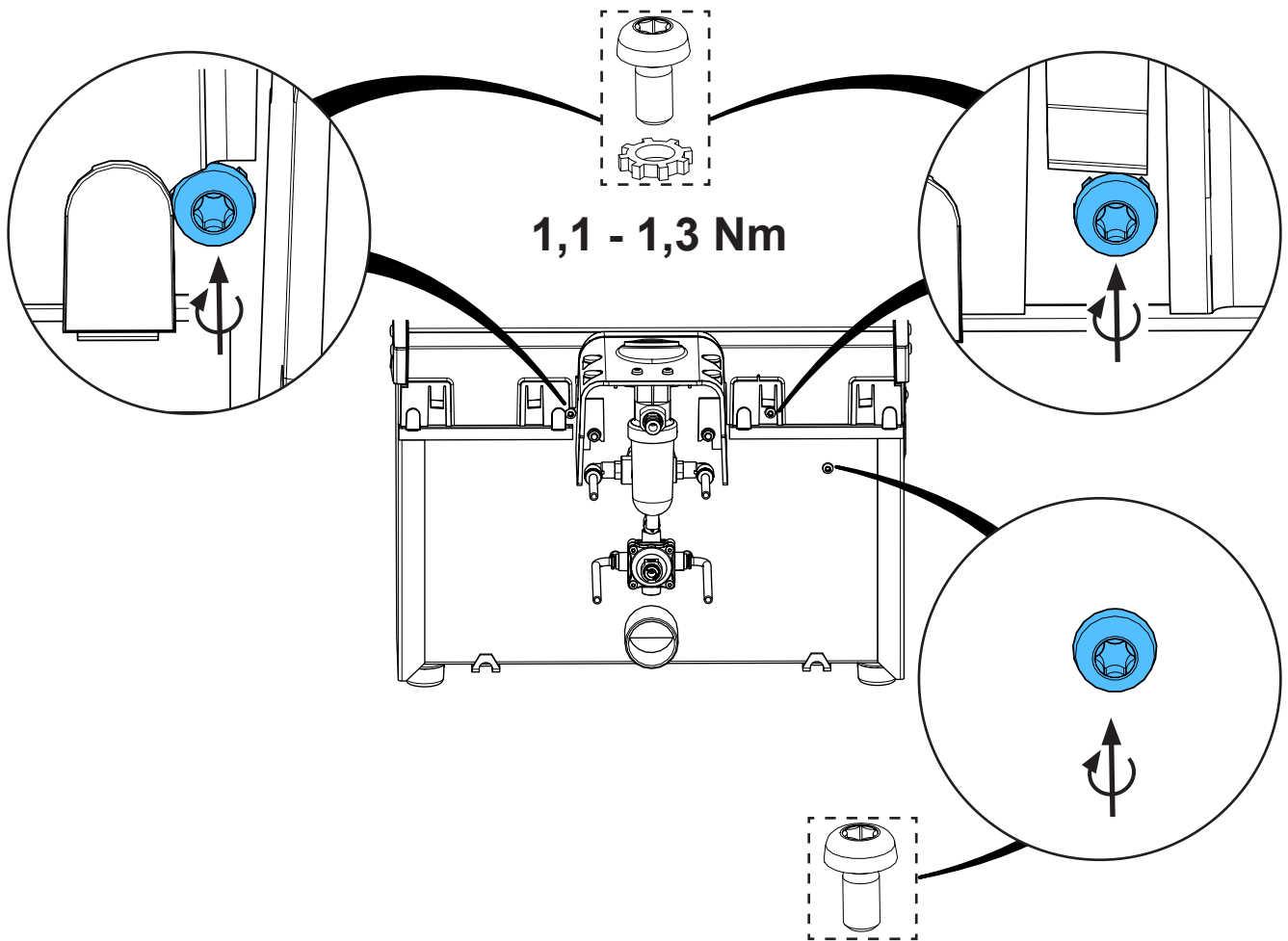






10.

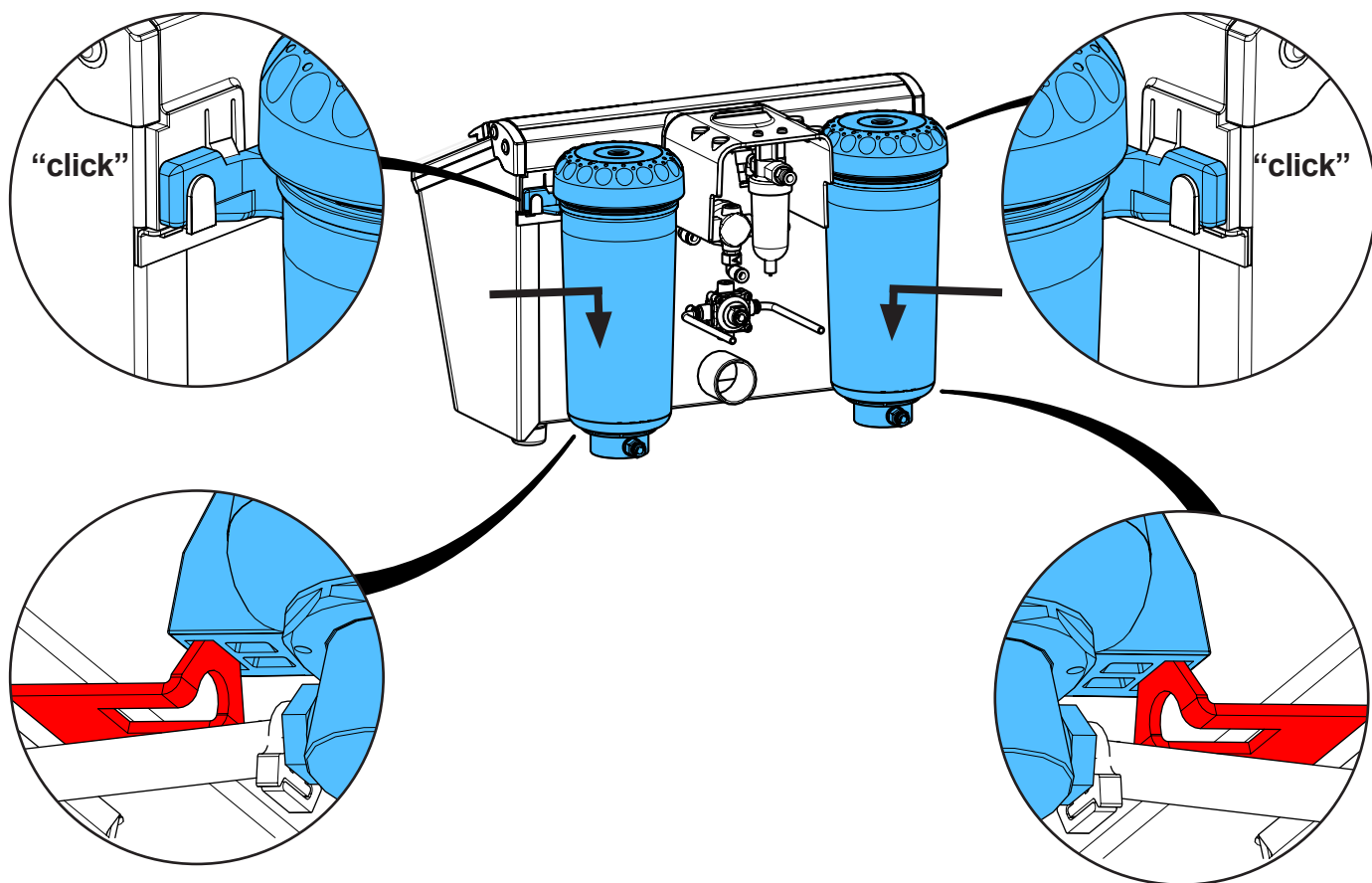

11.

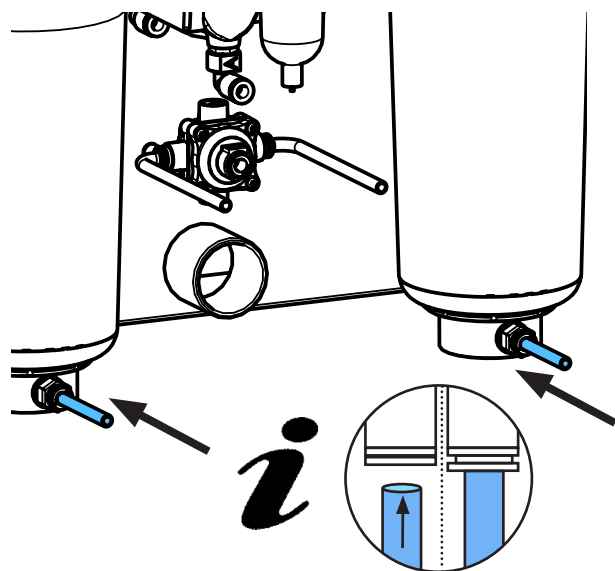

12.

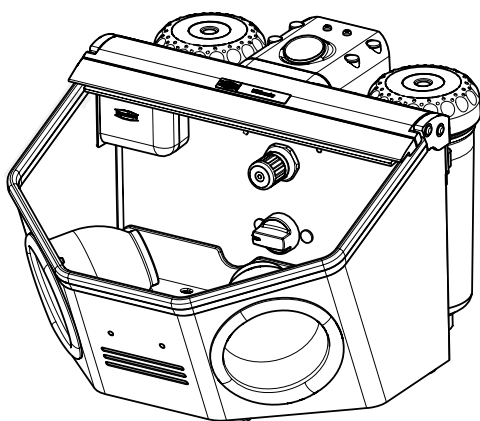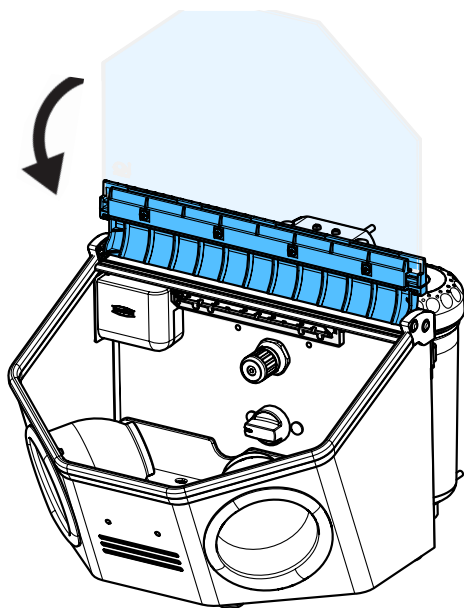





Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу



making work easy

90003 6659


Lampenabdeckung • Cover for lamp • Protection pour lampe • Protezione per la  
lampada Cobertura de lámpara • Крышка для светильника

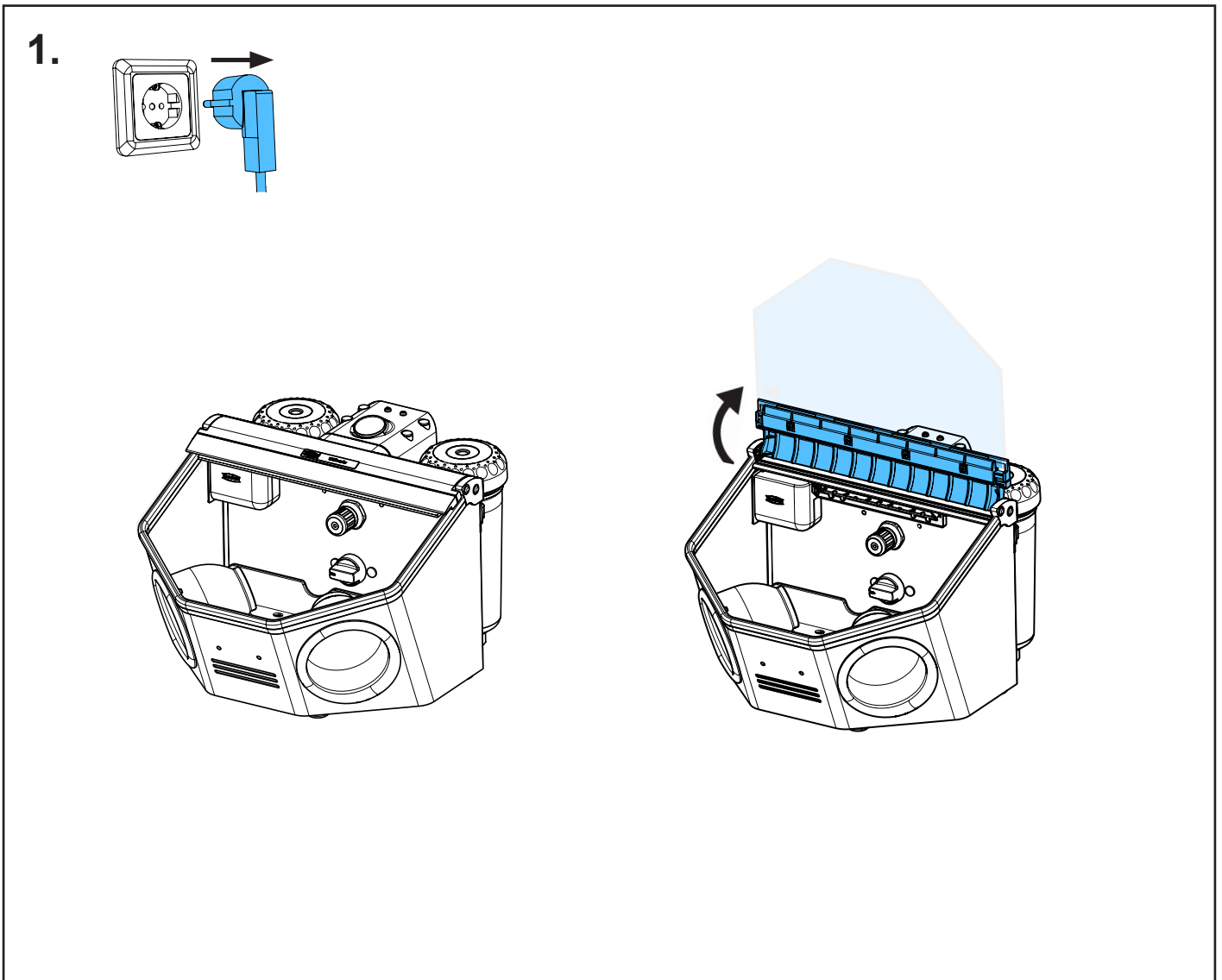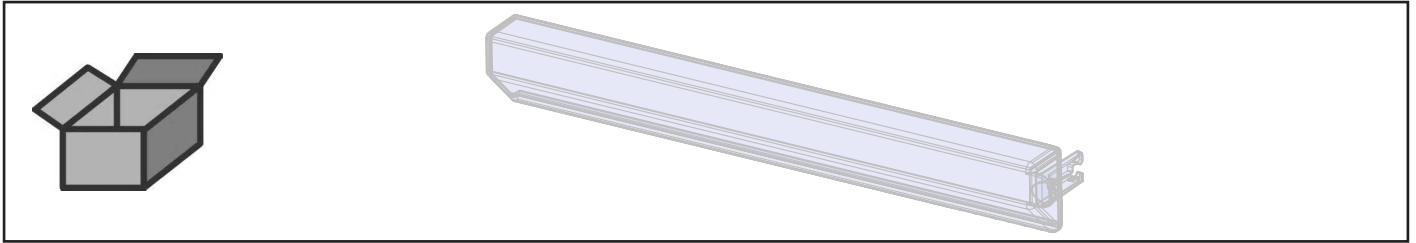


2.

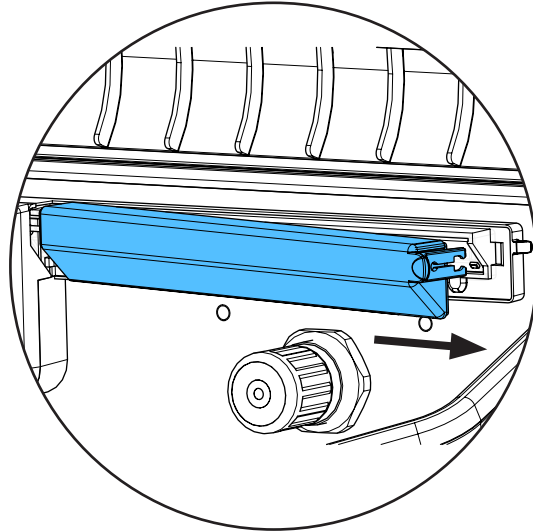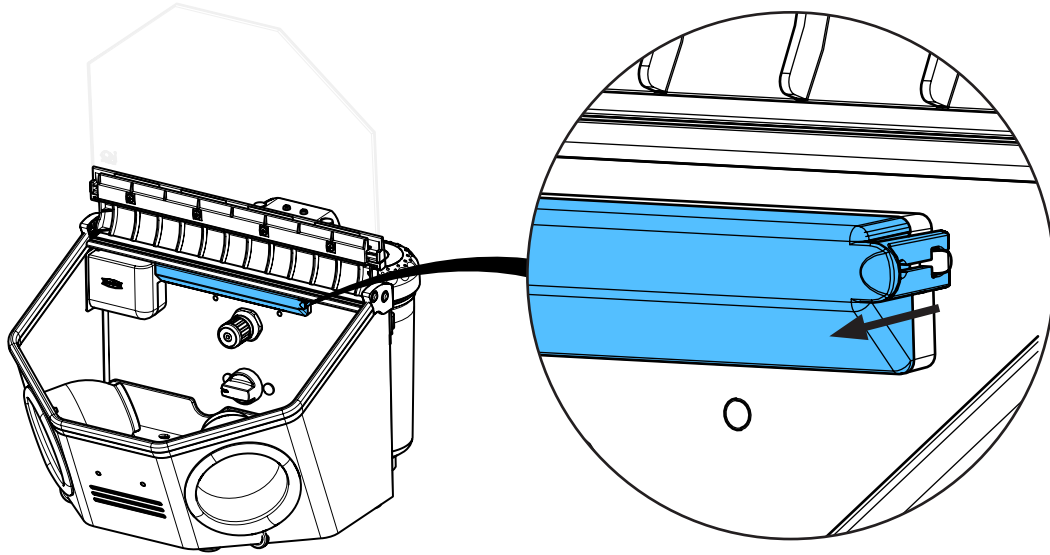

3.

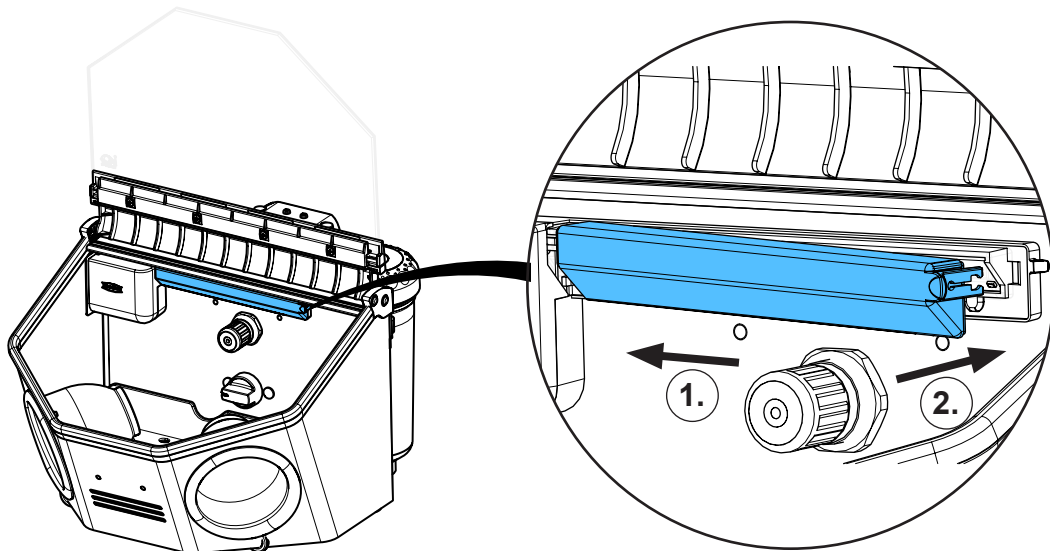

4.

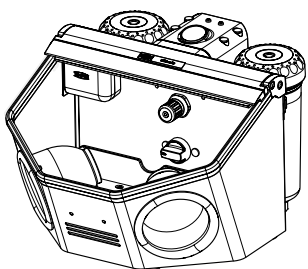






Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по монтажу

making work easy

**90003 6121/...6122/...5938/...2769/...6070/...2767/...2768**

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Netzkabel • Power cord • Cordon d'alimentation • Cavo di rete • Cable de red •  
Сетевой кабель

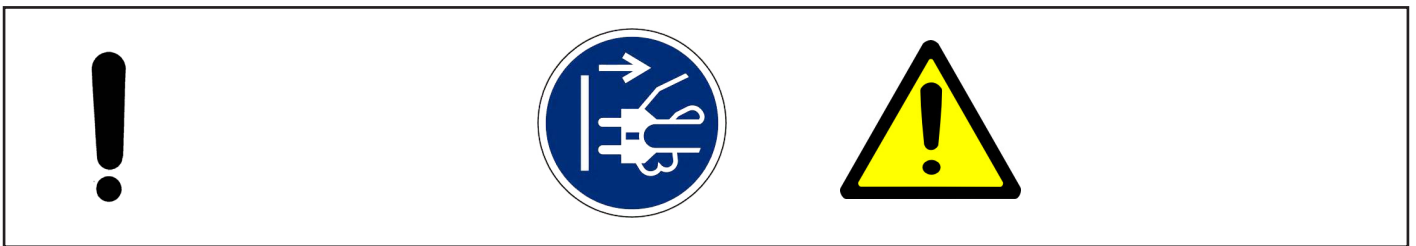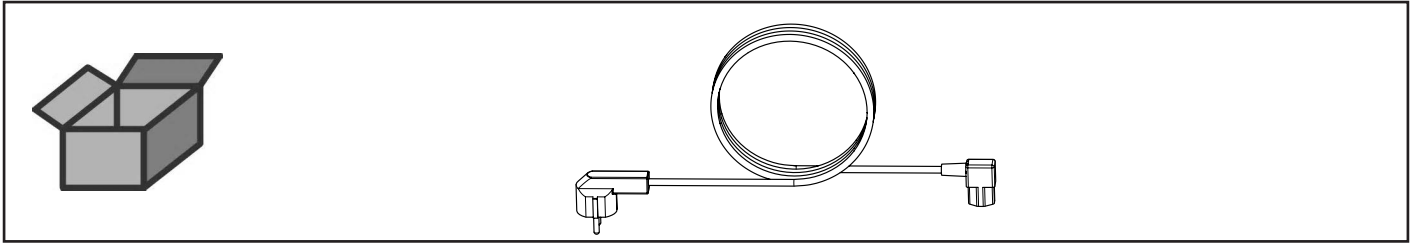







Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по монтажу


making work easy

**2958 0050 / 2958 0250**


Zusatztank • Additional tank • Silo supplémentaire • Serbatoio supplementare  
Depósito adicional • Дополнительные резервуары

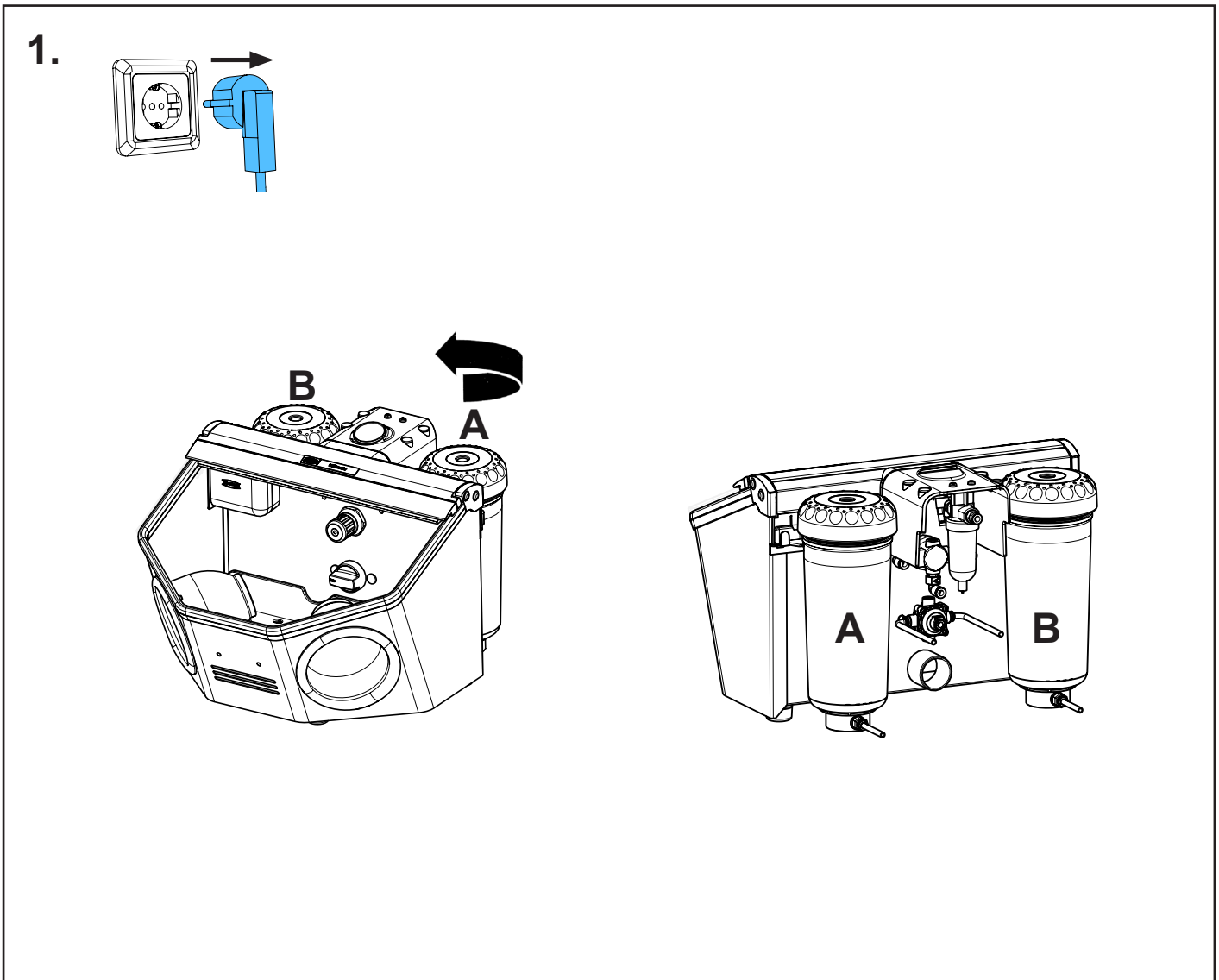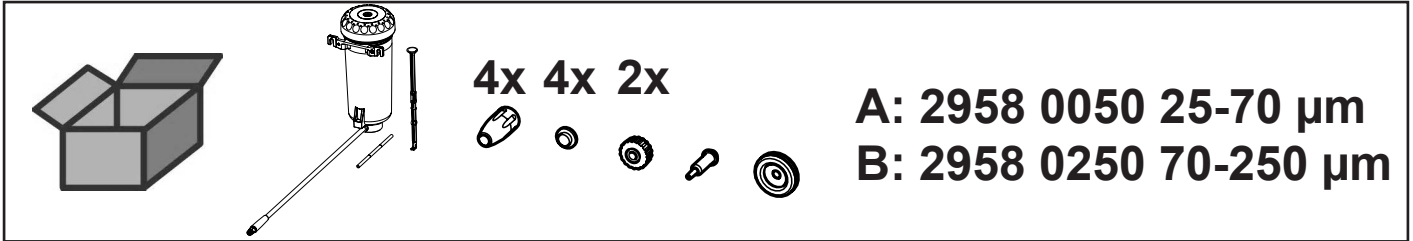


2.

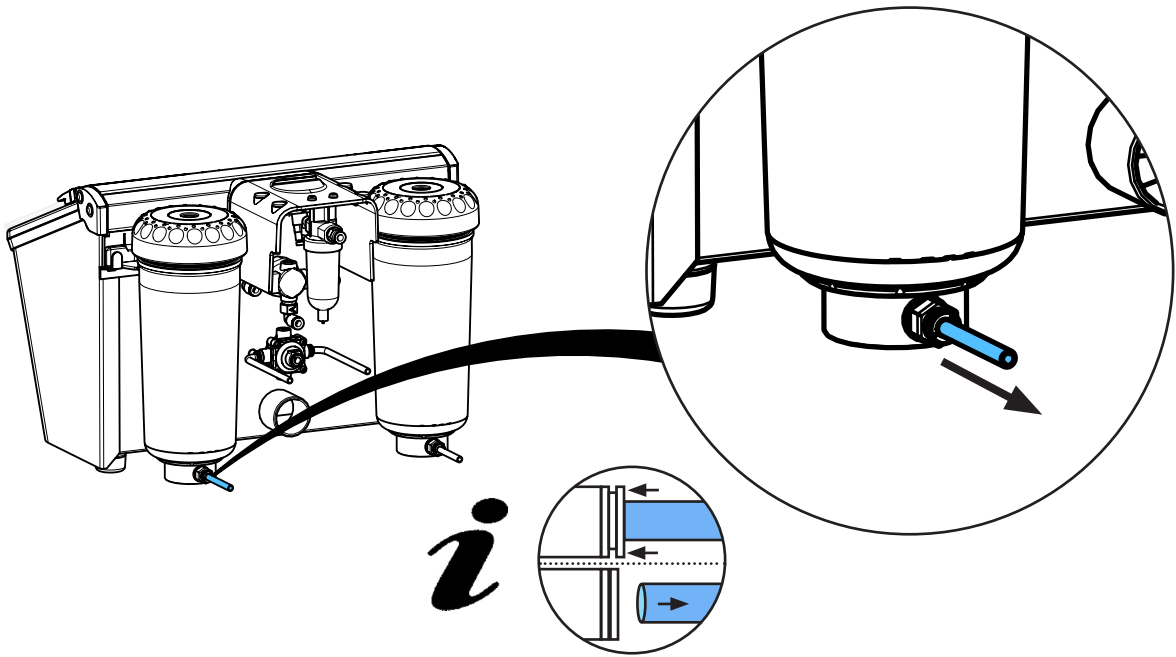

3.

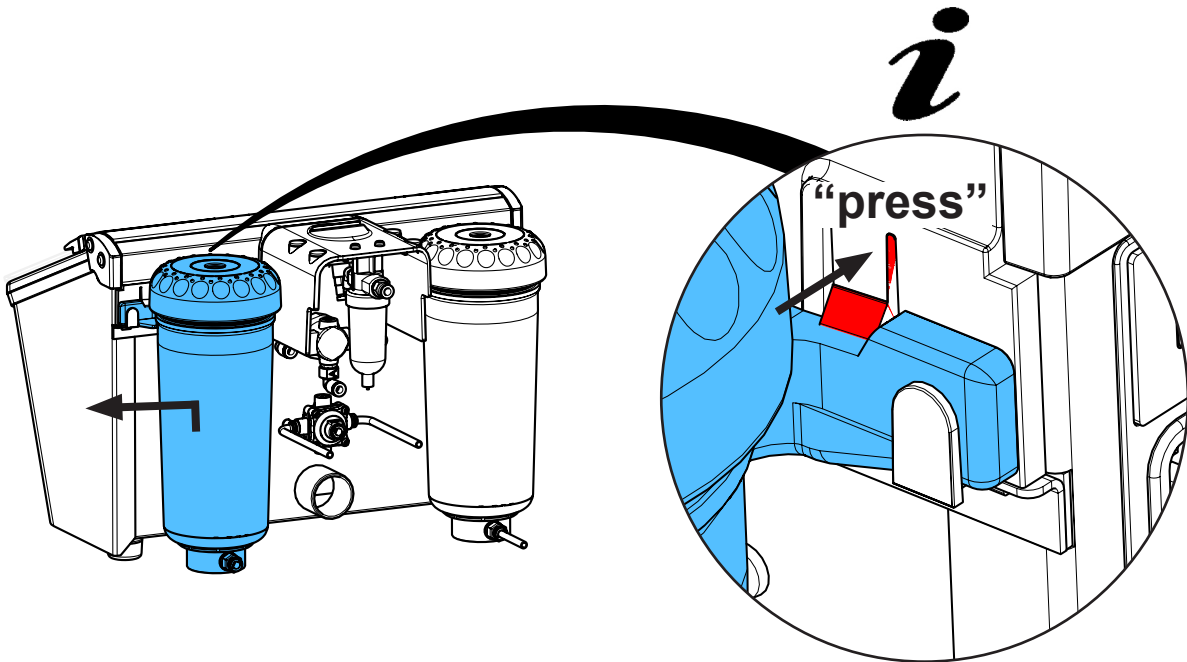


4.




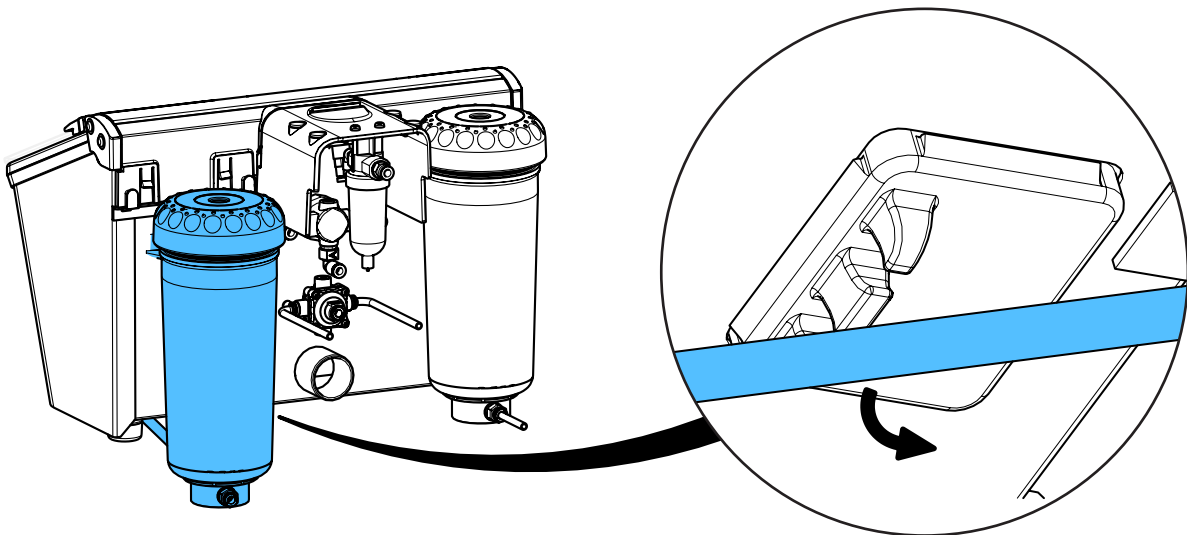




5.

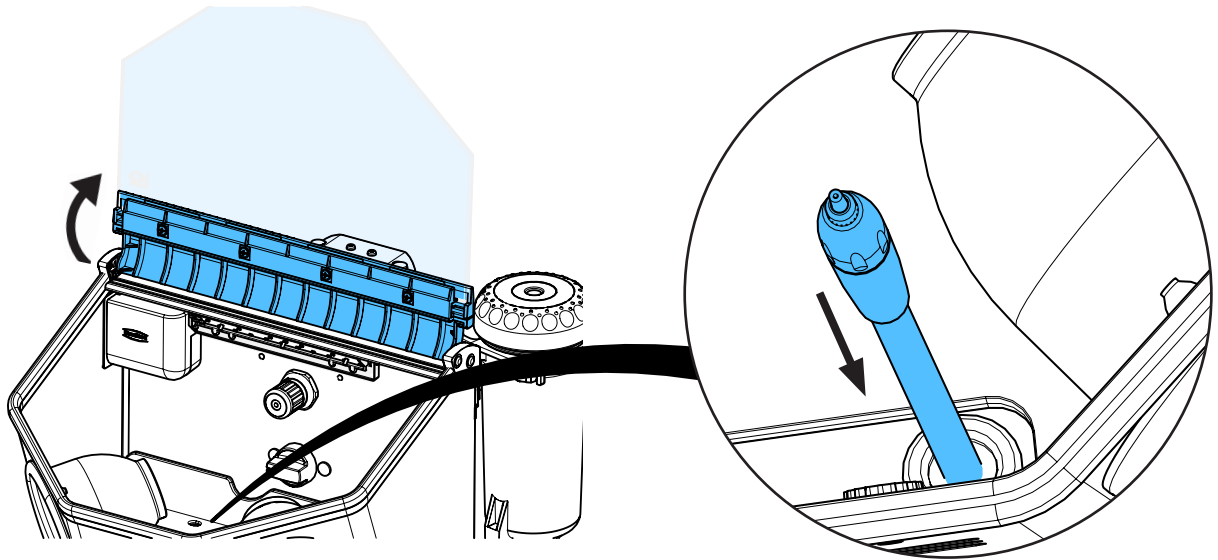


6.

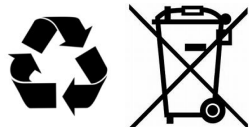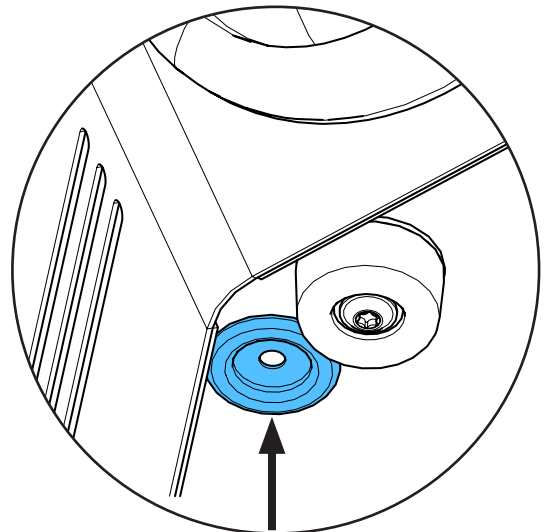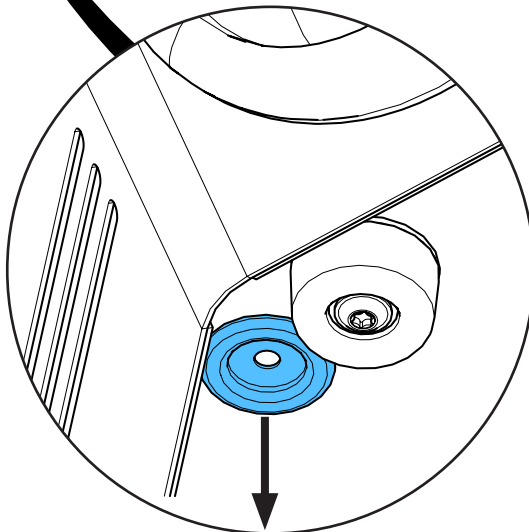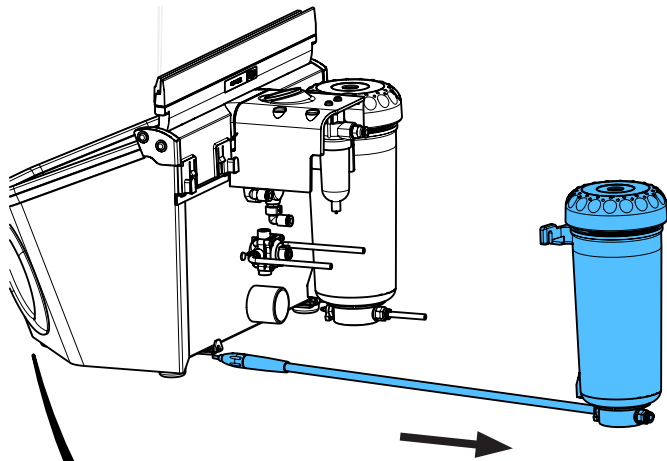


7.


yellow

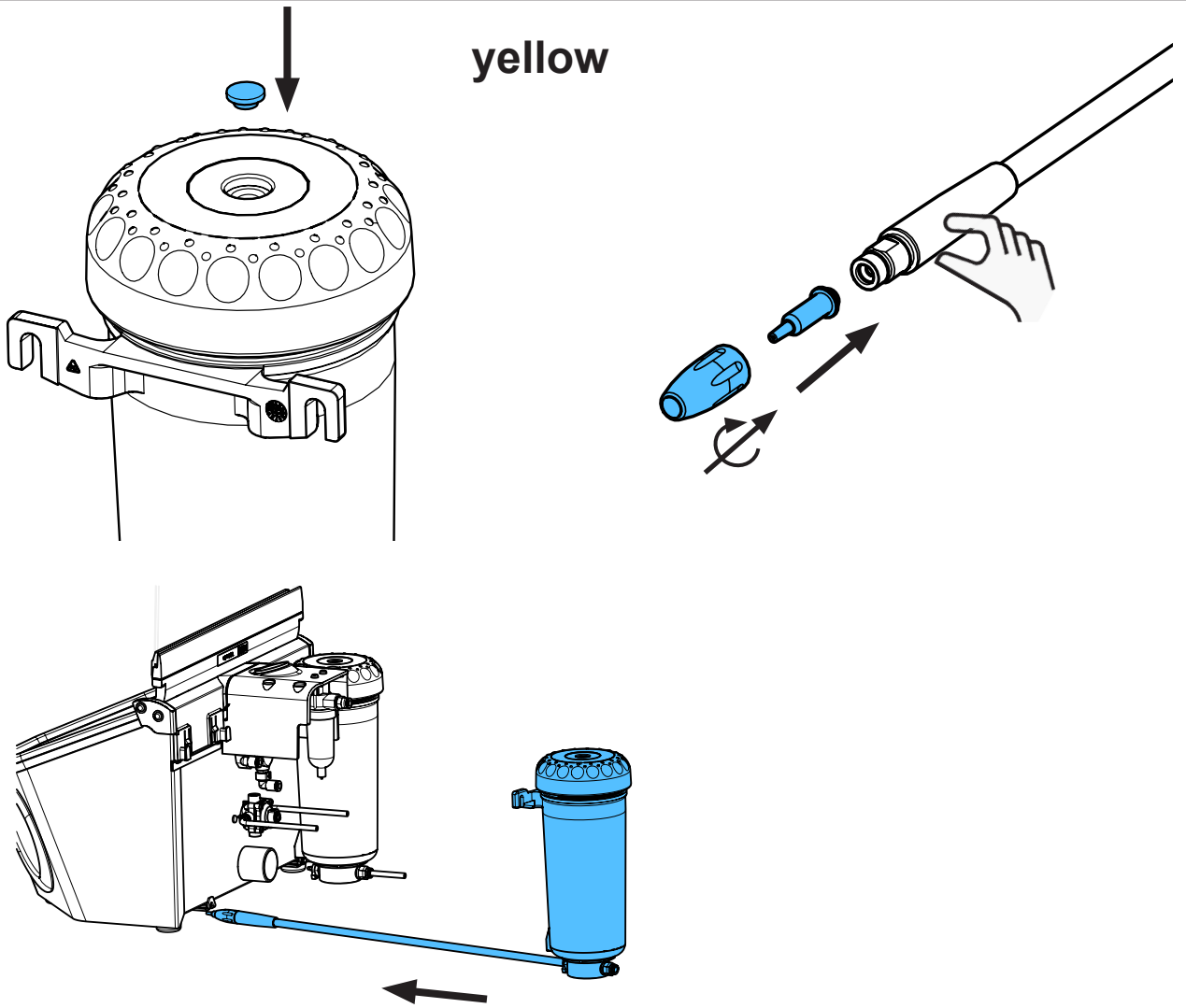


8.

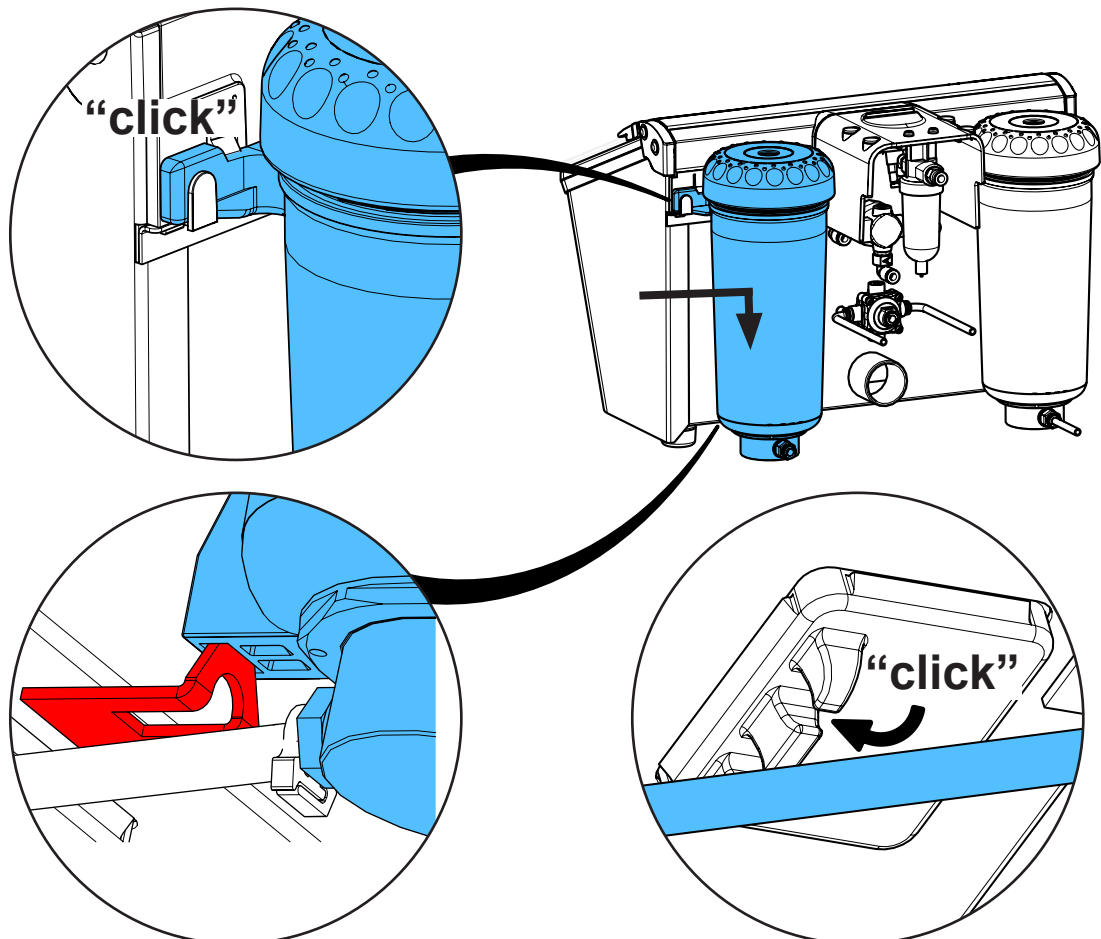


9.

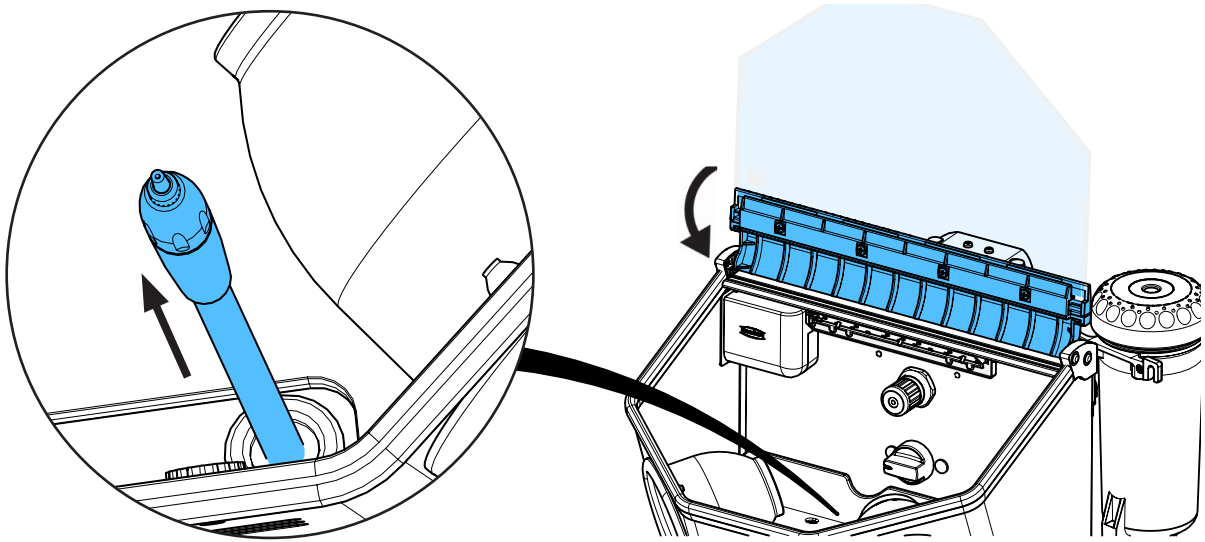


10.

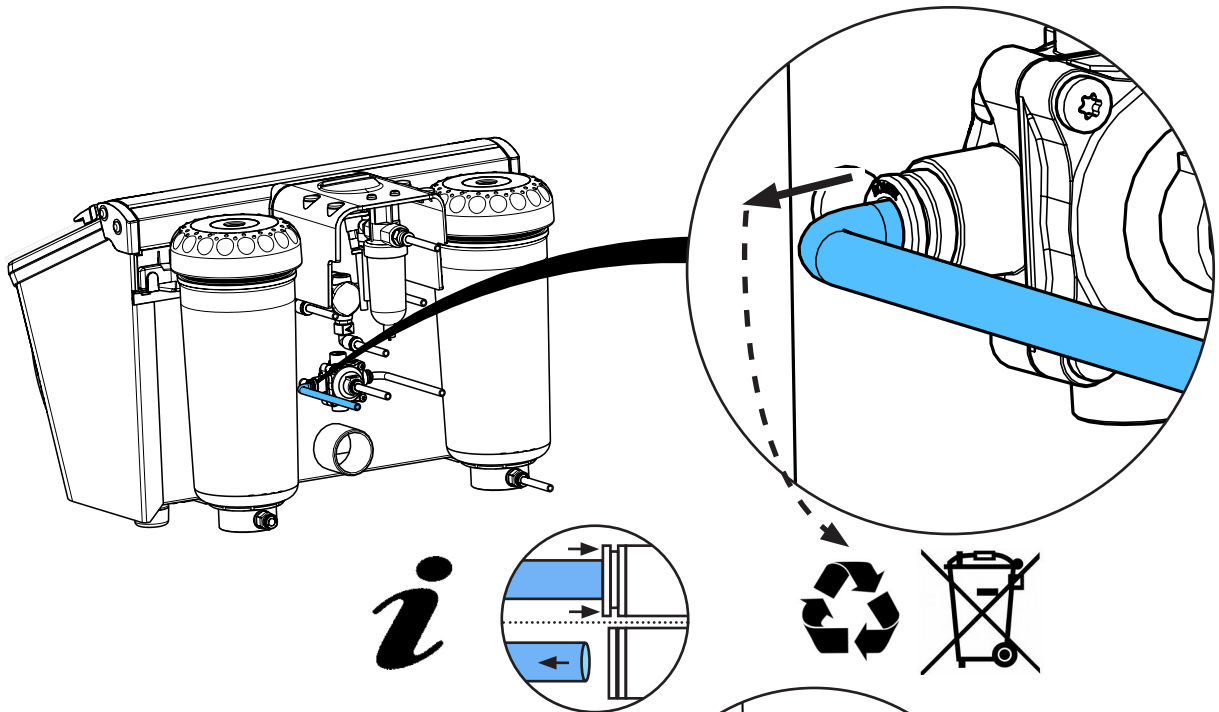

11.

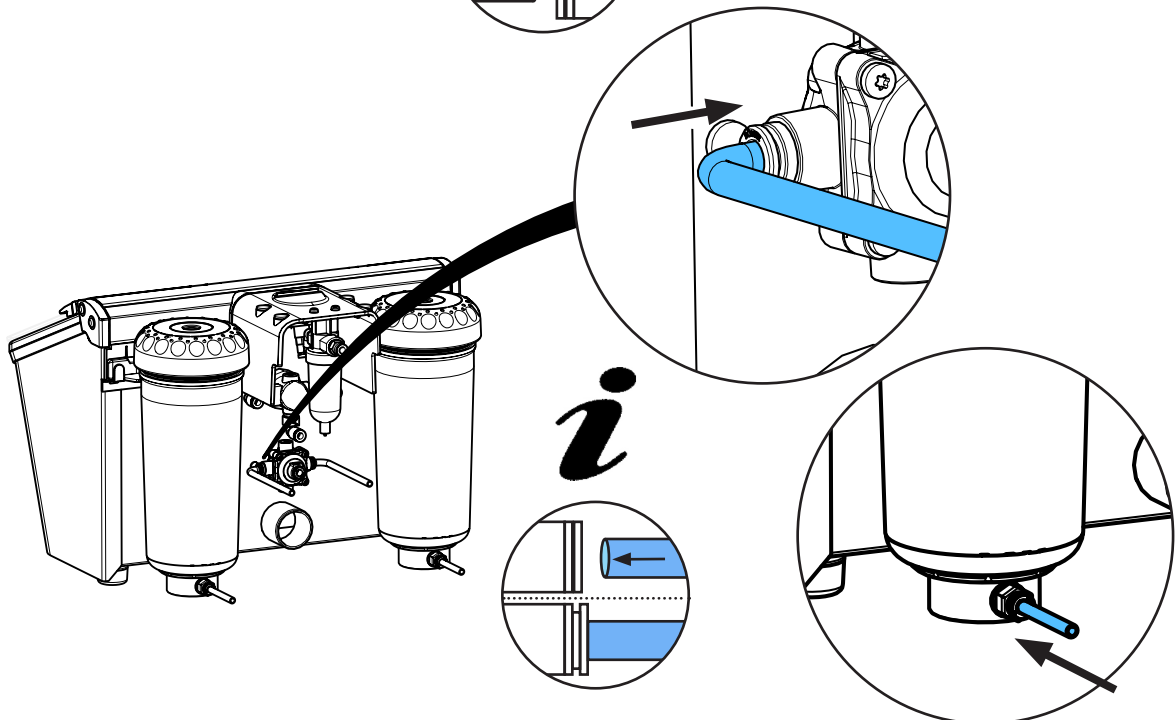

12.

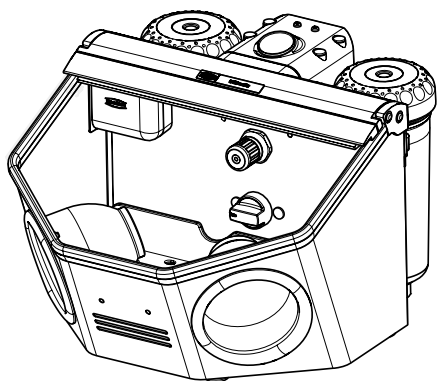





Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу


making work easy

90003 5801


Ersatztankdeckel • Lid f.replacement tank • Couvercle p. silo rec. • Coperchio  
p.dep. ric. • Tapa p.depósito de rec. • Крышка запасного бачка

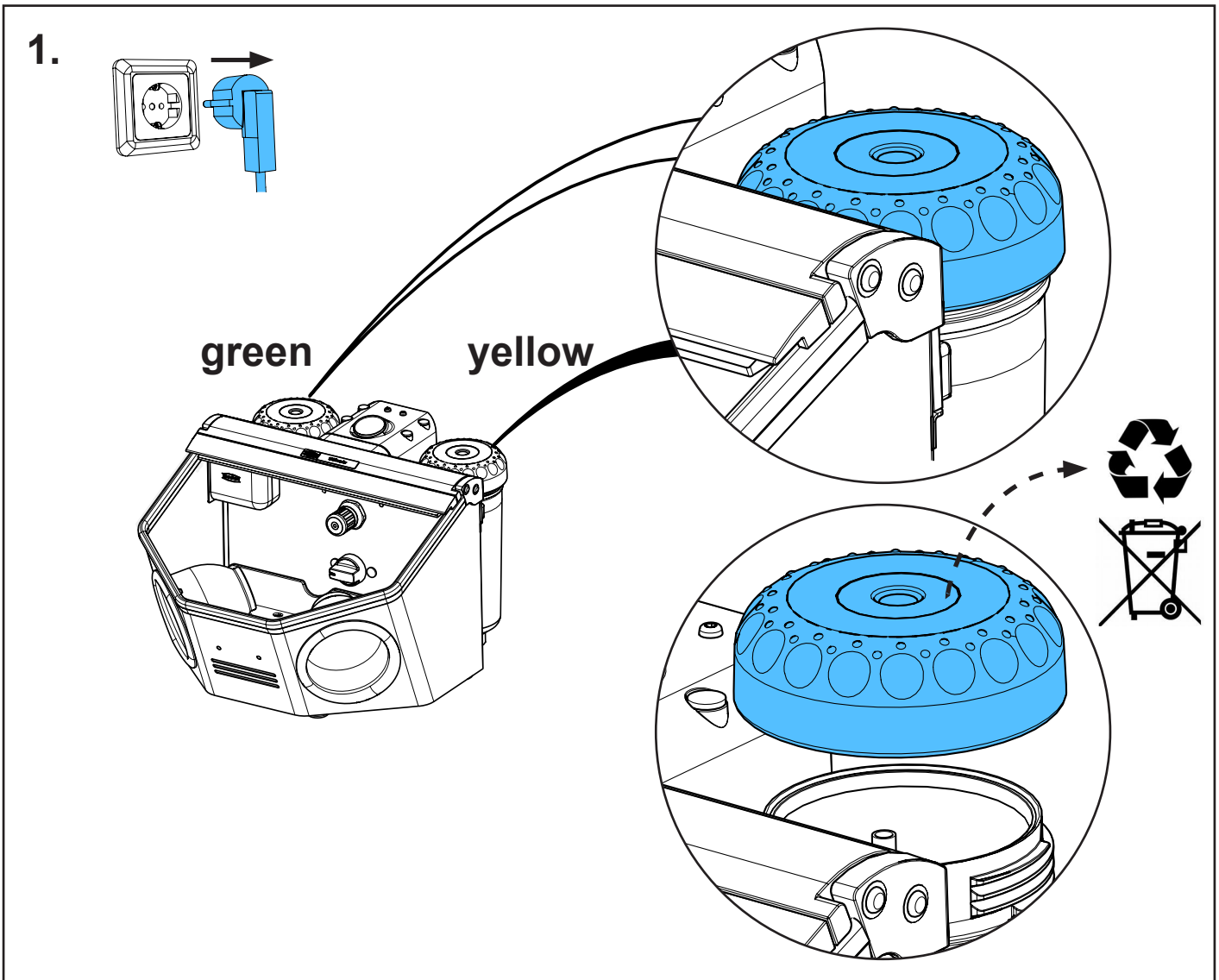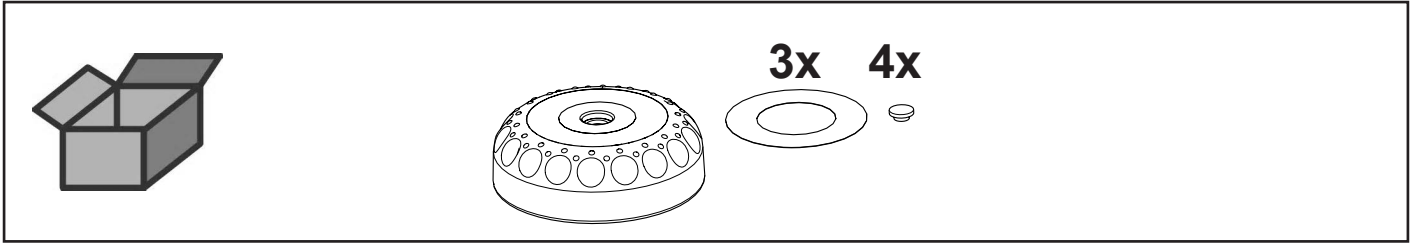


2.

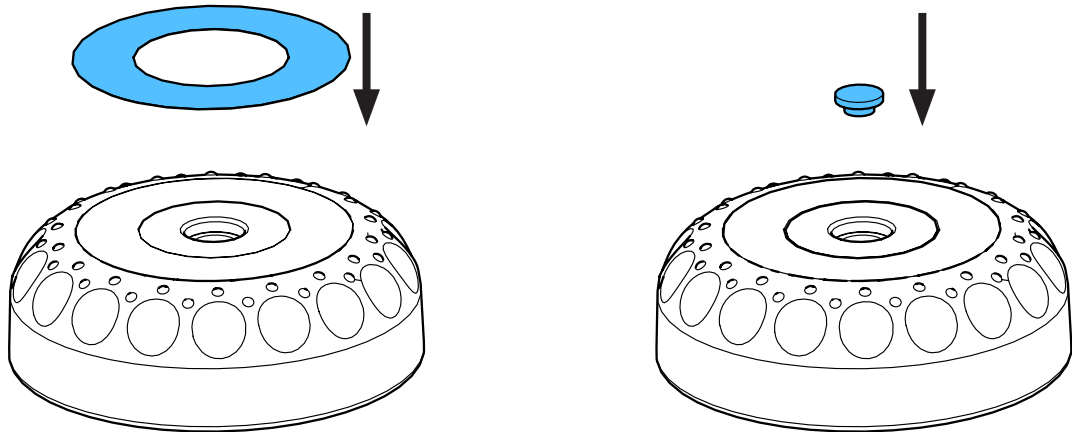

3.

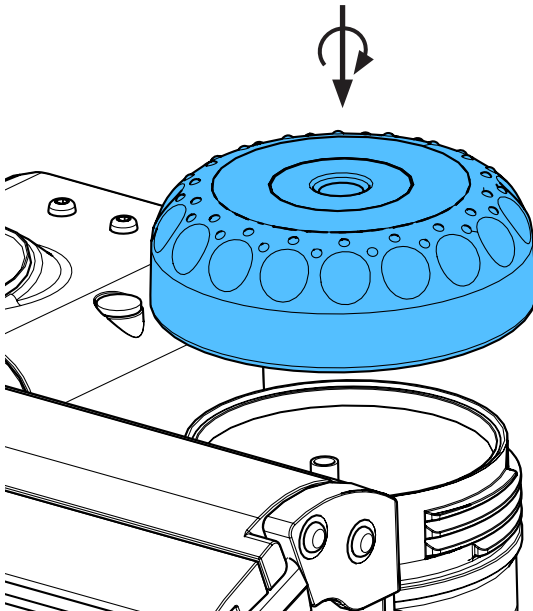

4.



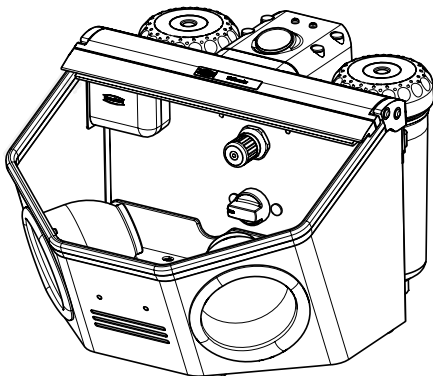





making work easy

## 90002 1431


Filterpatrone • Filter cartridge • Cartouche de filtres • Cartuccia dei filtri con  
raccordo • Cartucho de filtración • Патрон фильтра

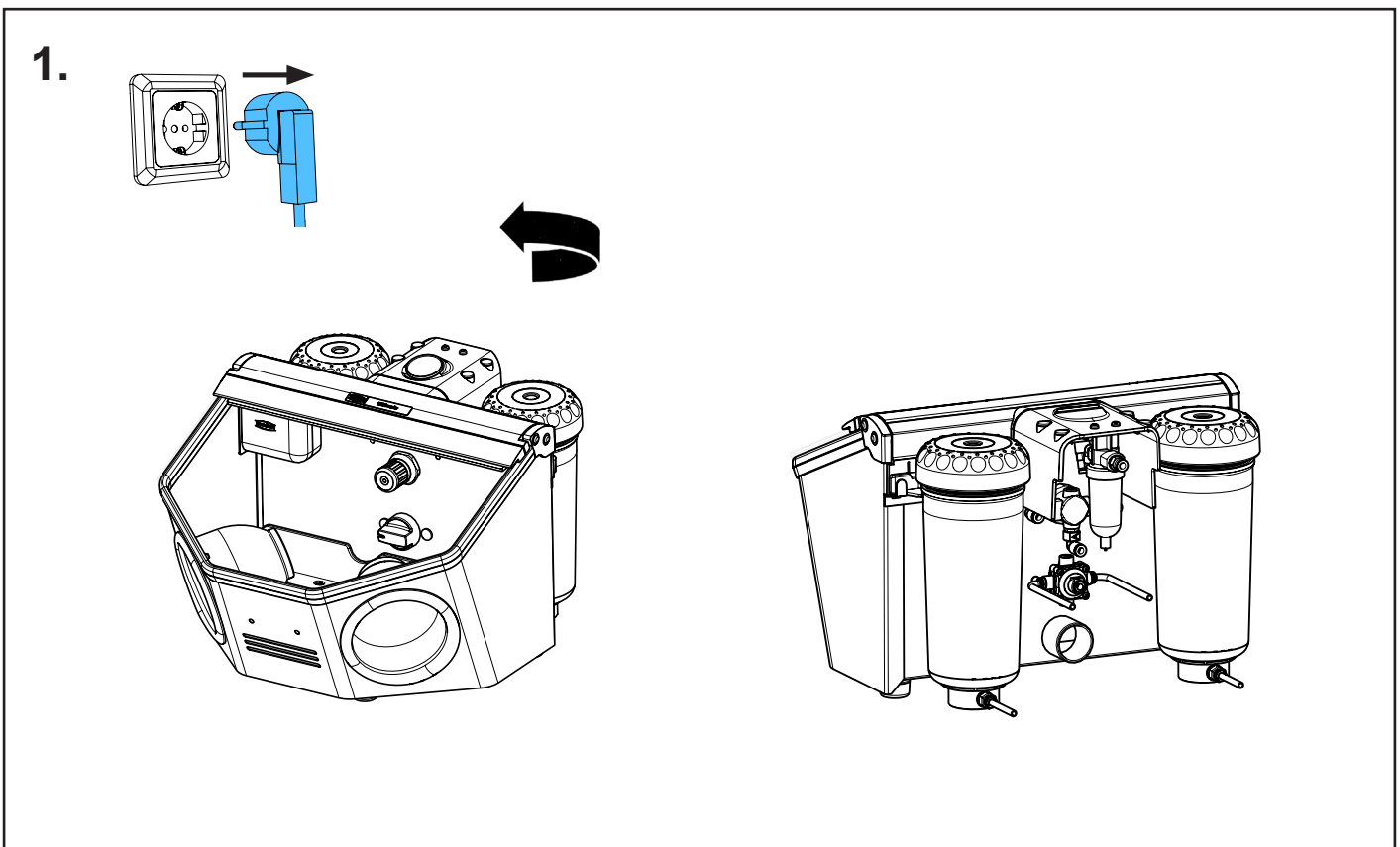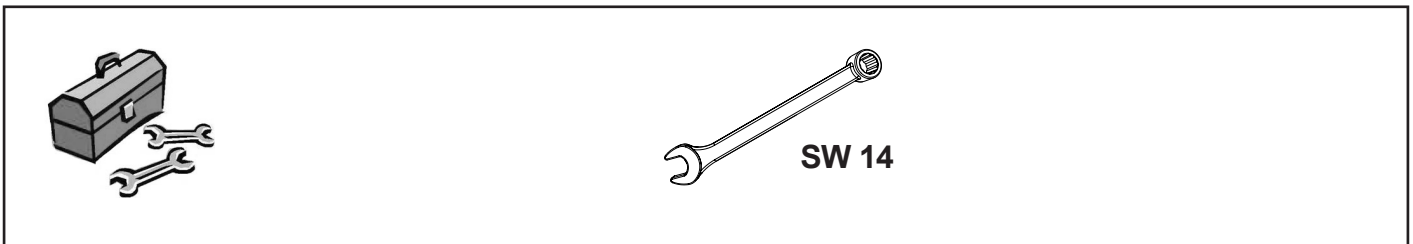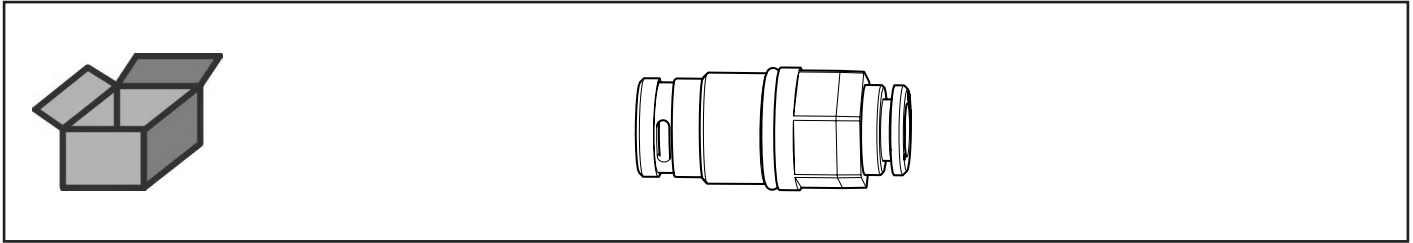


2.

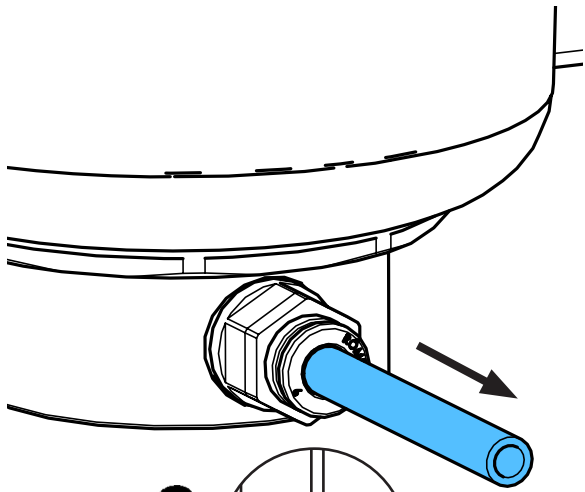


**i**

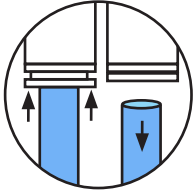


3.

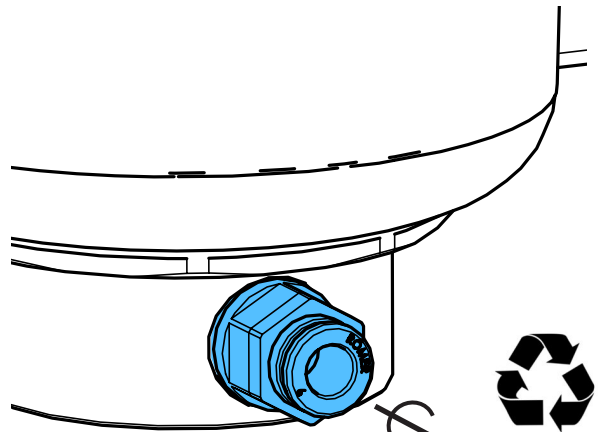



4.

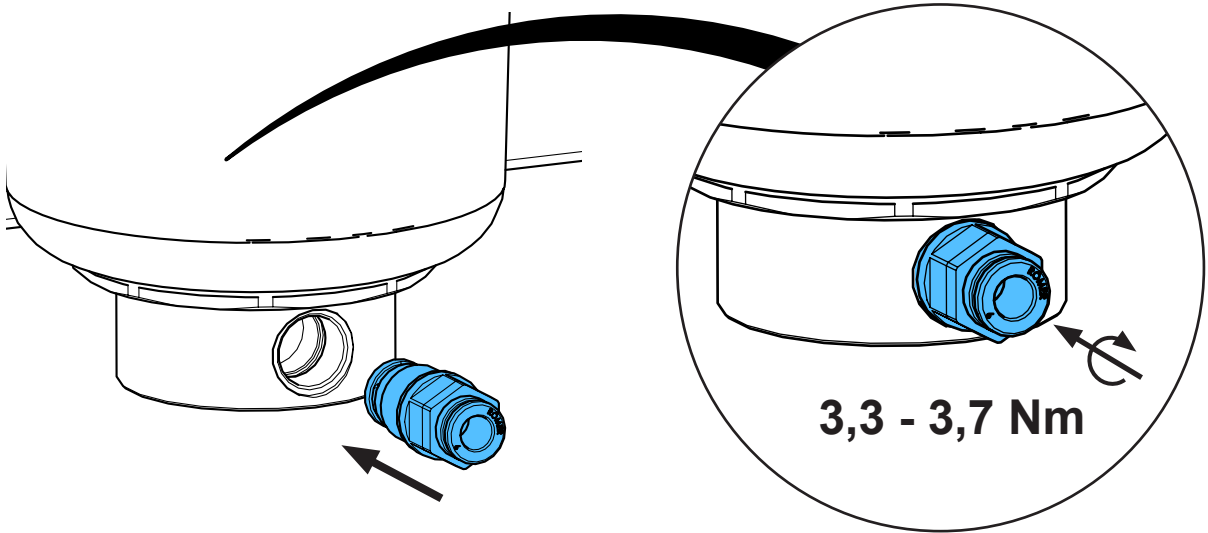

5.

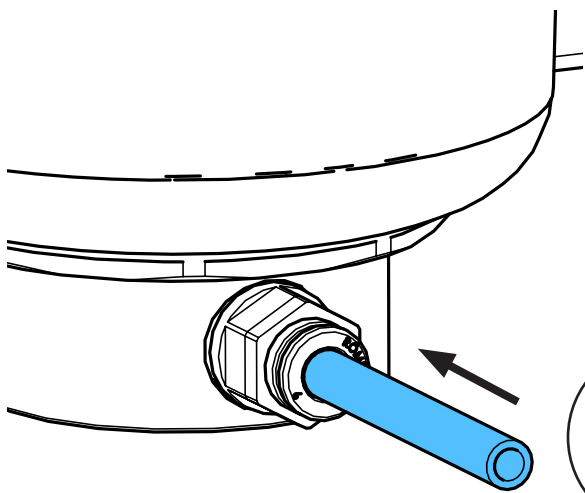

**i**

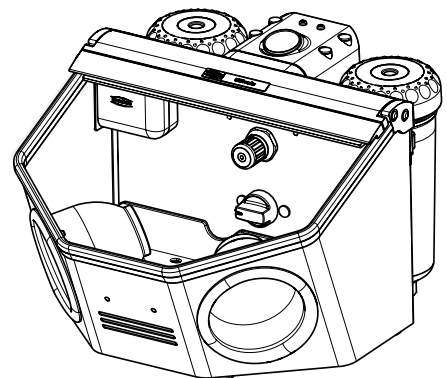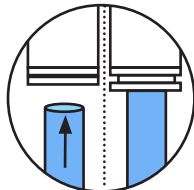









making work easy

## 90003 3035



O-Ring DIN 3771 82x4 NBR 55 • O-ring DIN 3771 82x4 NBR 55 • Joint thorique  
DIN 3771 82x4 NBR 55 • Guarnizione DIN 3771 82x4 NBR 55 • Anillo en O DIN  
3771 82x4 NBR 55 • Кольцо круглого сечения DIN 3771, 82x4 NBR 55

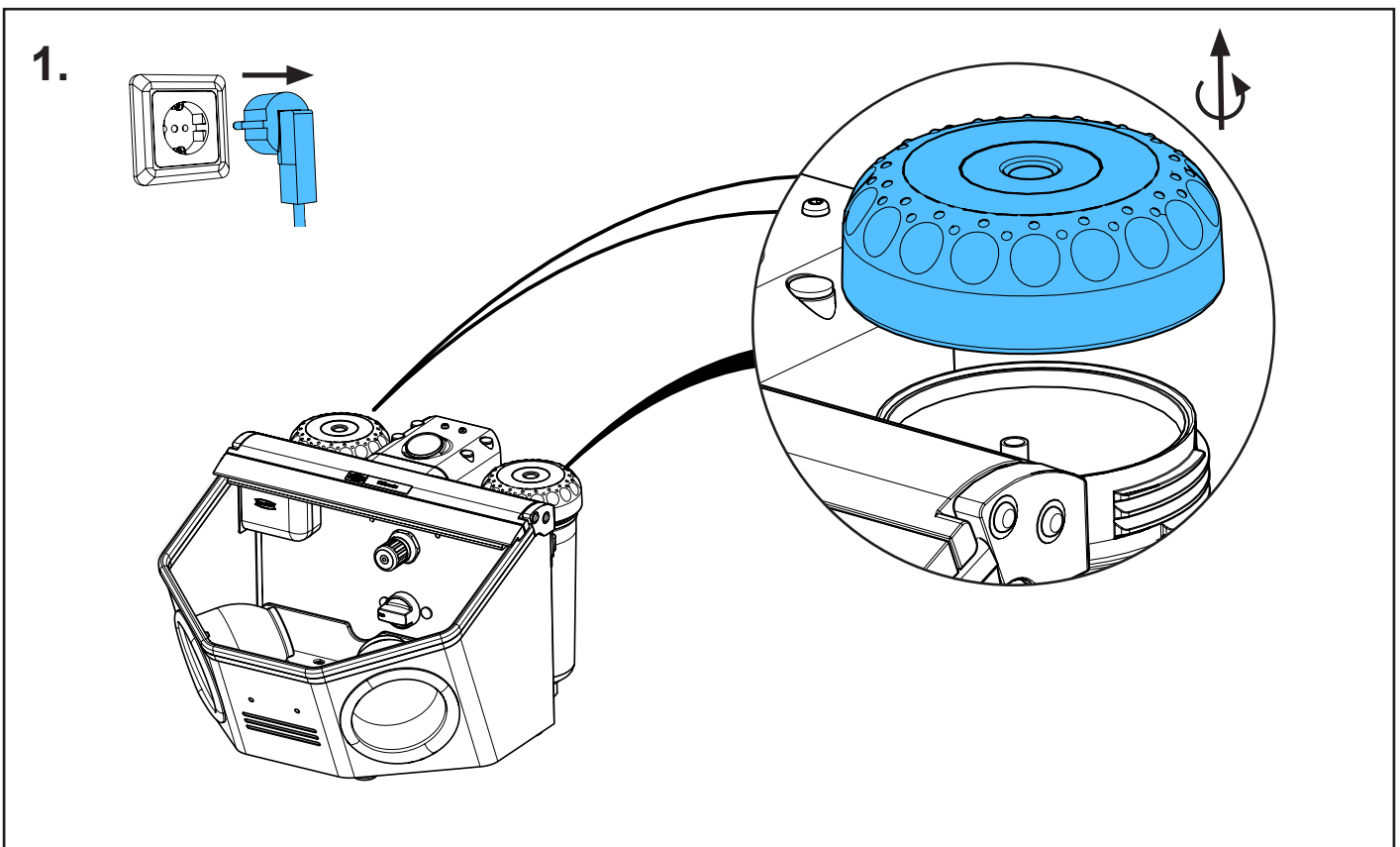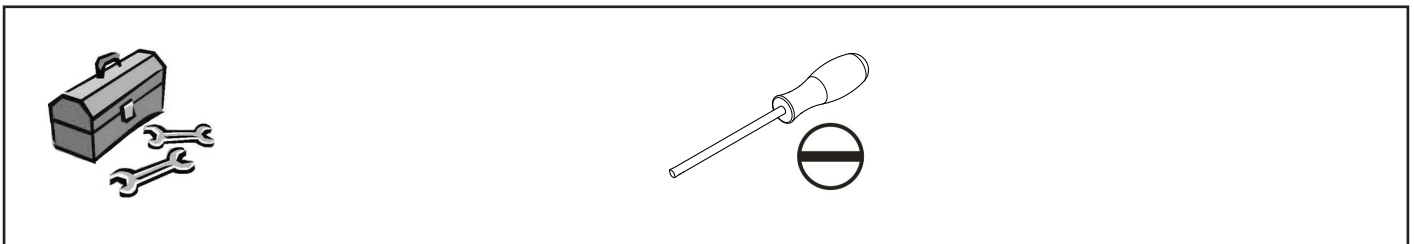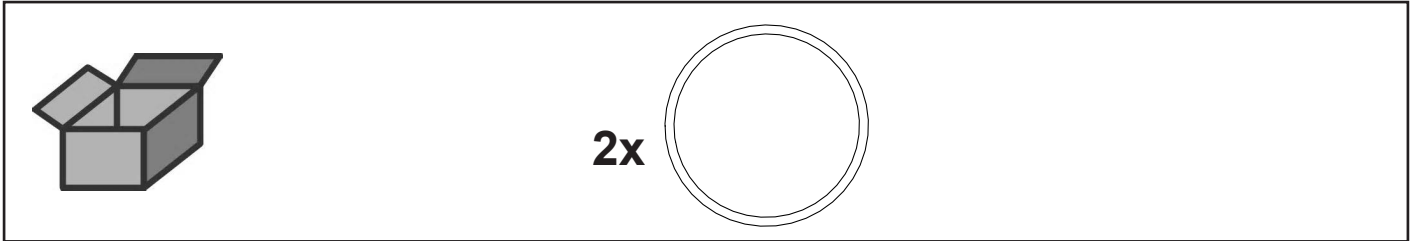



2.

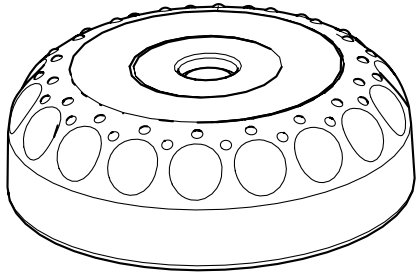



3.

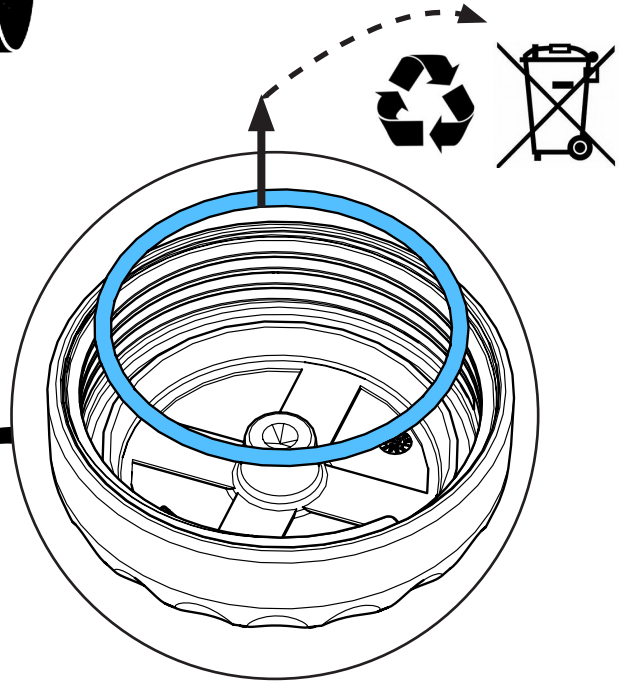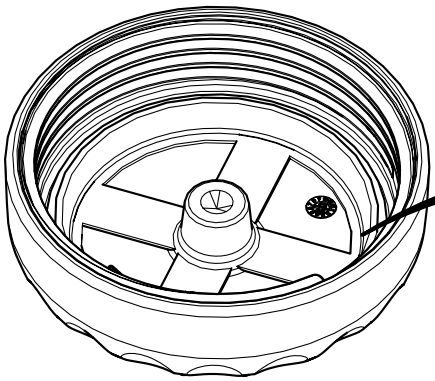

4.

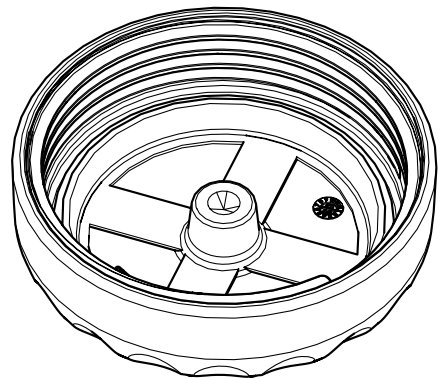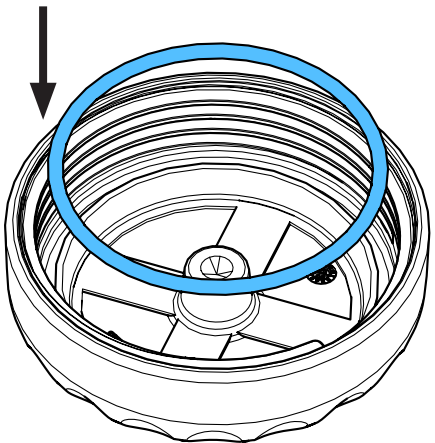

5.

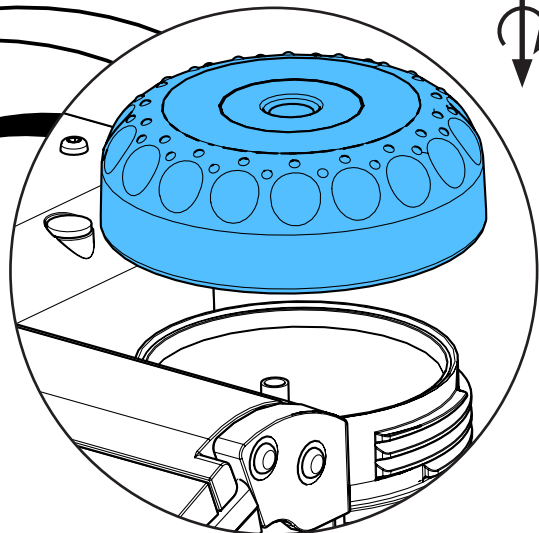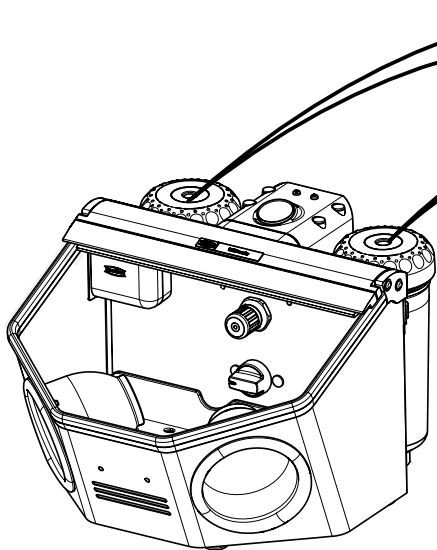





Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по монтажу


making work easy

**92947 0002**


Kleinteilesset Basic • Hardware Basic • Petit materiel d.montage Basic • Kit di minuteria Basic • Juego de piezas pequeñas Basic • Комплект мелких деталей

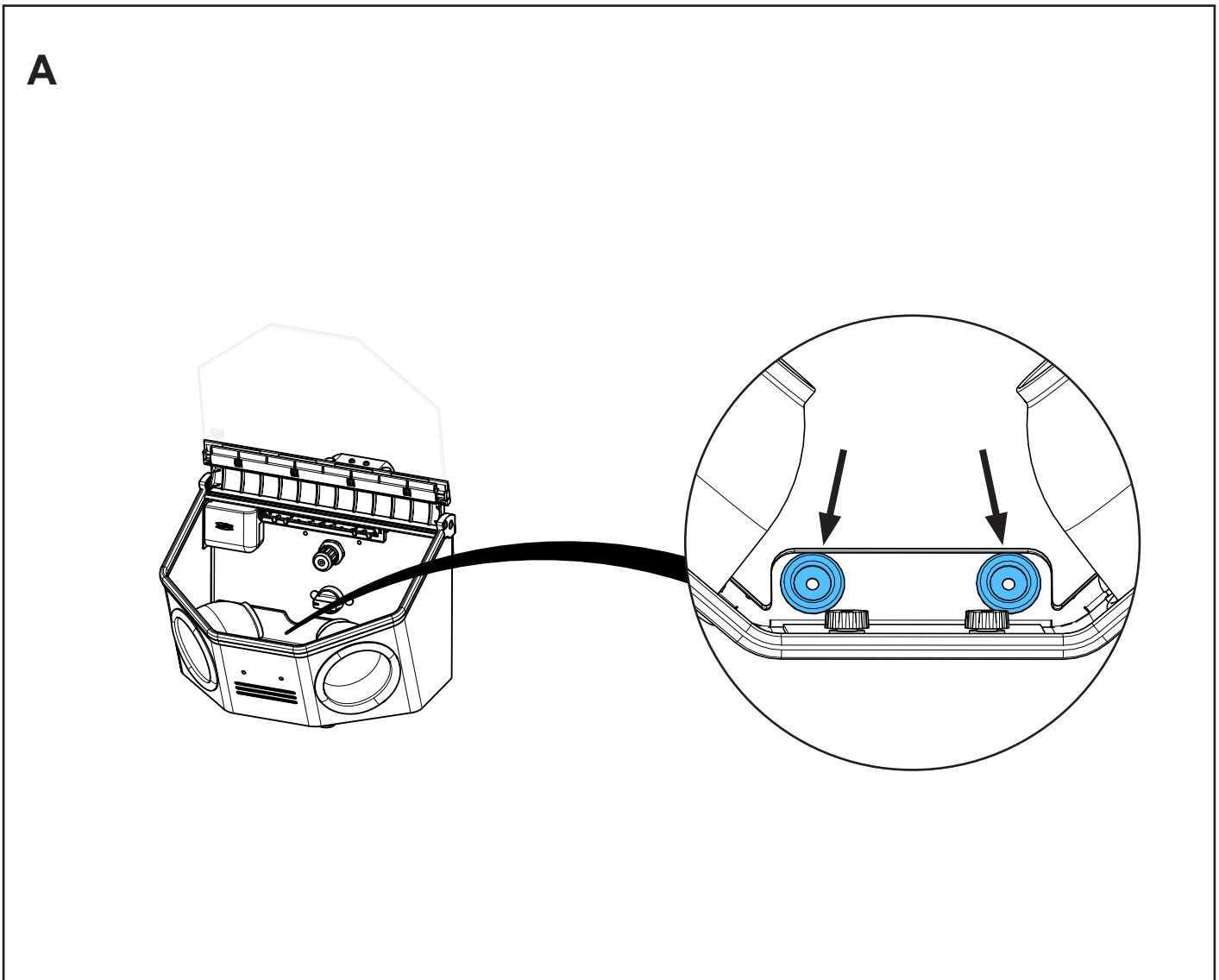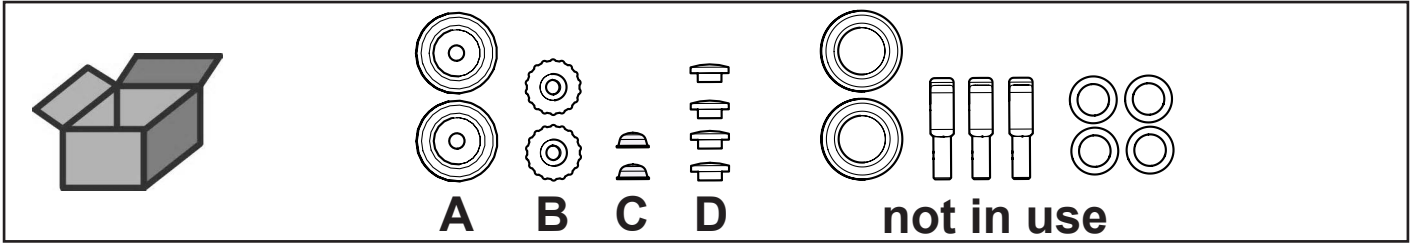


**B**

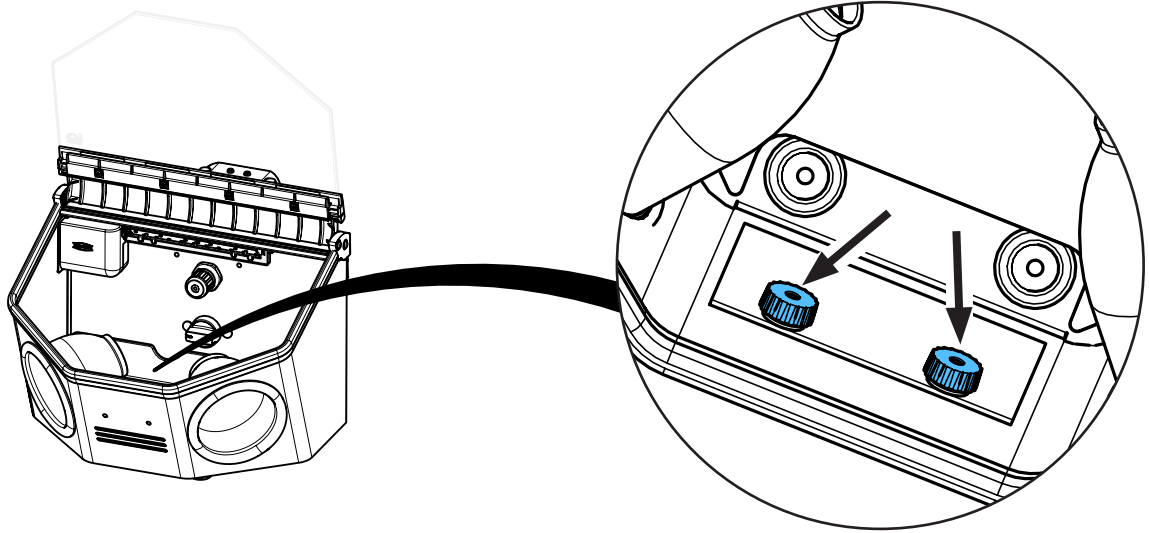

**C**

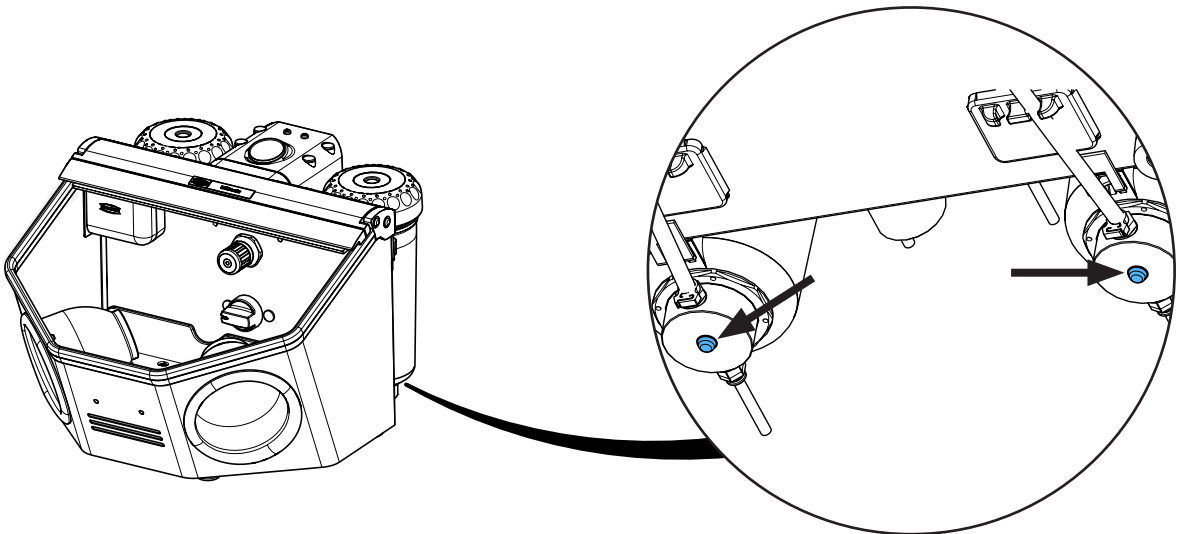

**D**

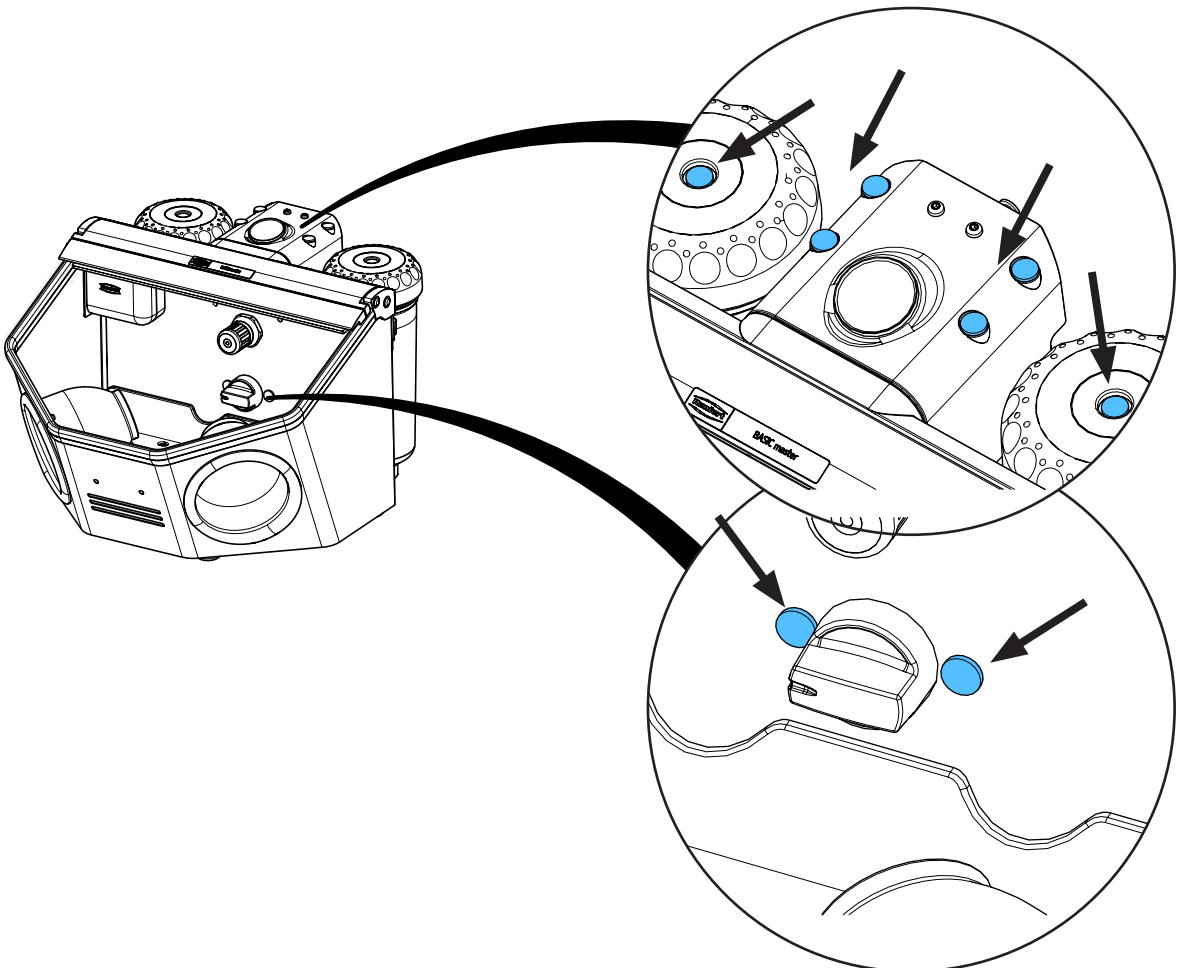





Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу


making work easy

**90003 6330**


Bodengitter • Base grid • Grille de fond • Griglia di fondo • Rejilla de fondo •  
Решетка для дна

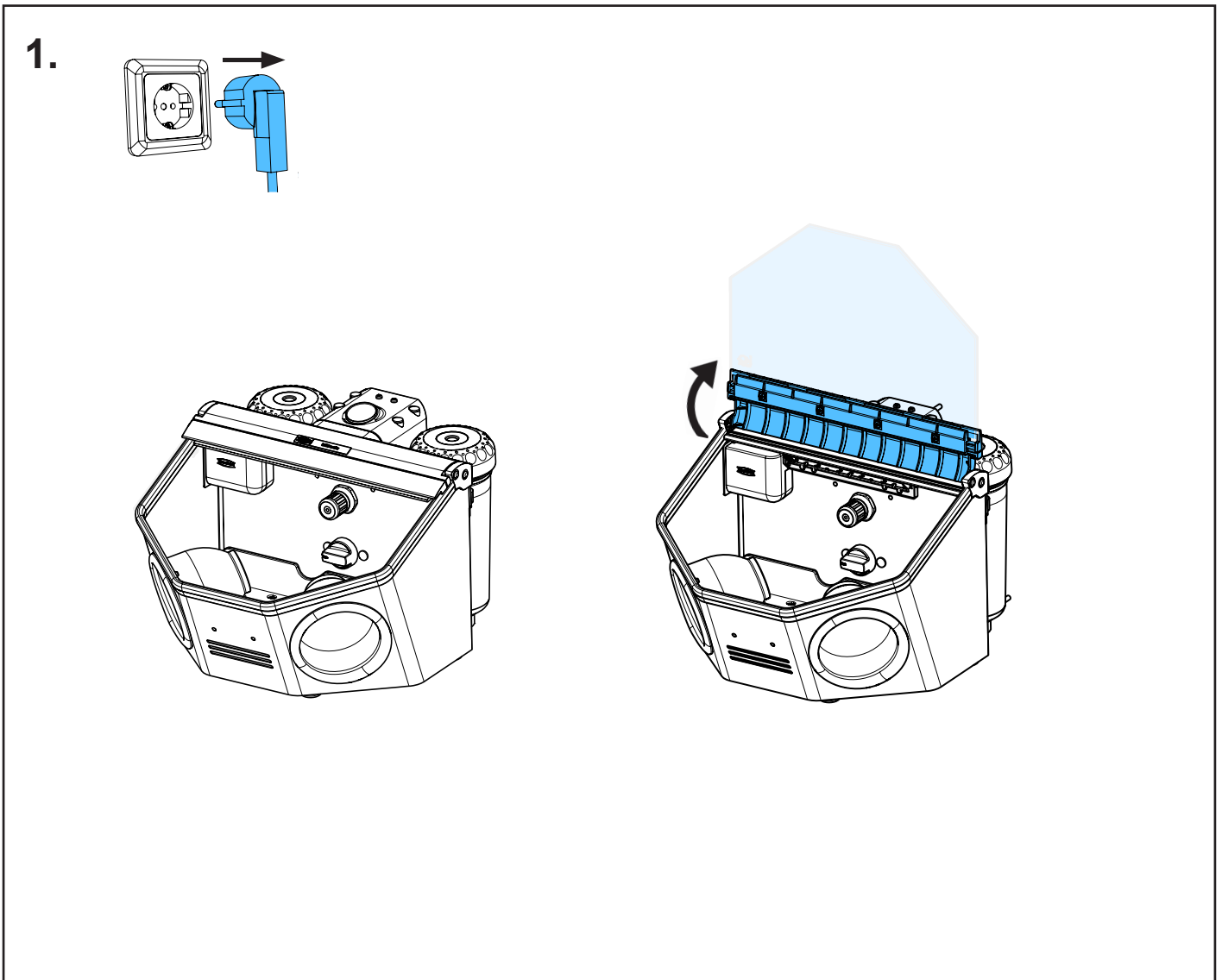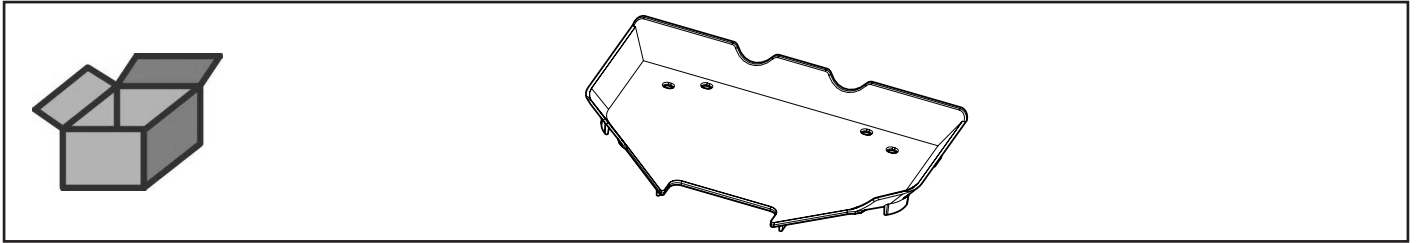


2.

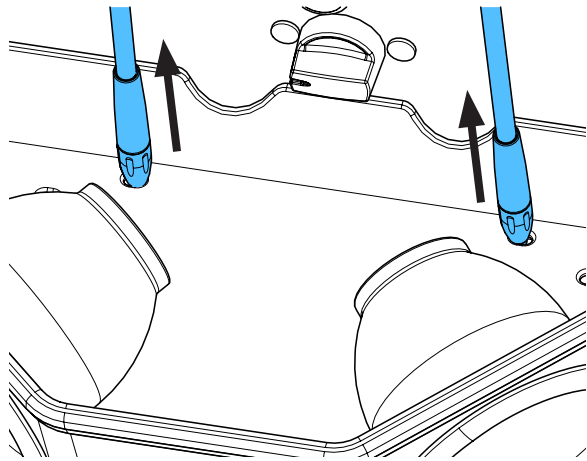


3.

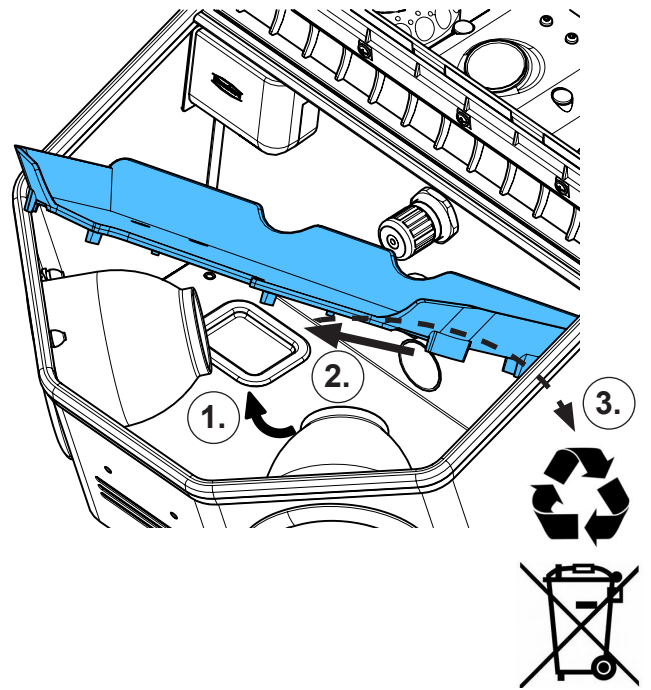


4.

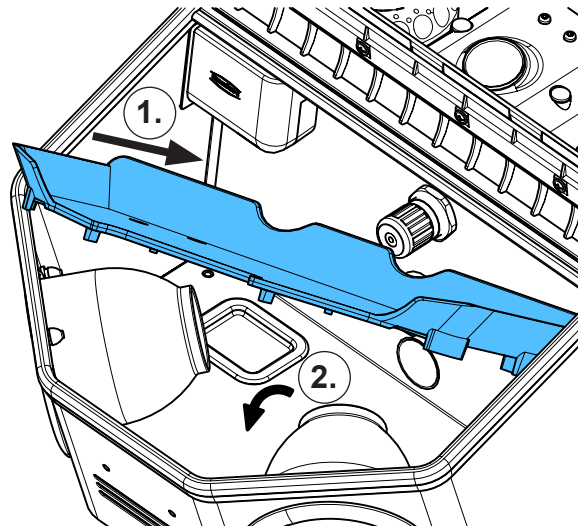

5.

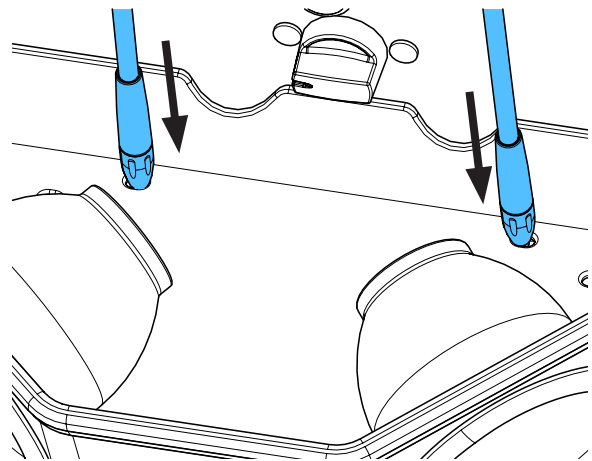

6.

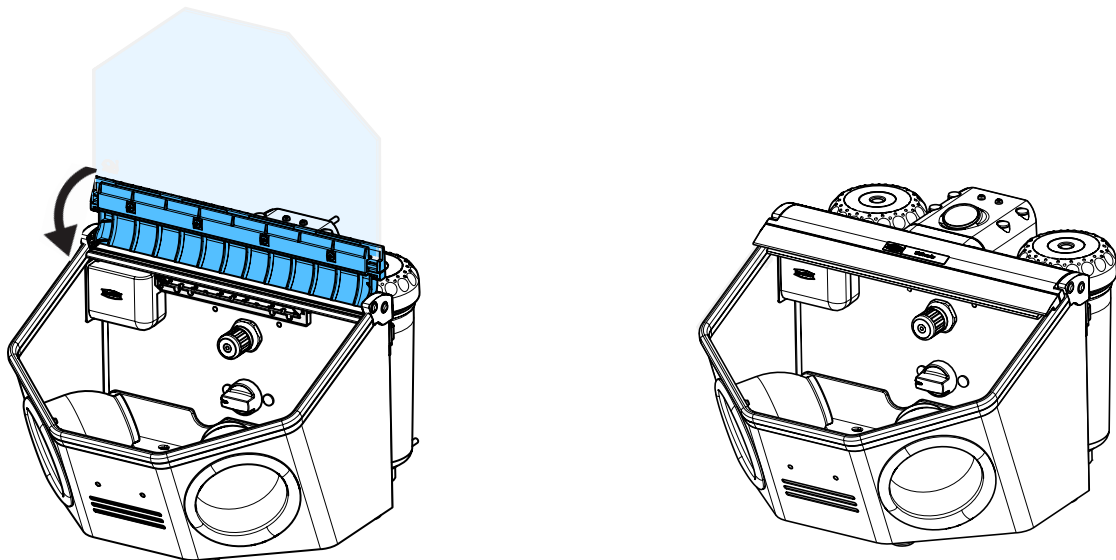





Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу


making work easy

90003 4751


Handlochmanschetten • Hand sleeves • Manchettes • Manicotto • Puños •  
Манжета серая

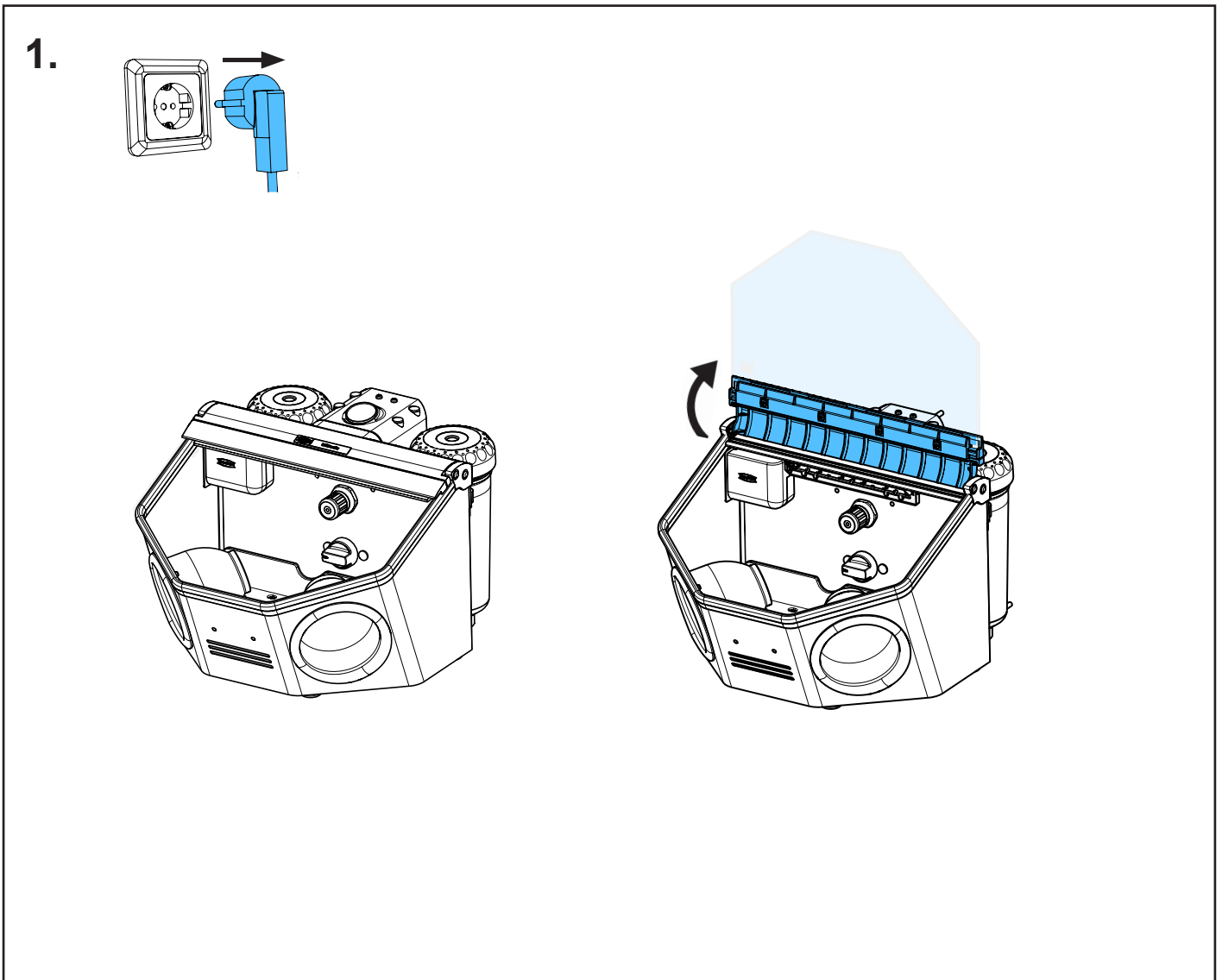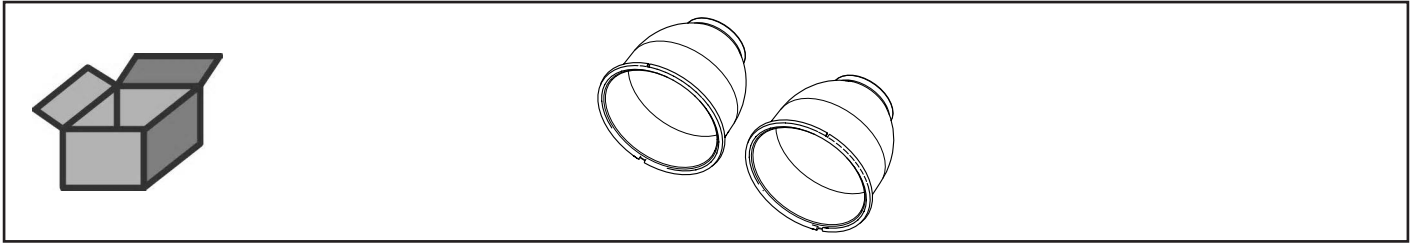


2.

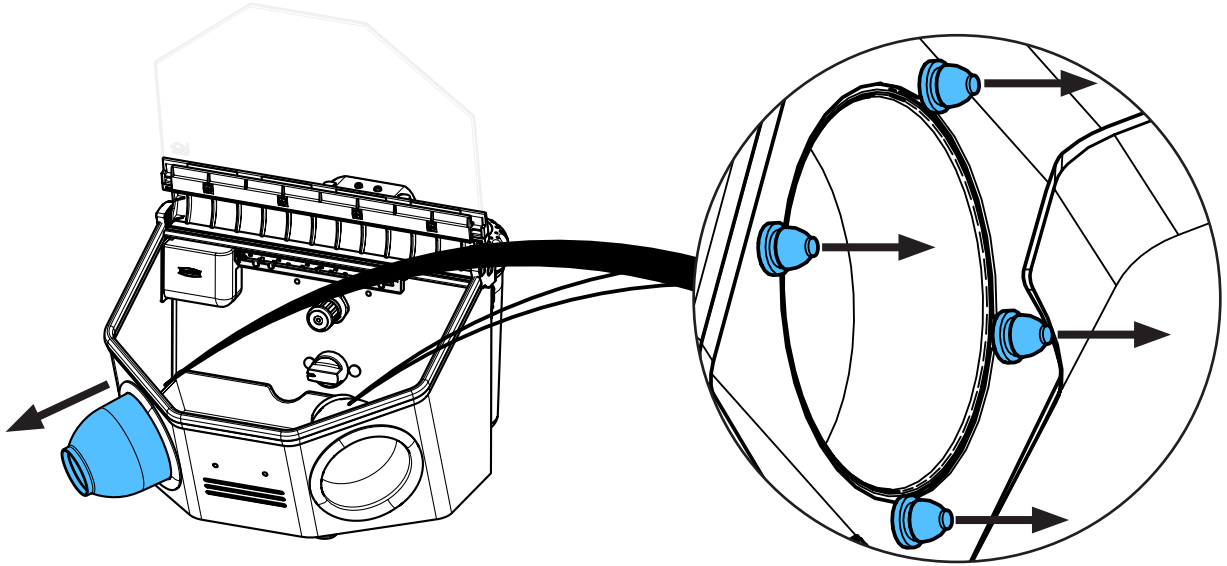

3.

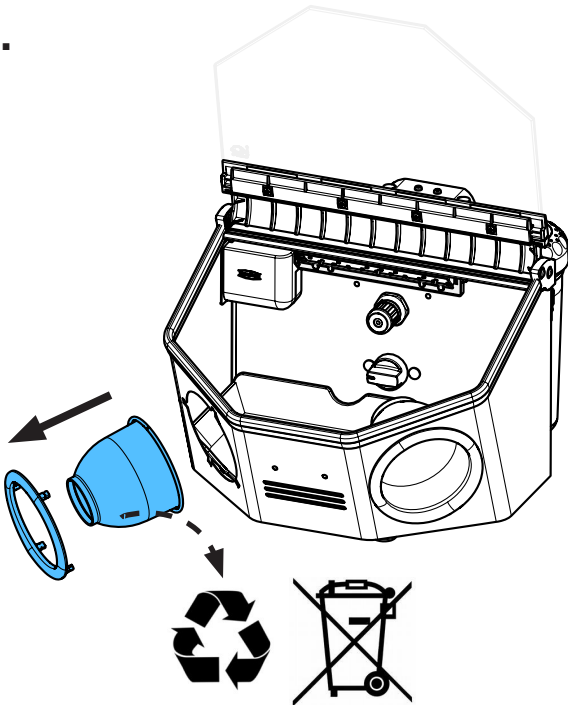


4.

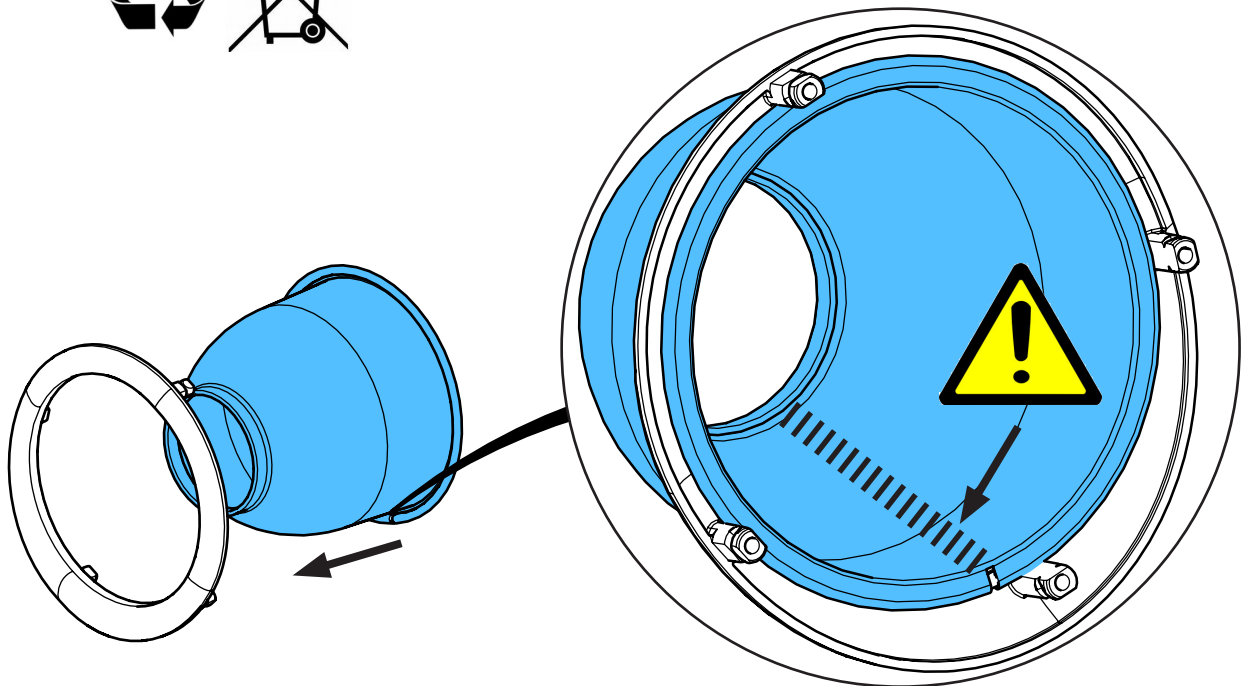





5.

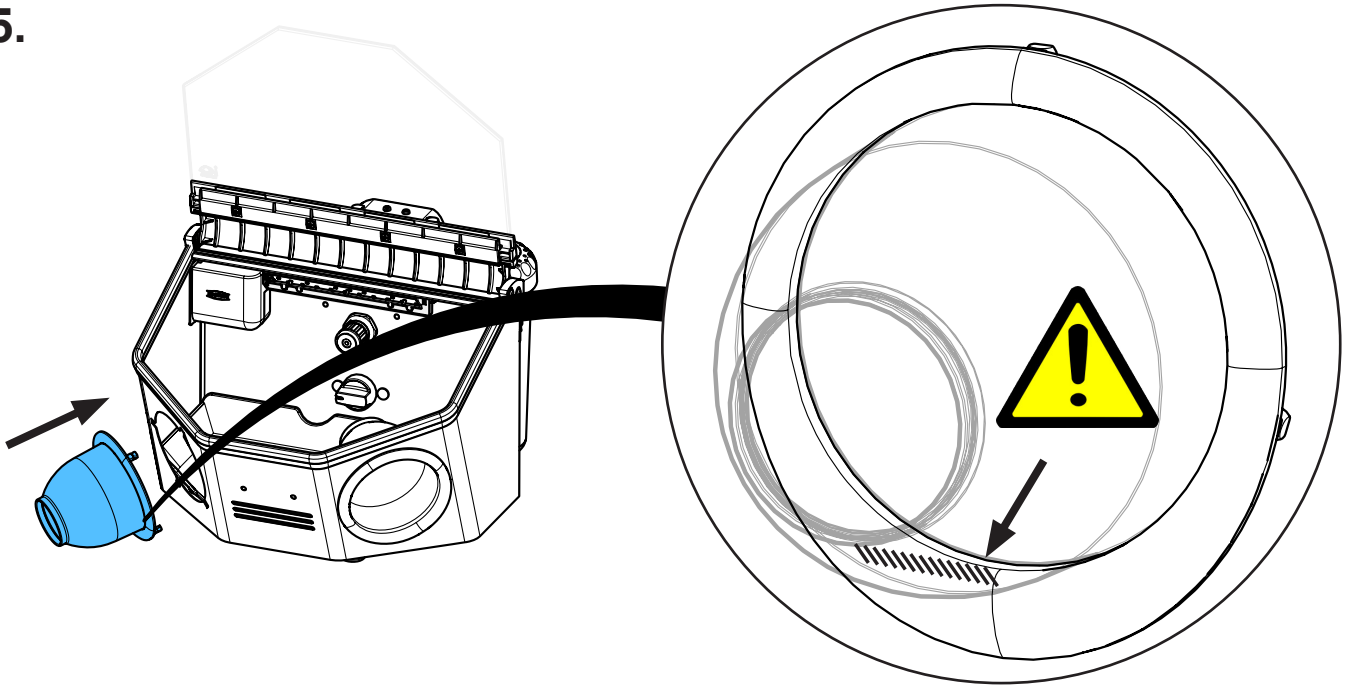

6.

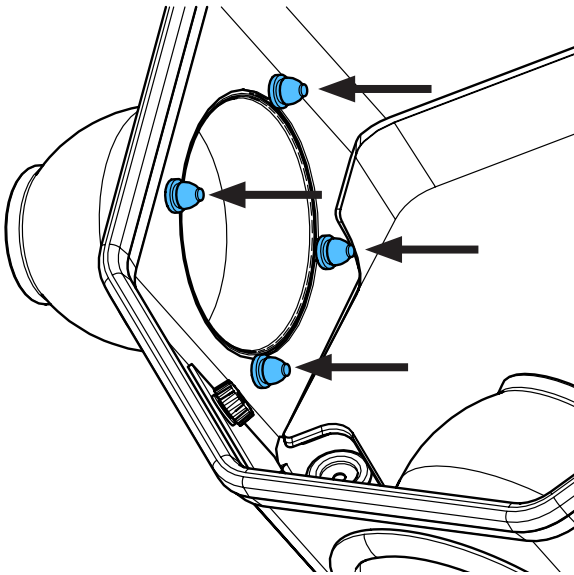

7.

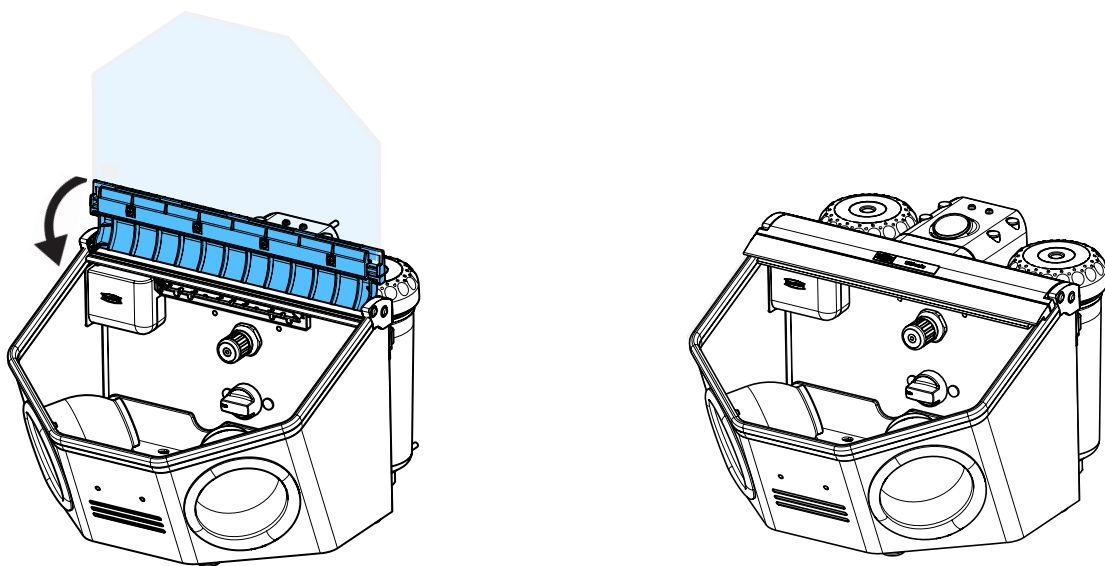







Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу


making work easy

**90003 6339**


Handstulpenring • Ring for hand sleeves • Anneau pour gants • Ghiera per mani-  
cotto • Anillo para puños • Кольцо манжеты

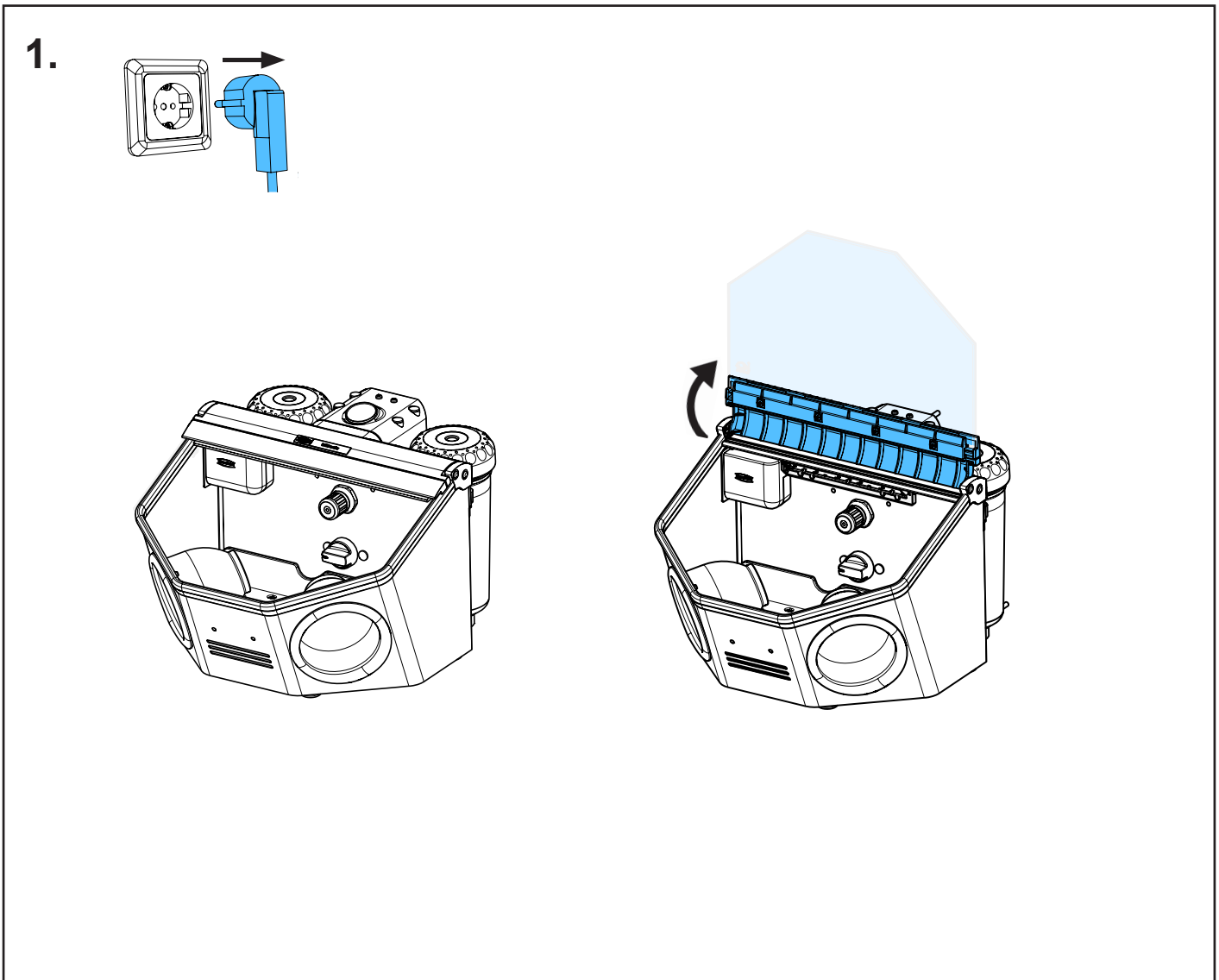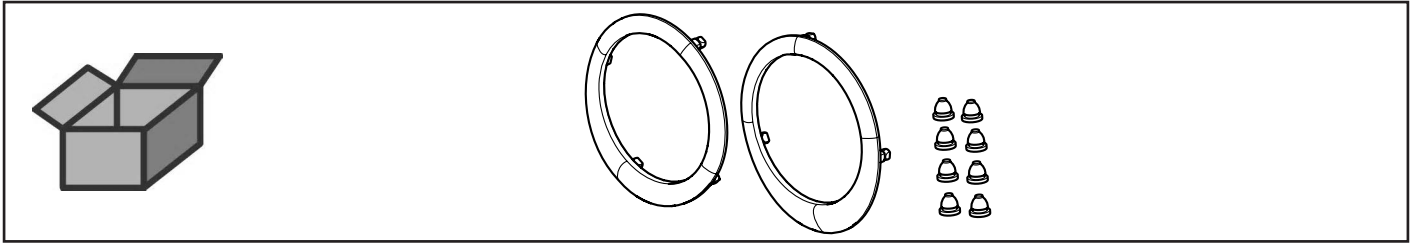


2.

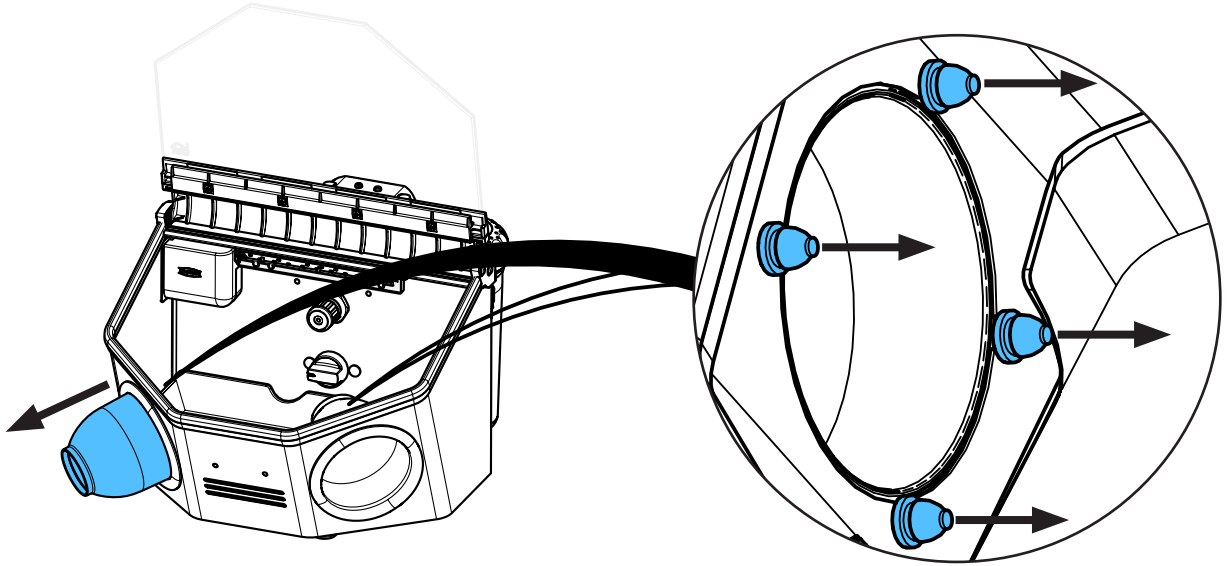


3.

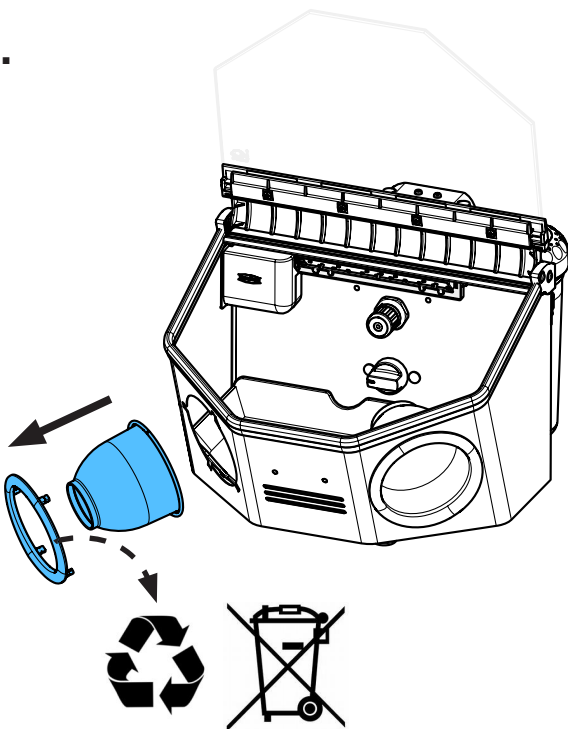


4.

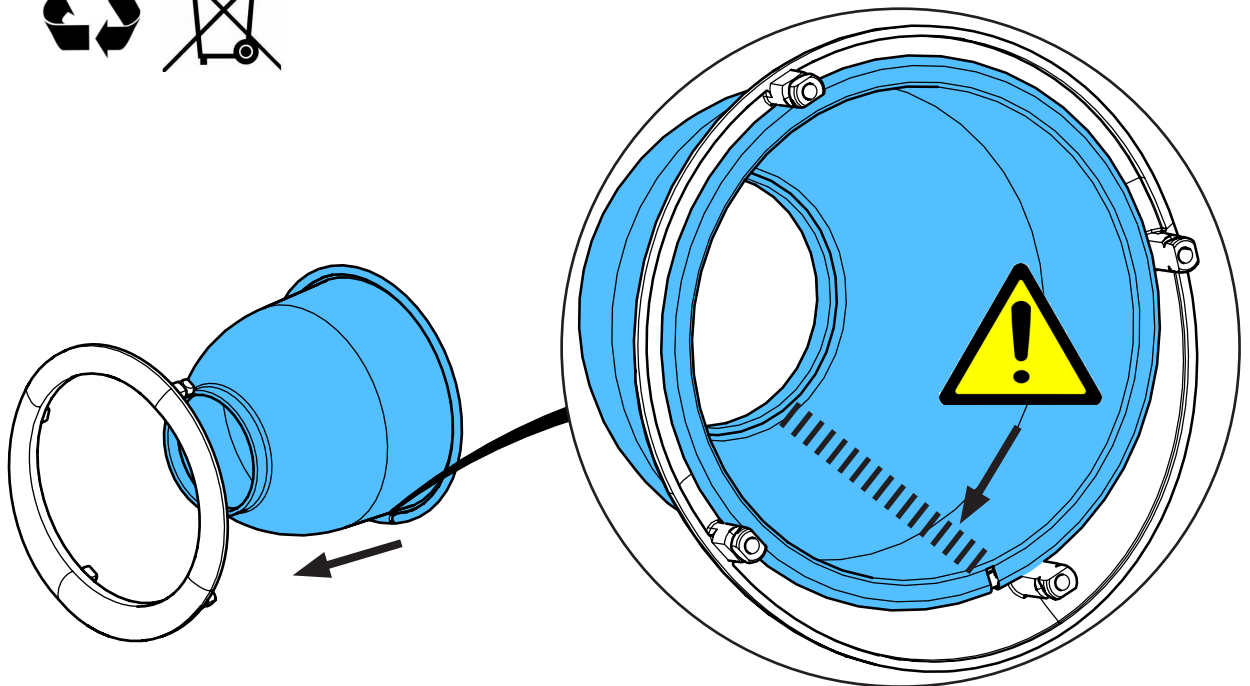


5.

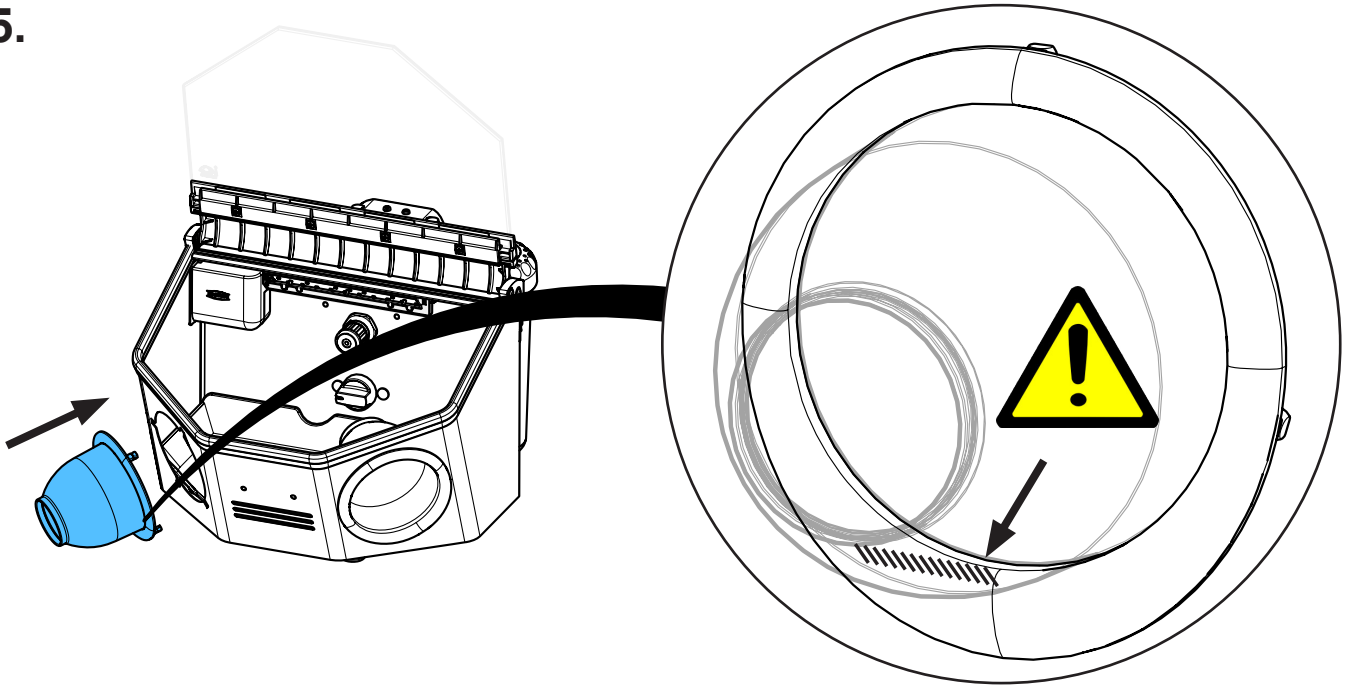

6.

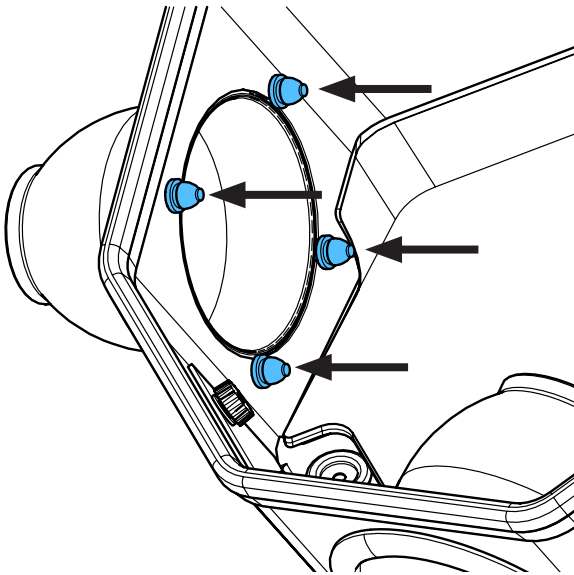

7.

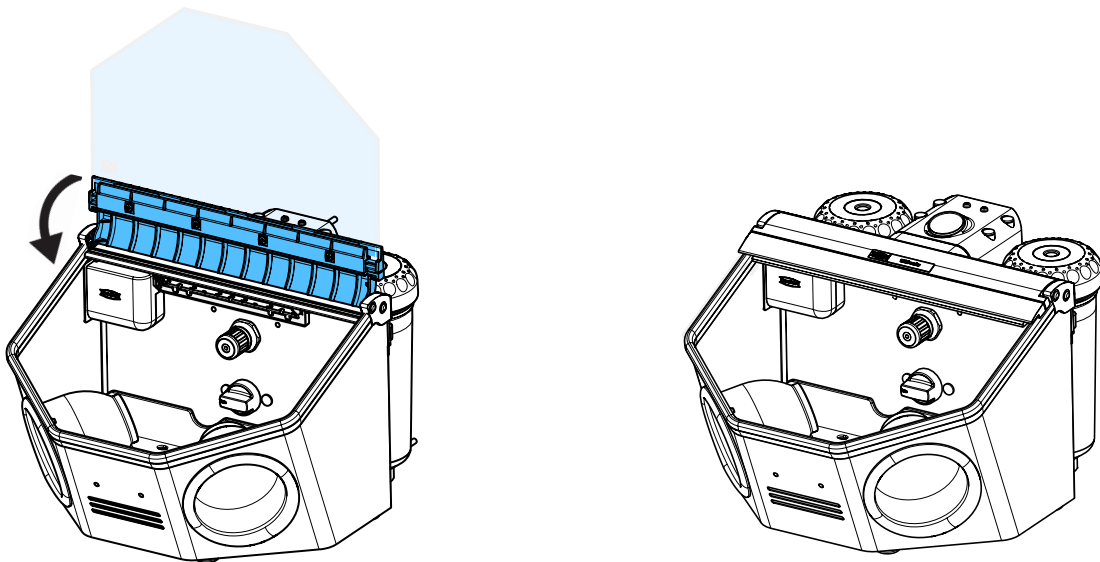









making work easy

90002 1435


Glasscheibe beschichtet • Coated glass screen • Vitre revêtu • Vetro rivestito •  
Cristal con revestimiento • Запасное стекло с покрытием

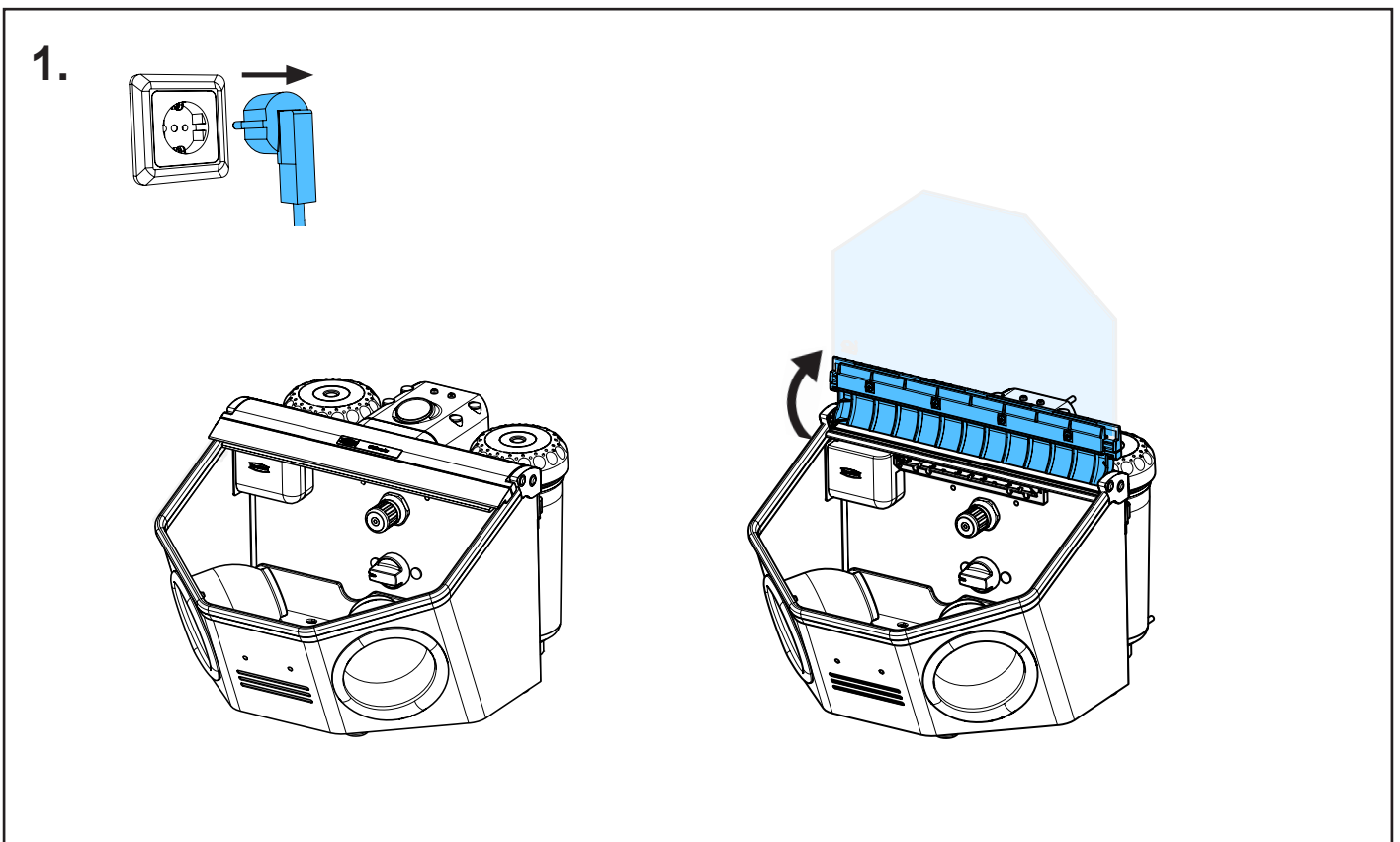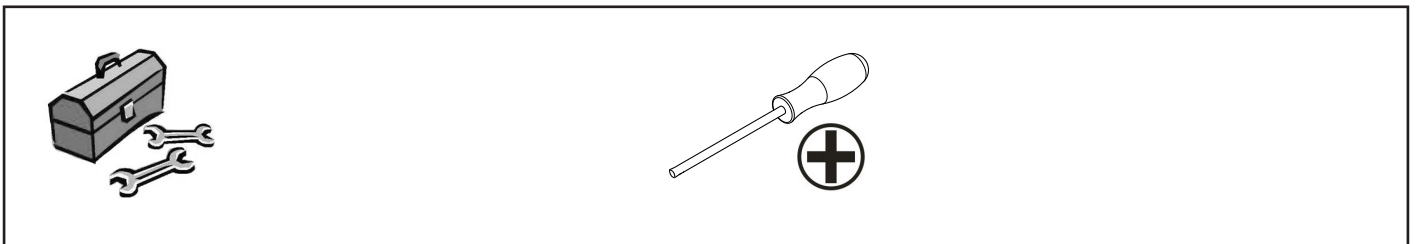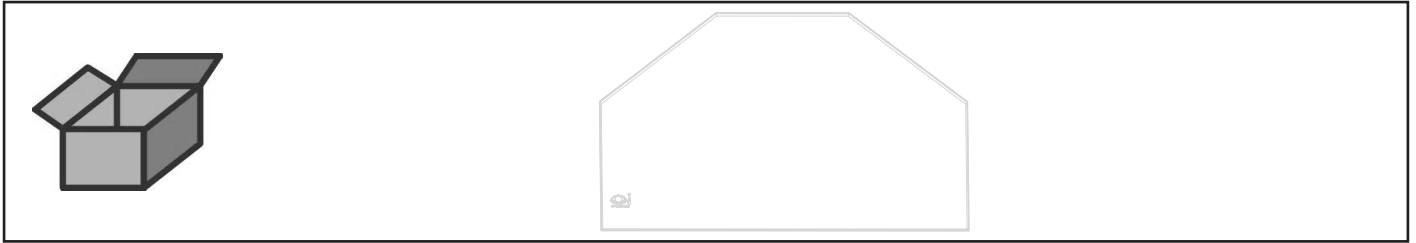


2.

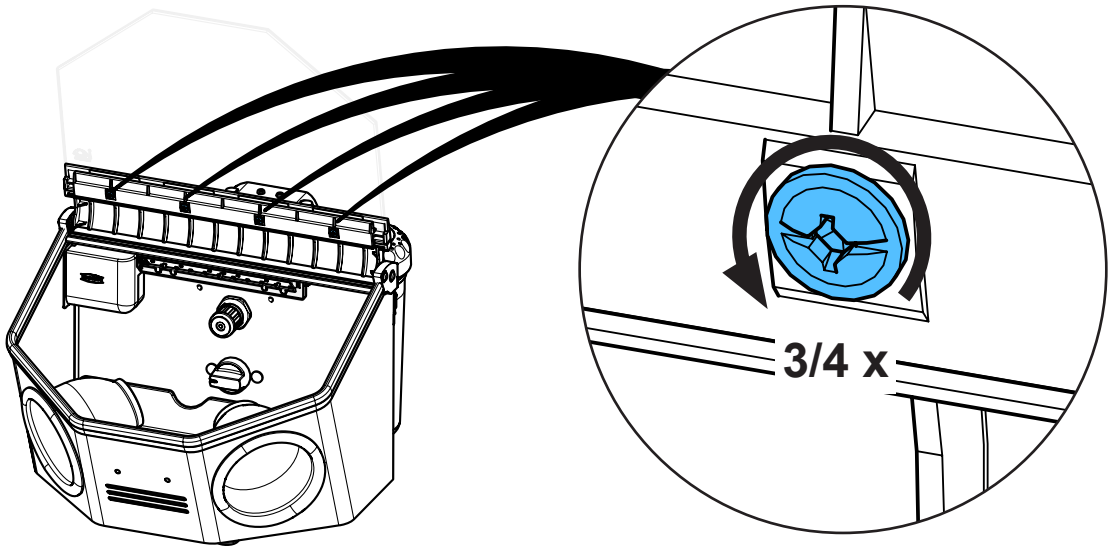

3.

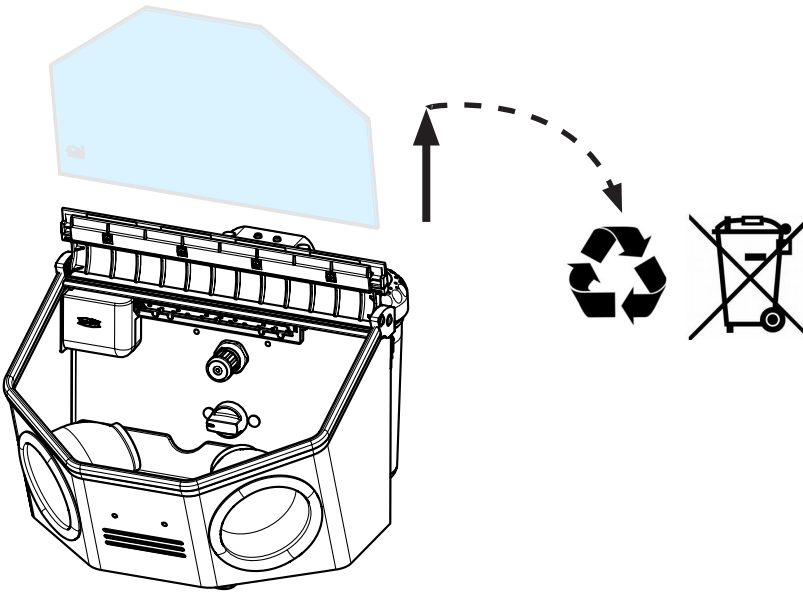


4.

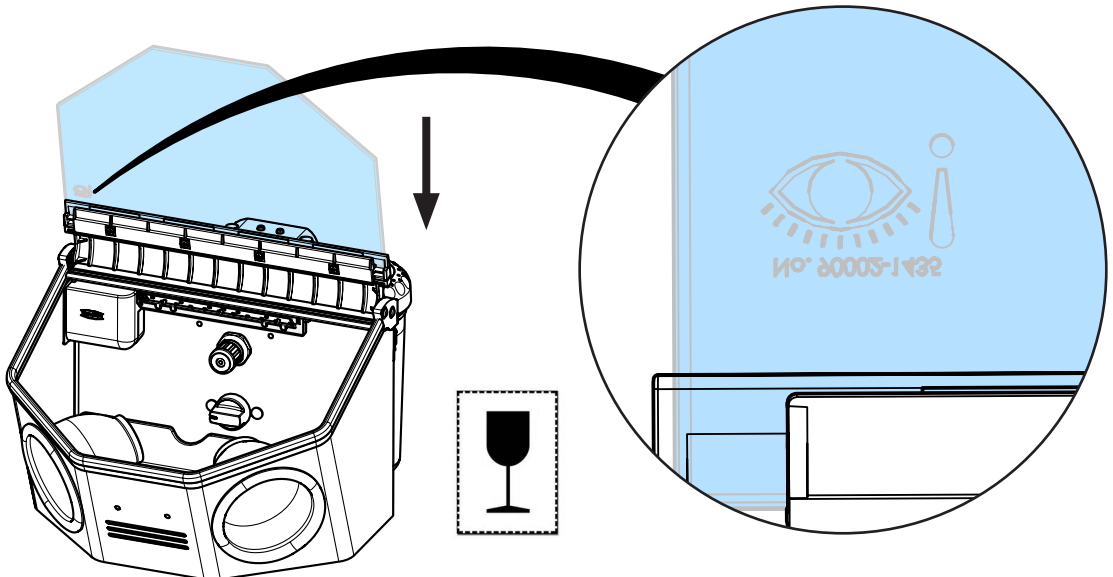




5.


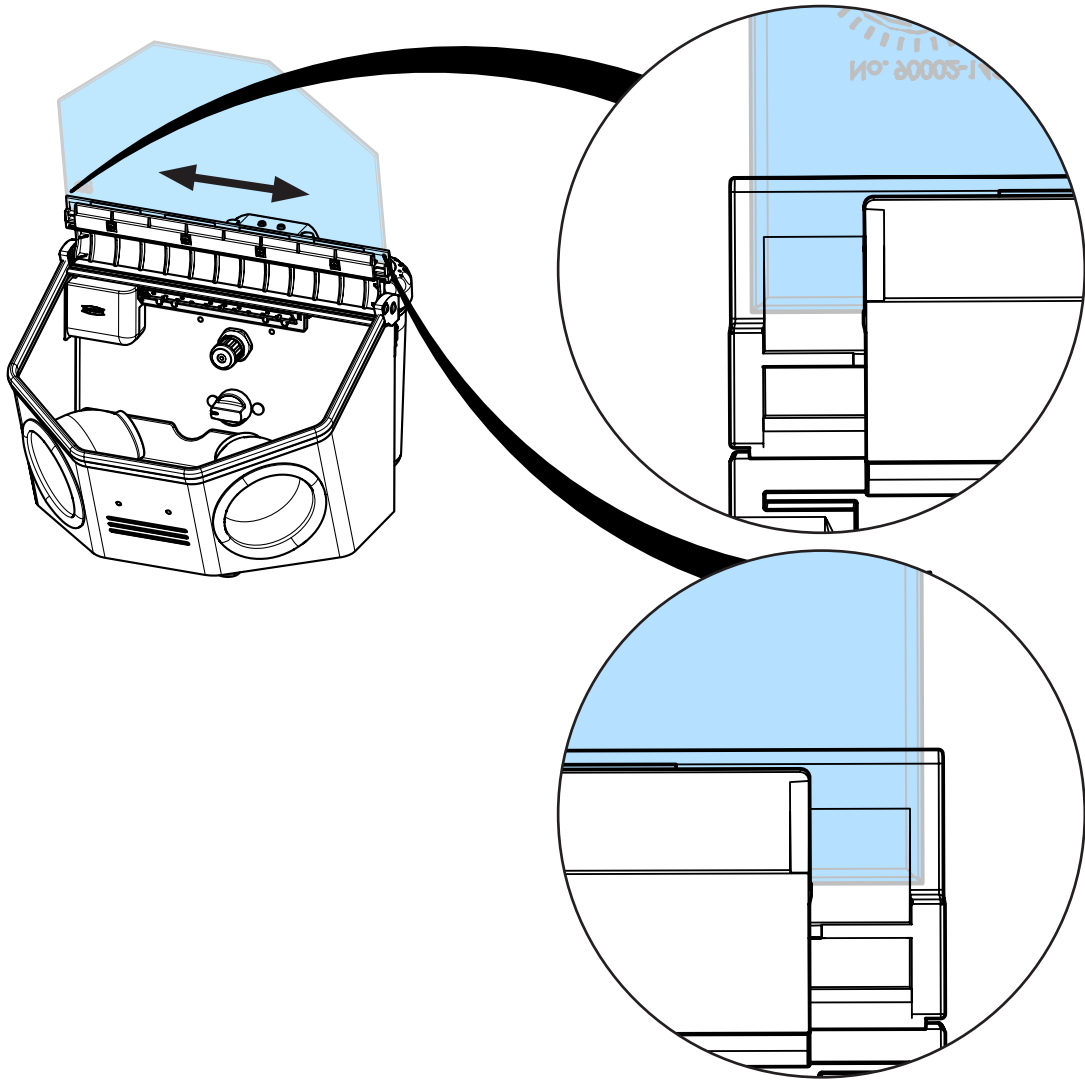


6.

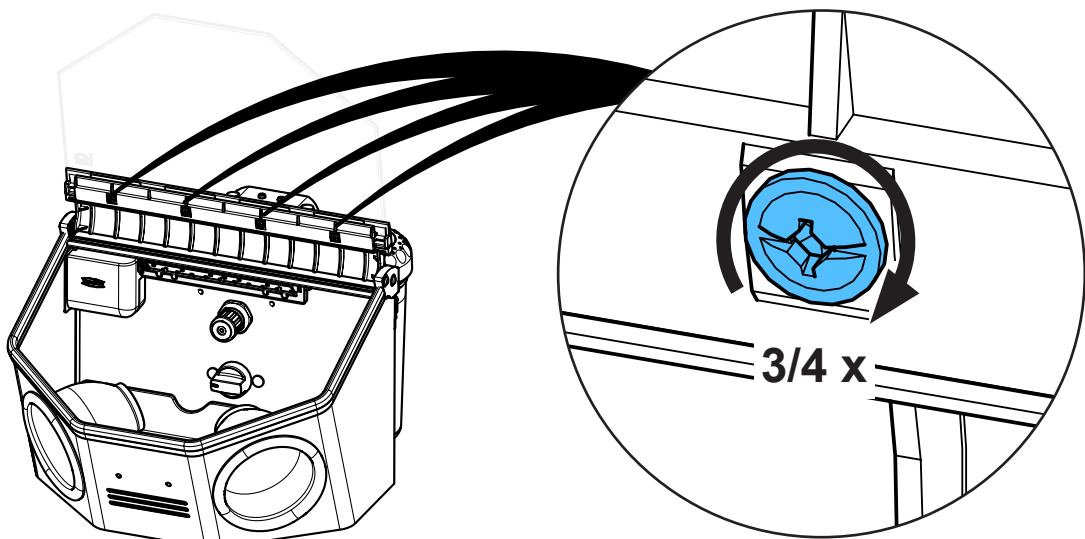

***i*** 1,2 - 1,4 Nm

7.


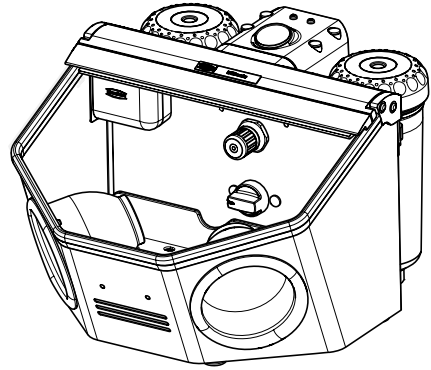
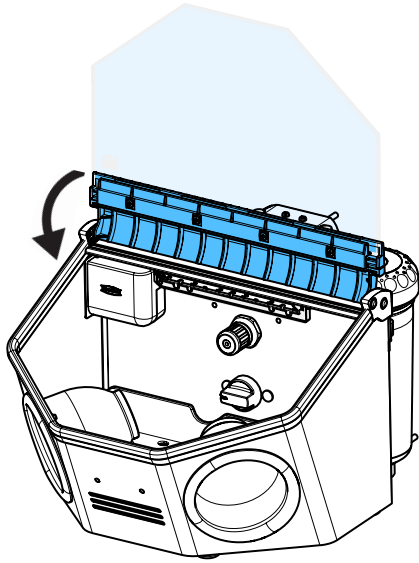





making work easy

90003 6311


Glasscheibe unbeschichtet • Uncoated glass screen • Vitre non revêtu • Vetro  
sen-za rivestimento • Cristal sin revestimiento • Запасное стекло без покрытия

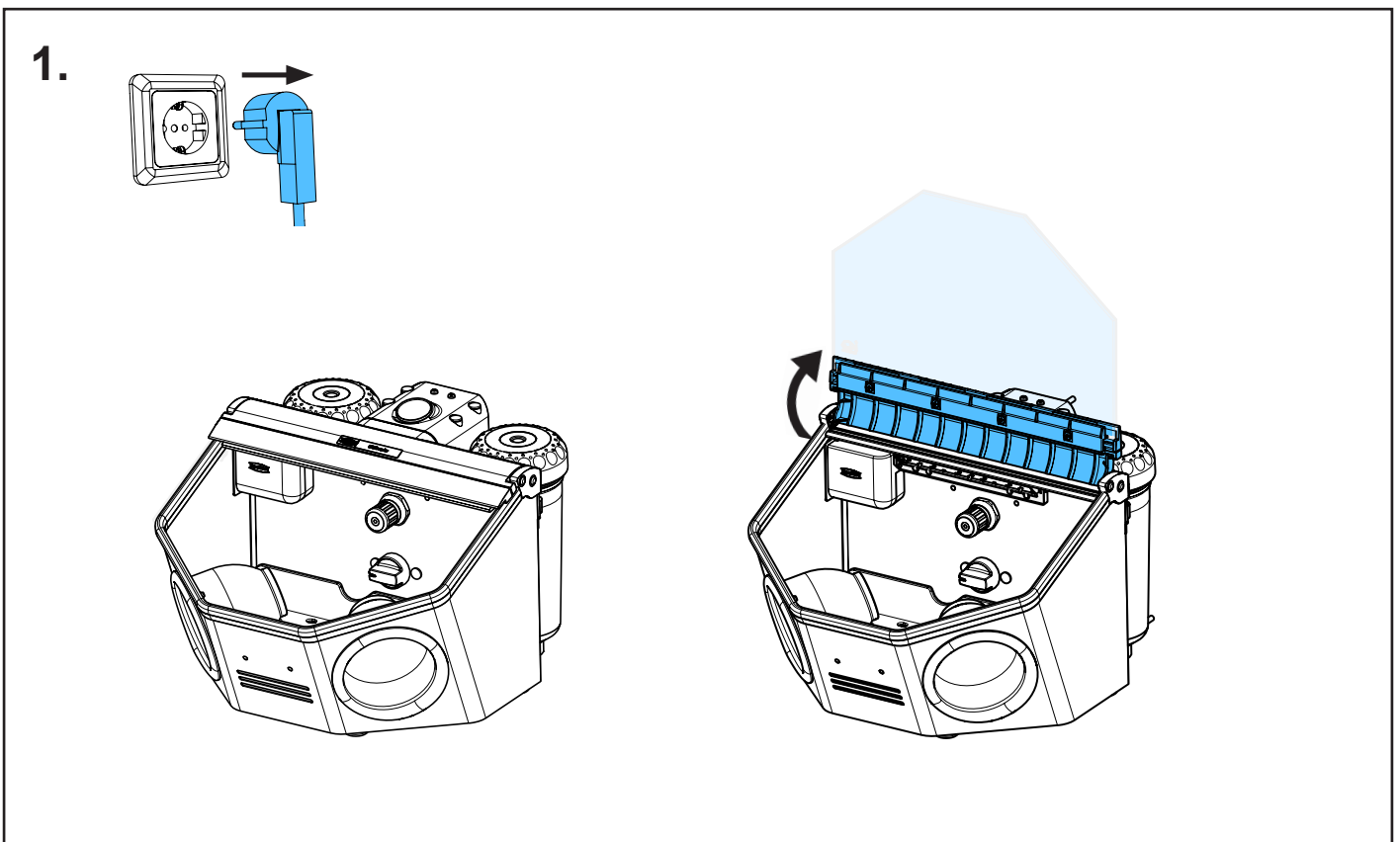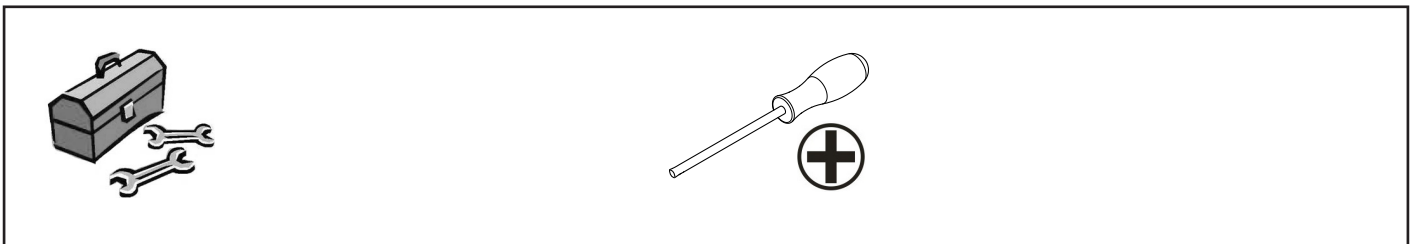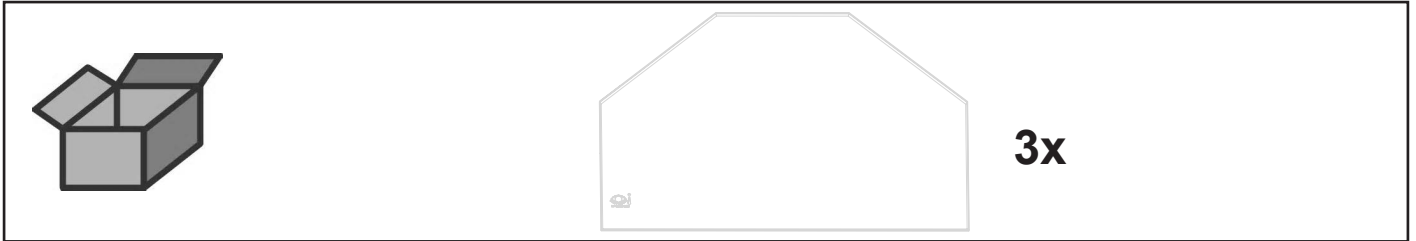


2.

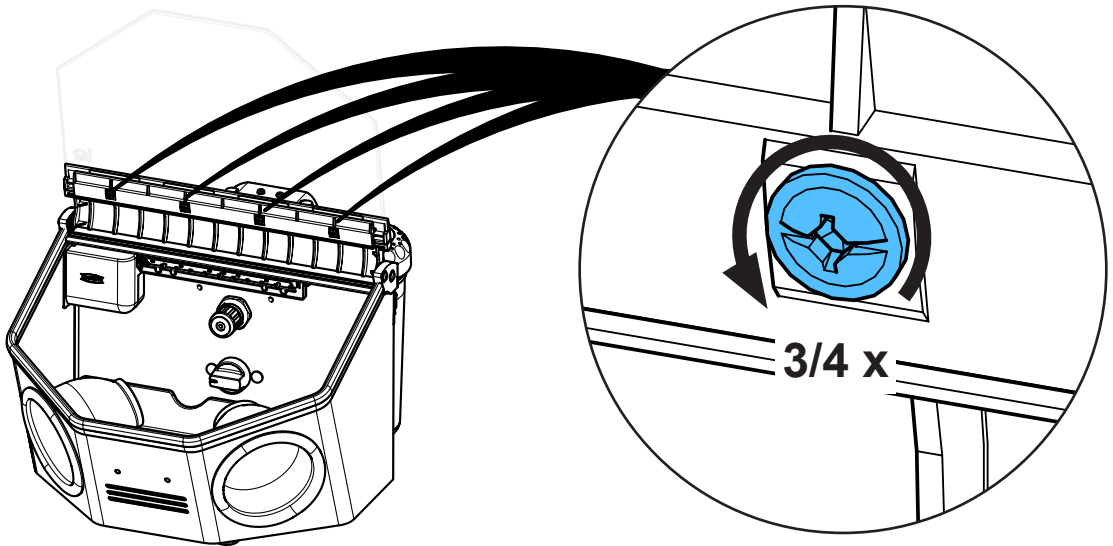


3.

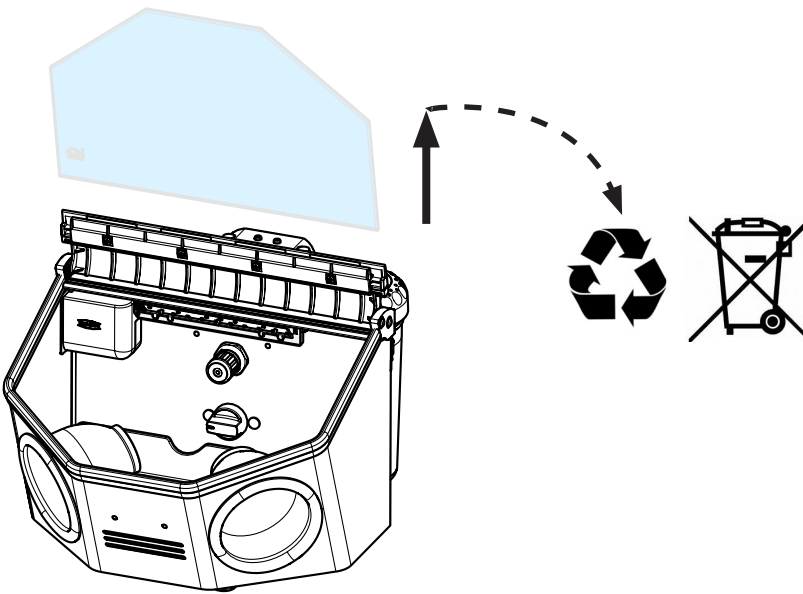


4.

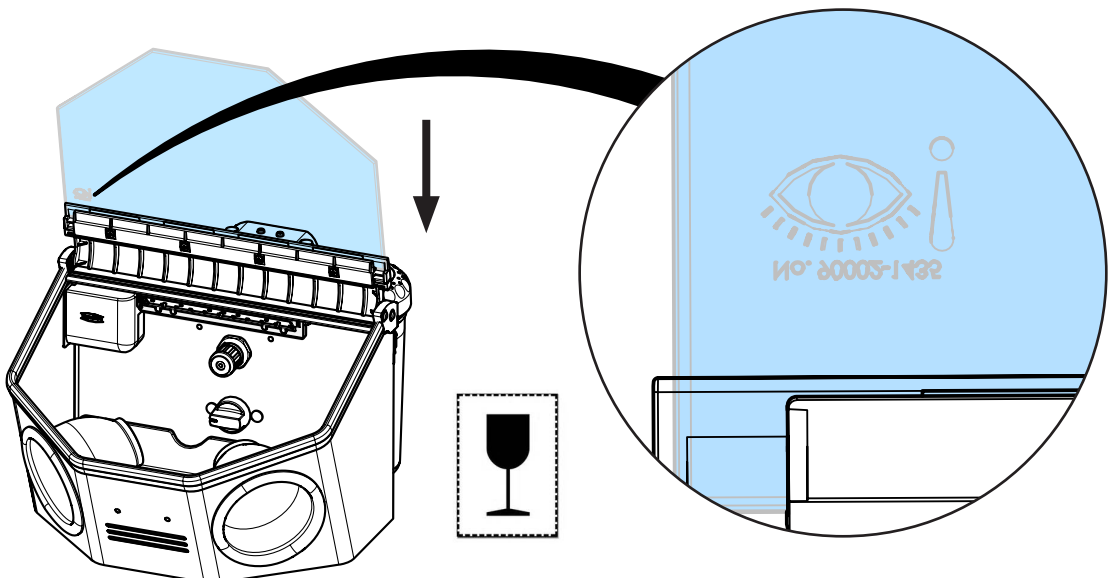

5.


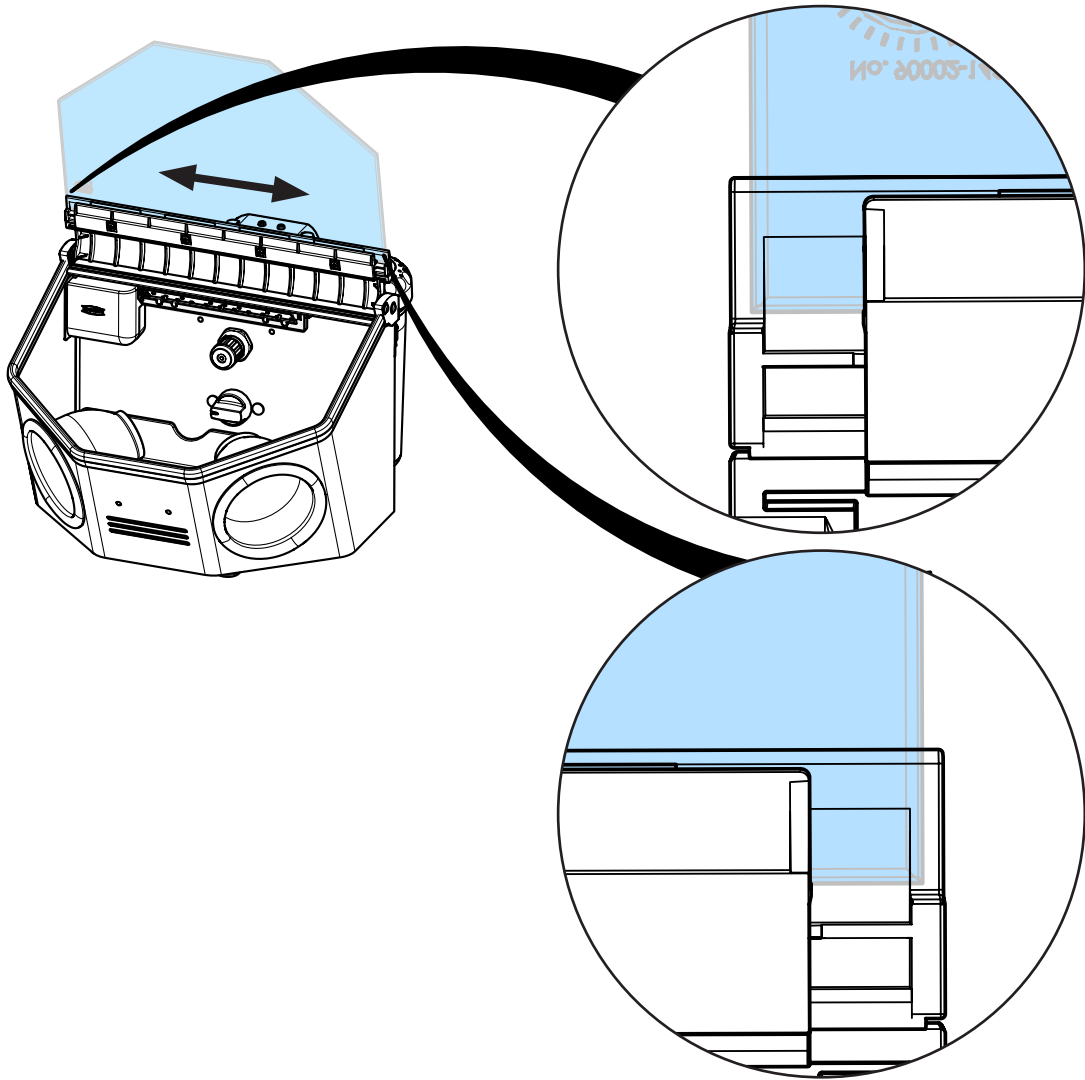


6.

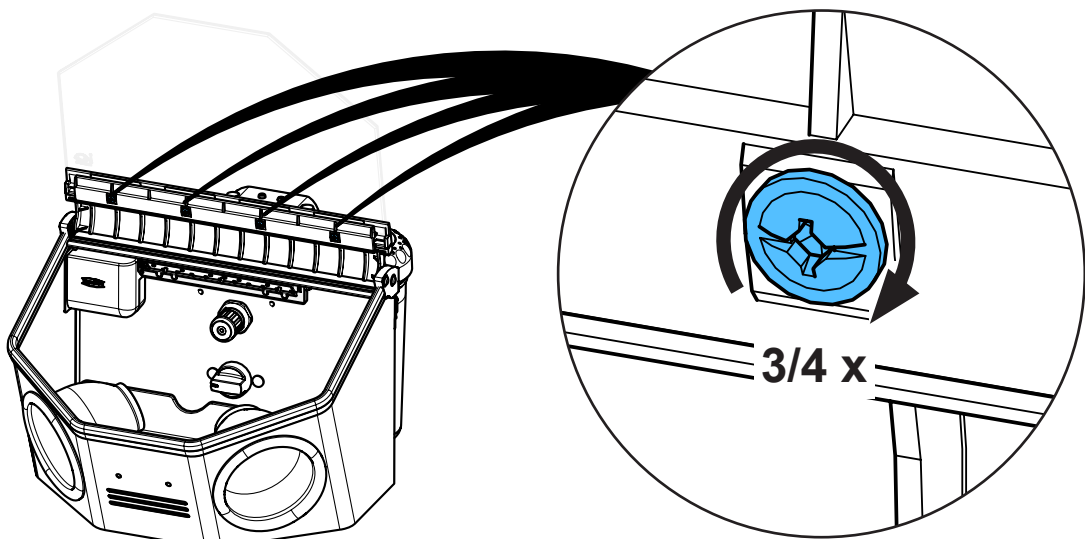

***i*** 1,2 - 1,4 Nm

7.

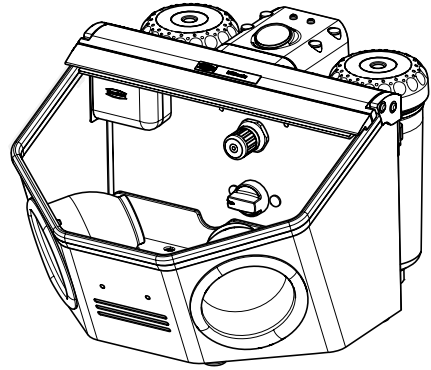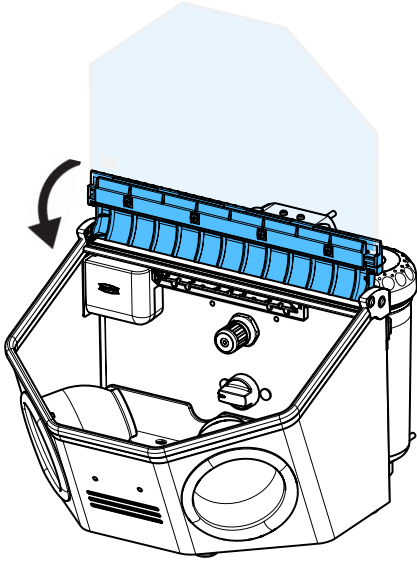





Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу


making work easy

92947 0001


Kantenschutzprofil • Edge protection • Bordure de protéction • Profilo protettivo  
Tira de protección • Краевой защитный профиль

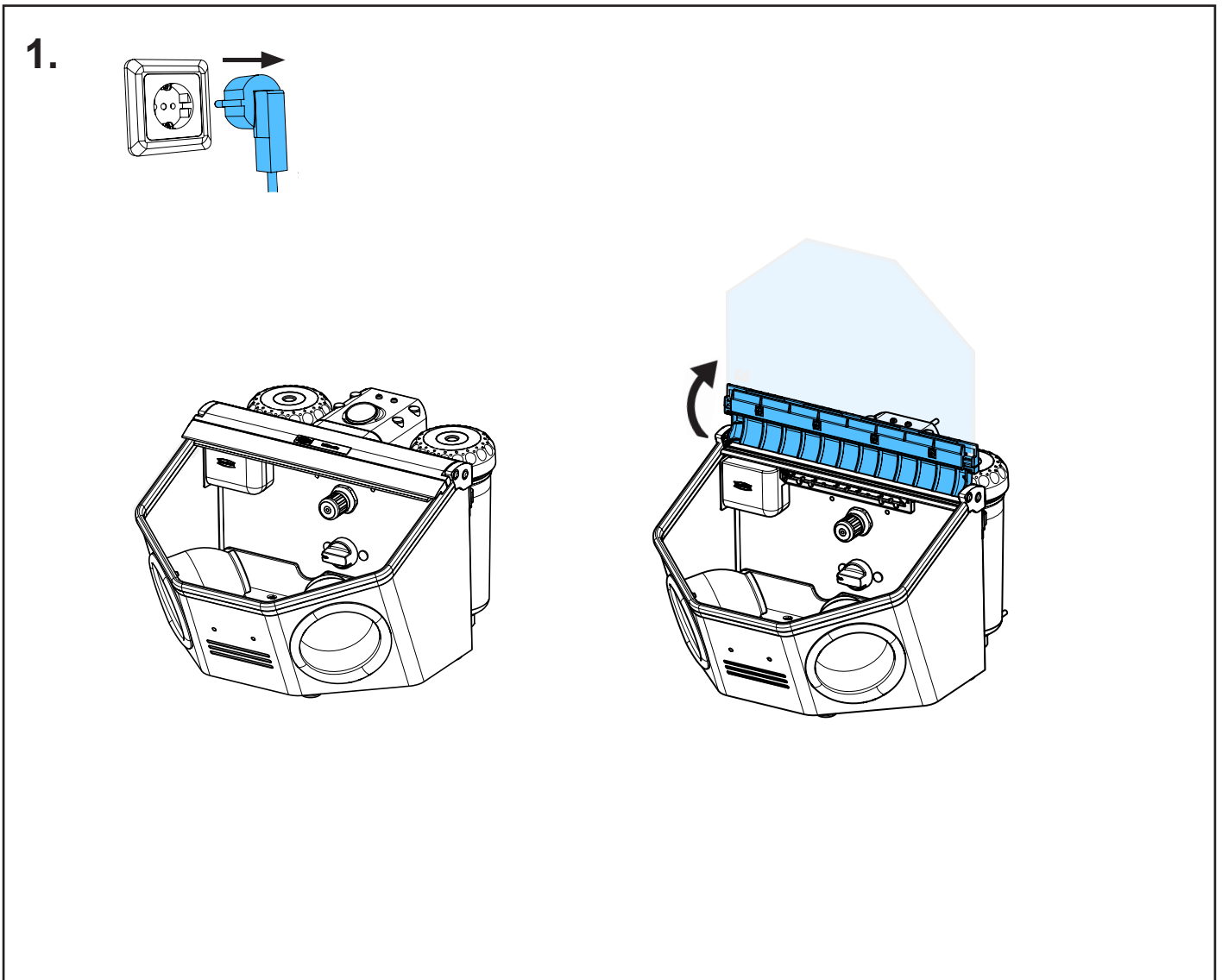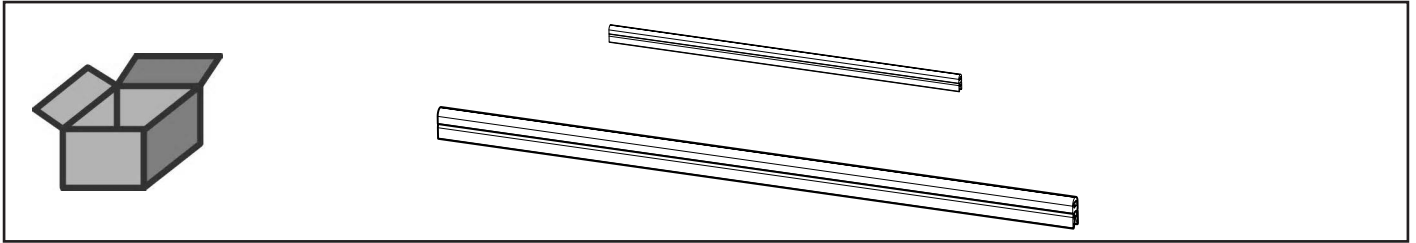


2.

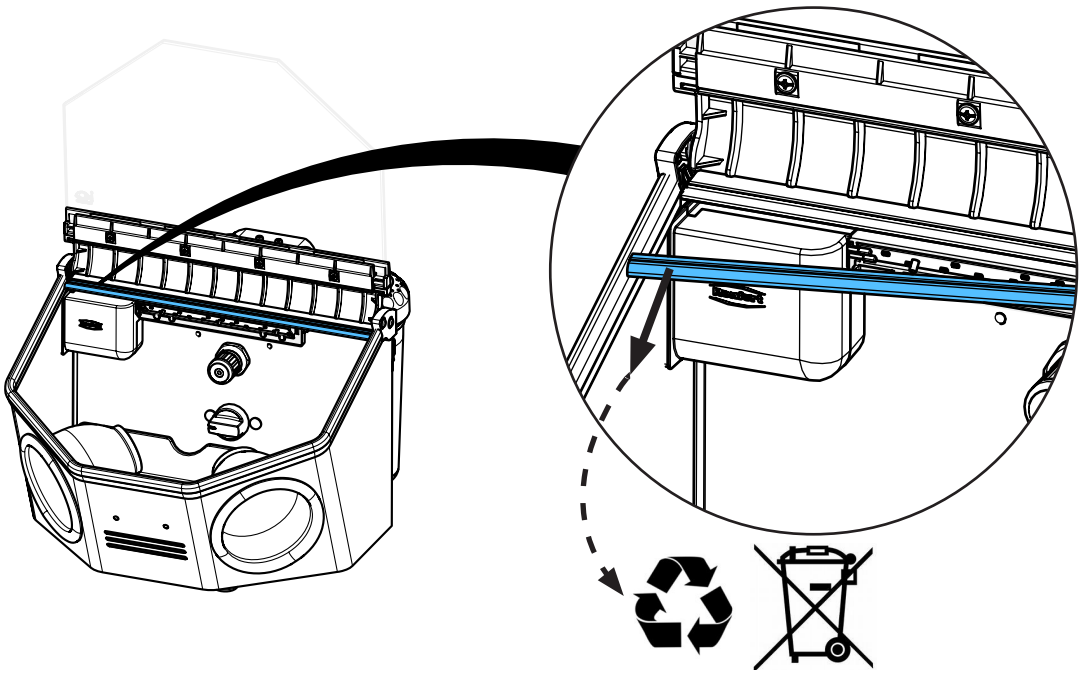

3.

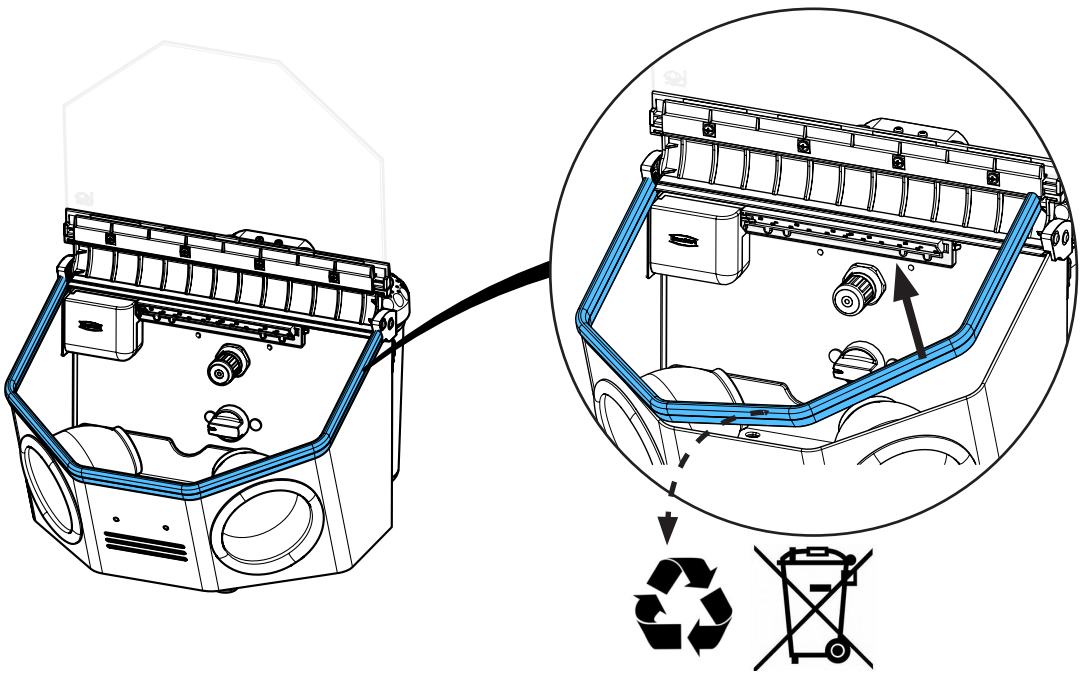


4.

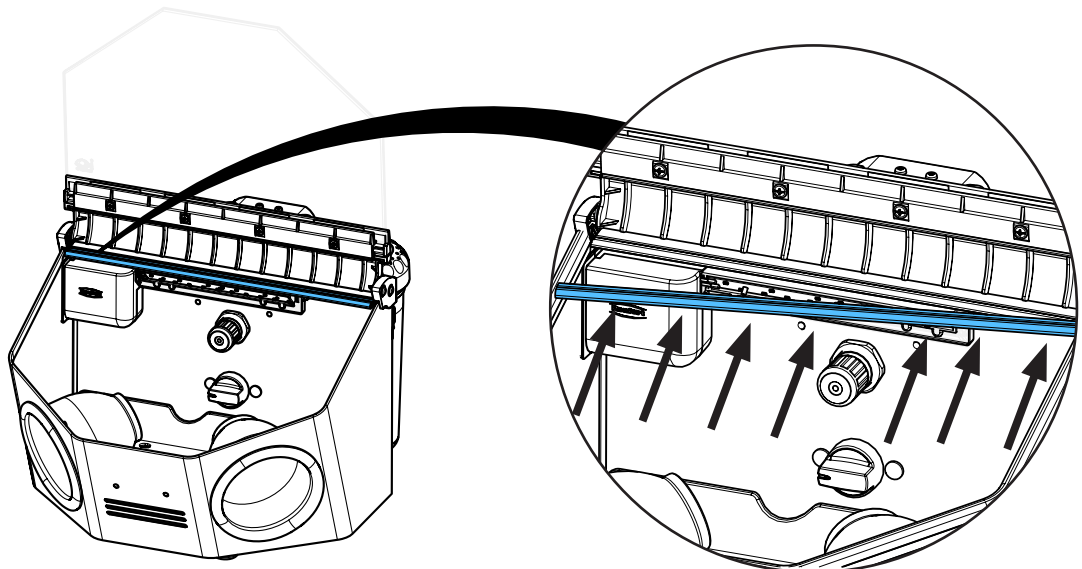






5.

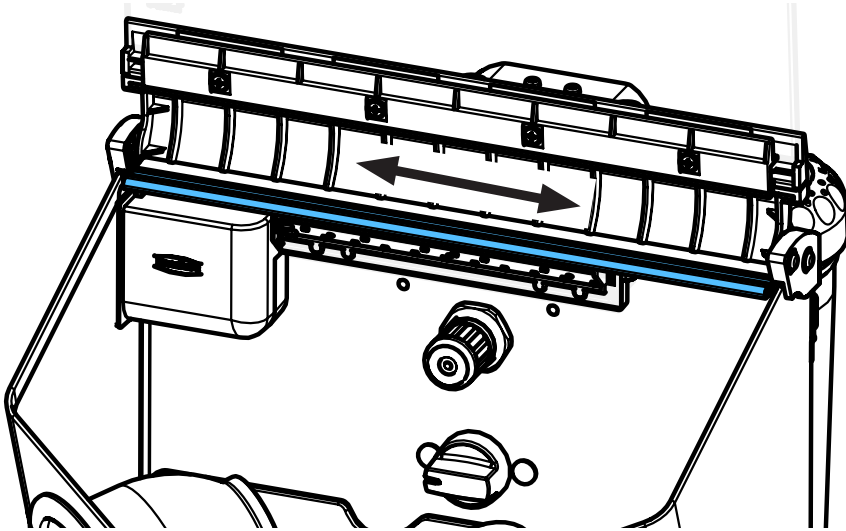

6.

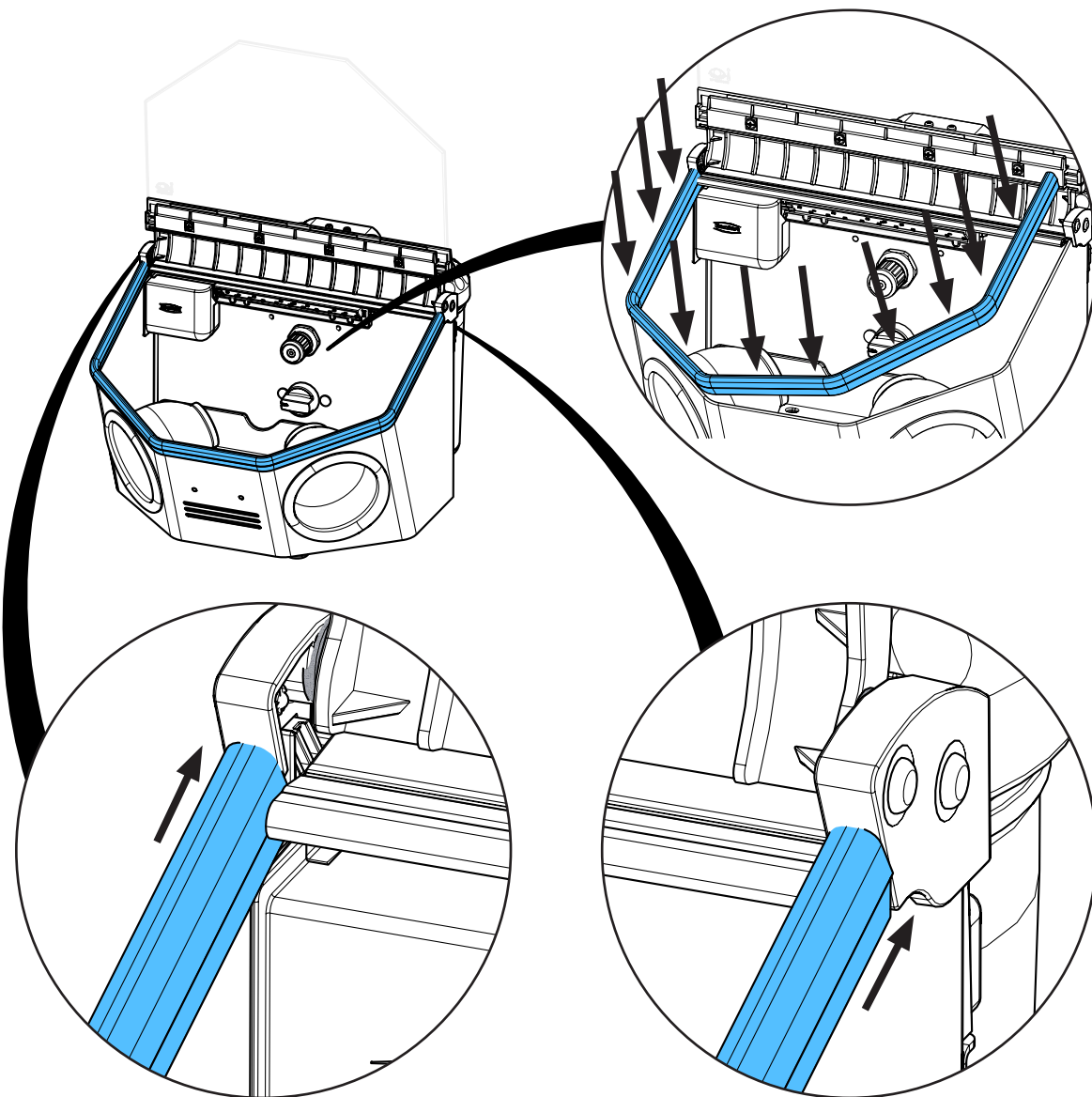

7.

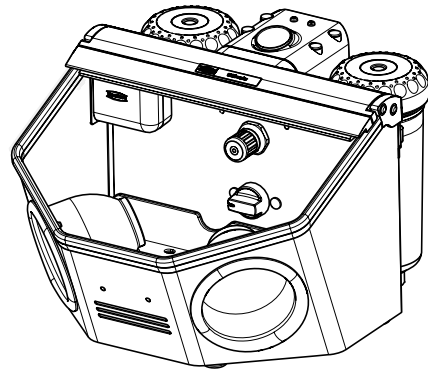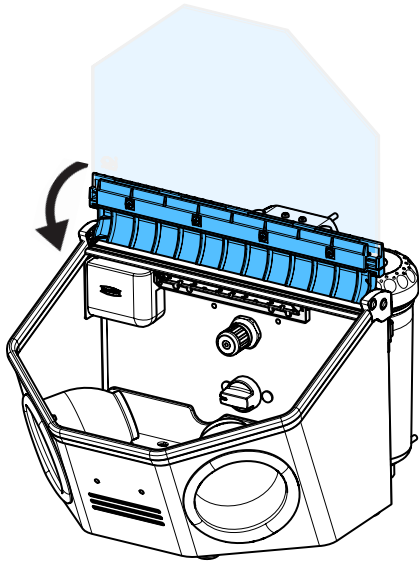





Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу



making work easy

90002 0921



Fußschalter pneumatisch • Pneumatic foot switch • Commande à pieds pneuma-  
tique • Comando a pedale pneumatico • Pedal interruptor neumático • Педаль  
пневматическая

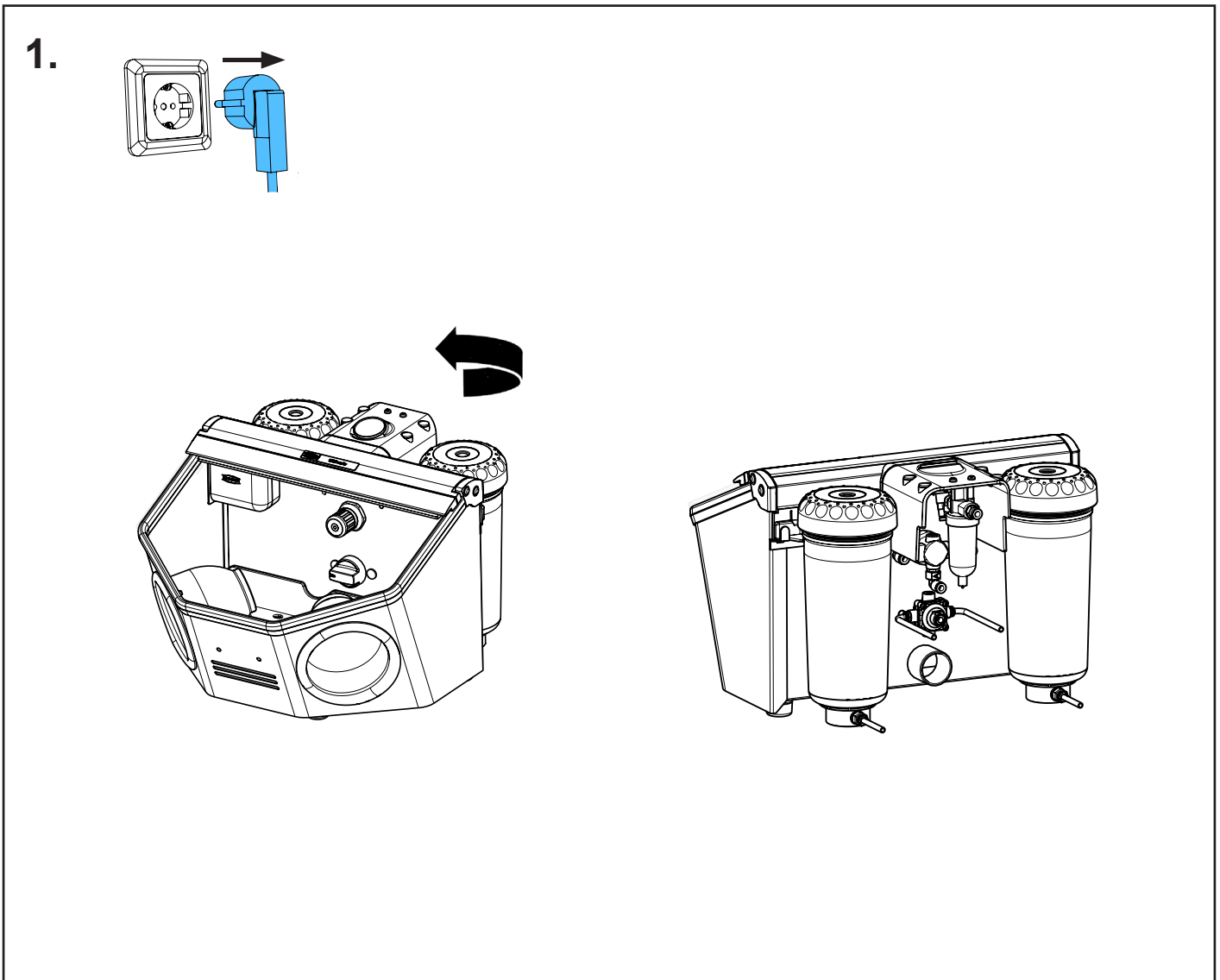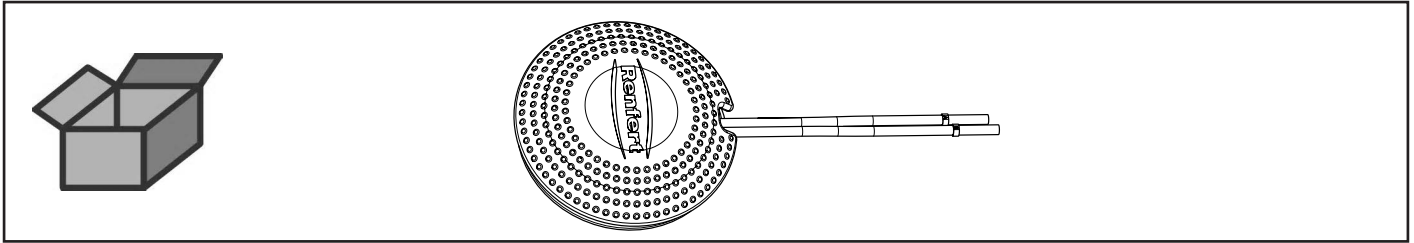



2.

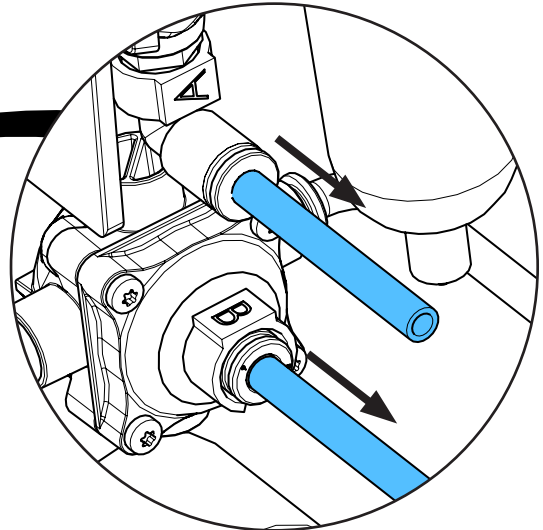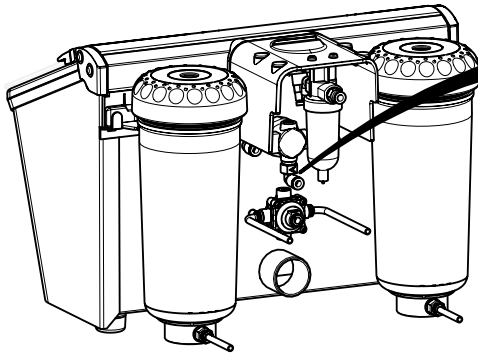


**i**

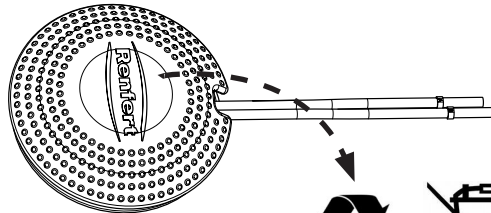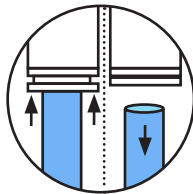

3.

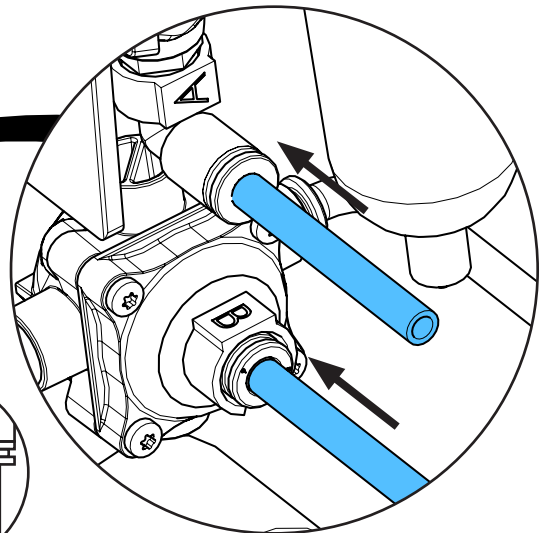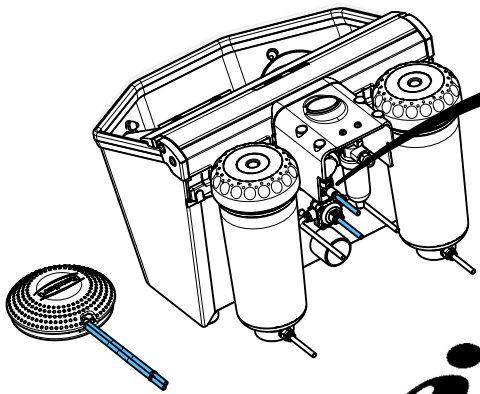


**i**

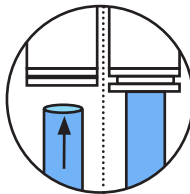

**!**

A - A  
B - B

4.

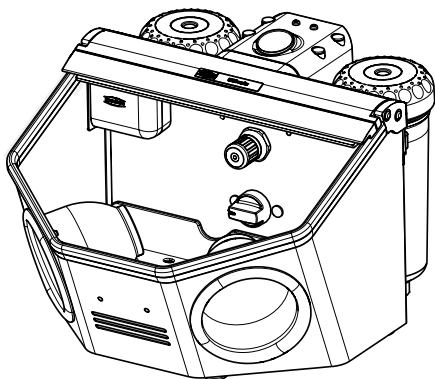






Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу


making work easy

90003 6598


EingangsfILTER • Inlet filter • Filtre d'entrée • Filtro d'entrata • Filtro de entrada  
входной фильТР

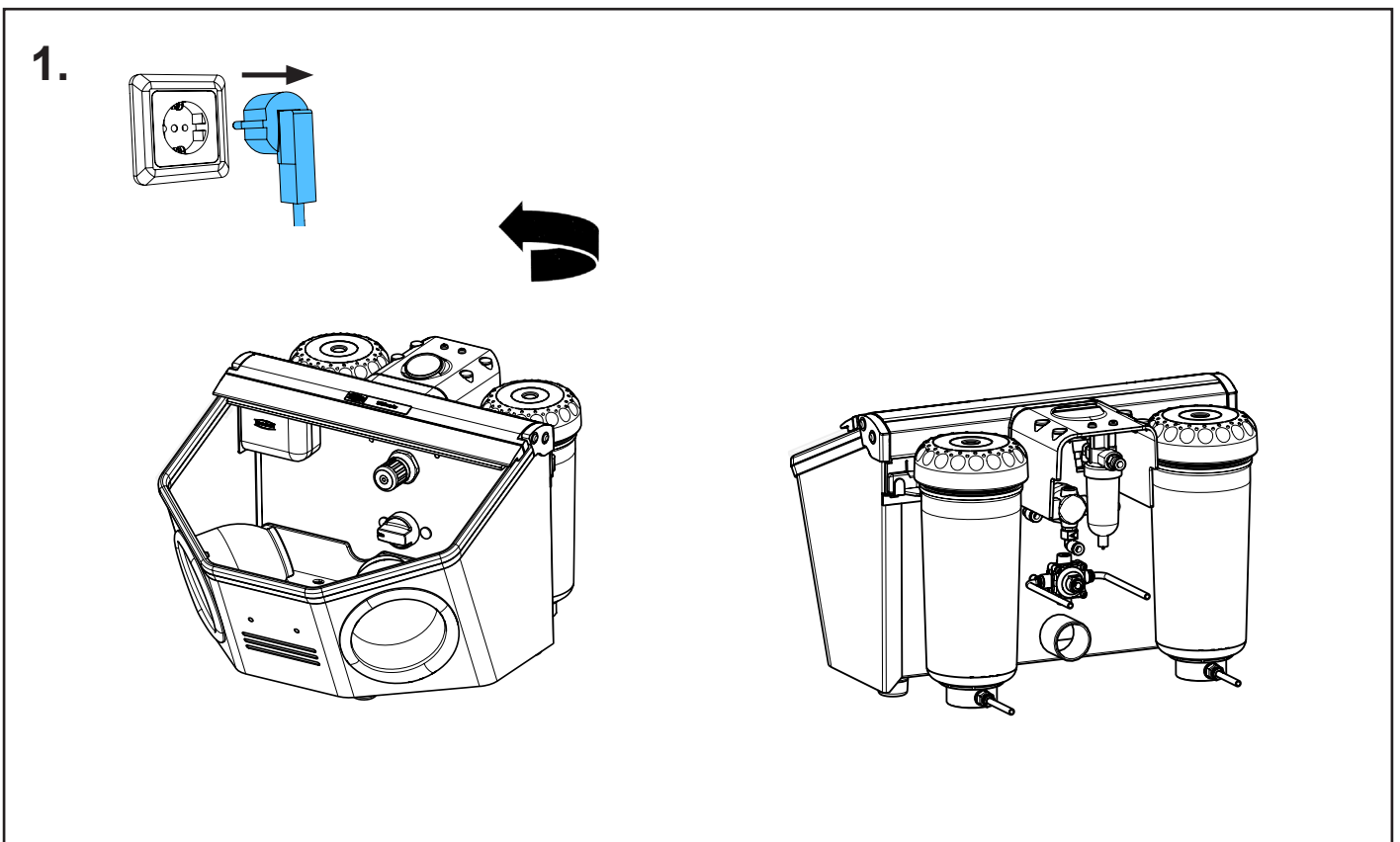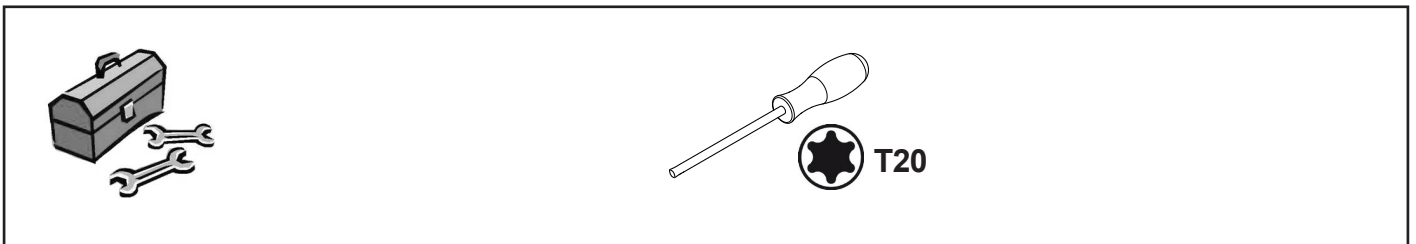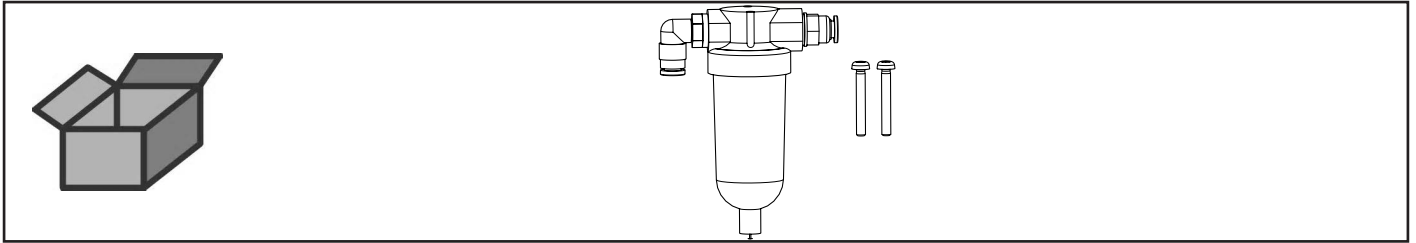




2.

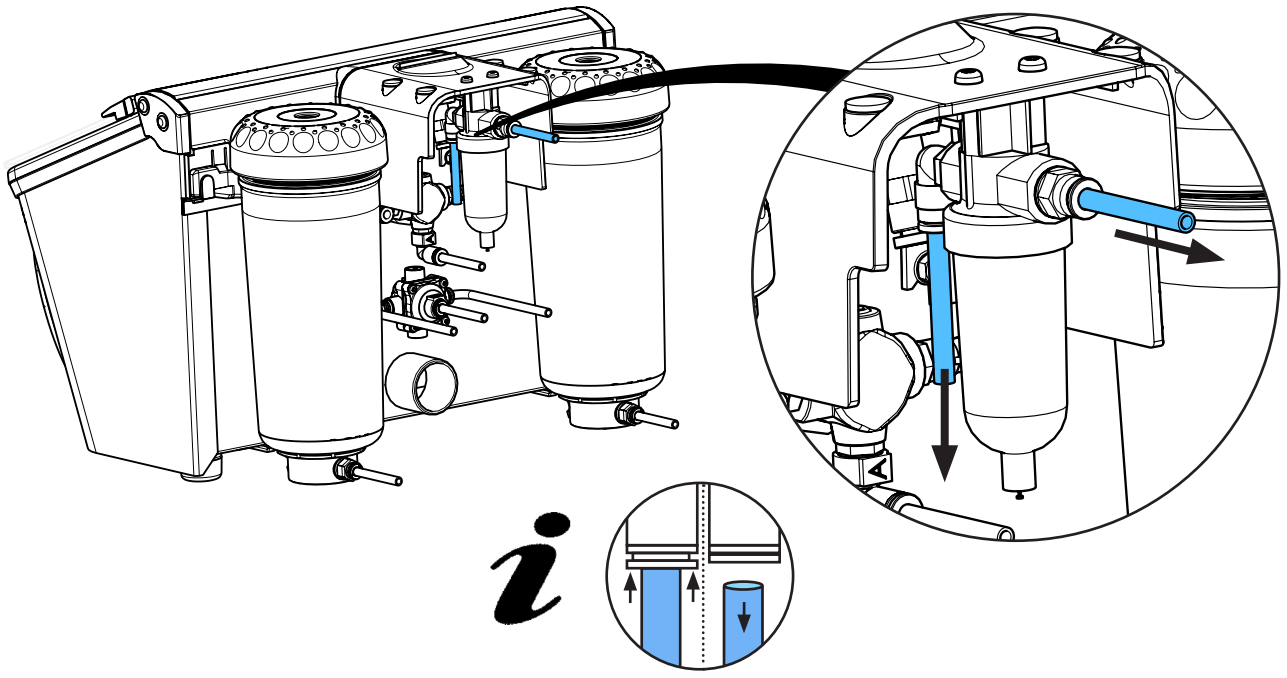

3.


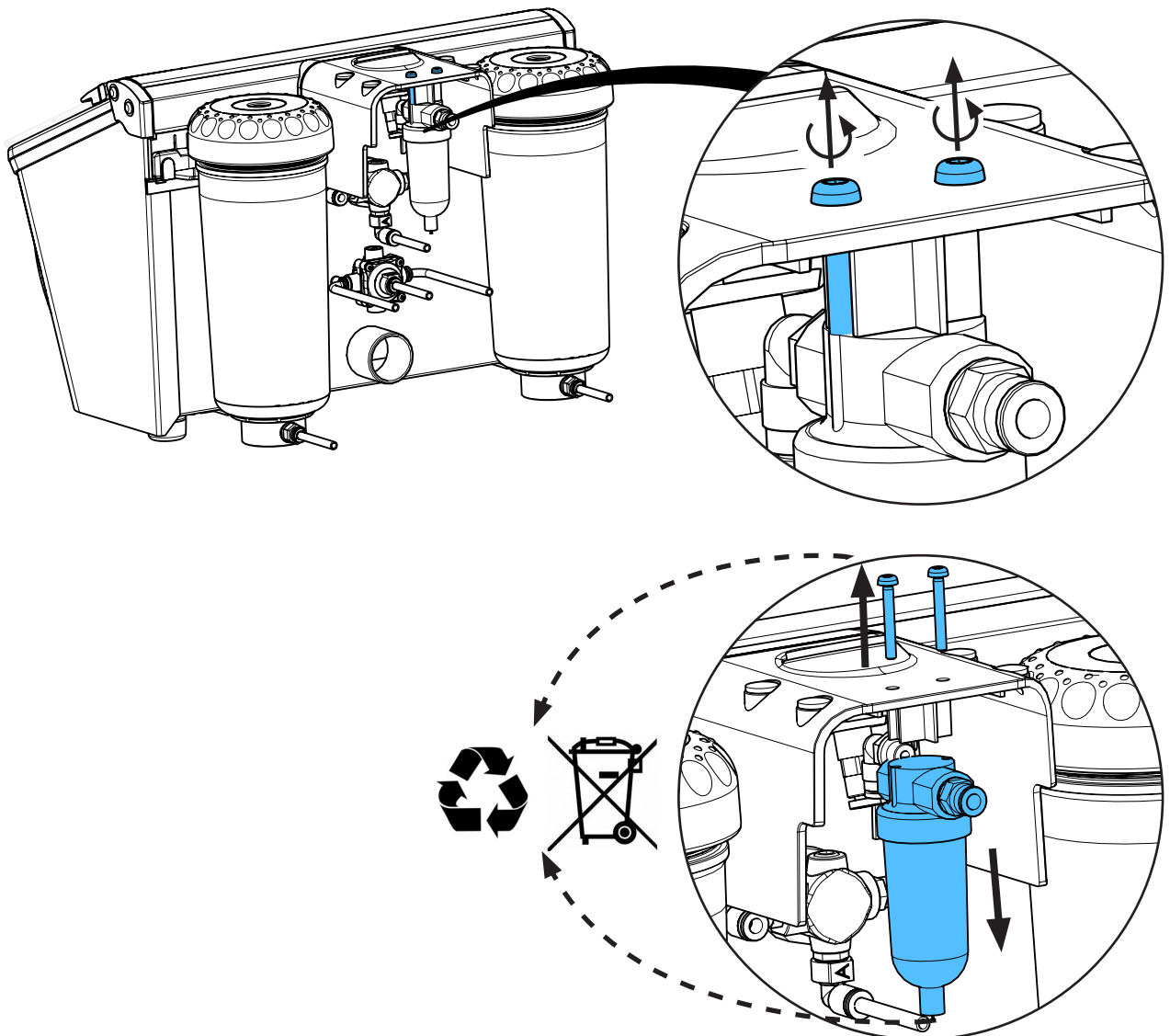


4.

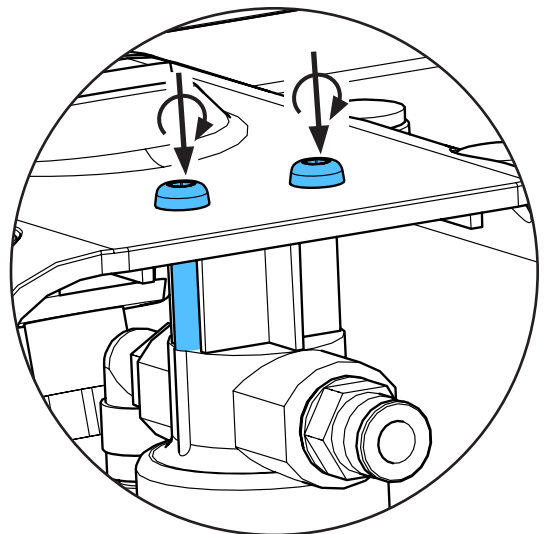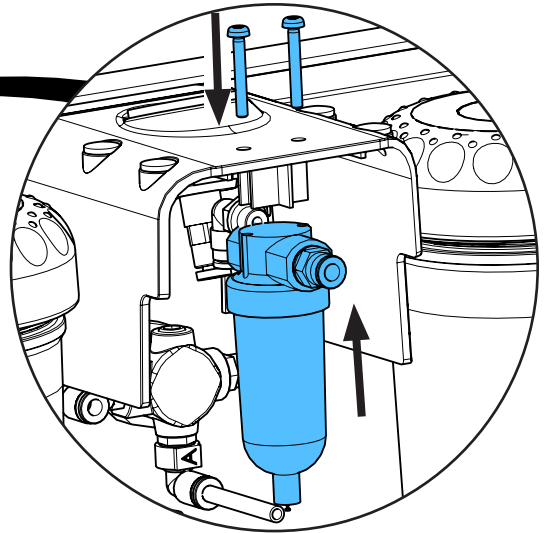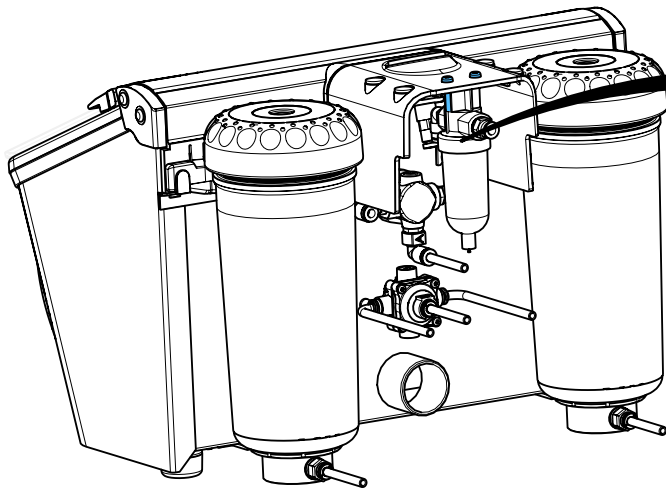


1,2 - 1,4 Nm


5.

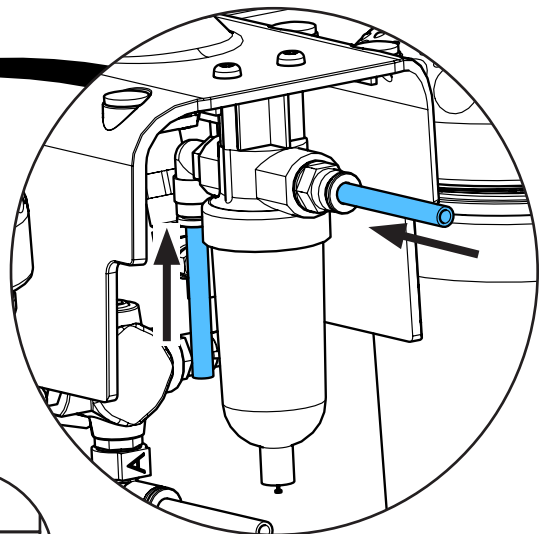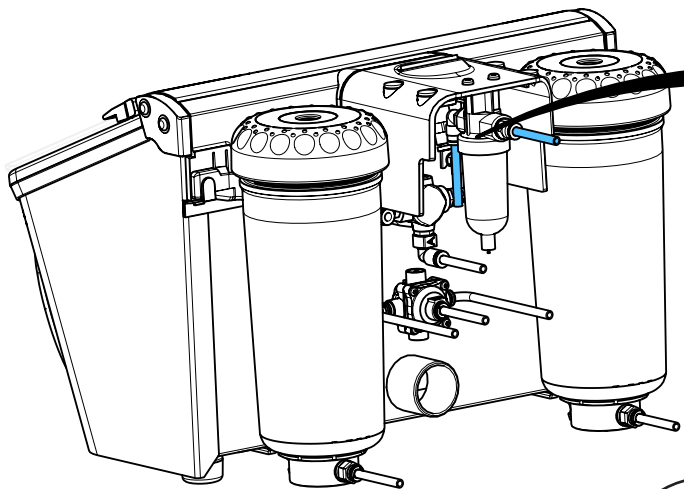

***i***

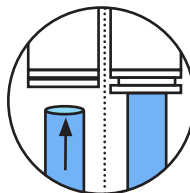

6.

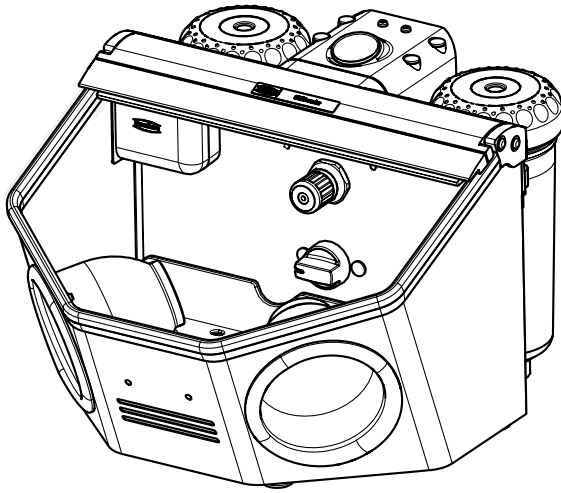







Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу


making work easy

**90003 4104**


Belüftungsklappe • Ventilating flap • Clapet de ventilation • Copertura per foro di  
aerazione • Chapaleta de ventilación • Вентиляционный клапан

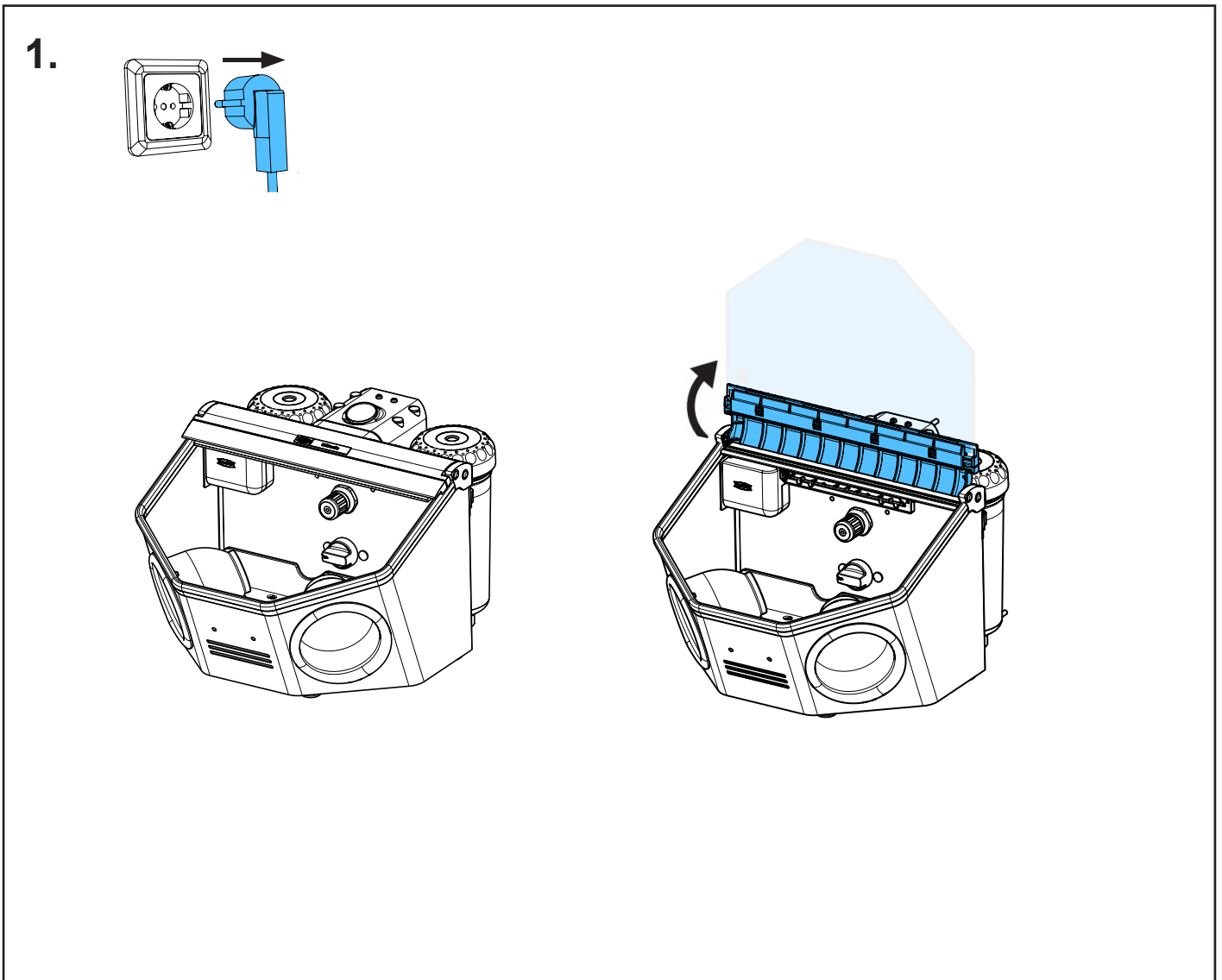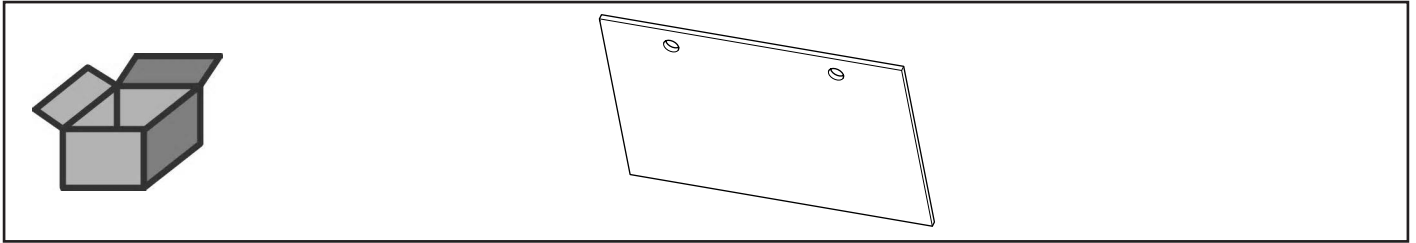


2.

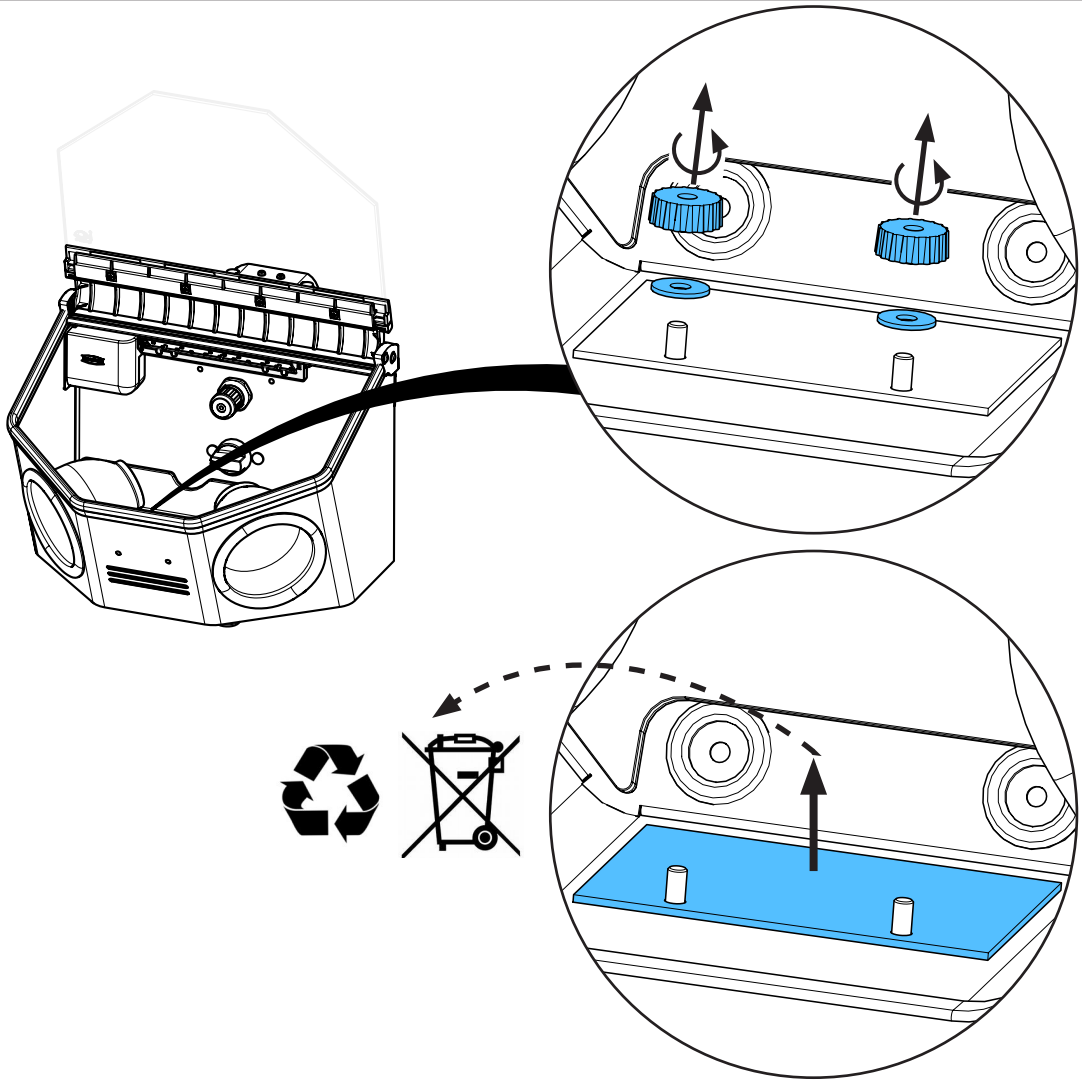

3.

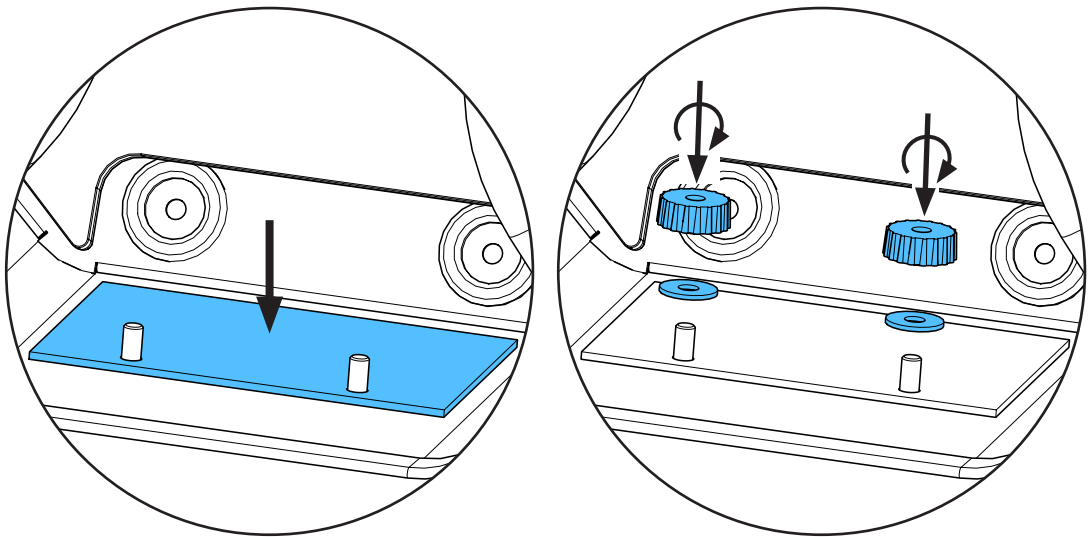

4.

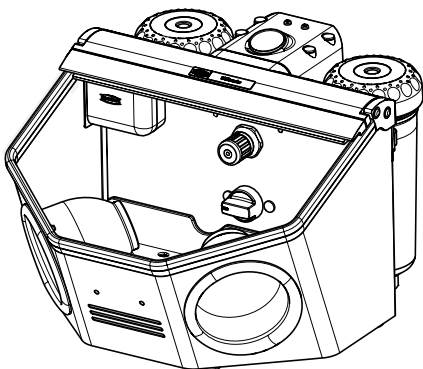






Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по монтажу


making work easy

## 92900 0050 / 92900 0500


IT Strahlschlauch inkl. Handstück • IT blasting hose incl. handpiece • IT tuyau de sablage avec pièce á main • IT tubo incl. manipolo • IT tubo incl. manipolo • Струйный шланг IT

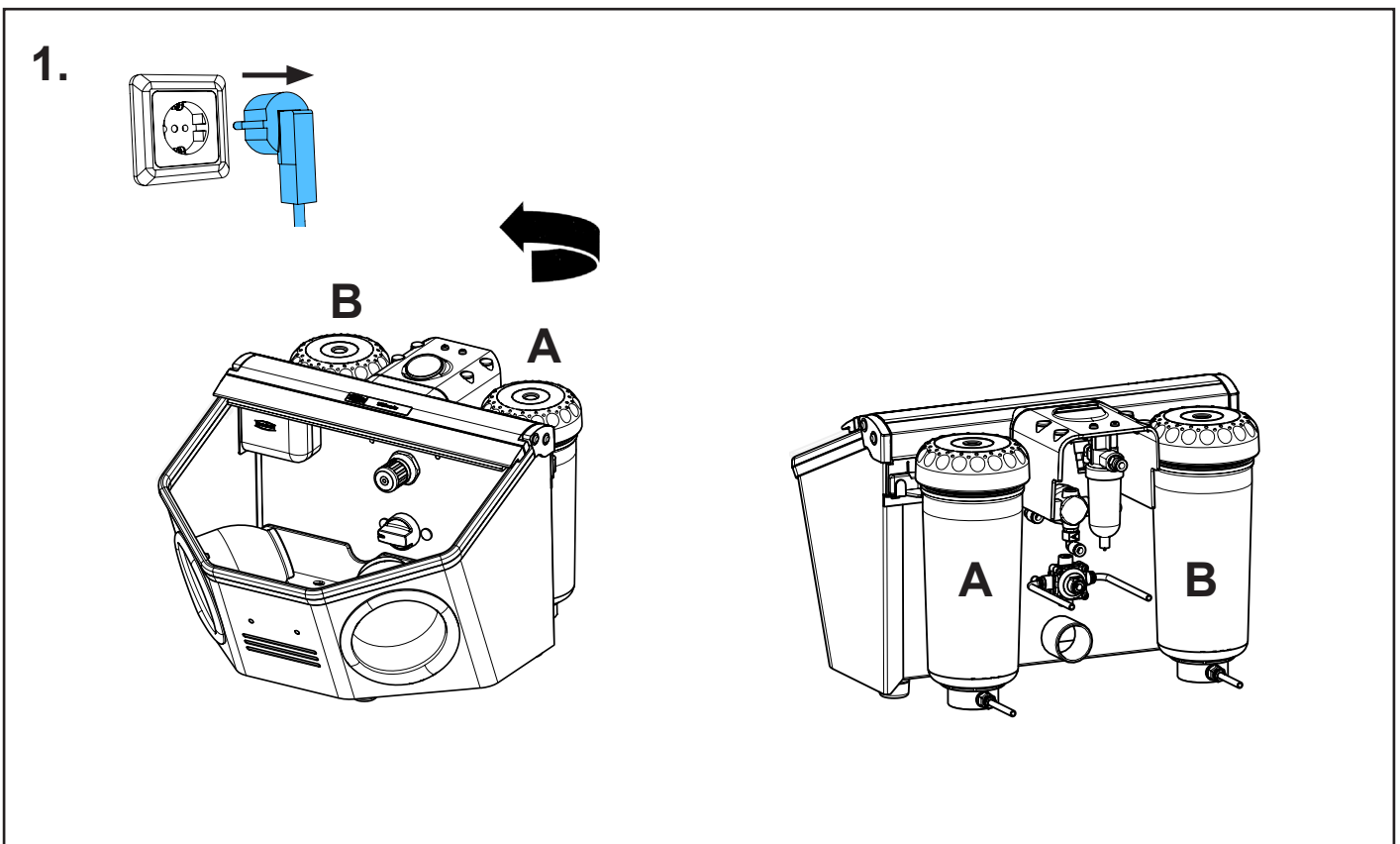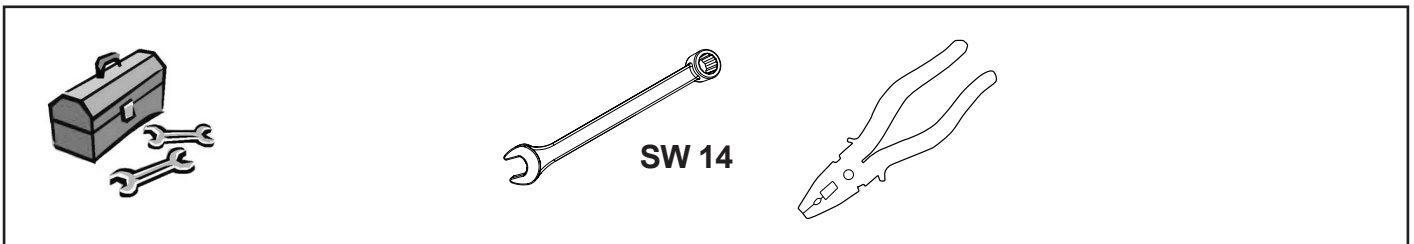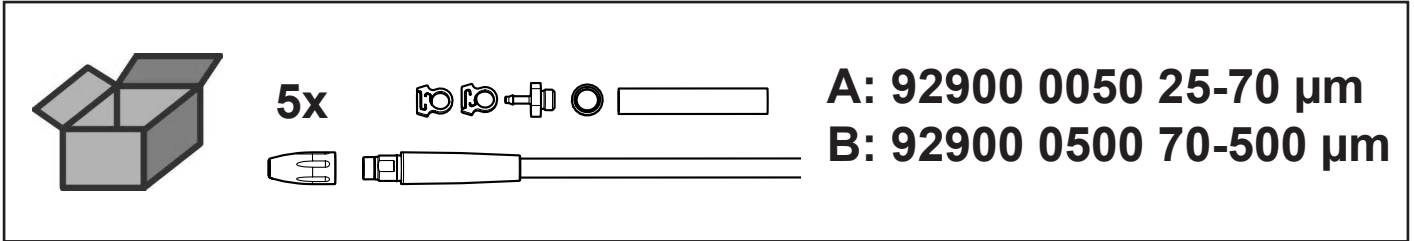


2.

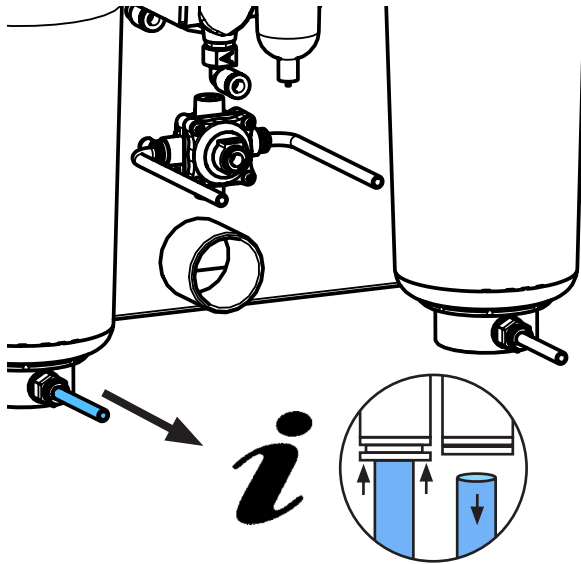


3.

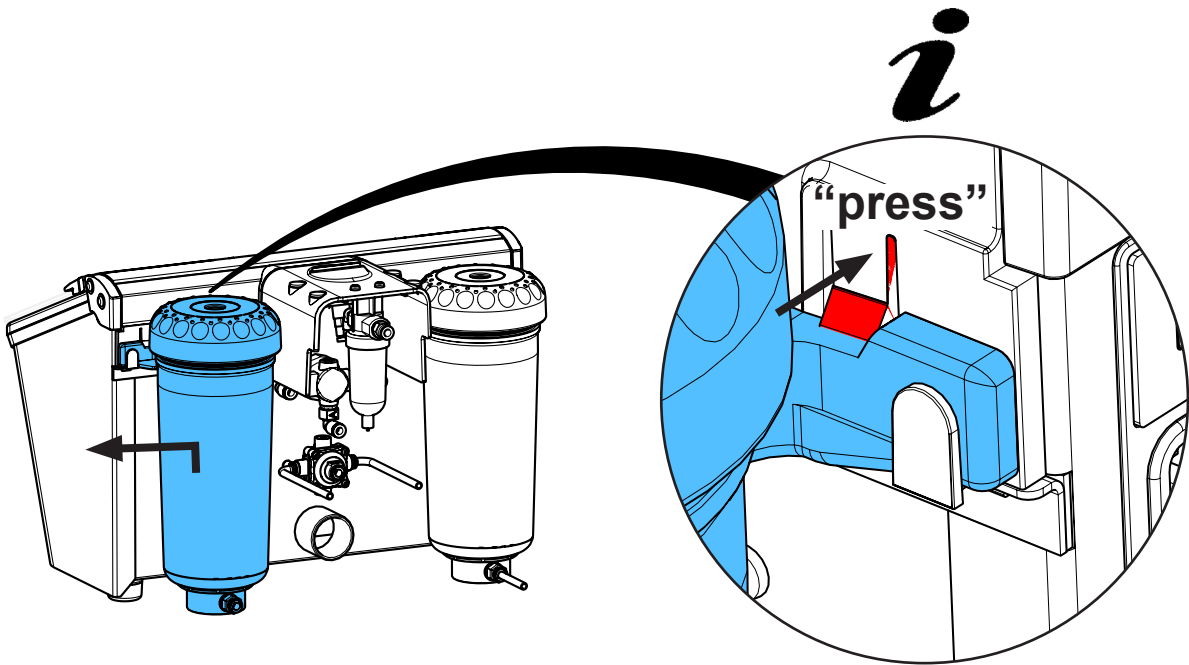


4.

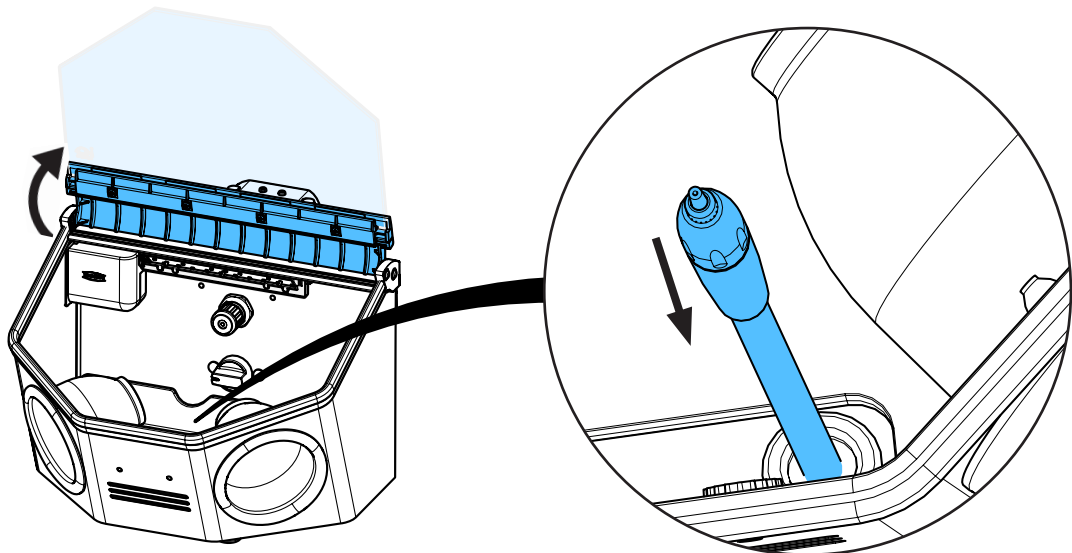


5.

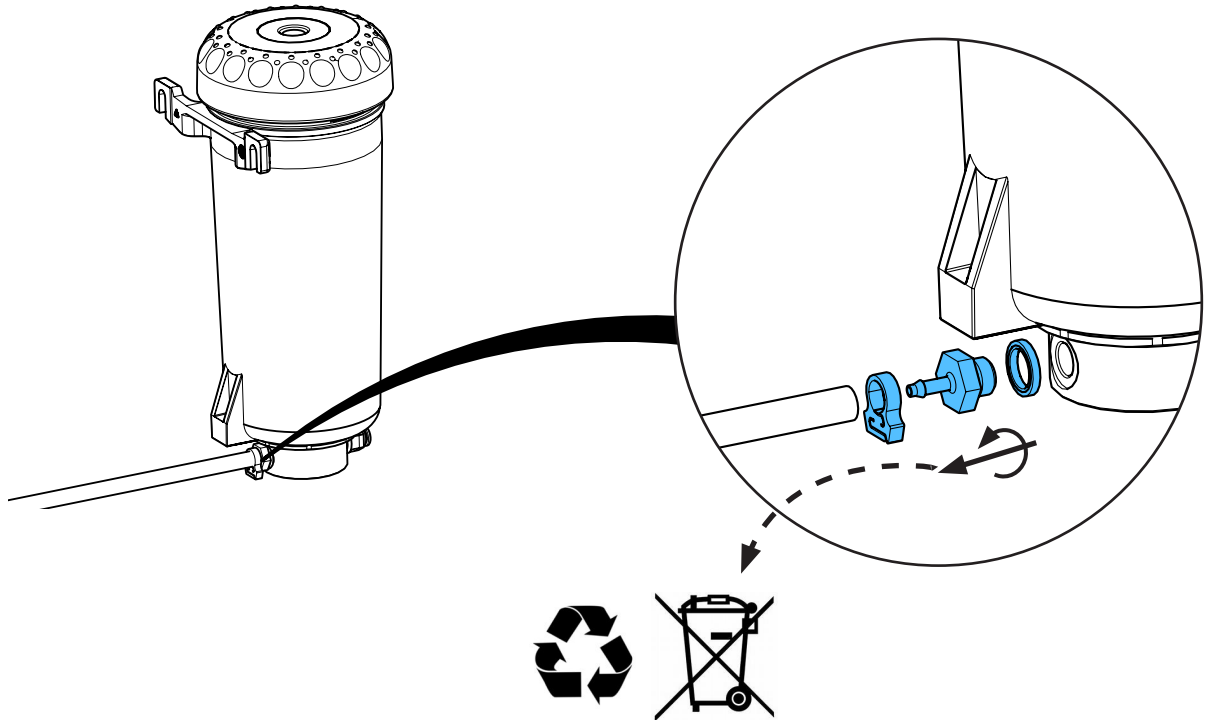


6.

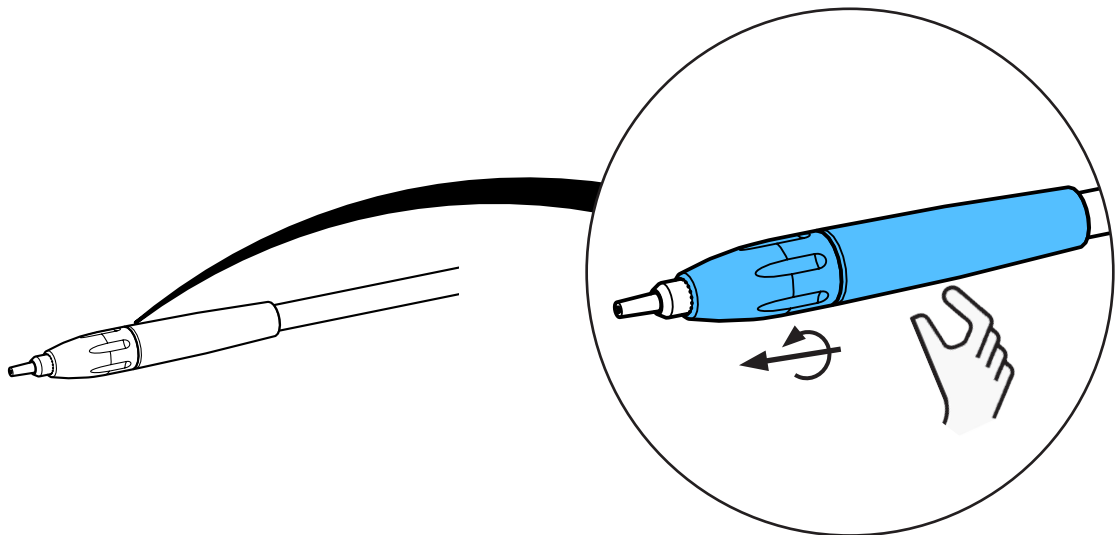


7.

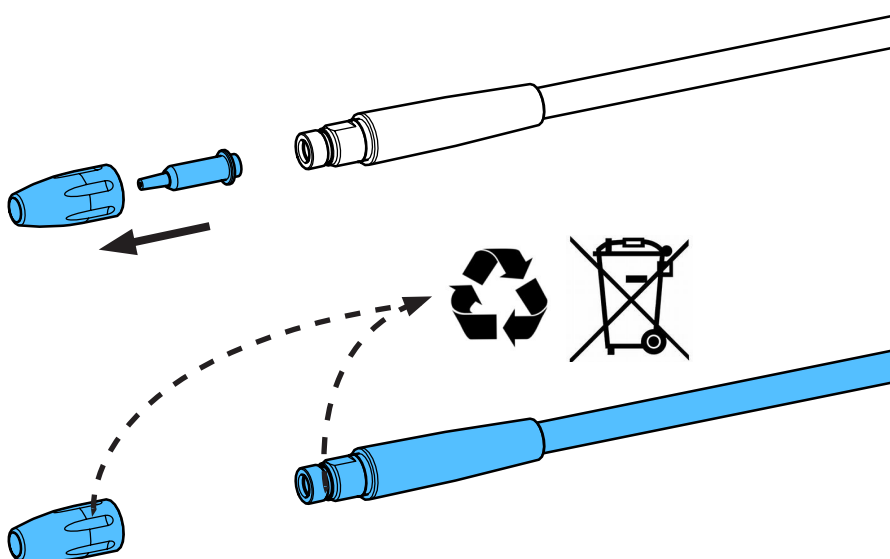

8.

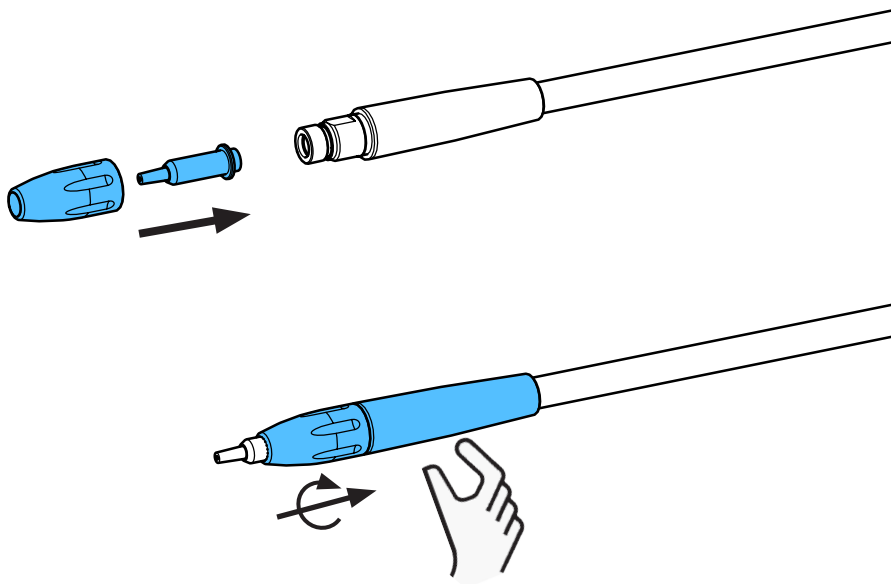


9.

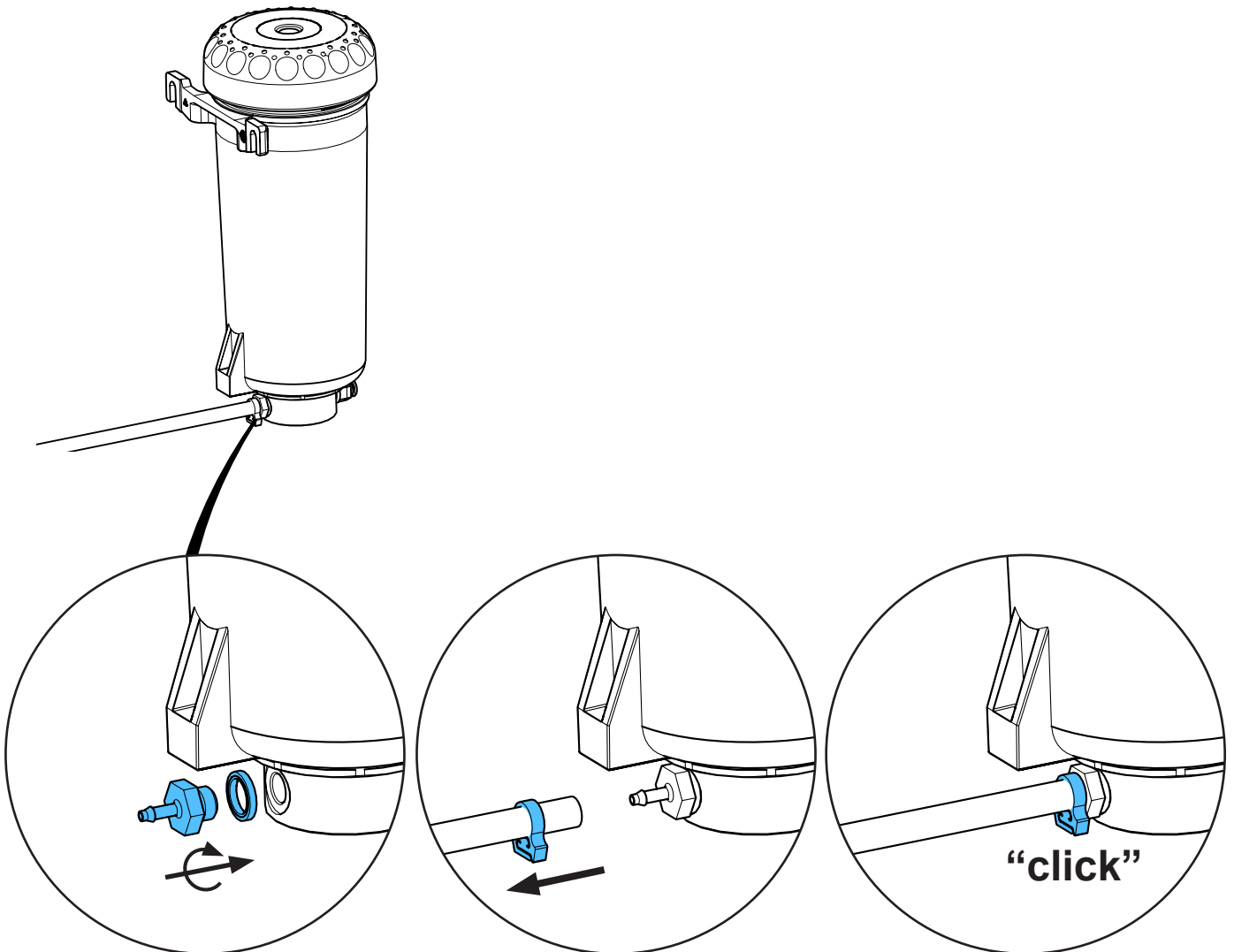


3,3 - 3,7 Nm


10.

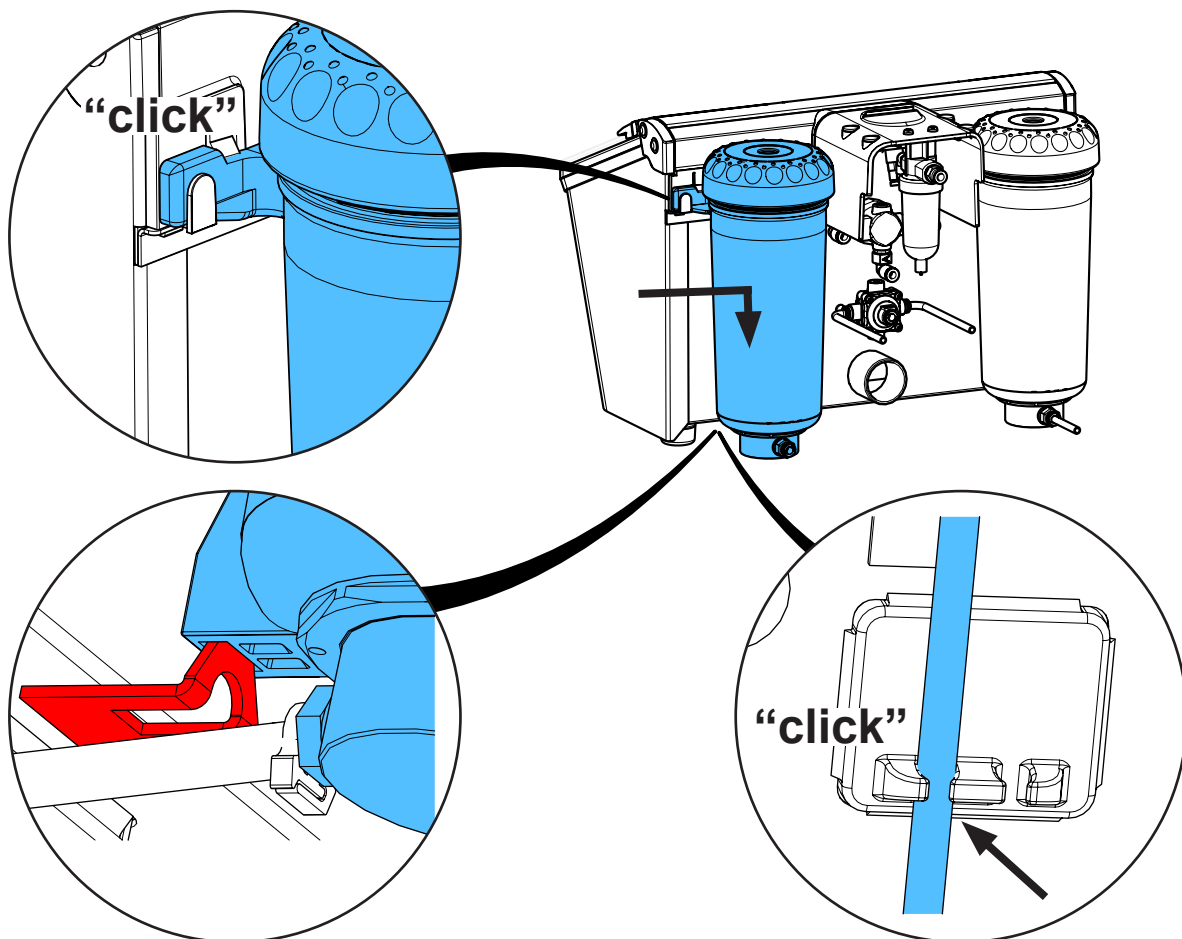


11.

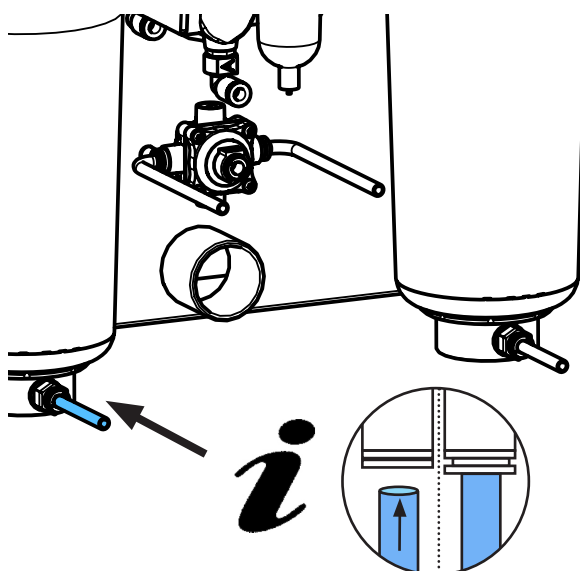

12.

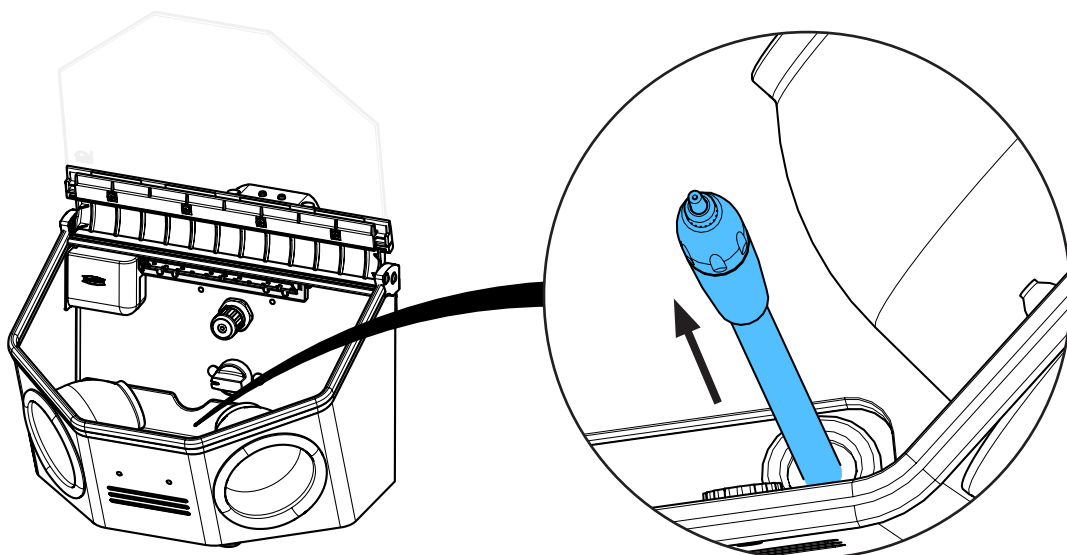

13.

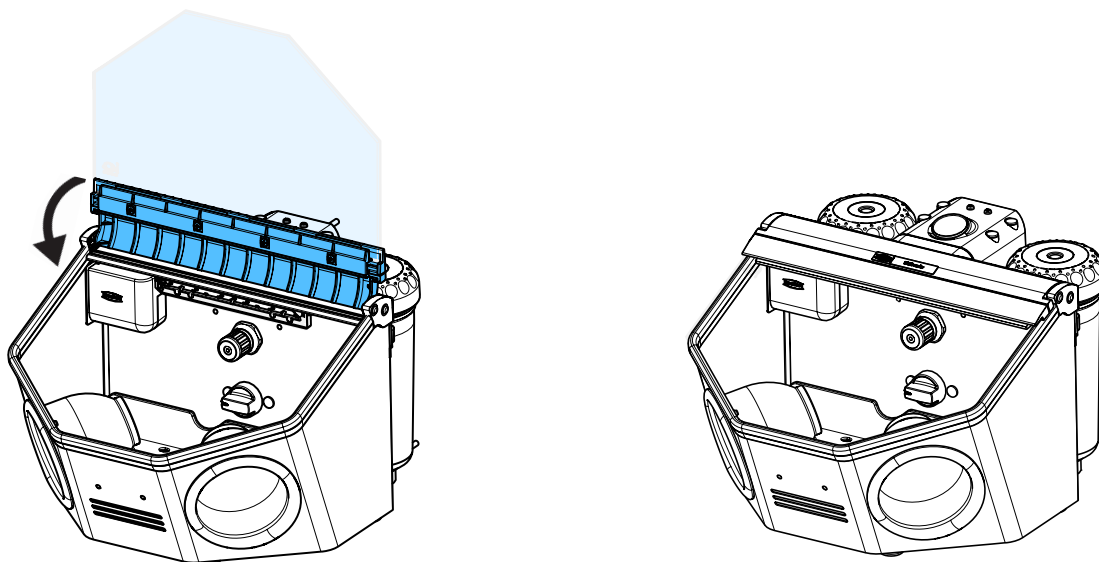









making work easy

### 90003 6353






Dosiernippel 25-70 µm steckbar • Plug-in dosing nipple 25-70 µm • Raccord de dosage 25-70 µm enfichable • Nipplo di dosaggio 25-70 µm innestabile • Boquilla de dosificación 25-70 µm enchufable • дозирующего сопла 25 мкм - 70 мкм

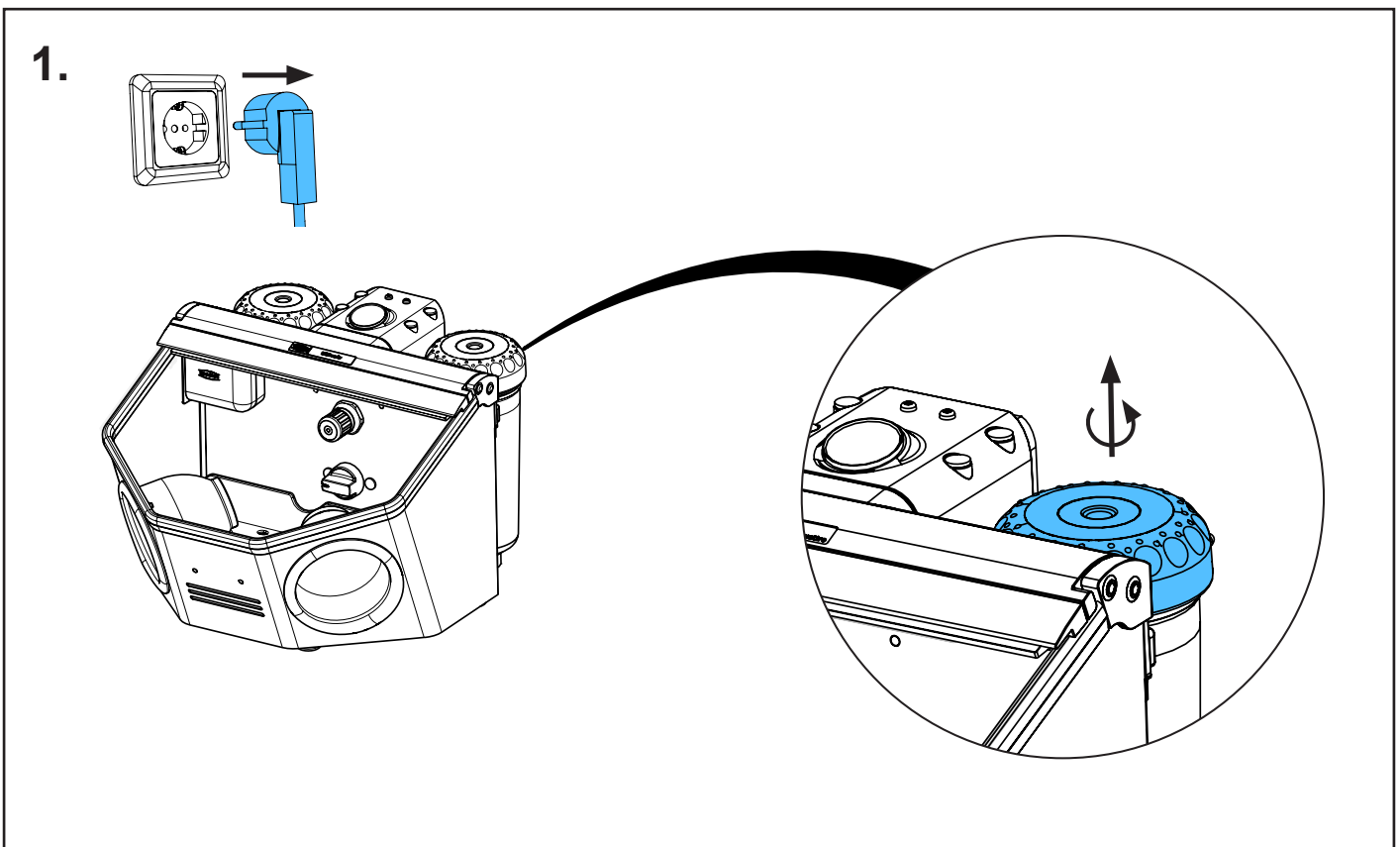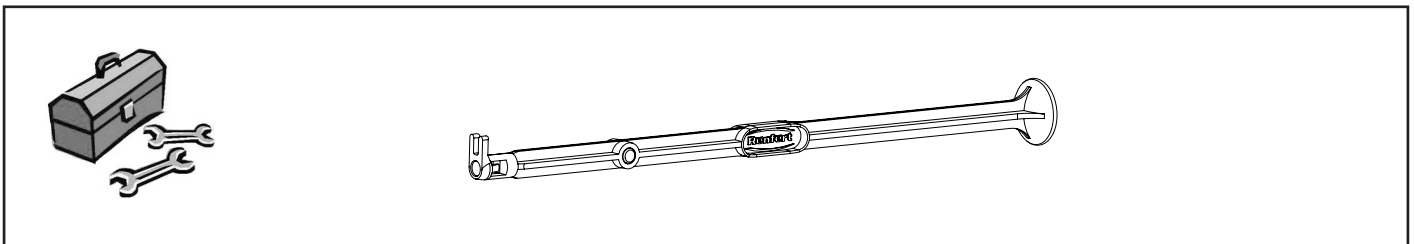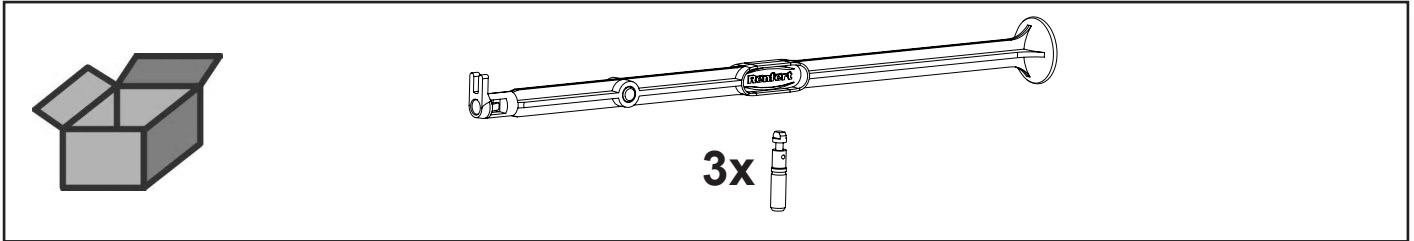


2.

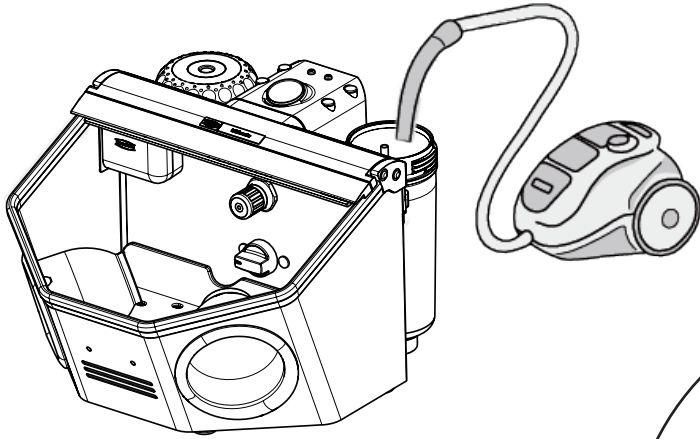


3.

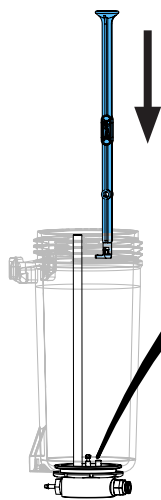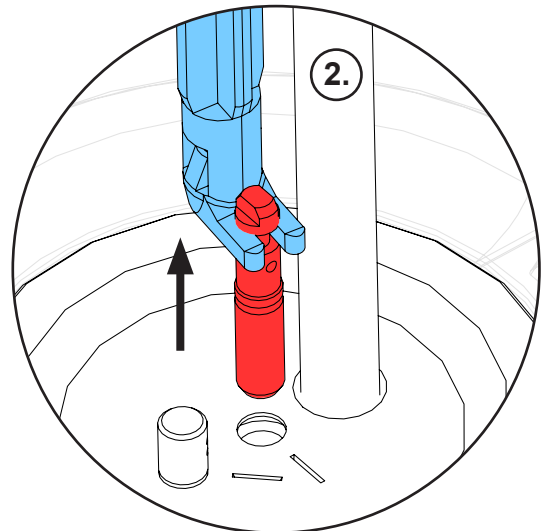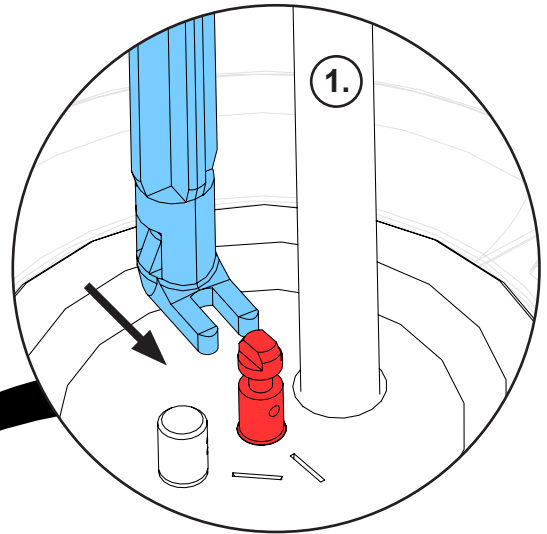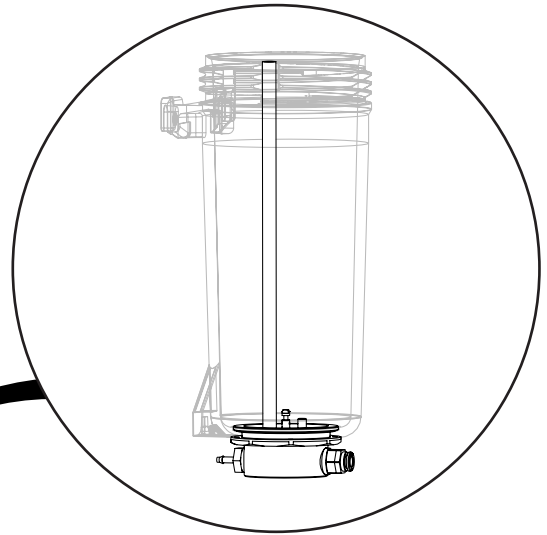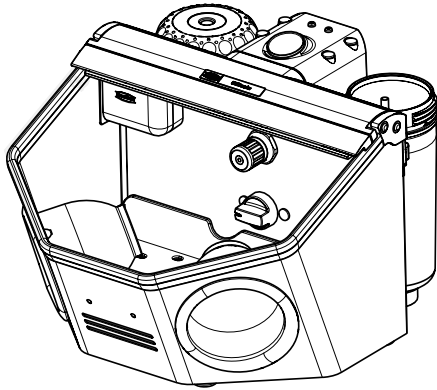


4.

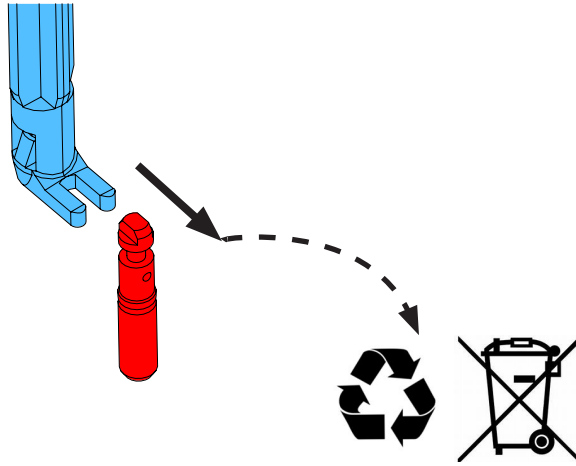




5.

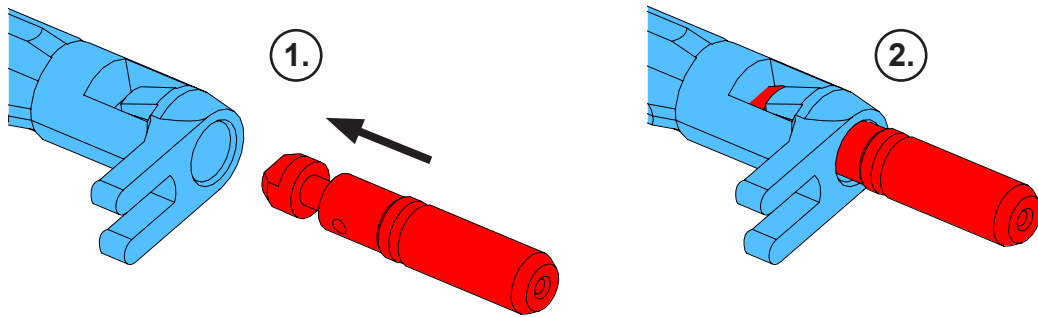


6.

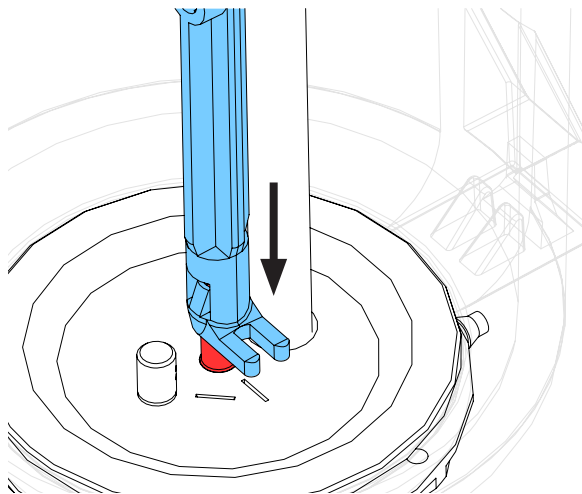


7.

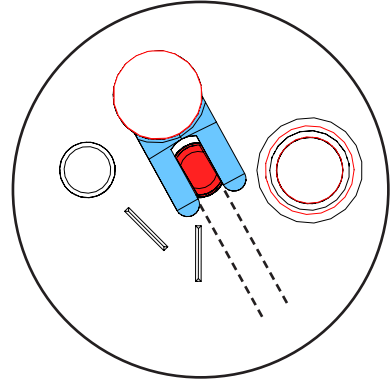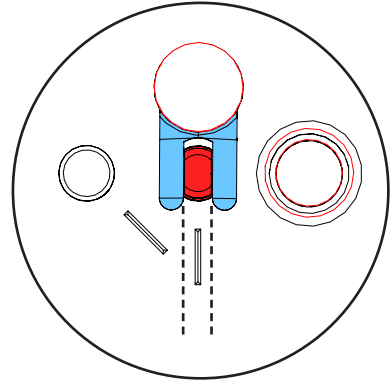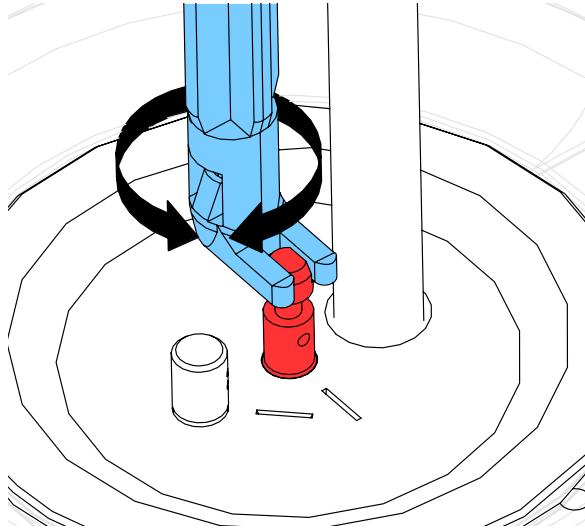

8.

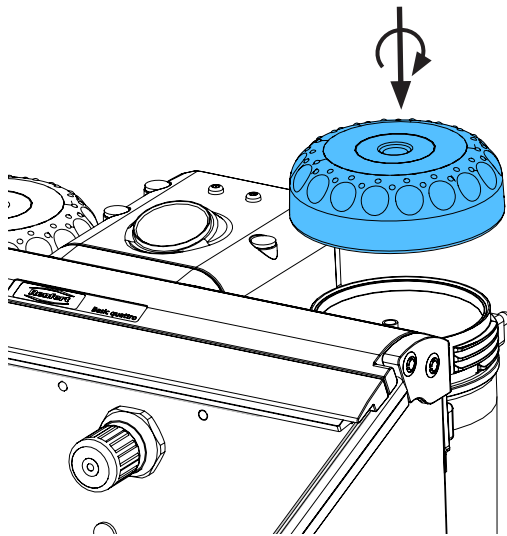

9.



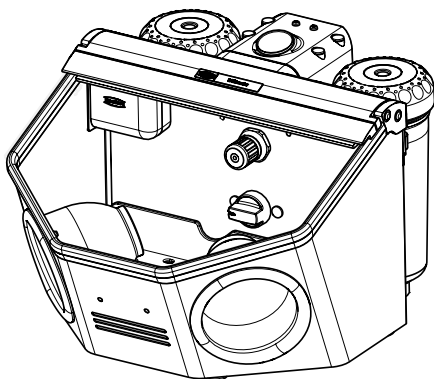





making work easy

**90003 6354**






Dosiernippel 70-250µm steckbar • Plug-in dosing nipple 70-250µm • Raccord de dosage 70-250 µm enfichable • Nipplo di dosaggio 70-250 µm innestabile • Boquilla de dosificación 70-250 µm enchufable • дозирующего сопла 70 мкм - 250 мкм

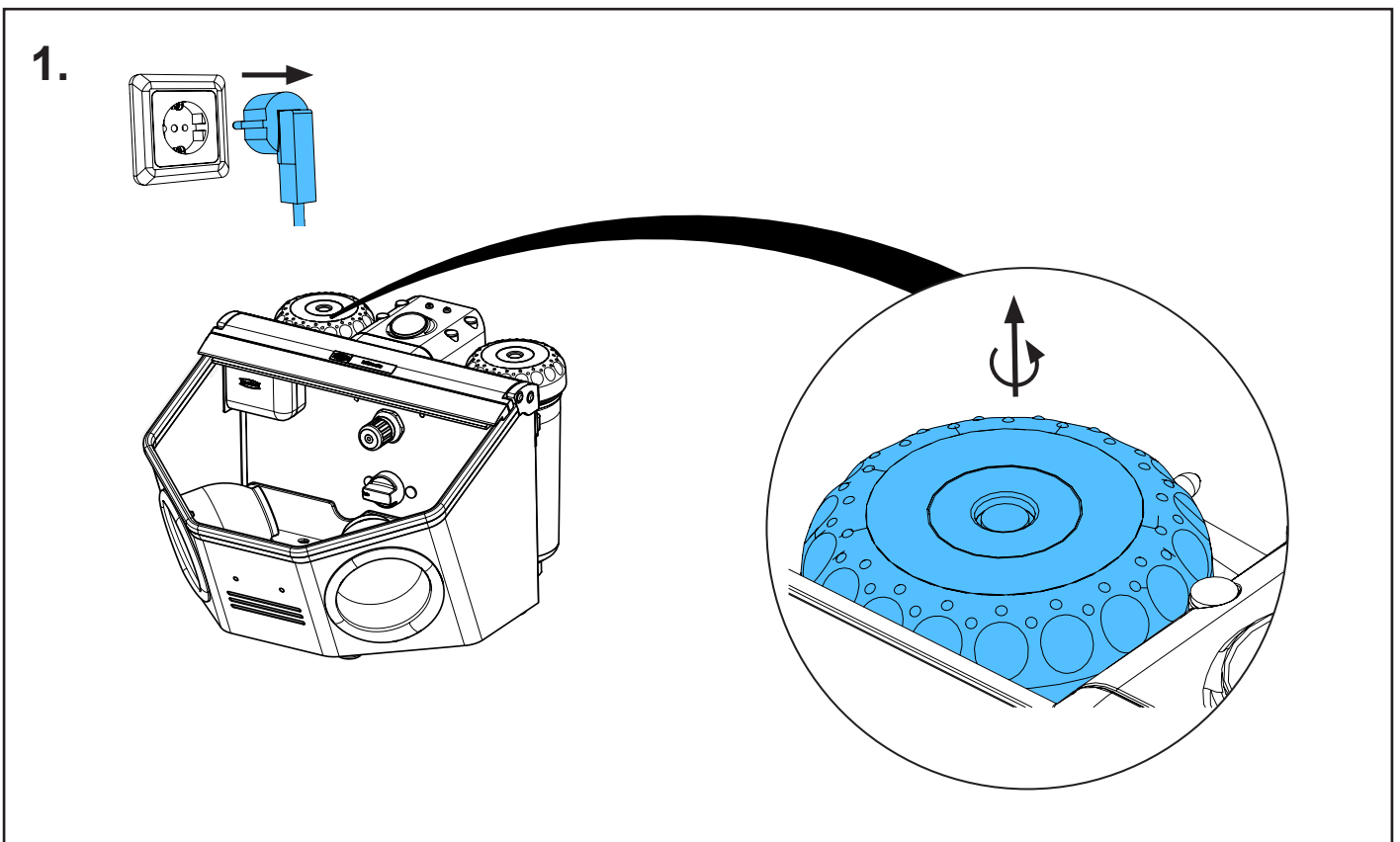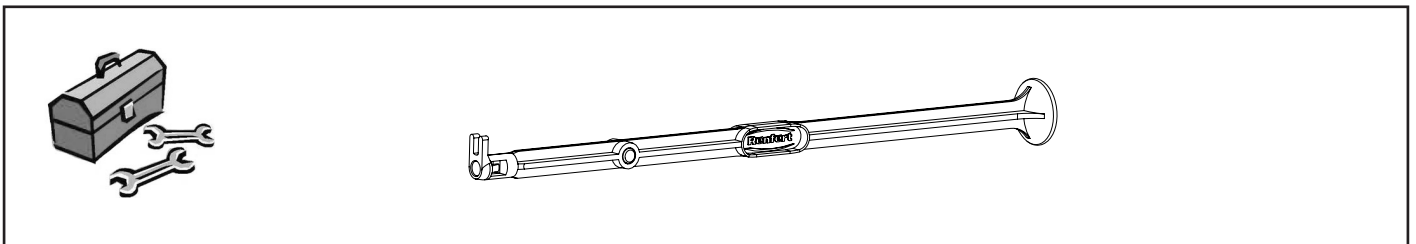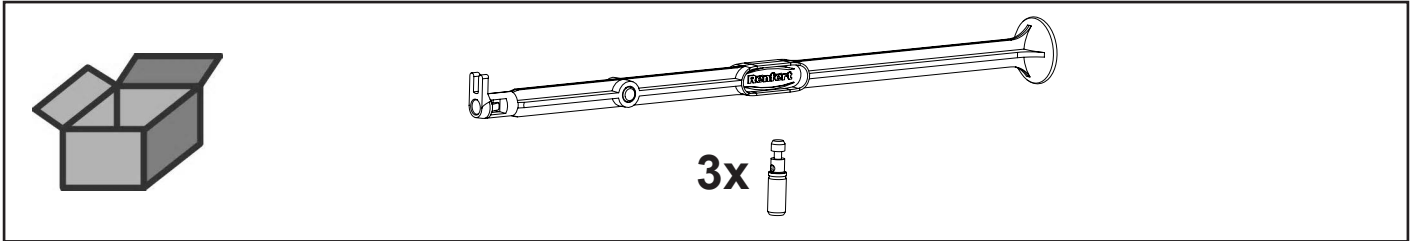


2.

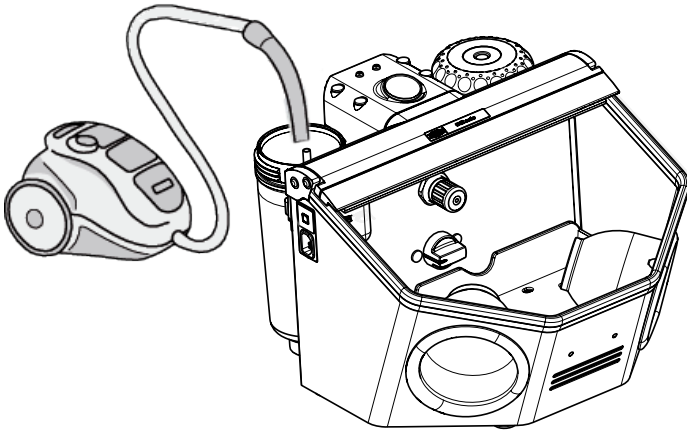


3.

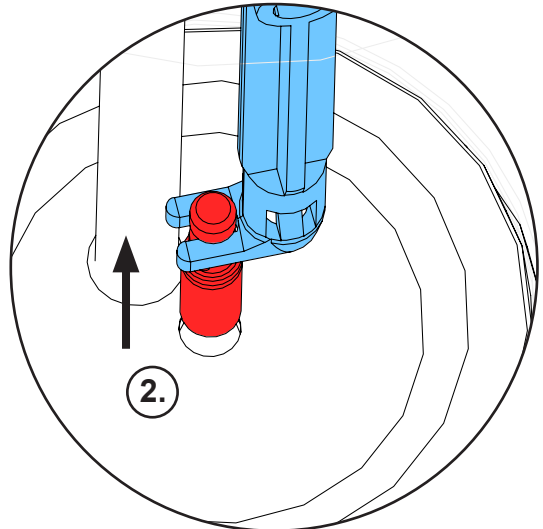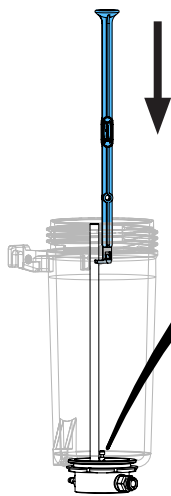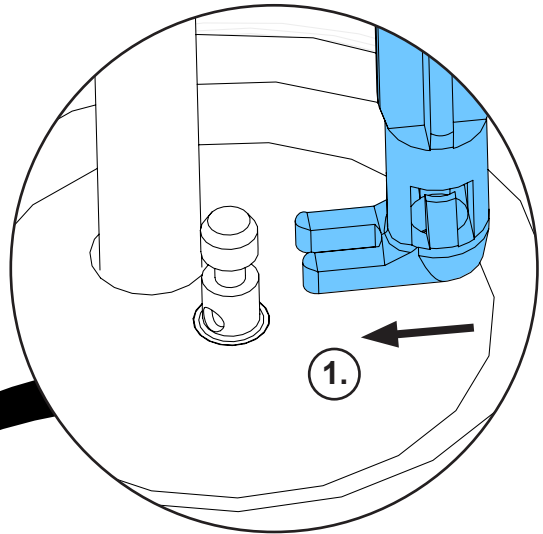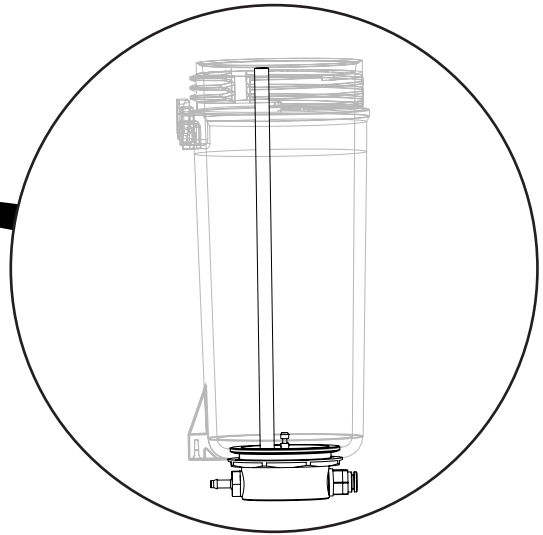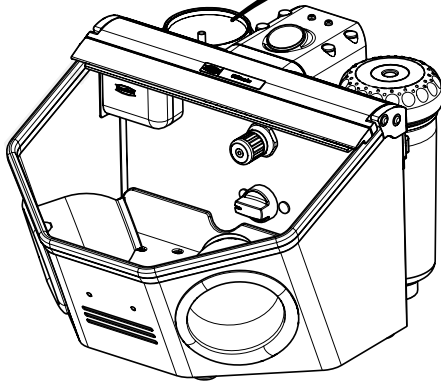


4.

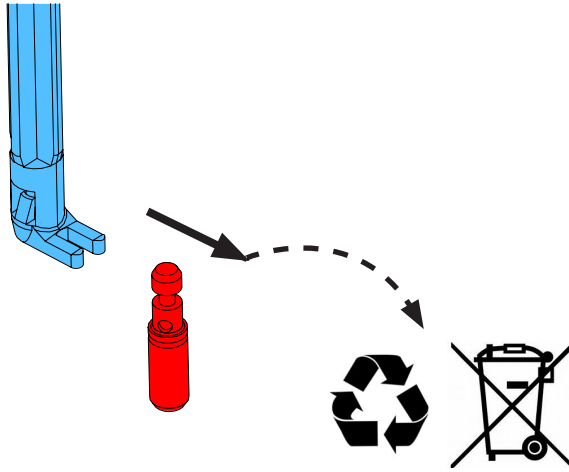


5.

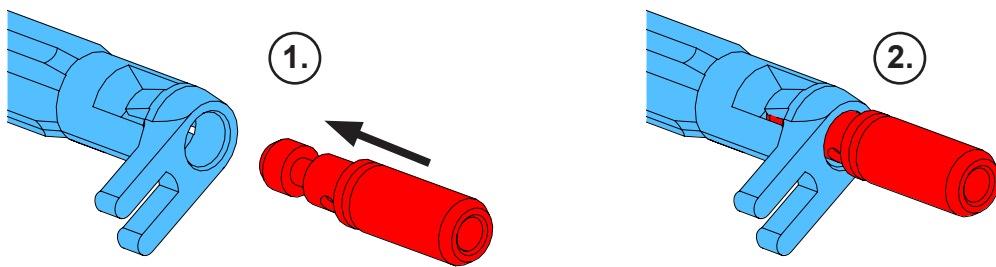


6.

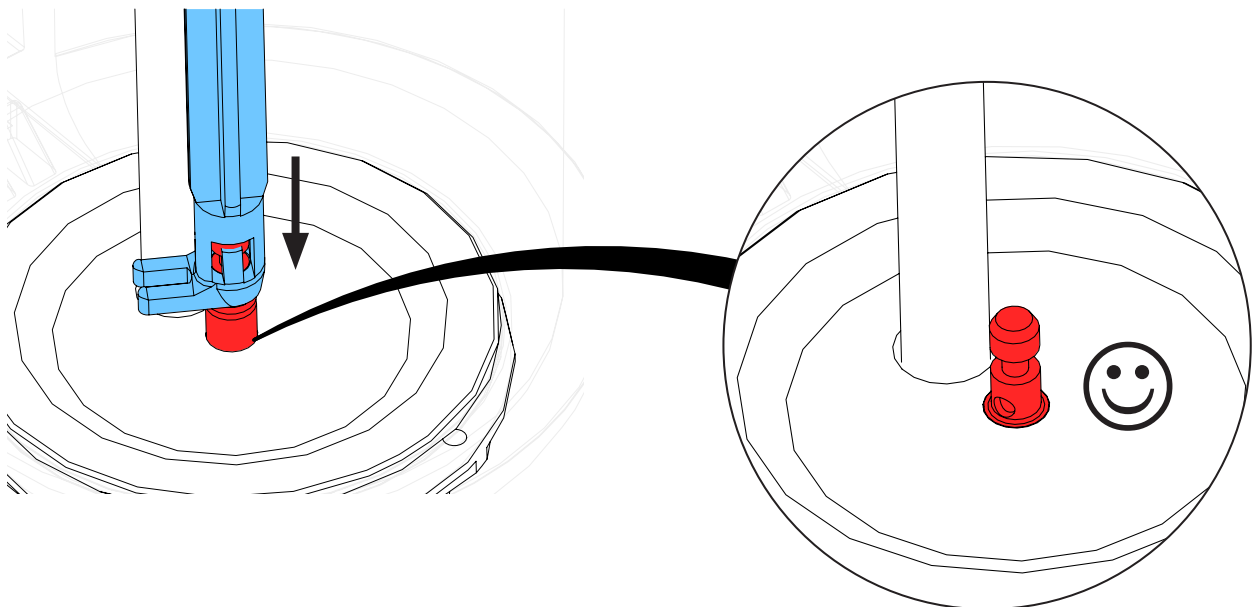

7.

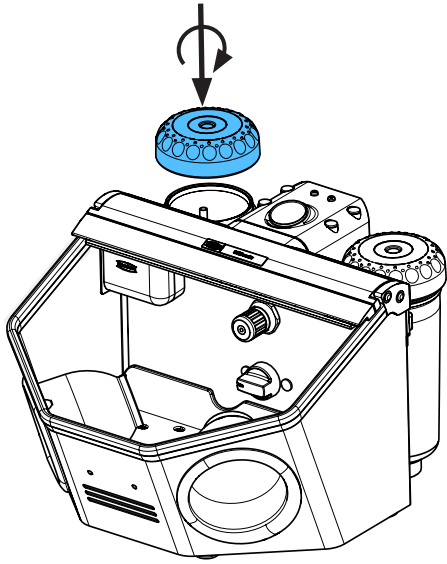

8.

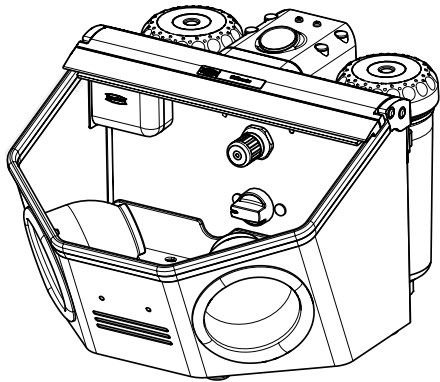







Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу


making work easy

92947 0003


Seitenteile Basic • Side panels Basic • Pièces latérales Basic • Parti laterali Basic  
Piezas laterales Basic • боковá я часть Basic

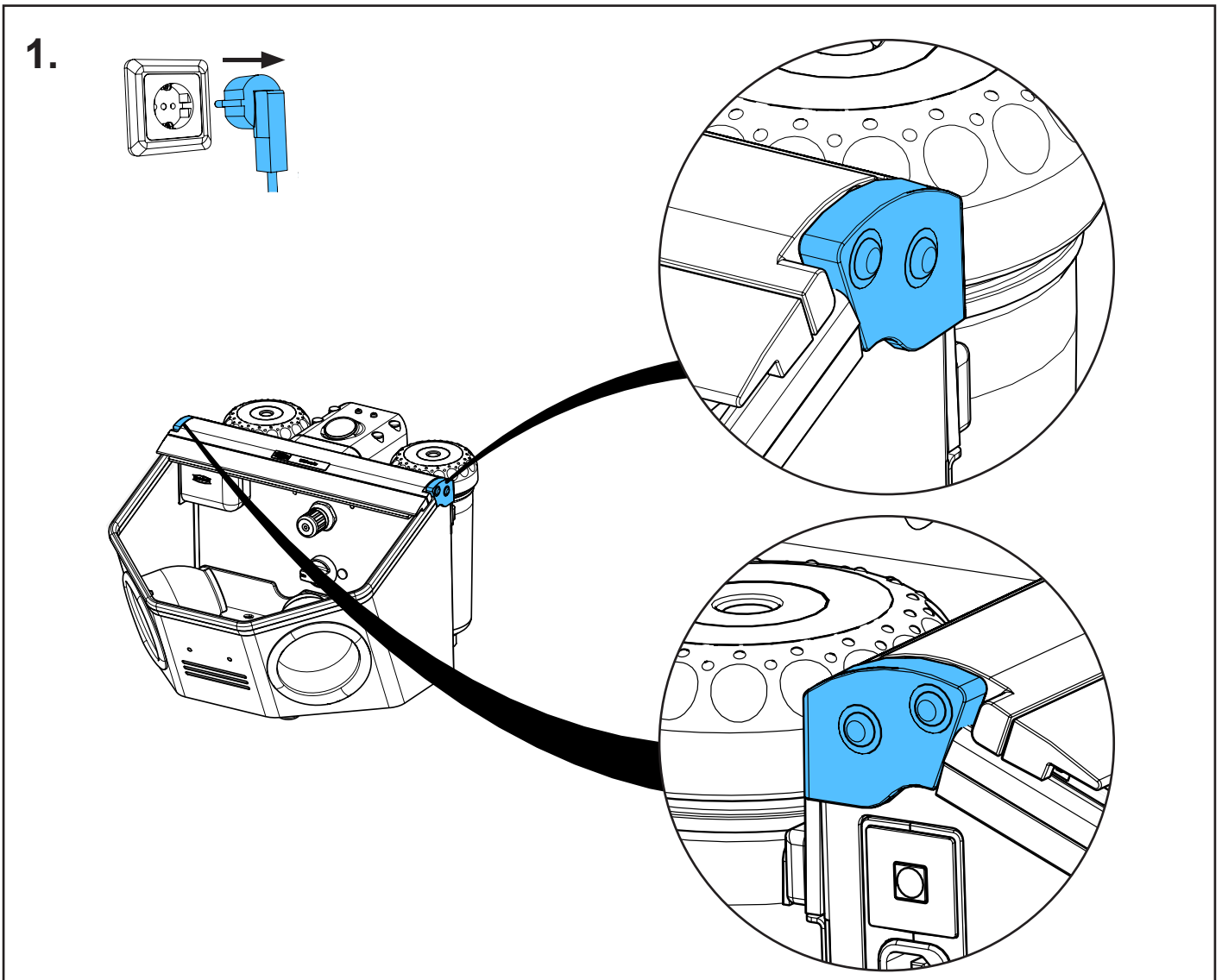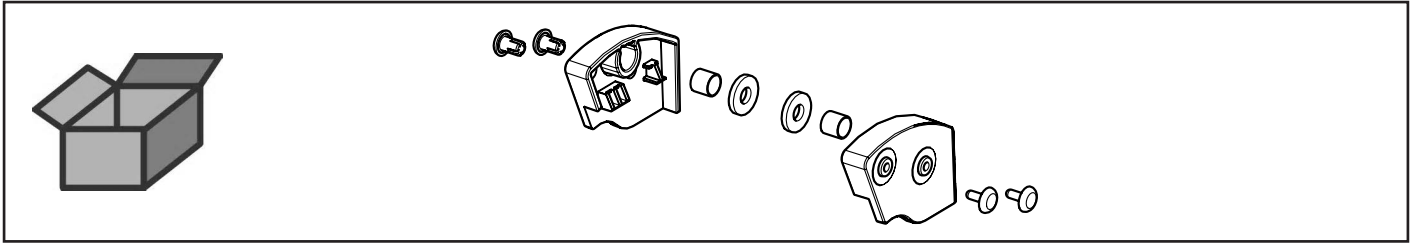


2.

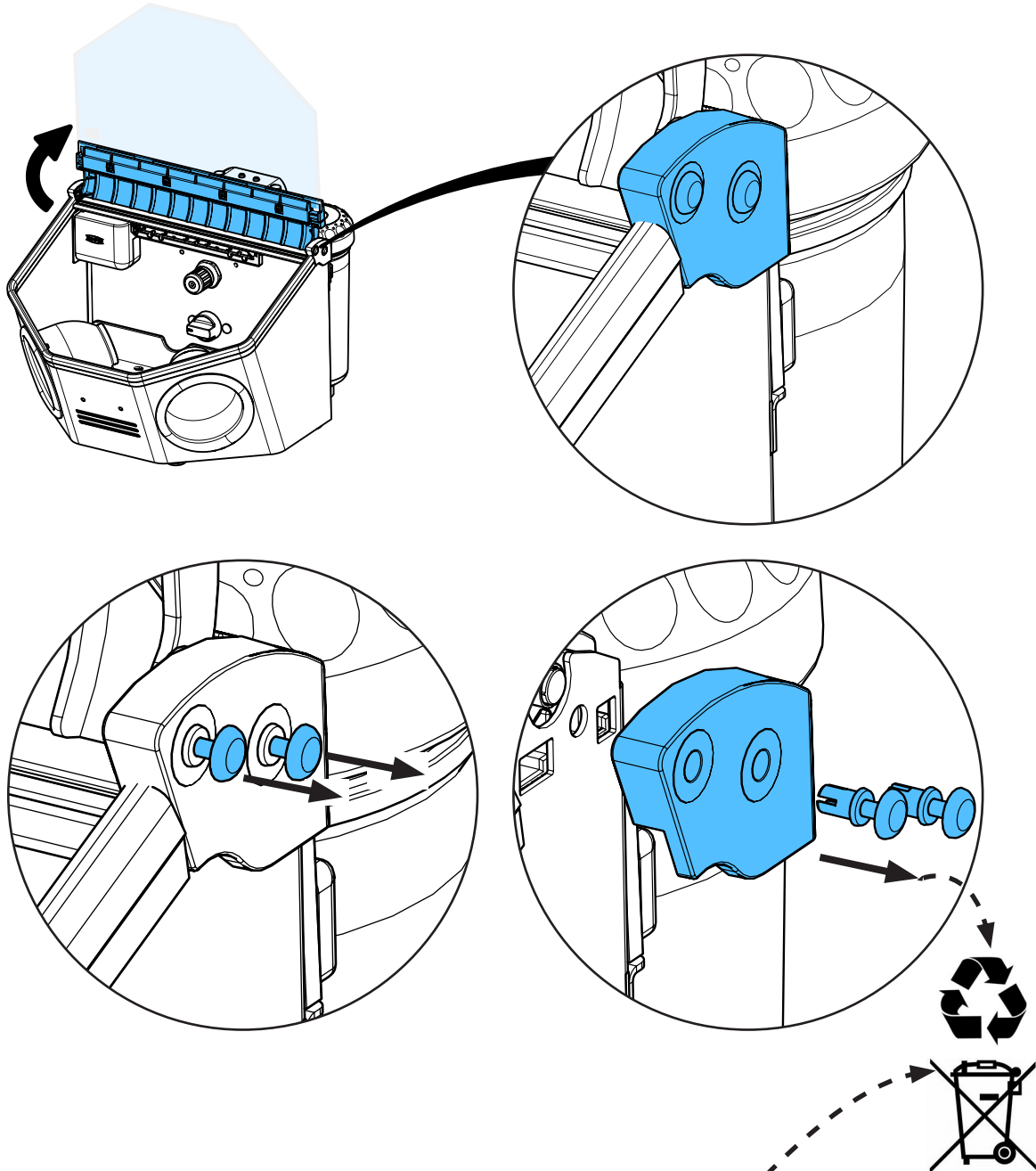


3.

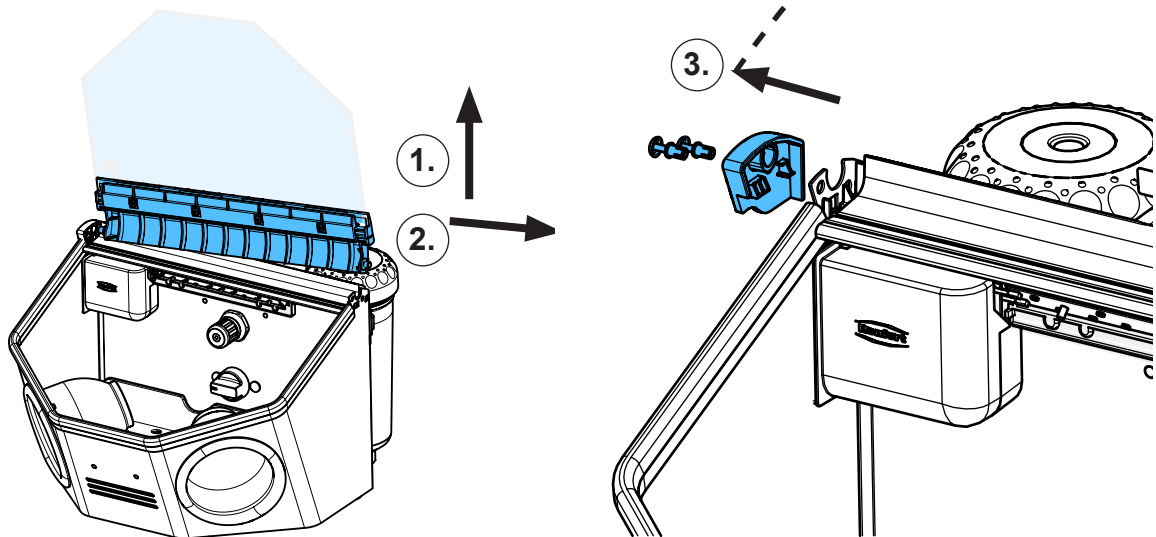


4.

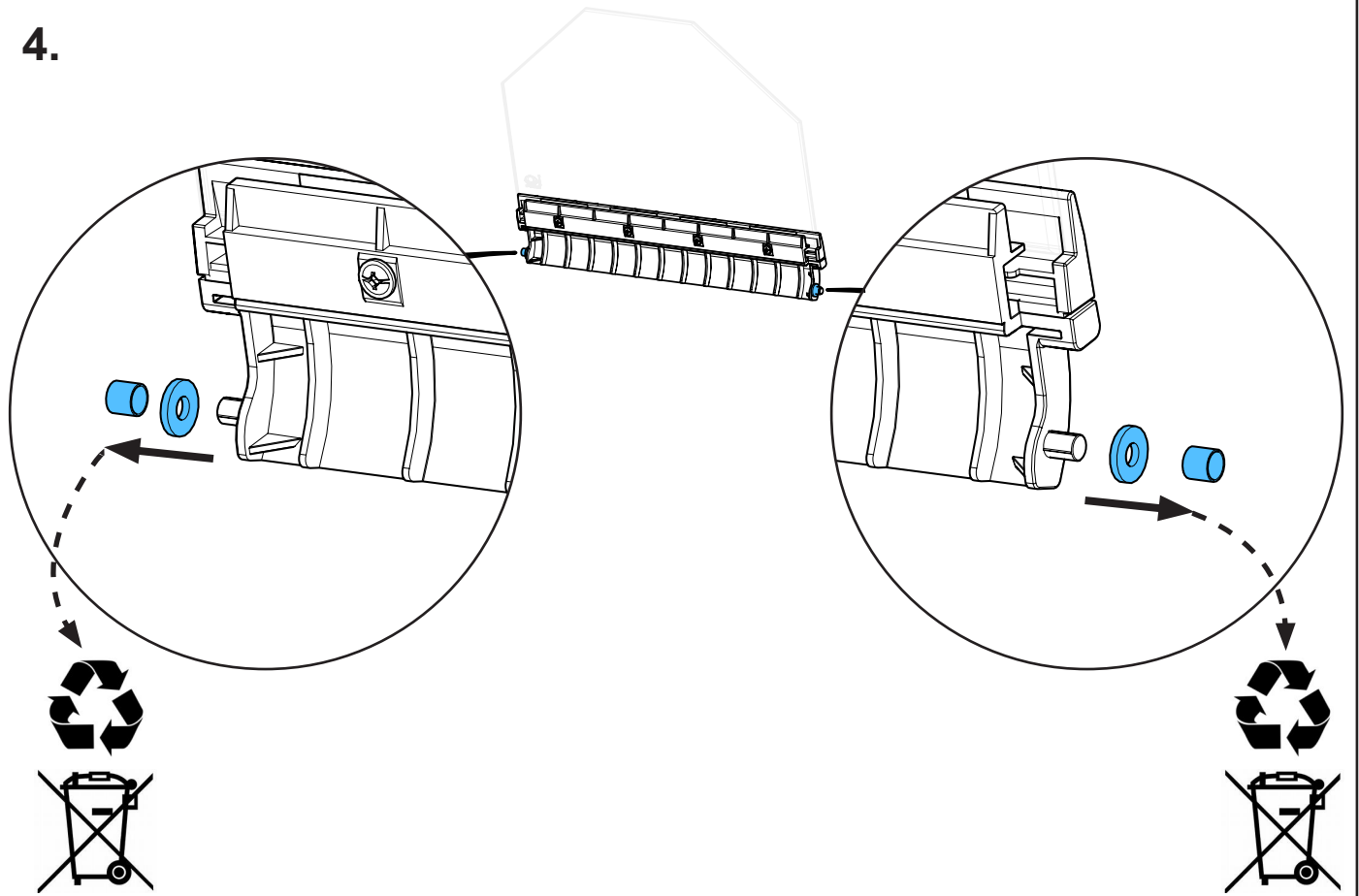


5.

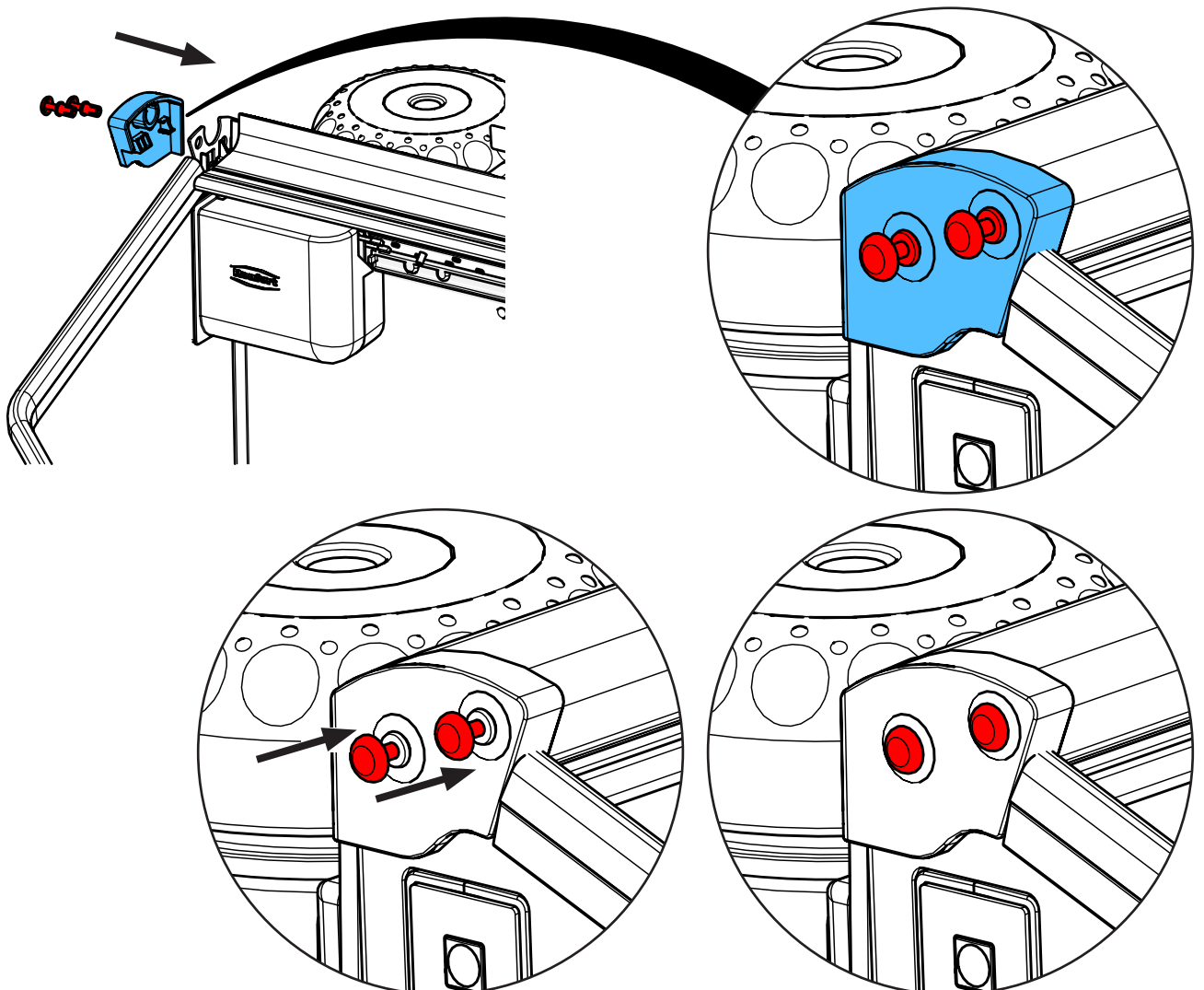


6.

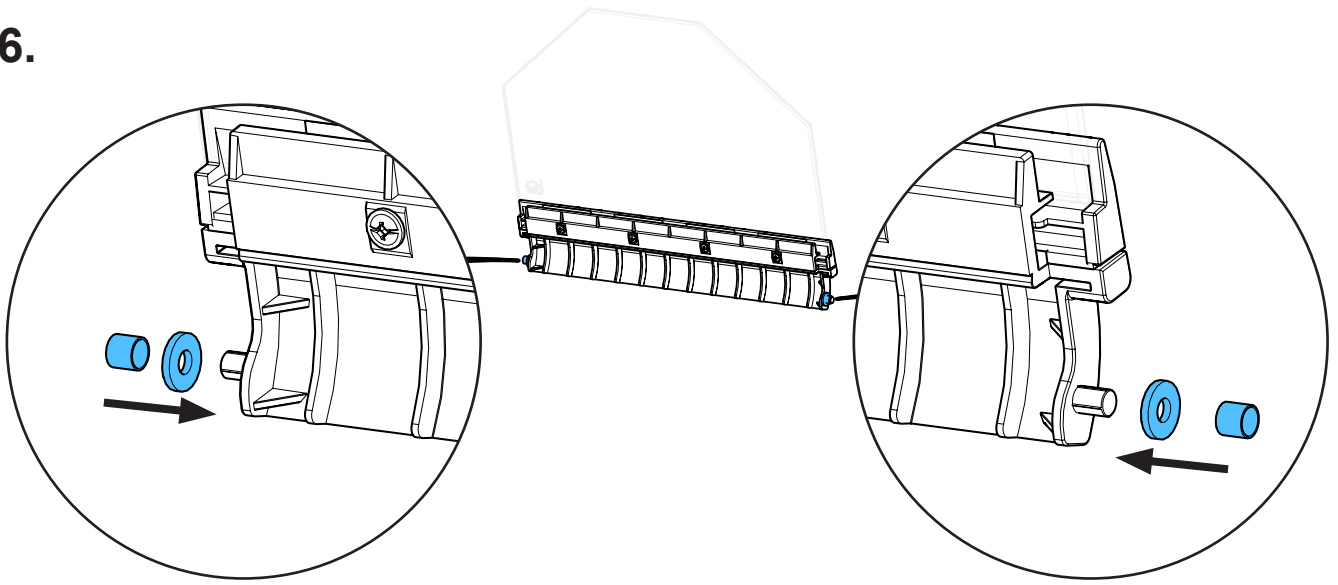

7.

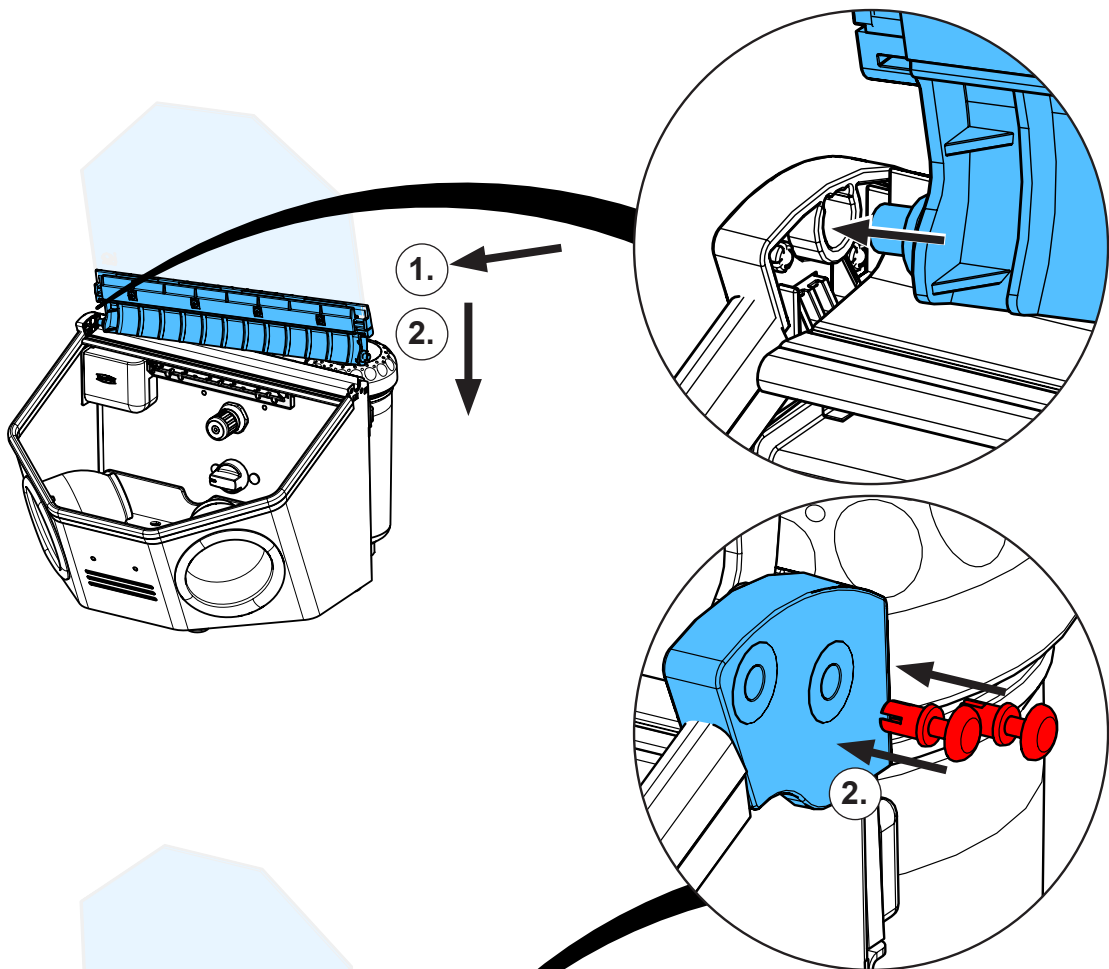

8.

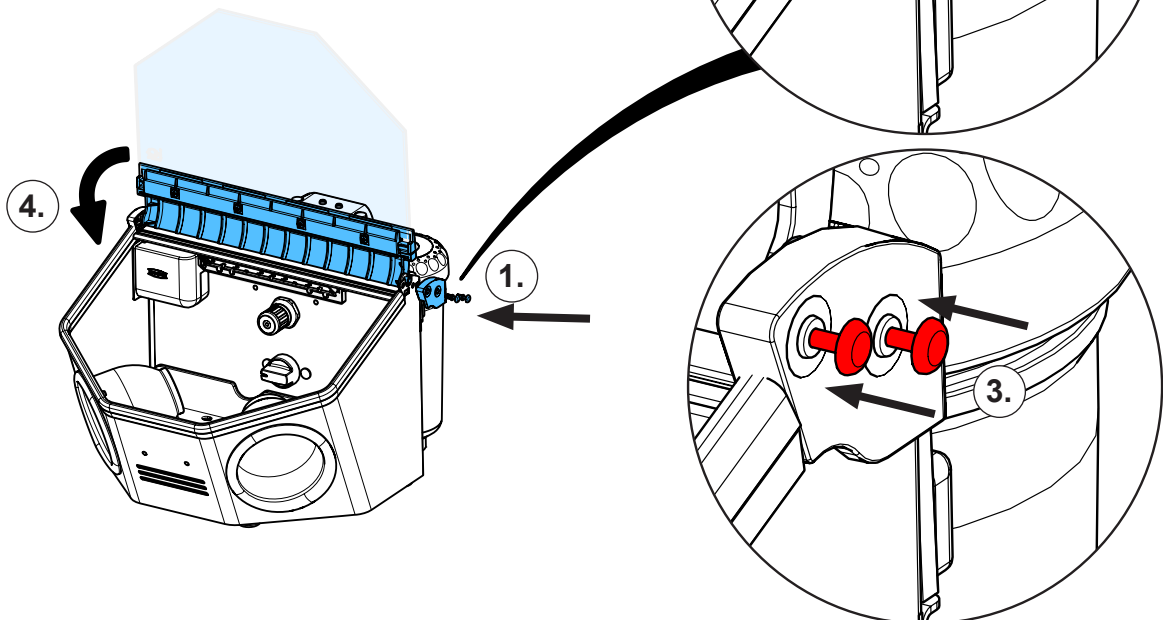





Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу


making work easy

90003 8600


Durchstrahlenschutz • Blast-through protection • Protection contre perforation • Pro-  
tezione contro perforazione • Protección contra perforación • Защита от сквозного  
прохождения струи песка

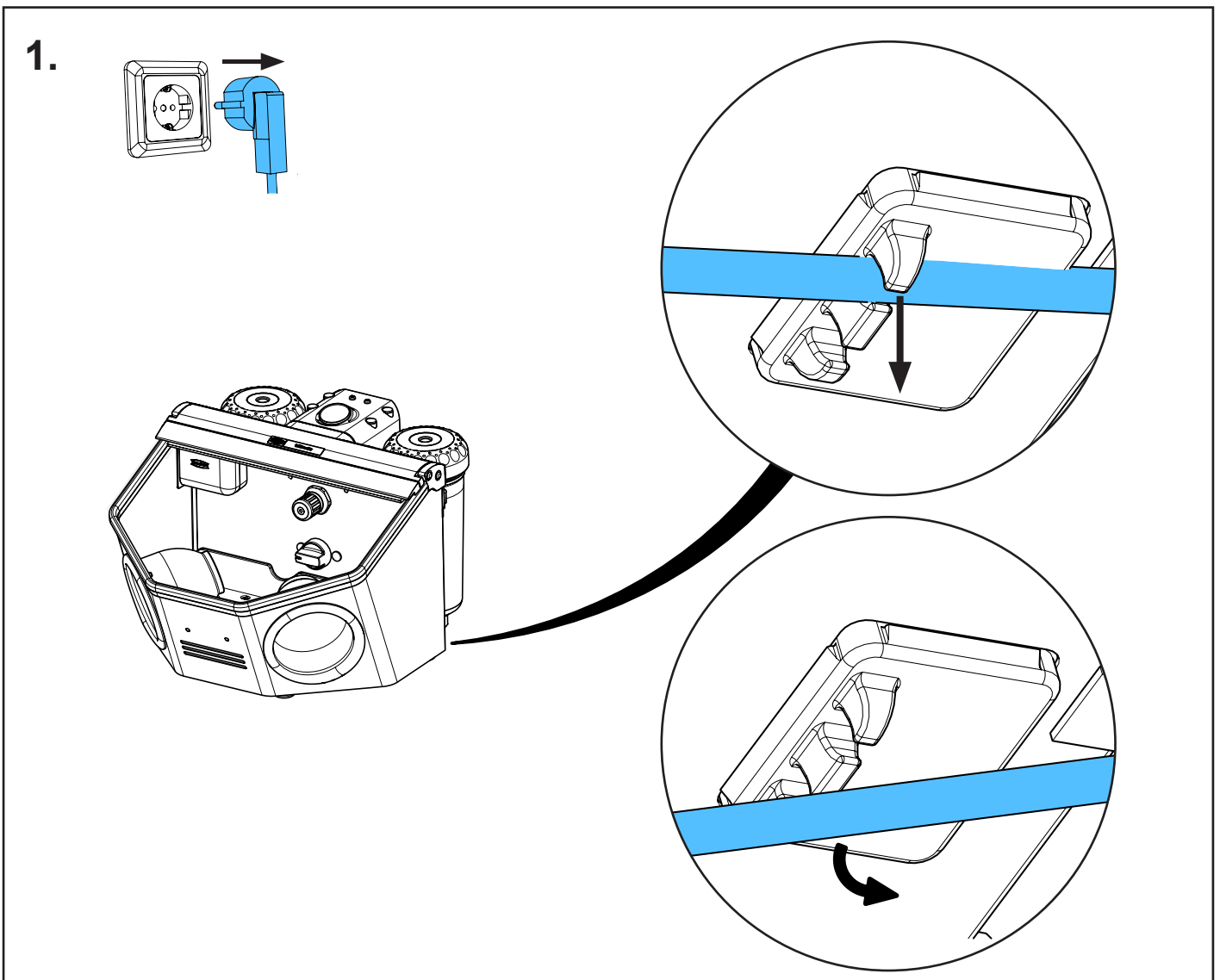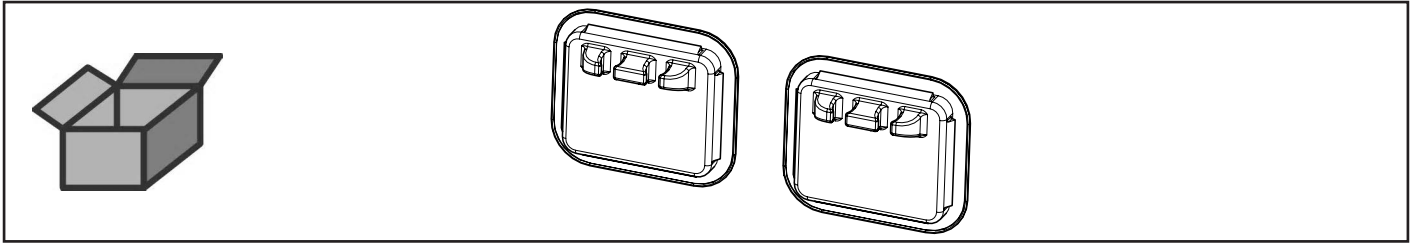


2.

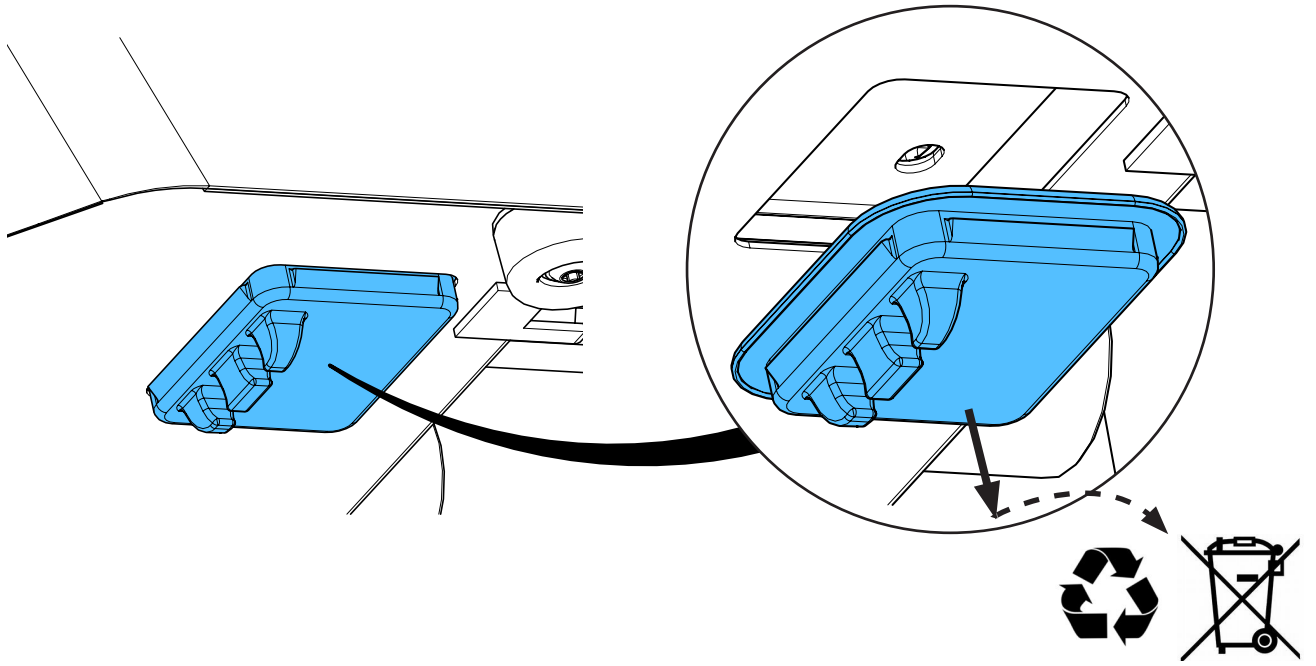

3.

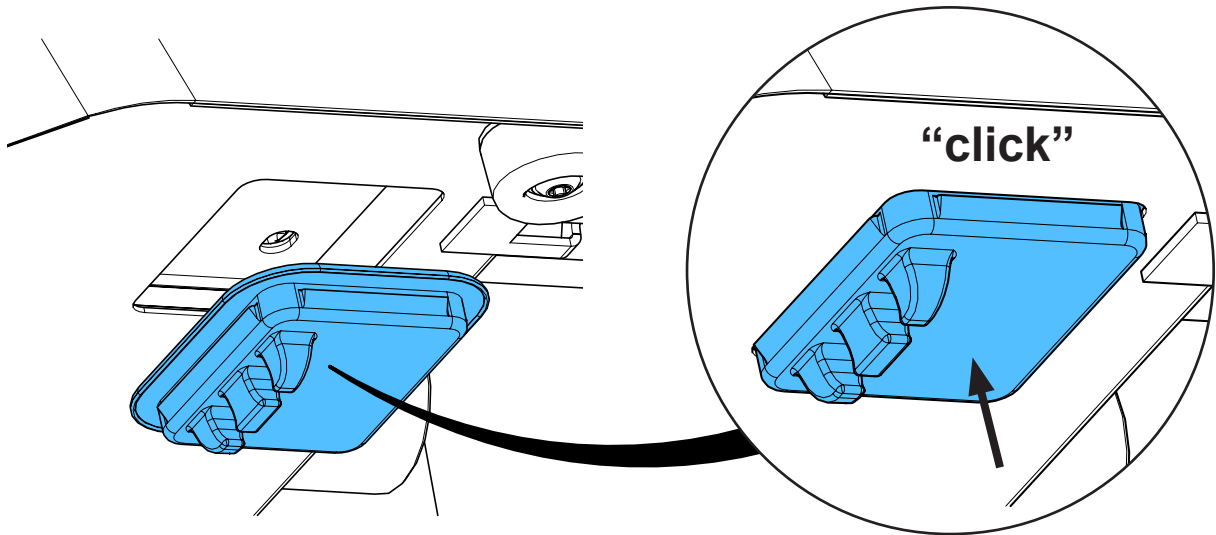

4.

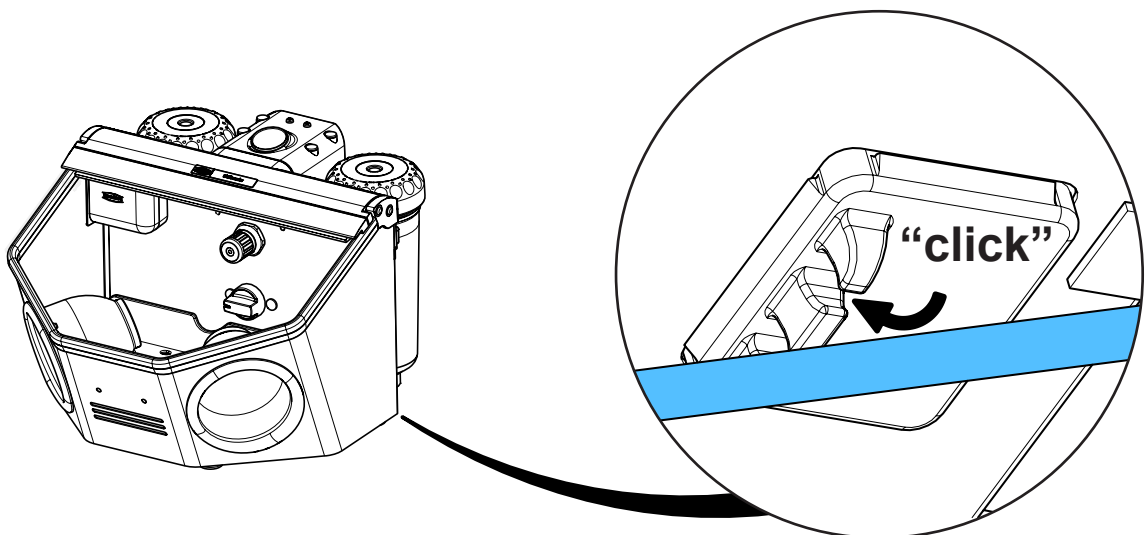





Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу

making work easy

**90015 0864**

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PU-Schlauch 6 x 3,9 mm • PU-hose 6 x 3.9 mm • Tuyau flexible en polyuréthane 6  
x 3,9 mm • Tubo flessibile in poliuretano 6 x 3,9 mm • Tubo flexible en poliuretano  
6 x 3,9 mm • PU-шланга 6 x 3,9 мм

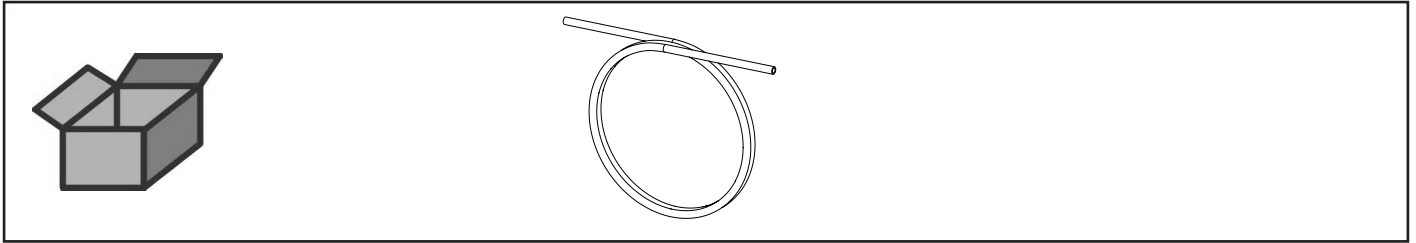











making work easy

**90003 5800**


Ersatztankmantel Basic • Replacement tank case Basic • Boîtier du silo Basic •  
Involucro deposito Basic • Envoltura depósito Basic • Корпус запасного бачка  
Basic

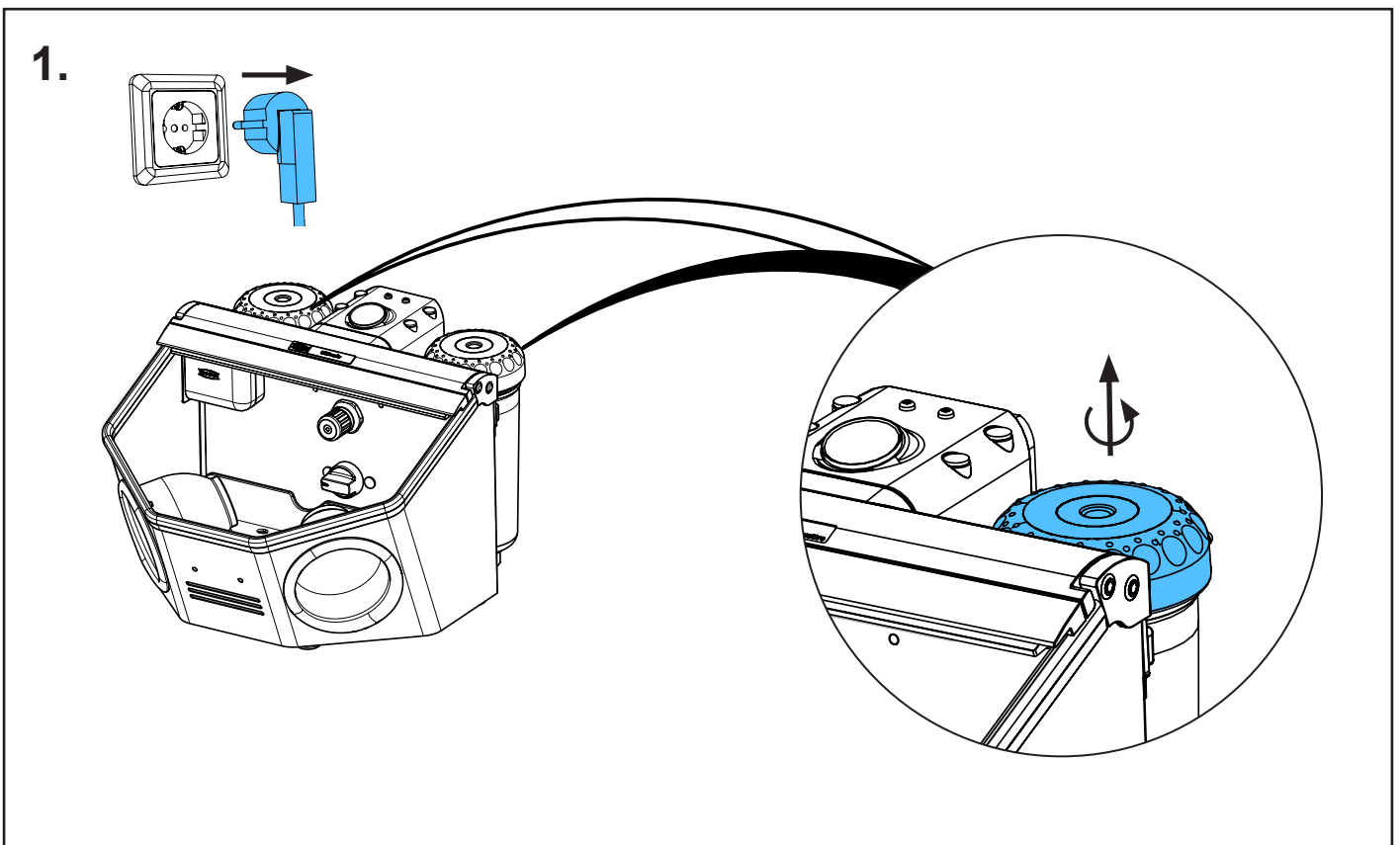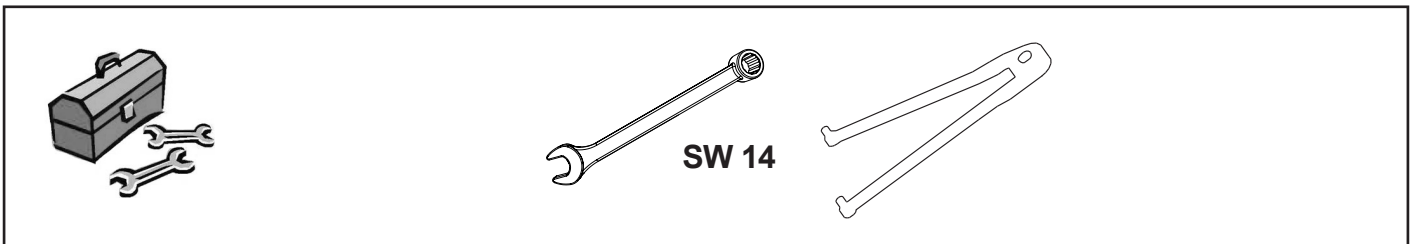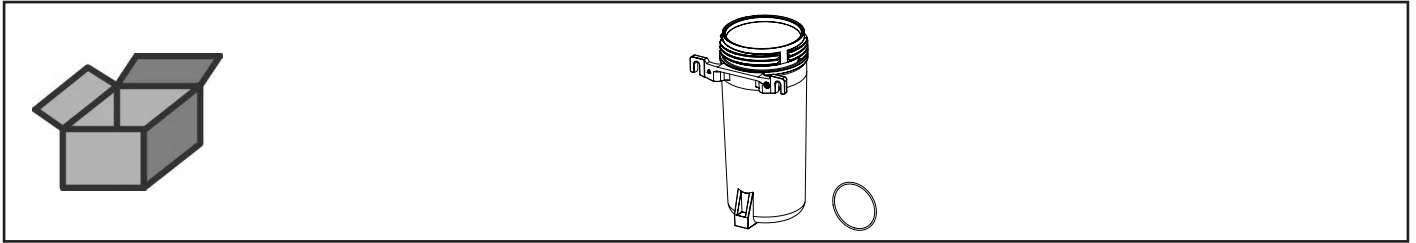


2.

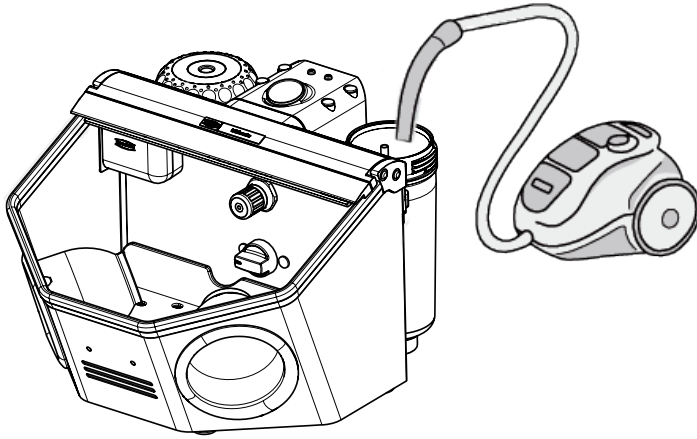


3.

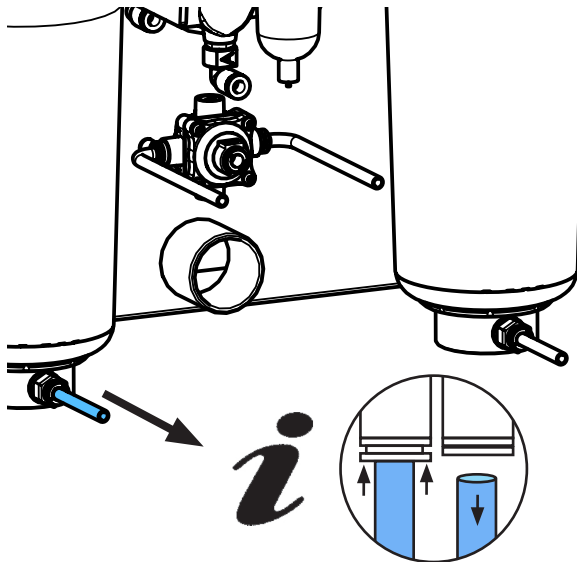


4.

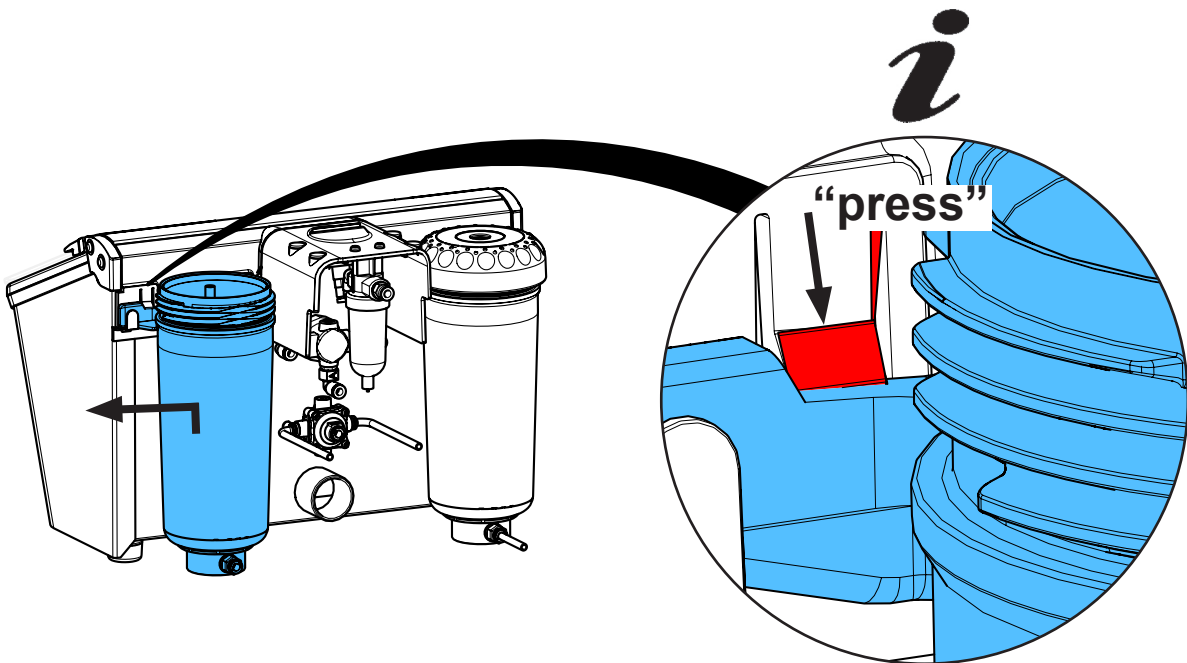


5.

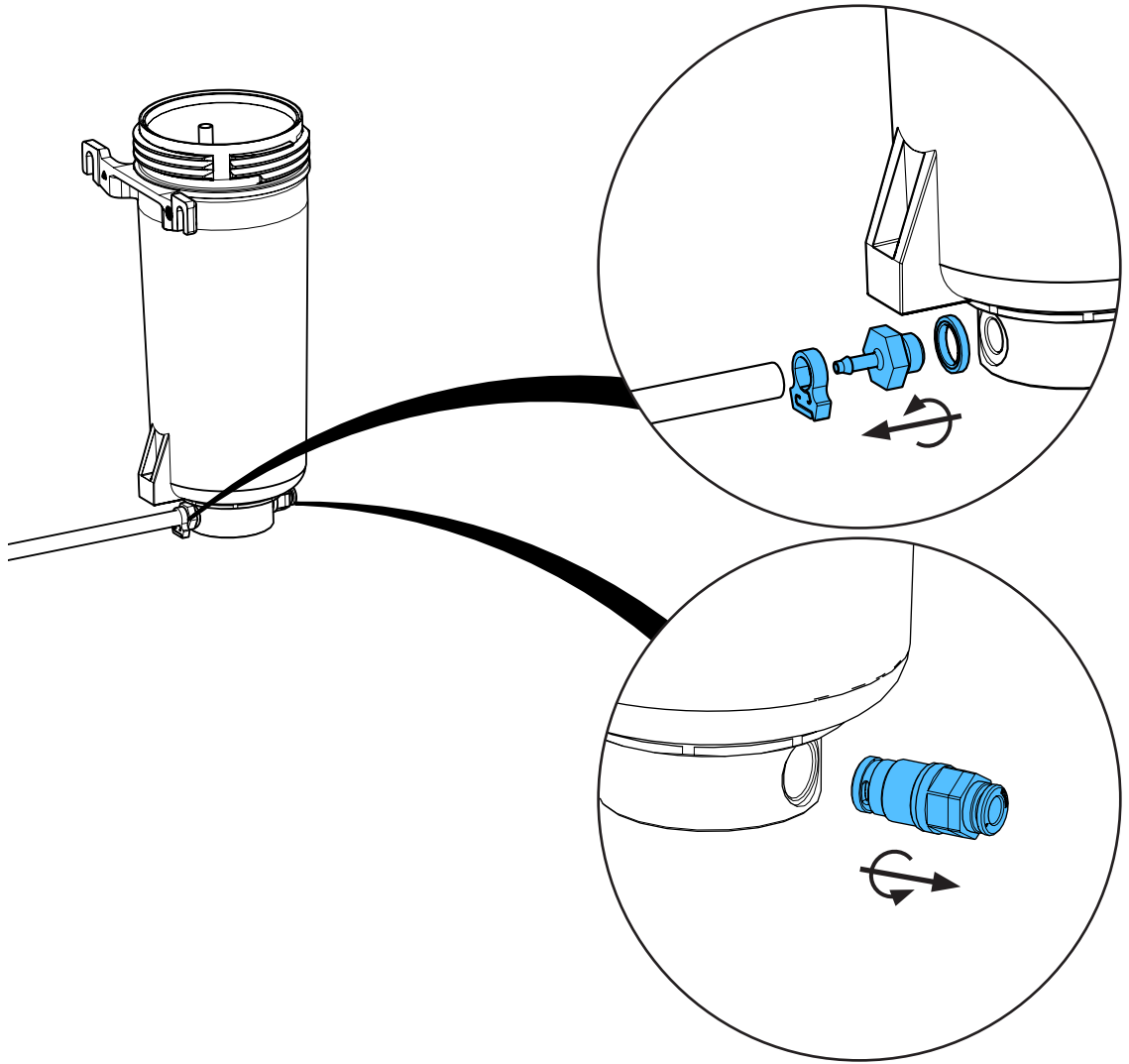


6.

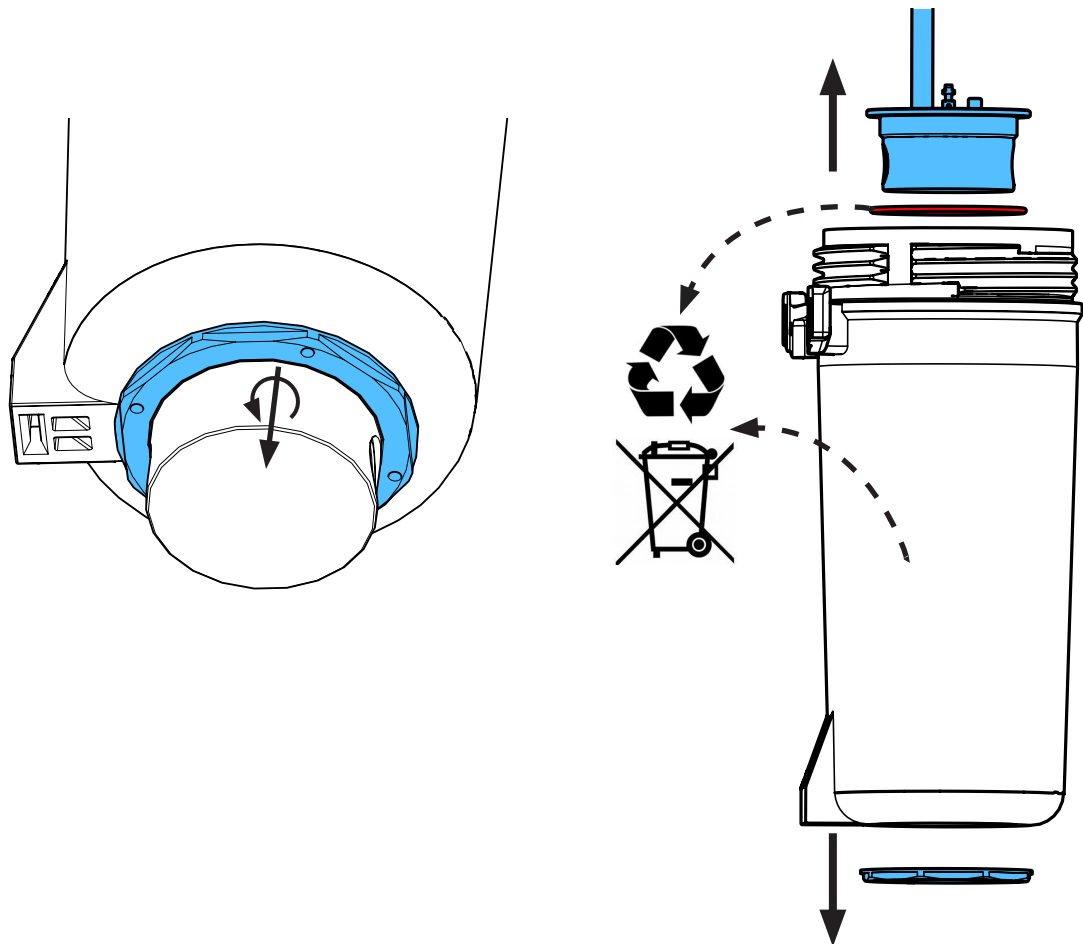


7.

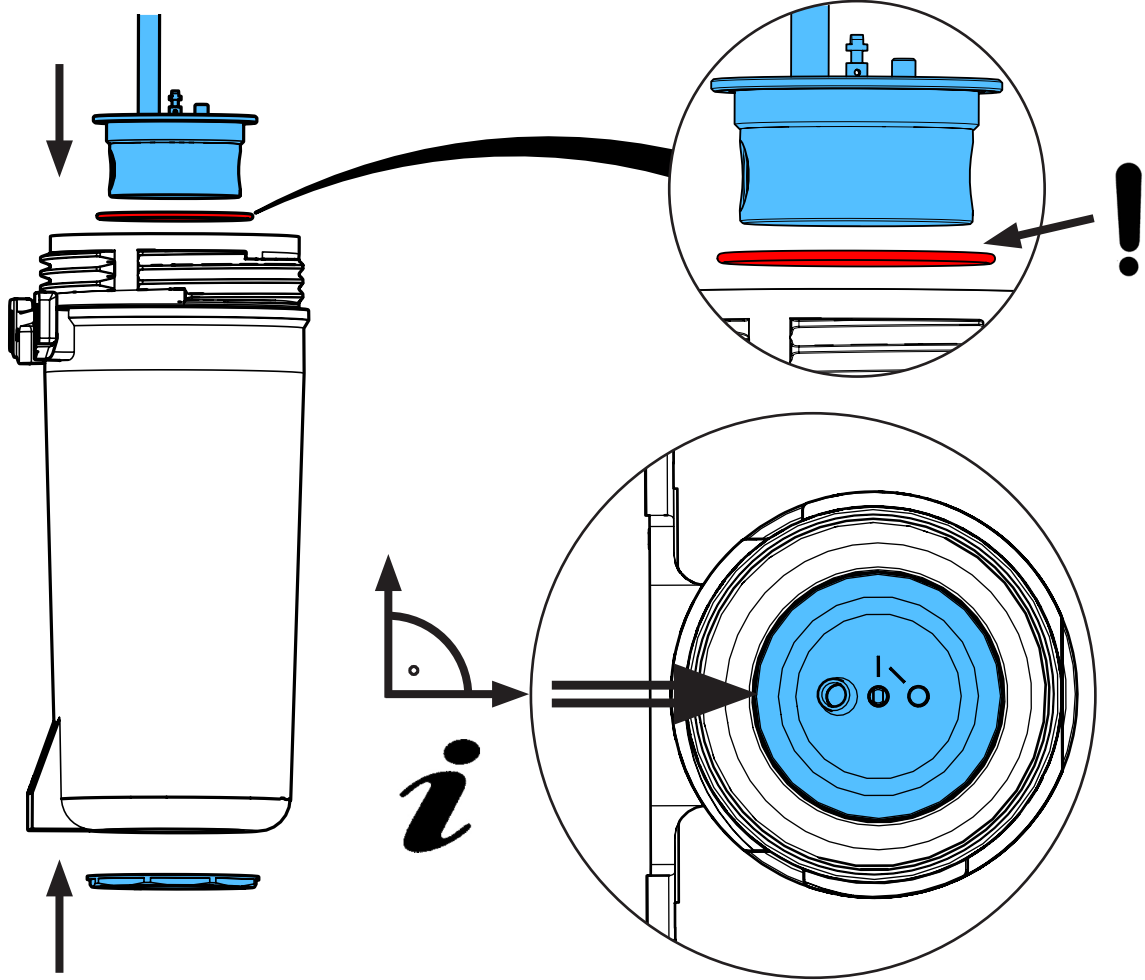


8.

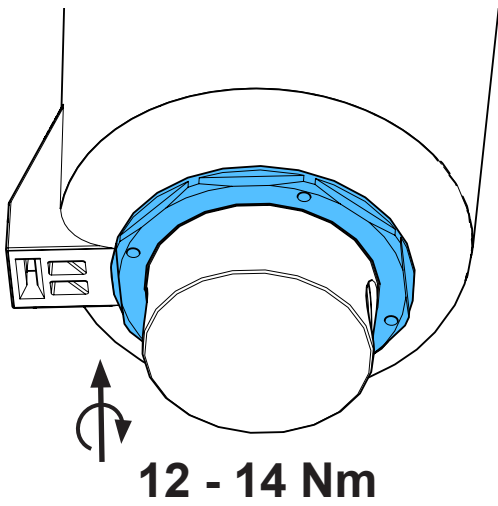


9.

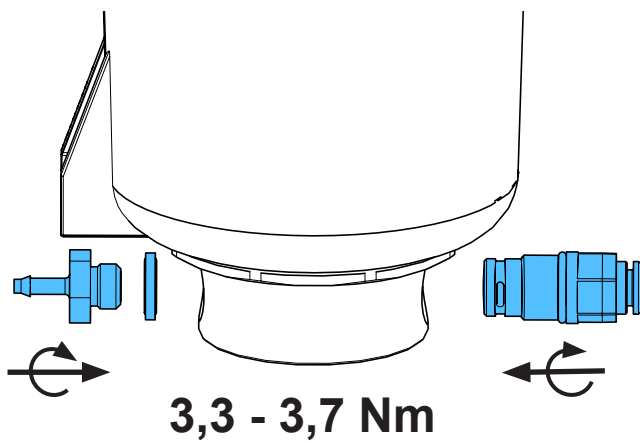


10.

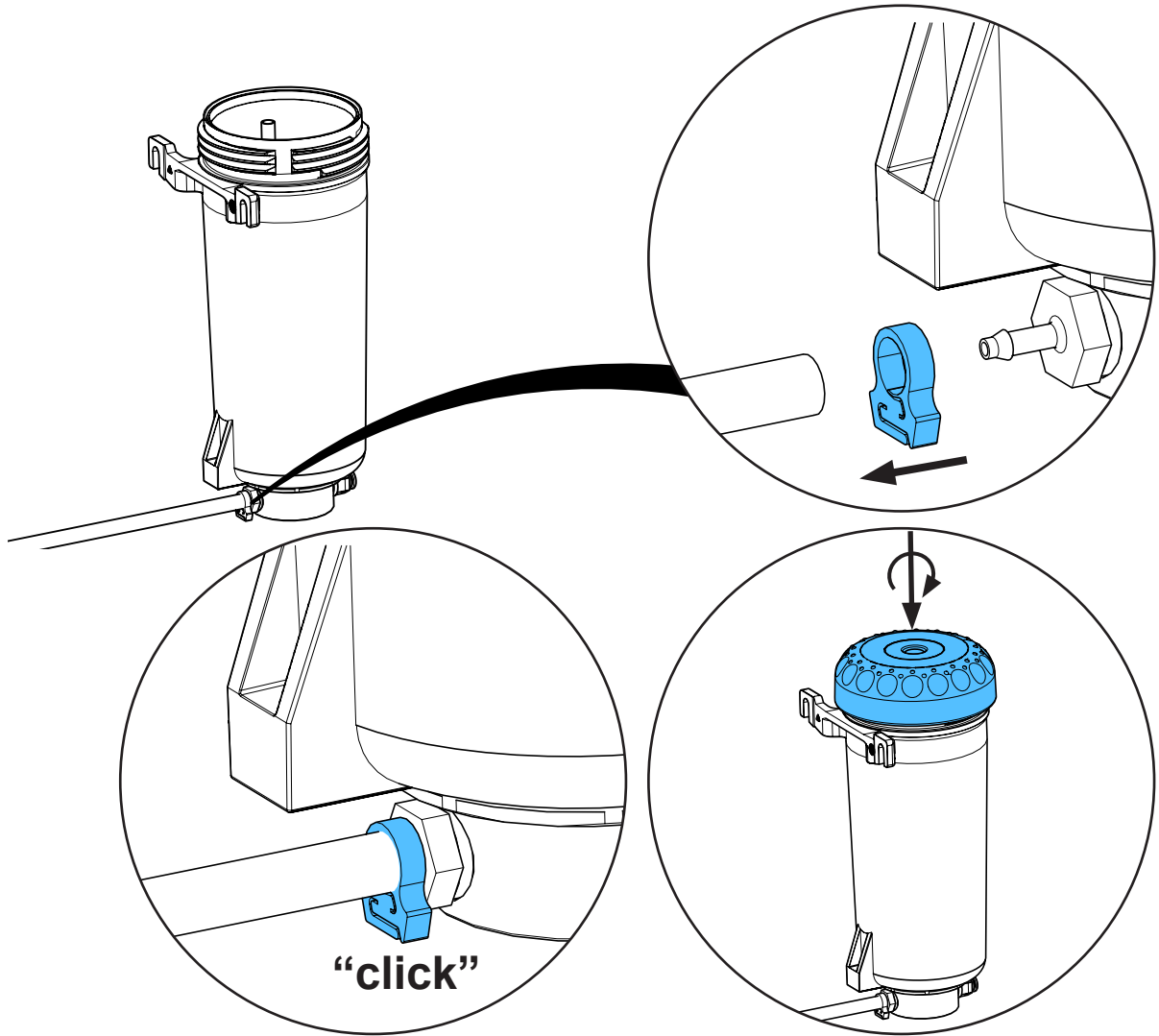


11.

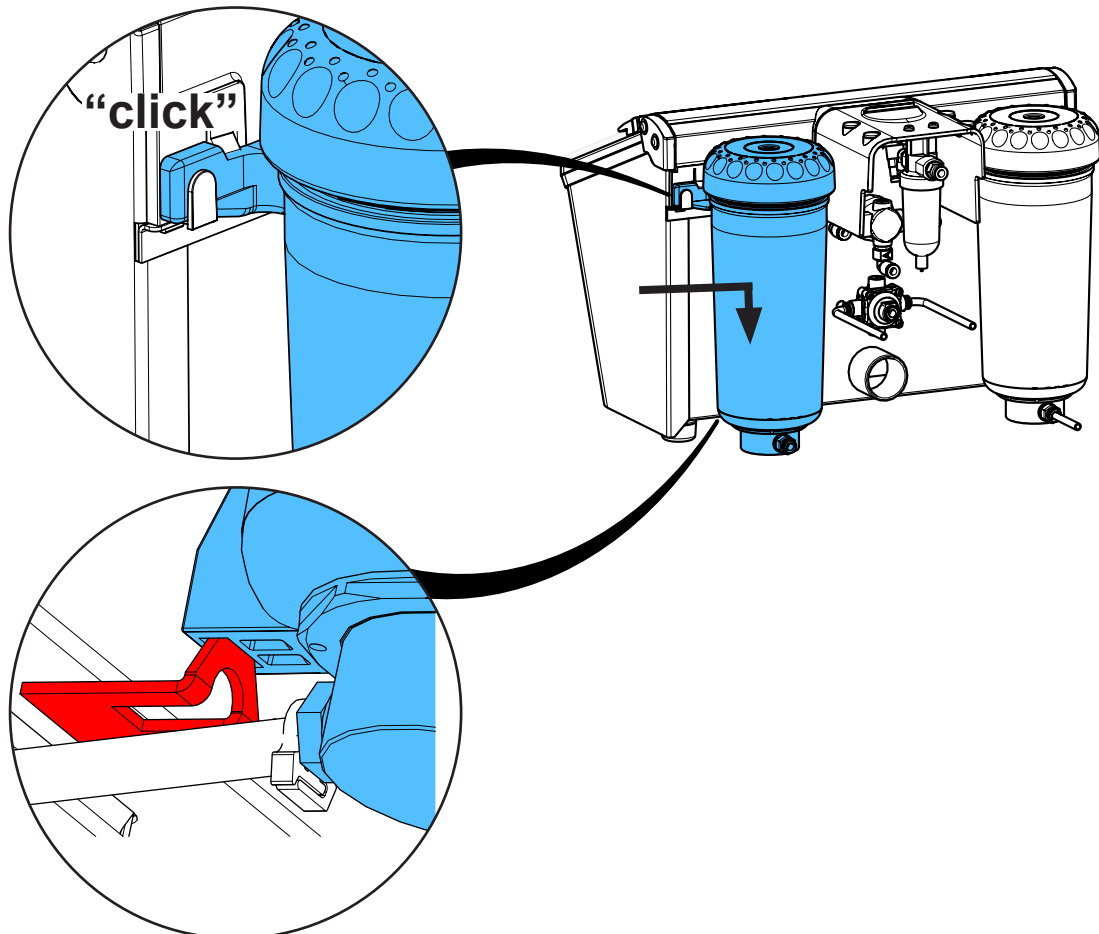

12.

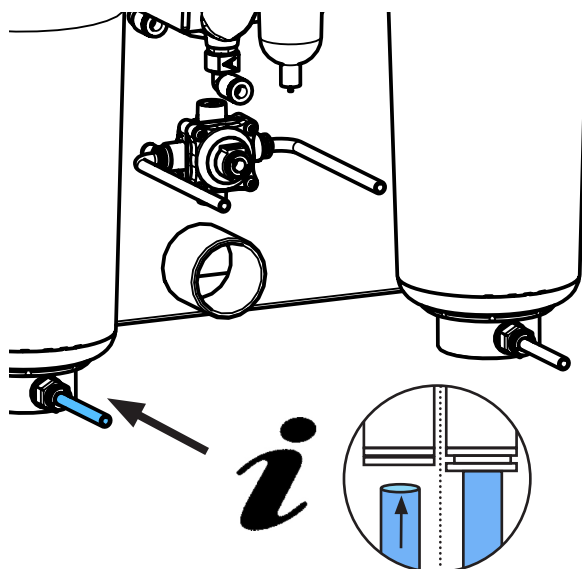

13.

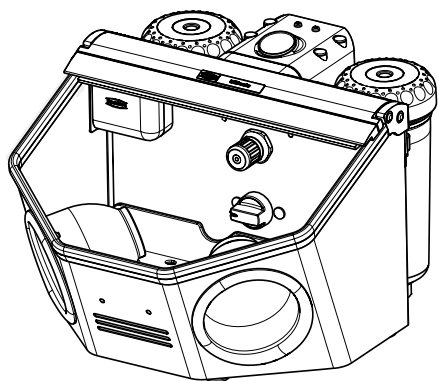






Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу


making work easy

**90003 6386**


Mischkammer • Mixing chamber • Chambre de malaxage • Camera di miscela •  
Camara de mezcla • Смесительная камера

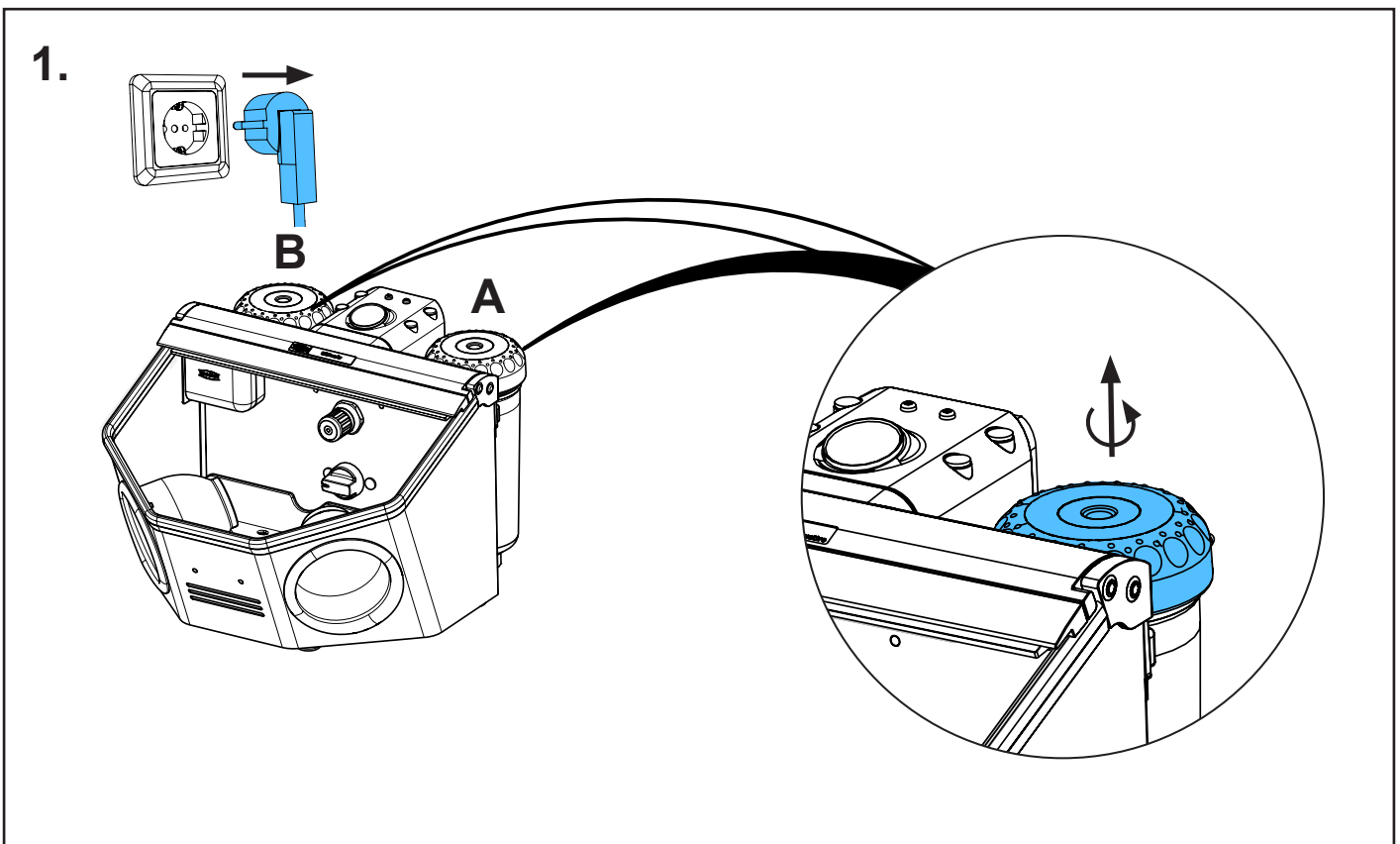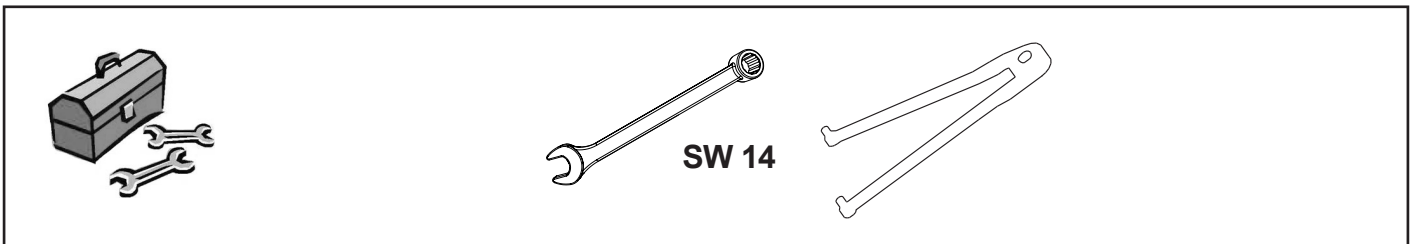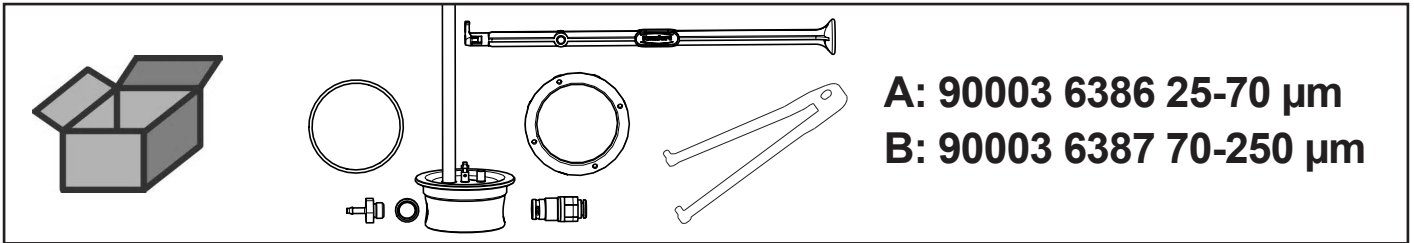


2.

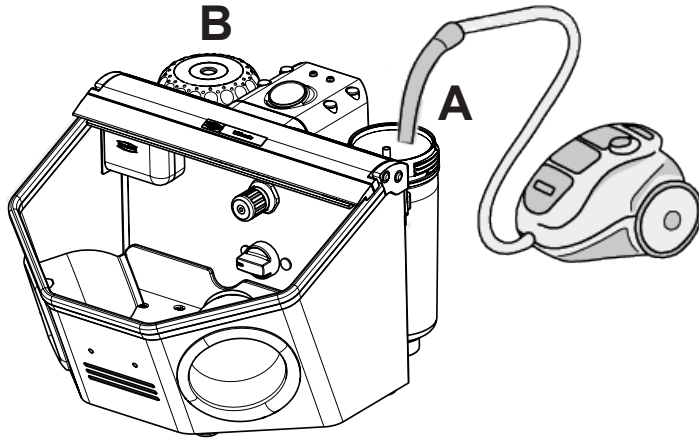

3.

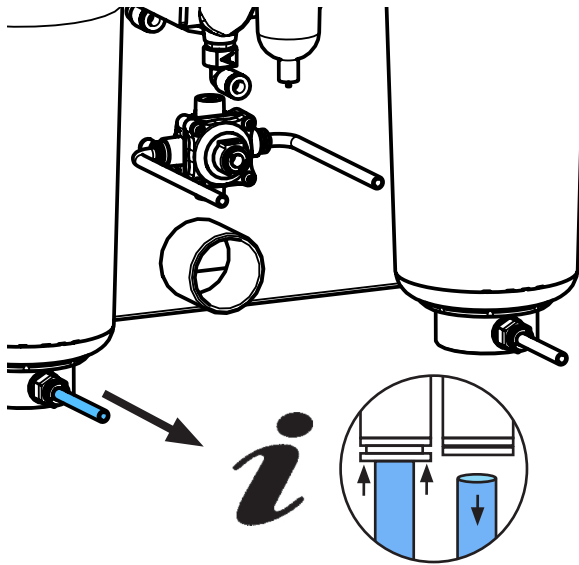


4.

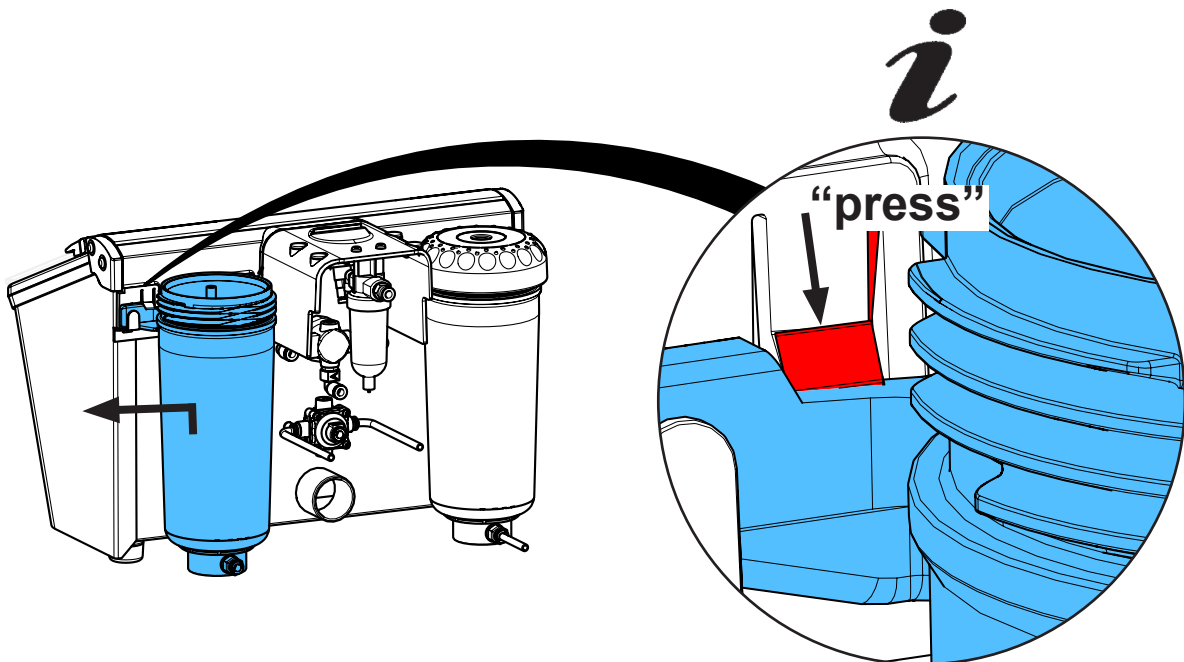





5.

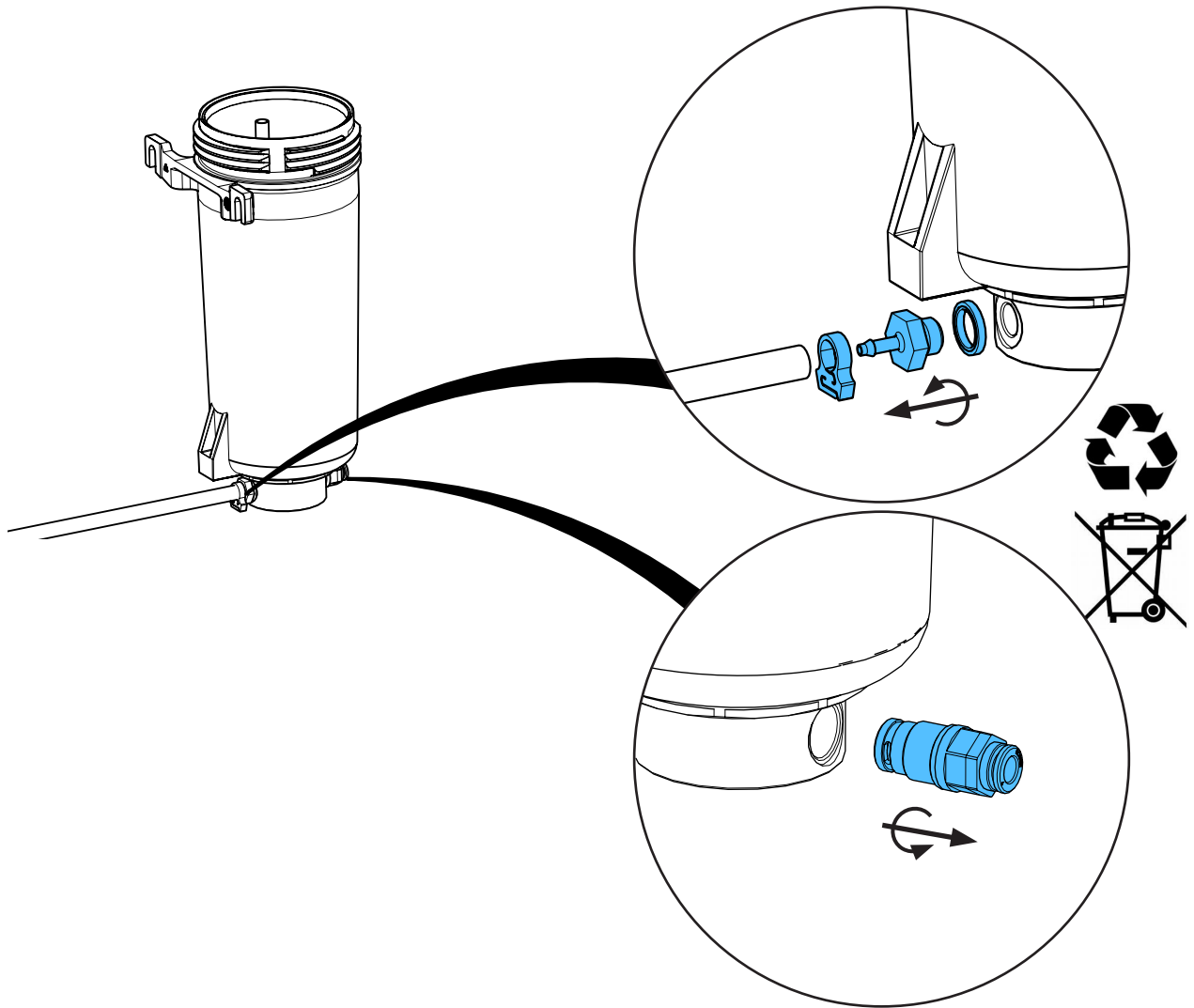


6.

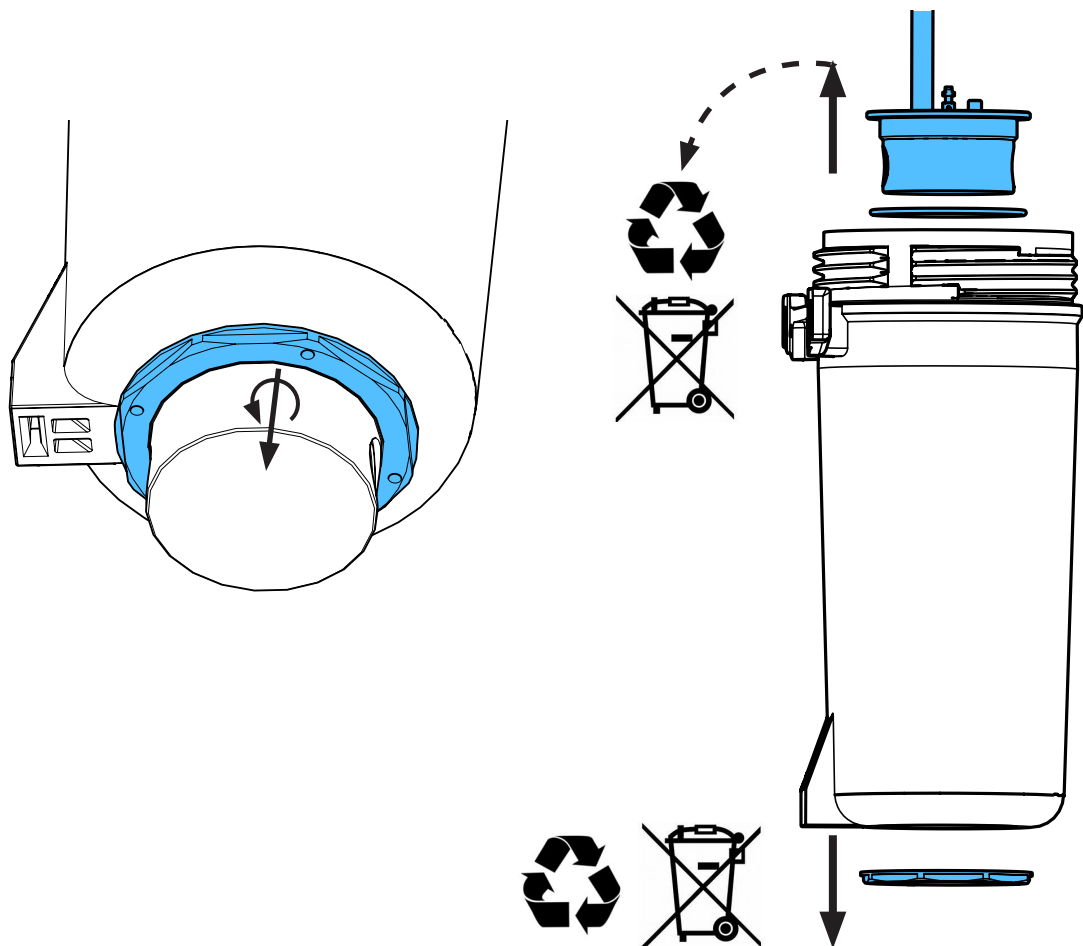


7.

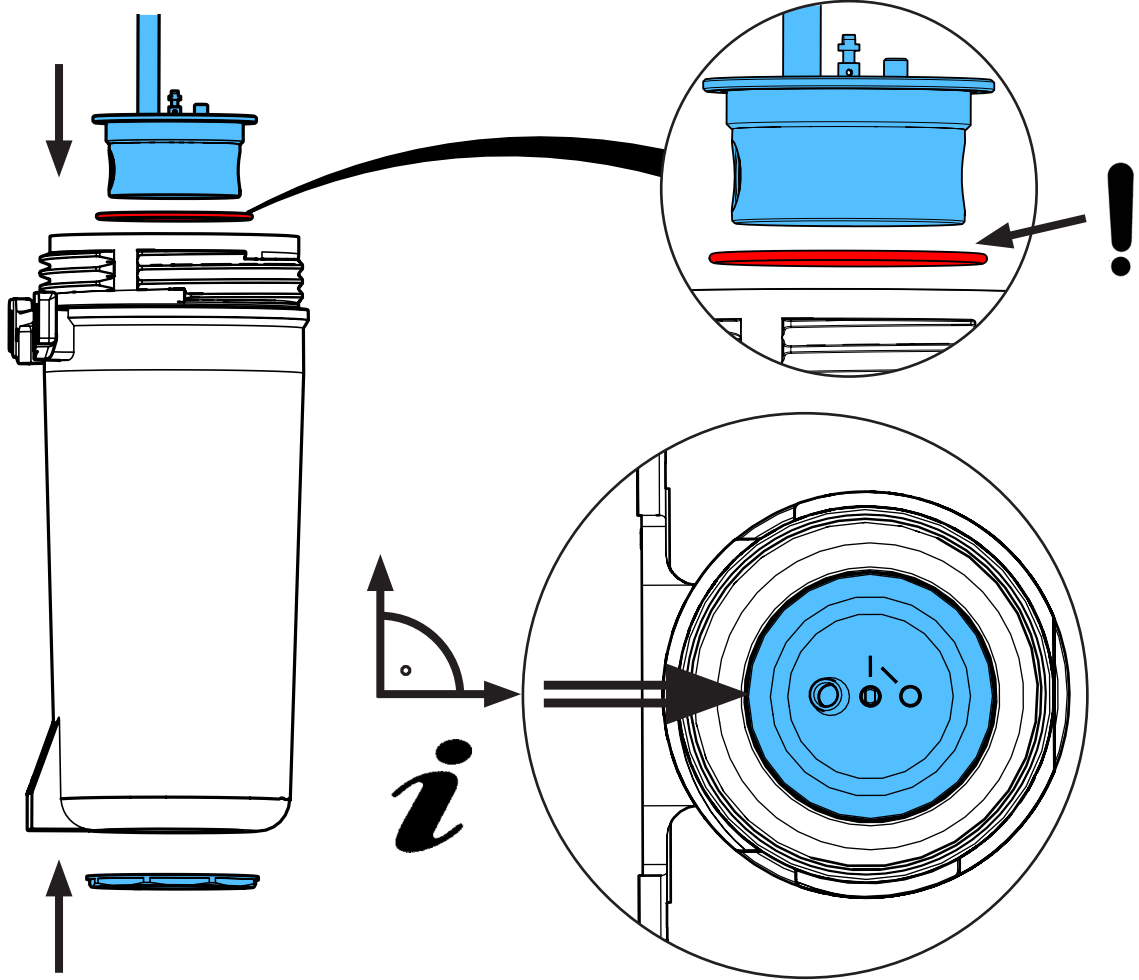


8.

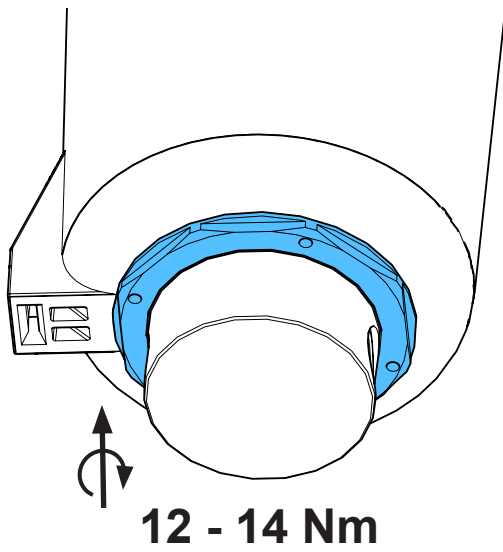


9.

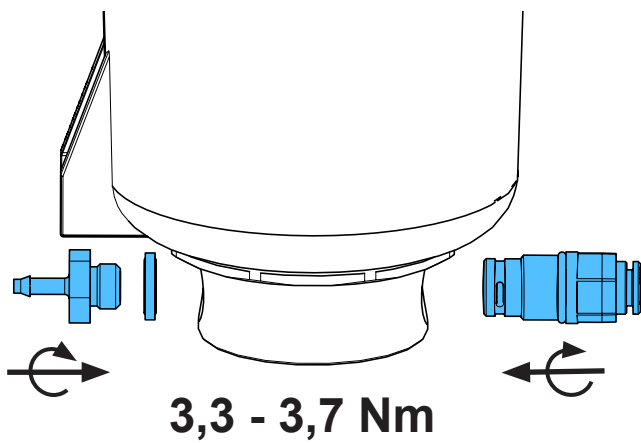


10.

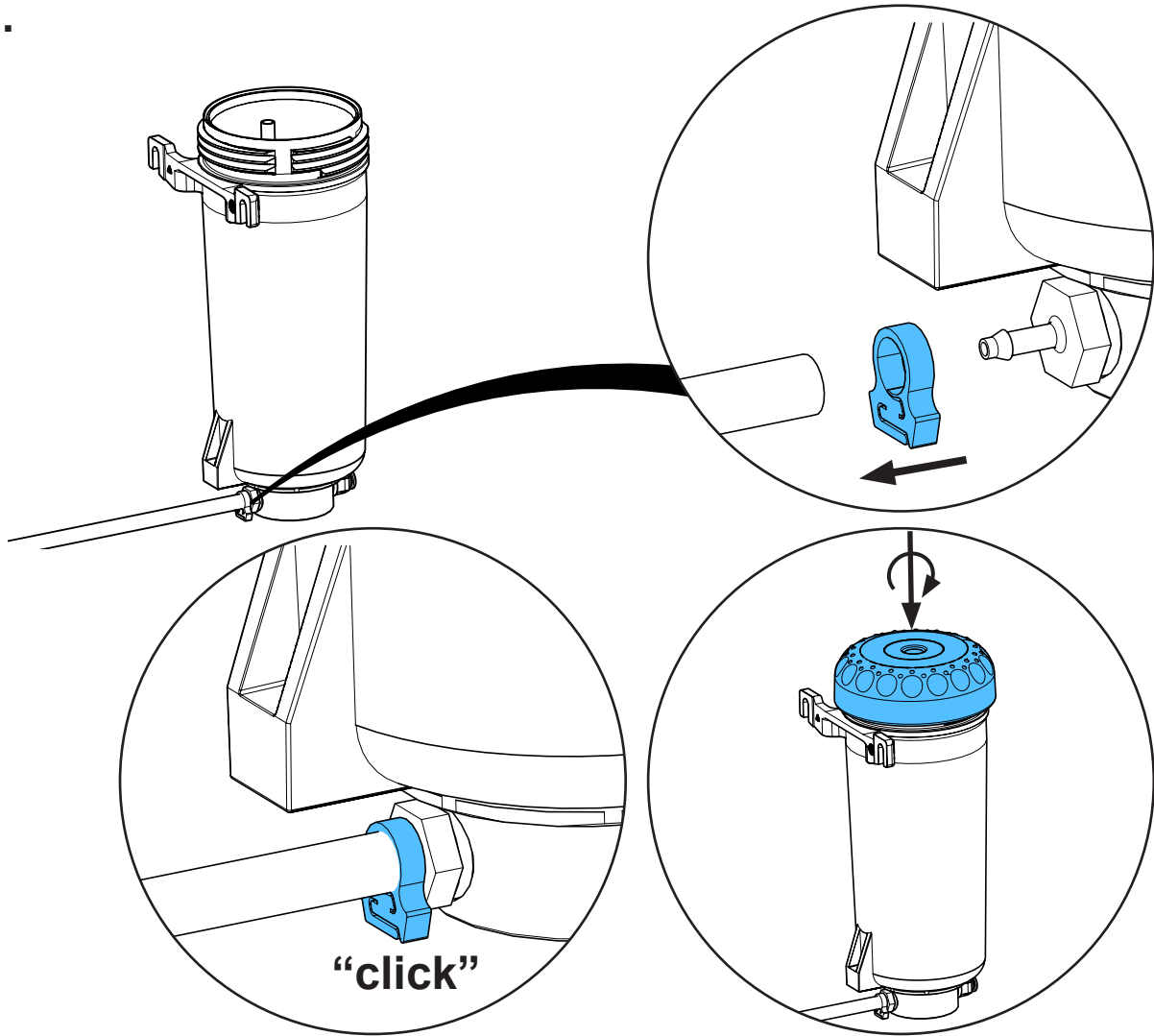


11.

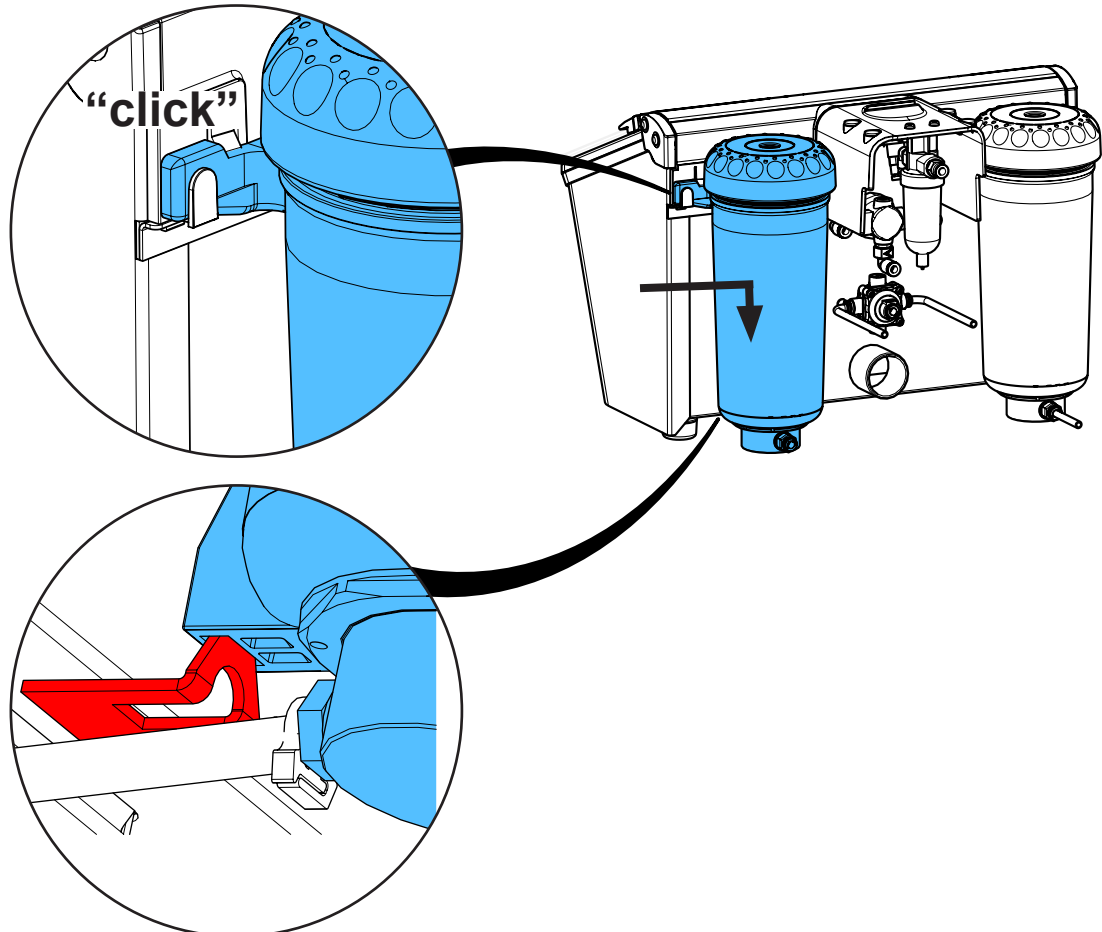

12.

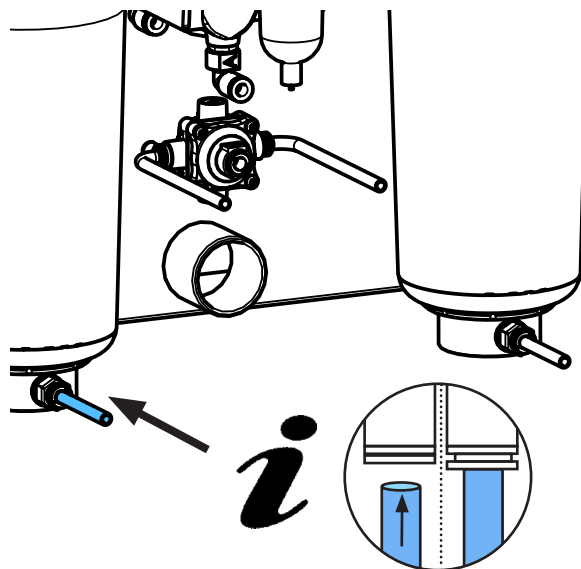

13.



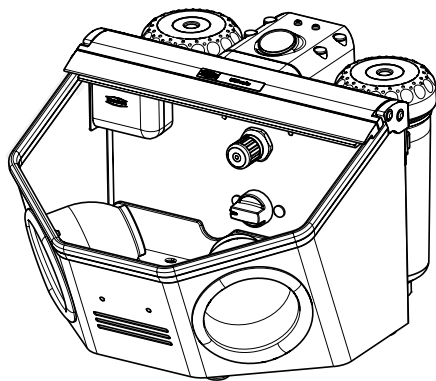





making work easy

## 2 1119


Manometer mit Verschraubung • Manometer with screw fitting • Manomètre avec  
raccord • Manometro con filettatura • Manómetro con atornilladura • Манометр с  
резьбовым соединением

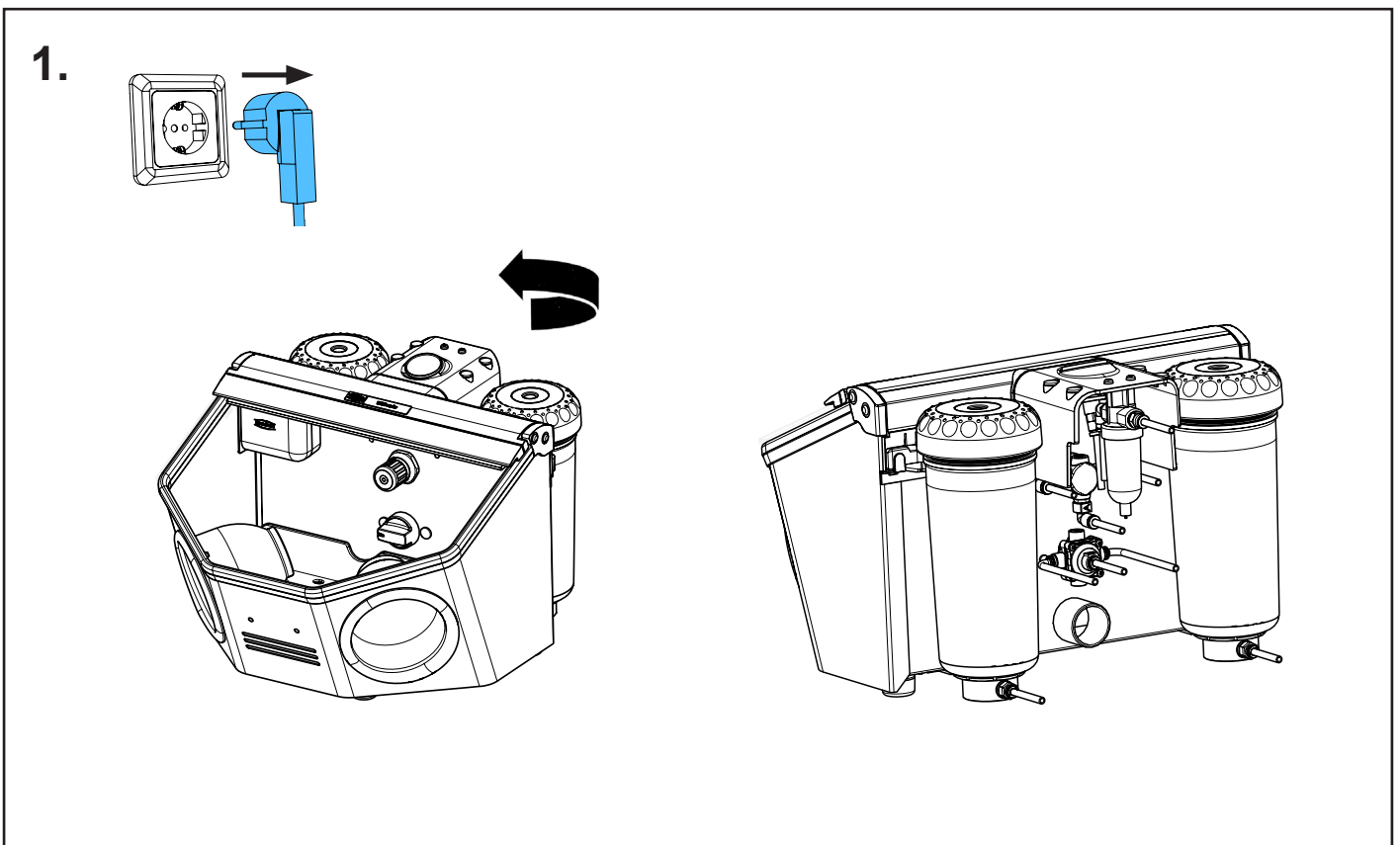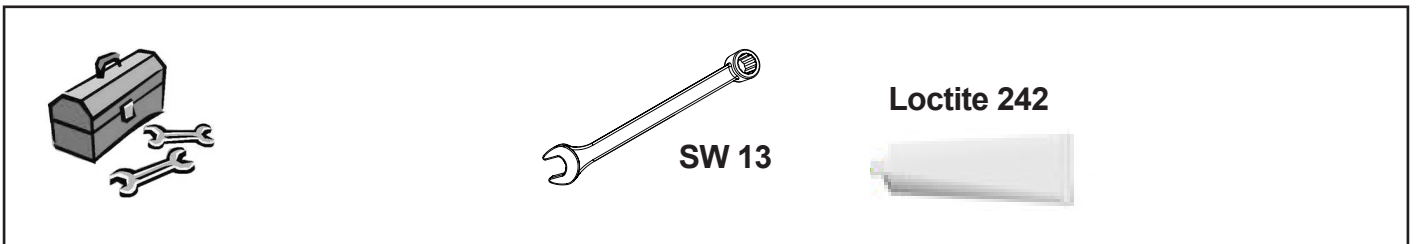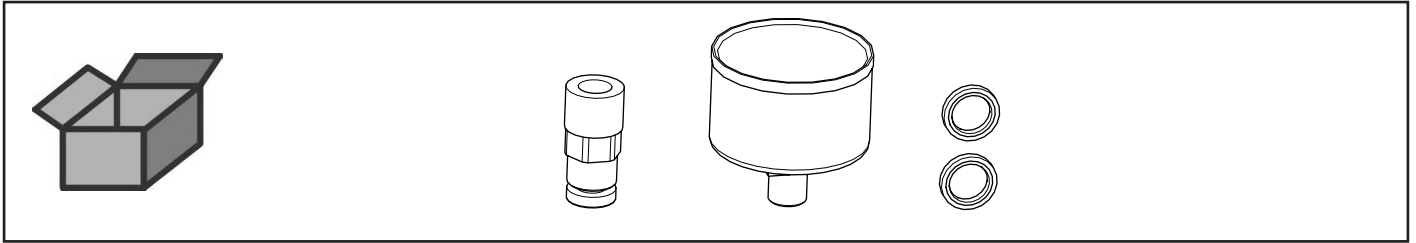


2.

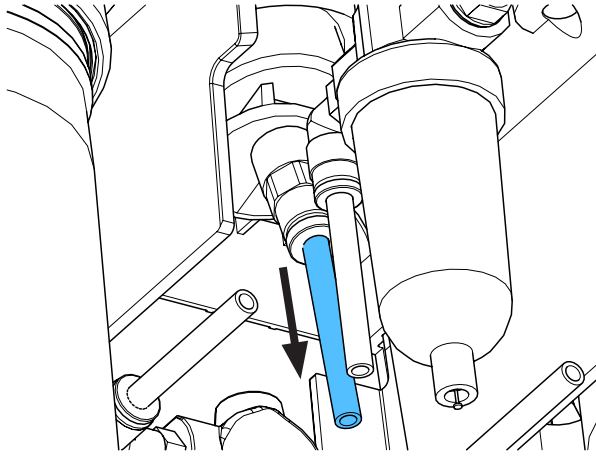



***i***

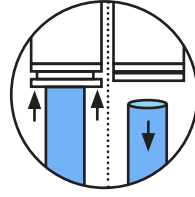

3.

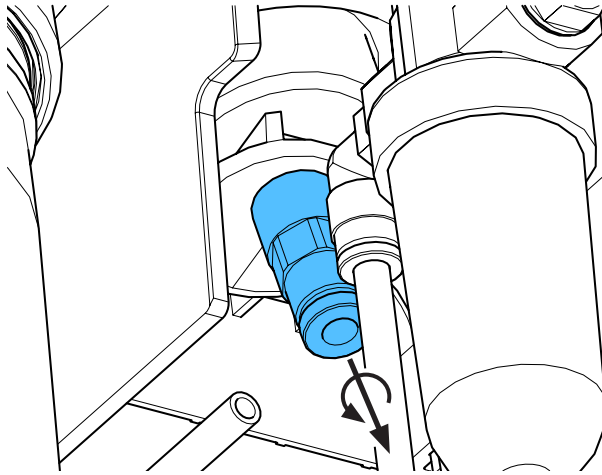




4.

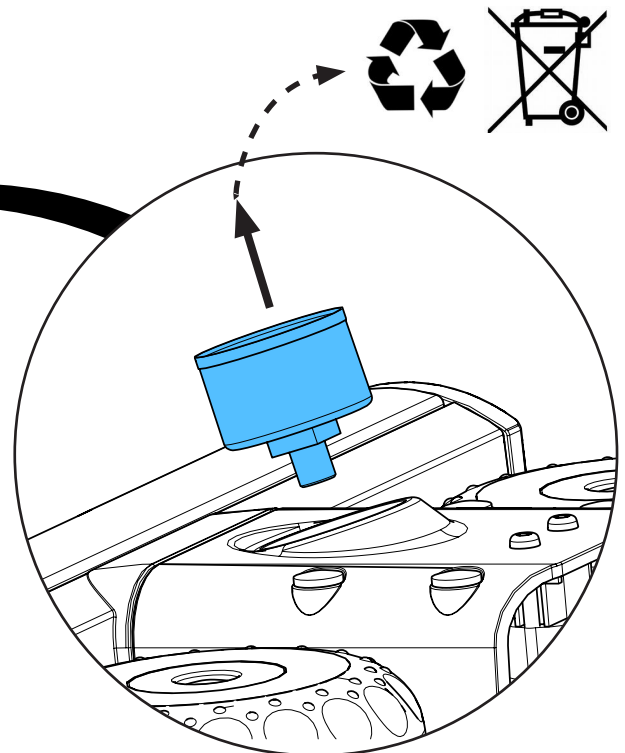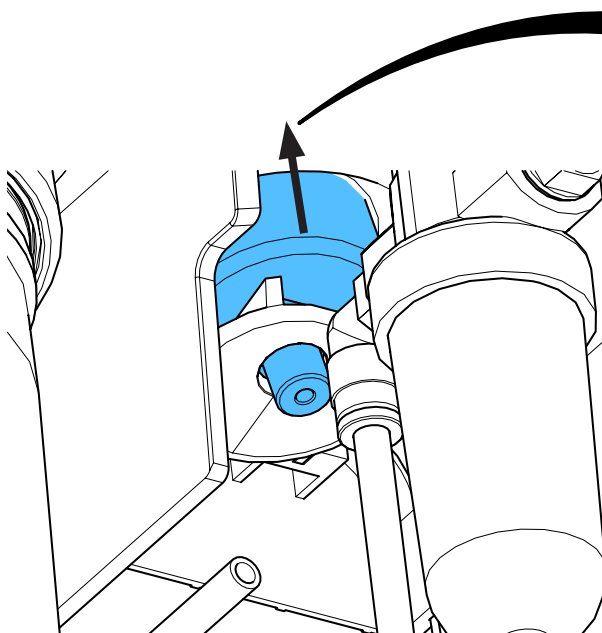


5.




6.

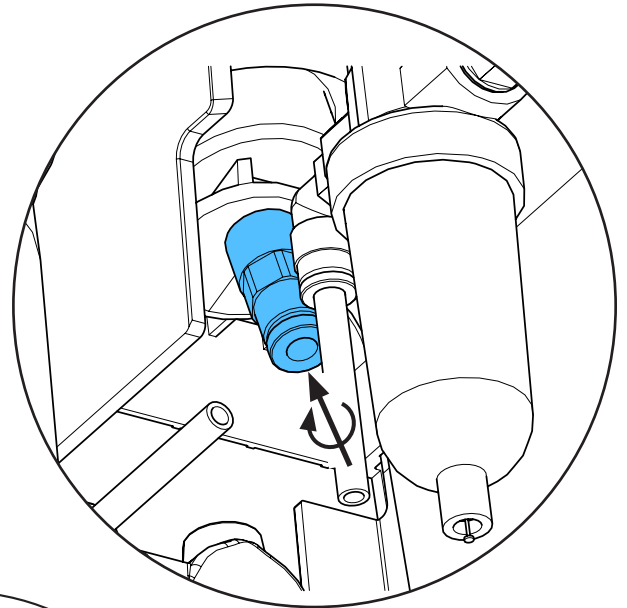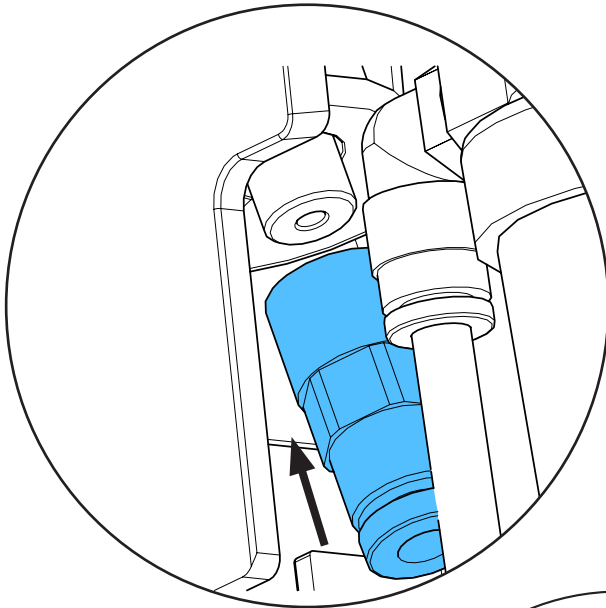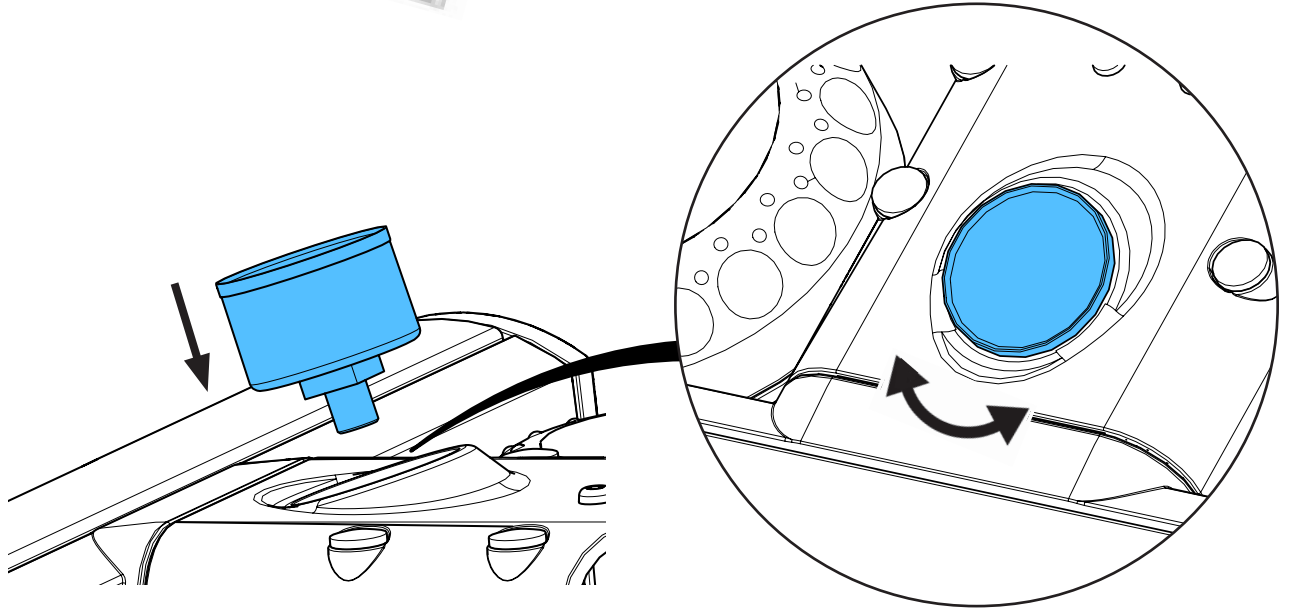


3,5 - 3,7 Nm

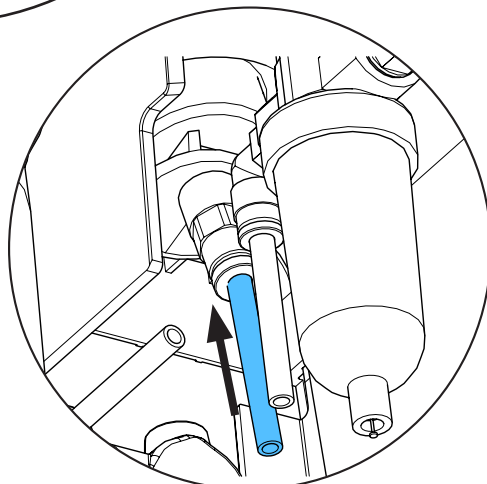

***i***

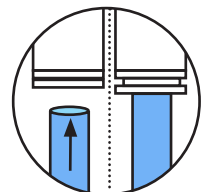

7.

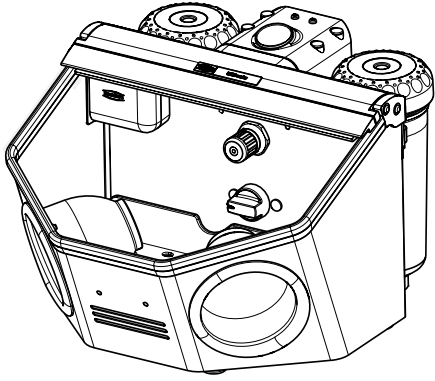








Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу


making work easy

**90002 1551**


Tankwahlschalter 2-fach • 2-fold tank selector switch • Commutateur de silo  
2-voies • Selettore di serbatoio 2-vie • Conmutador de depósito doble •  
Переключатель на 2 бачка

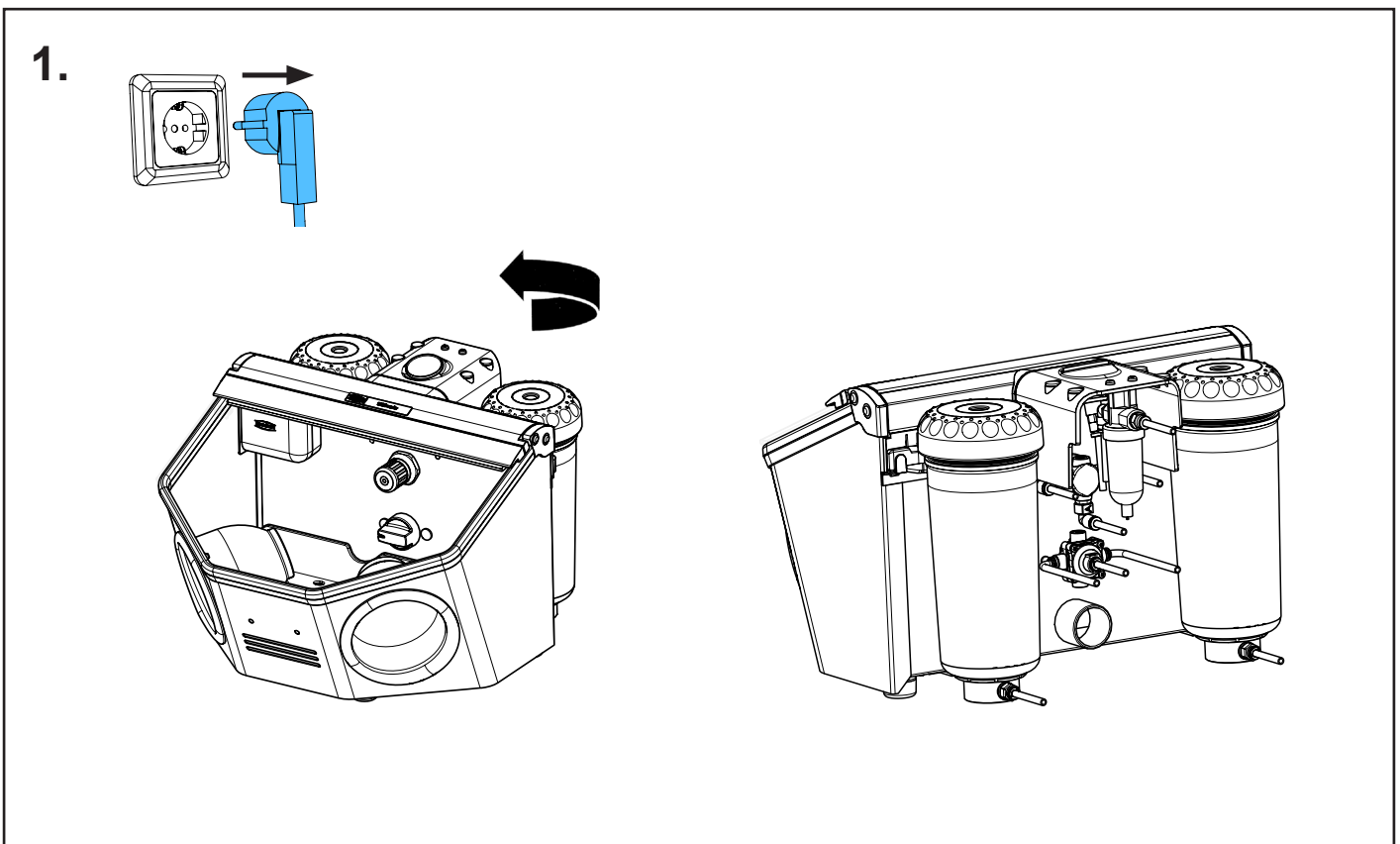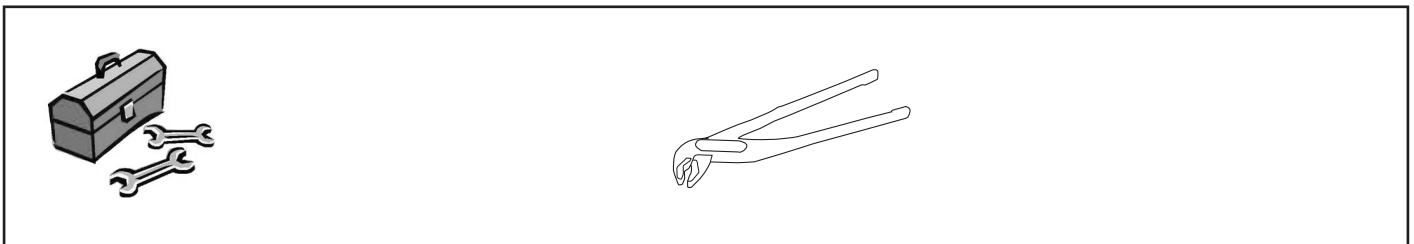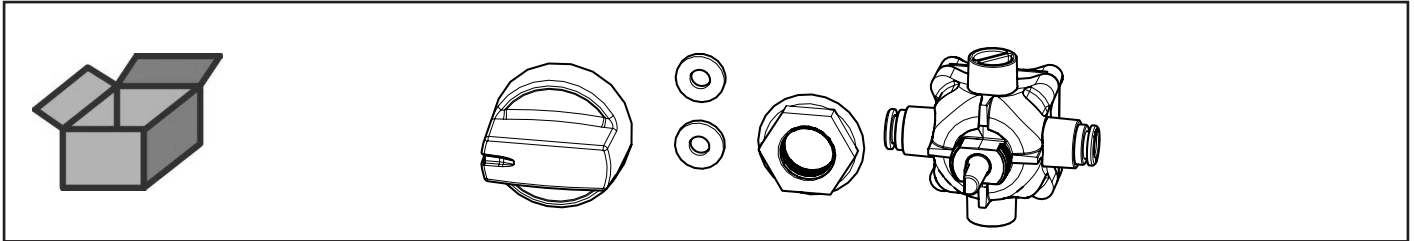


2.

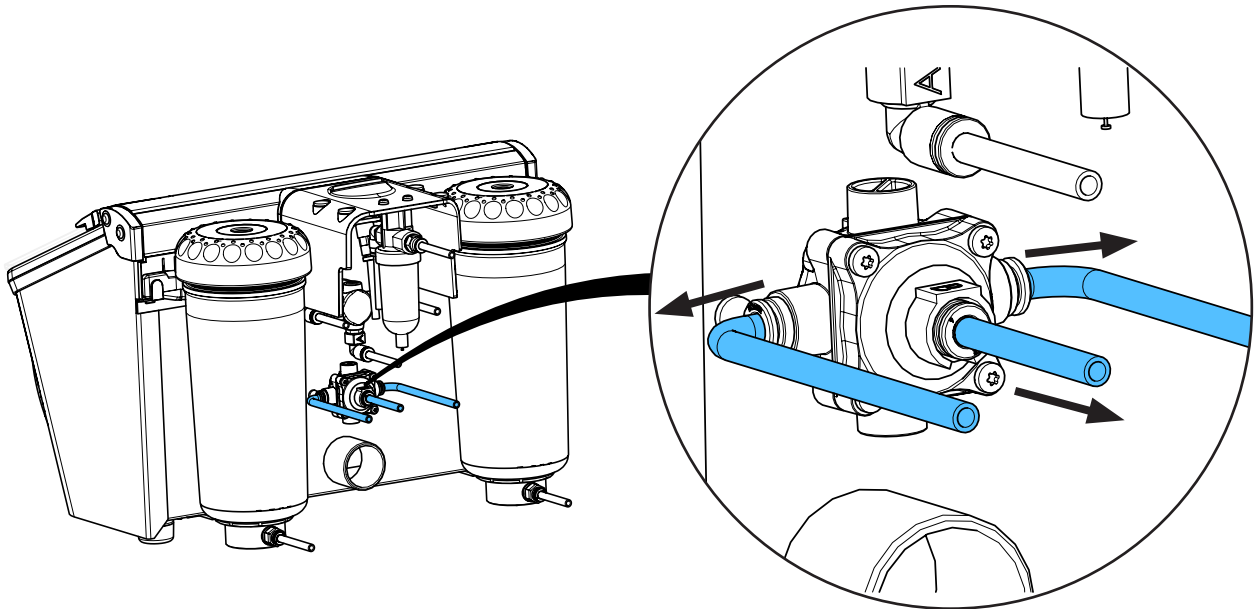




**i**

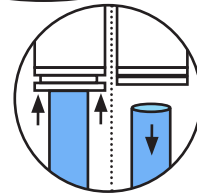



3.

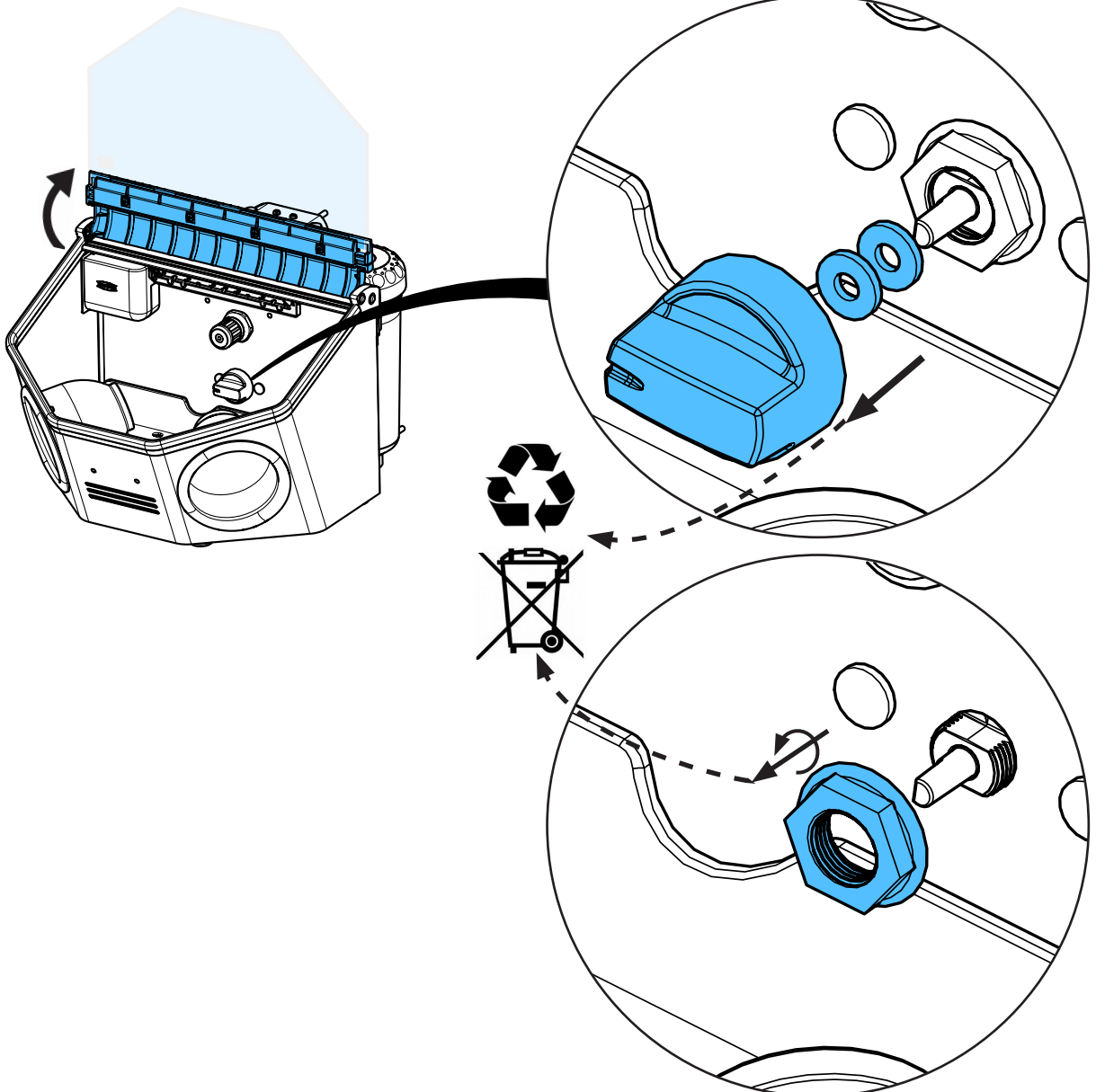




4.

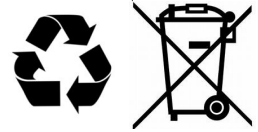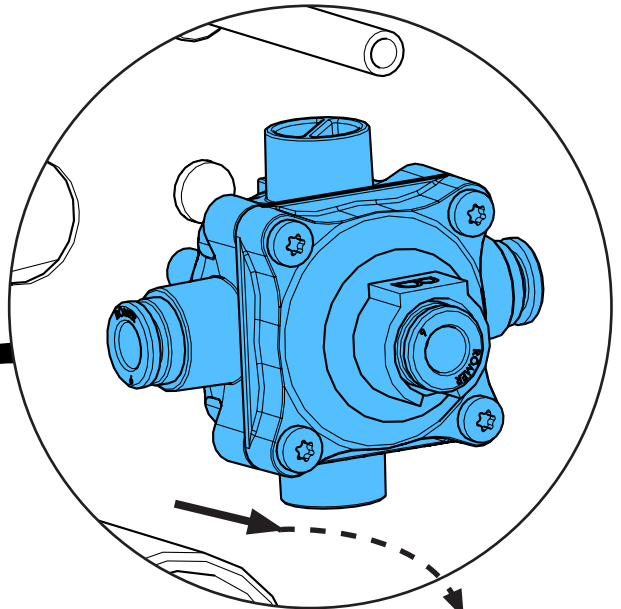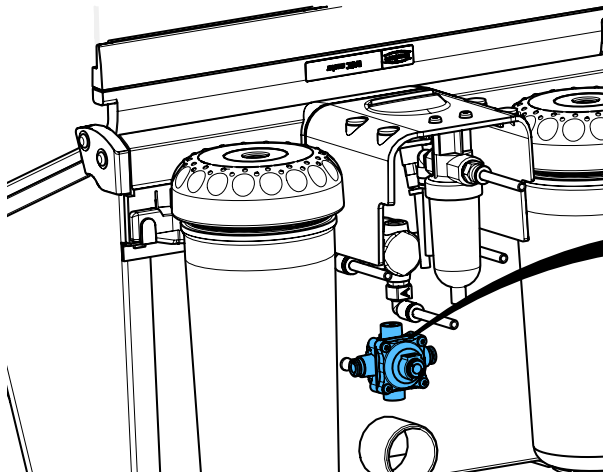


5.

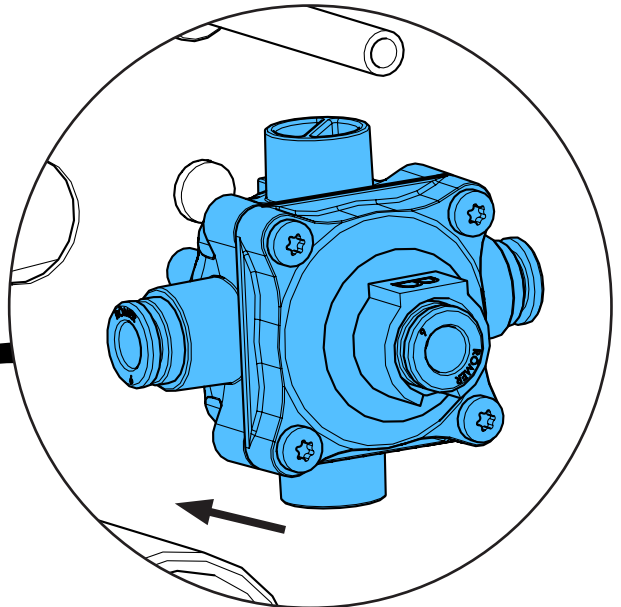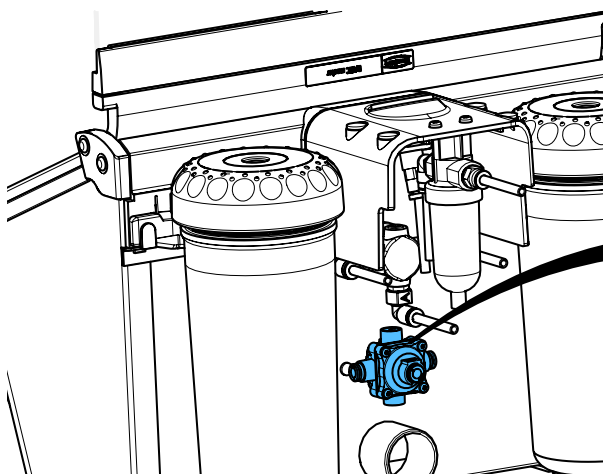


6.

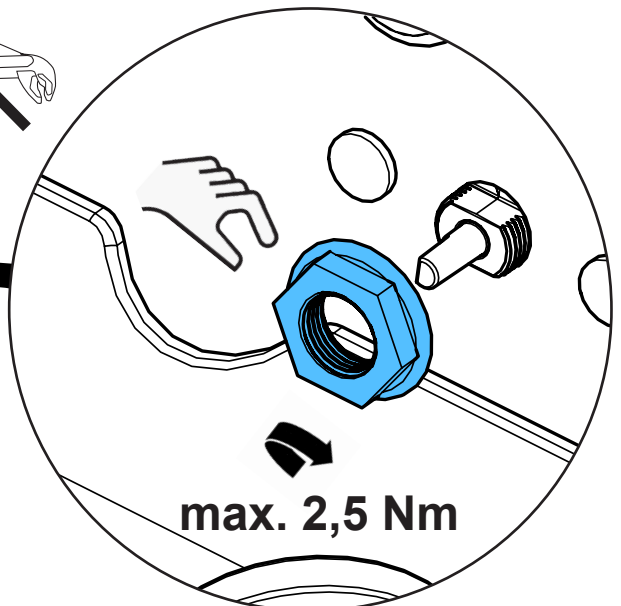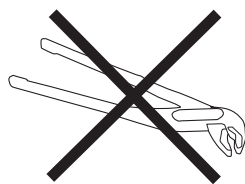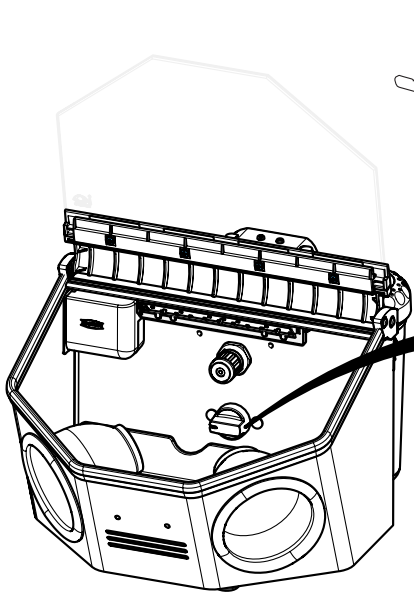


7.

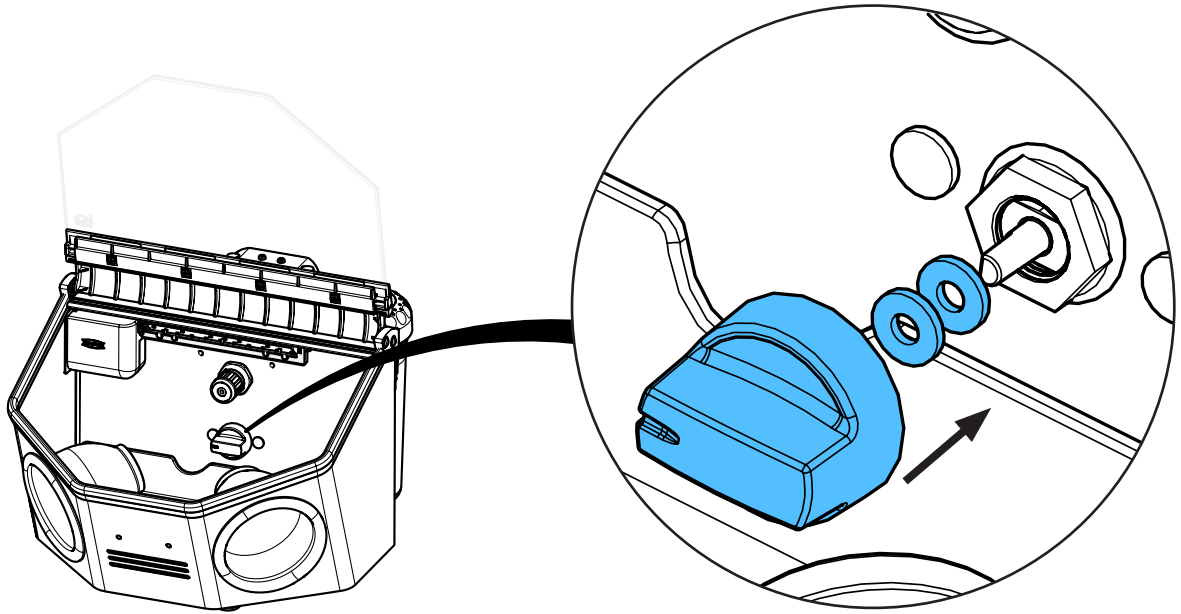


8.

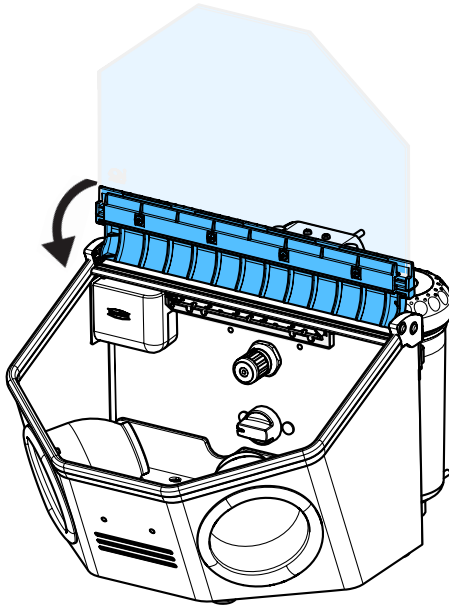


9.

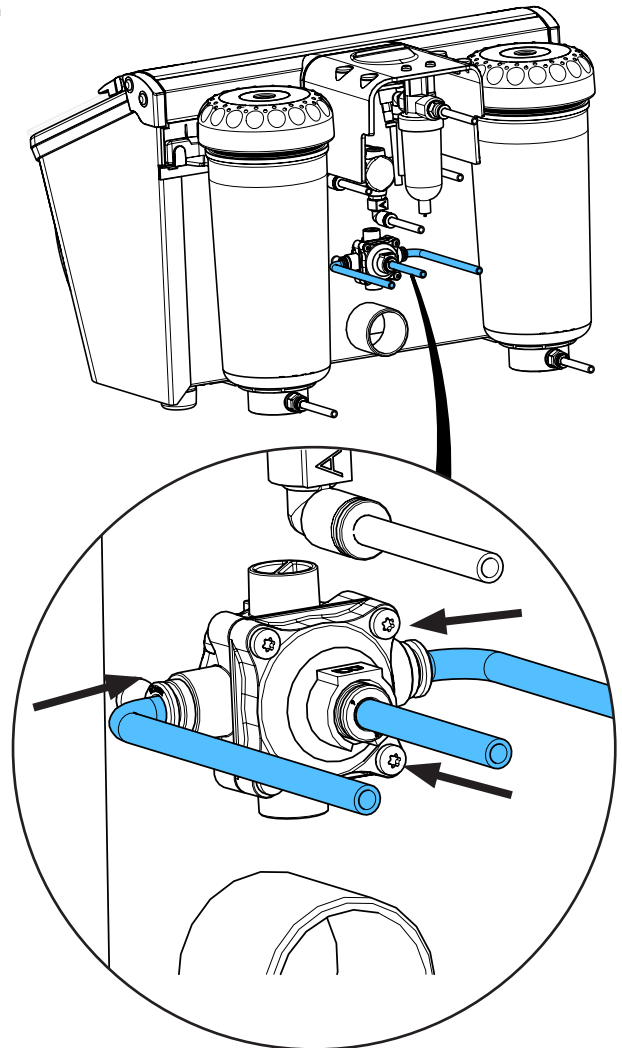

10.

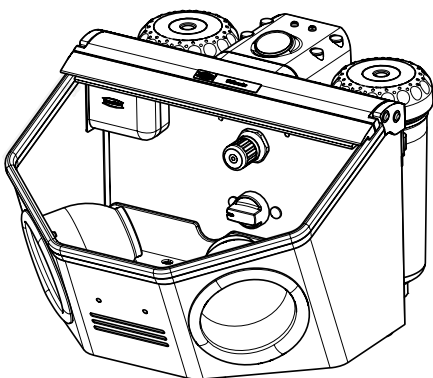

**i**

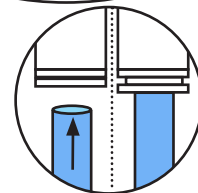







Montageanleitung • Assembly Instruction • Instruction d'installation  
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по  
монтажу


making work easy

**90003 6413**


Druckregler • Pressure regulator • Régulateur de pression • Regolatore di pressio-  
ne • Regulador de presión • Регулятор давление

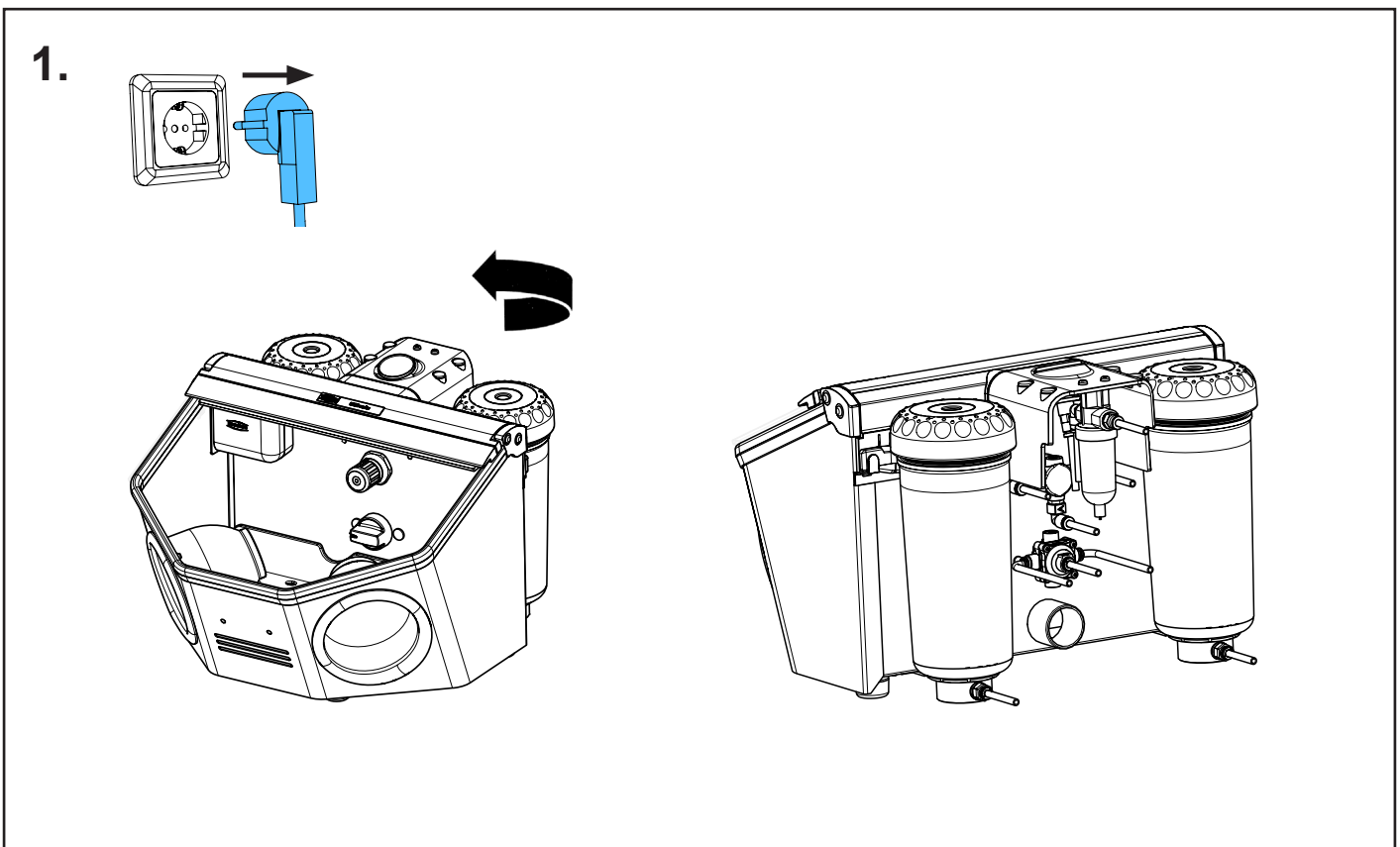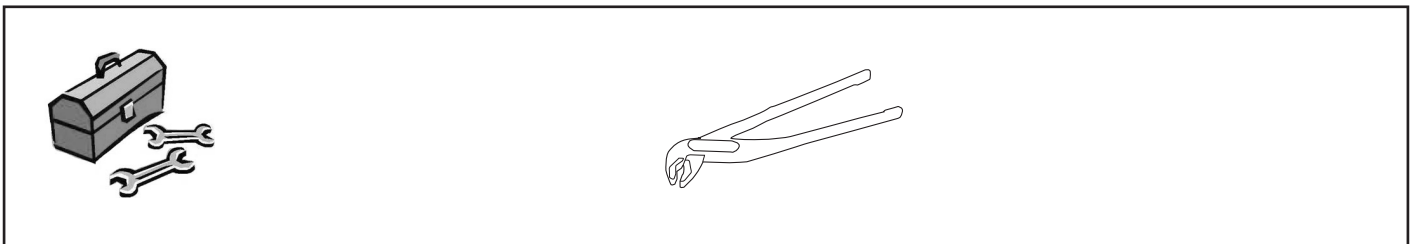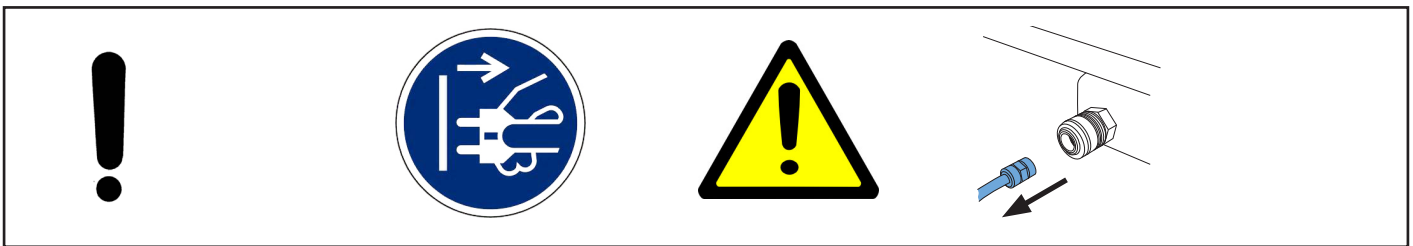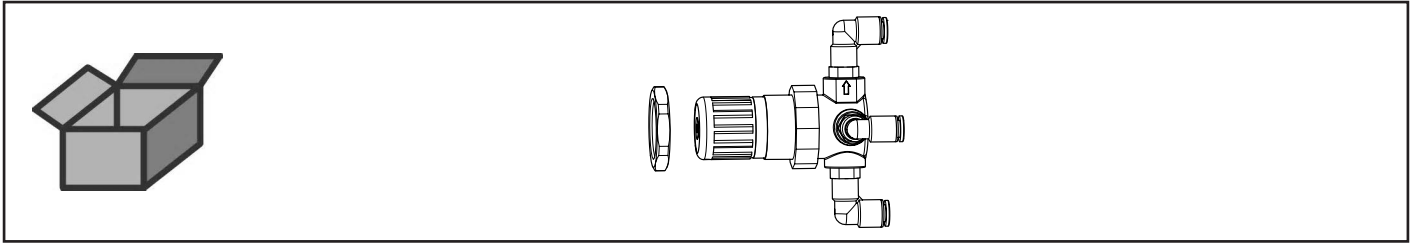


2.

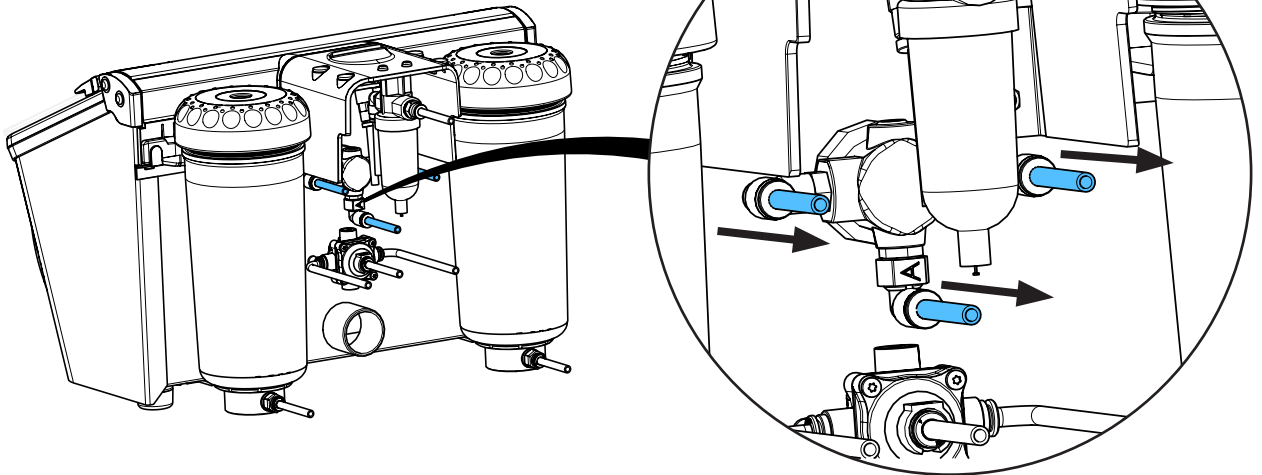


**i**

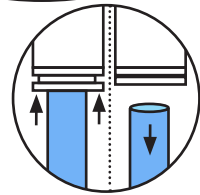


3.

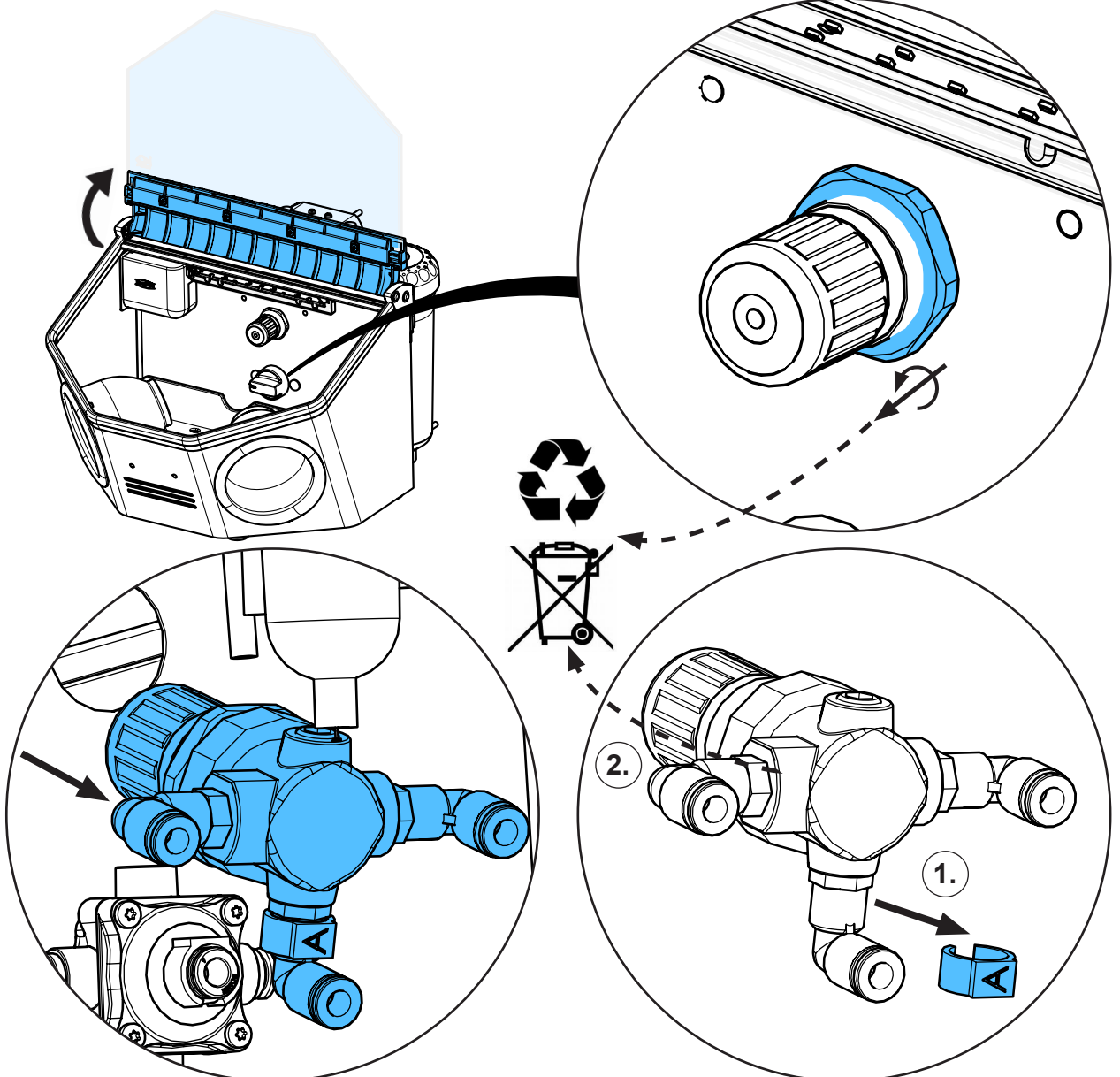


4.

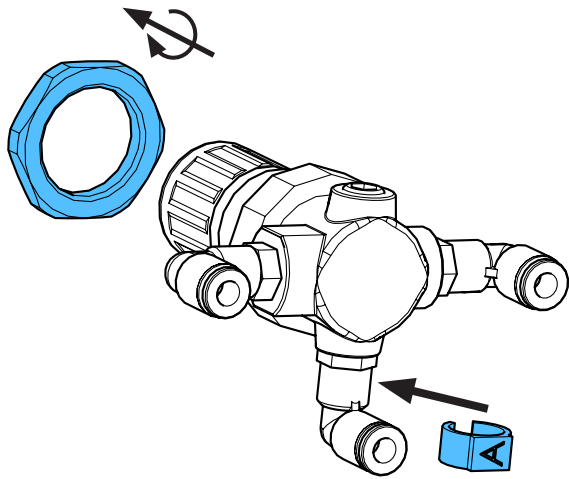


5.

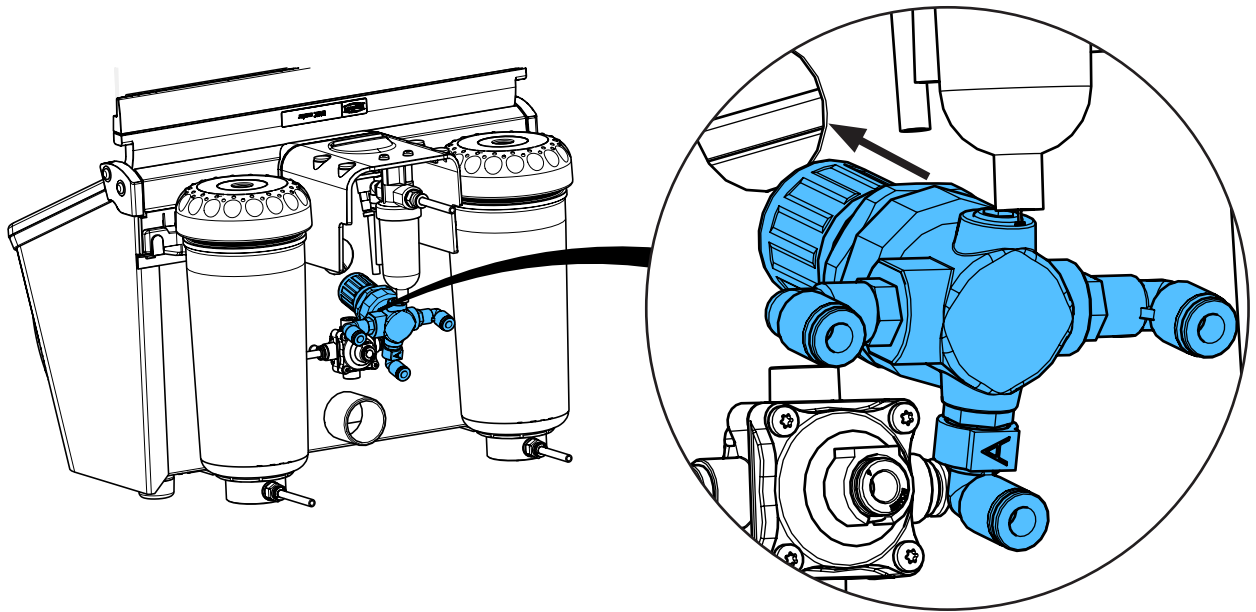


6.

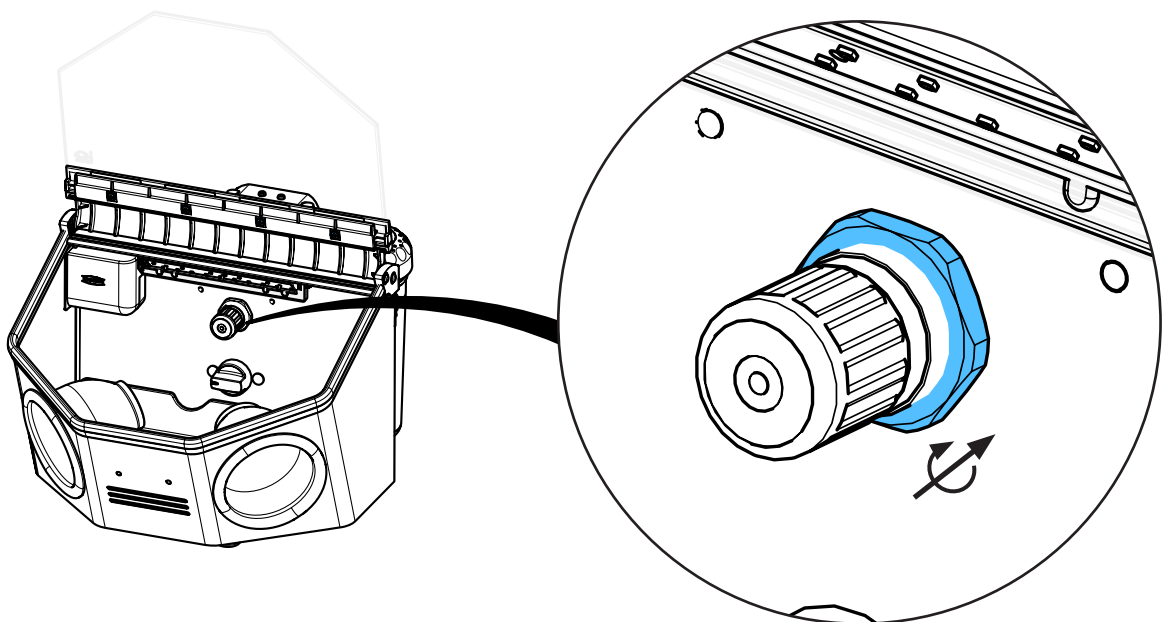


7.

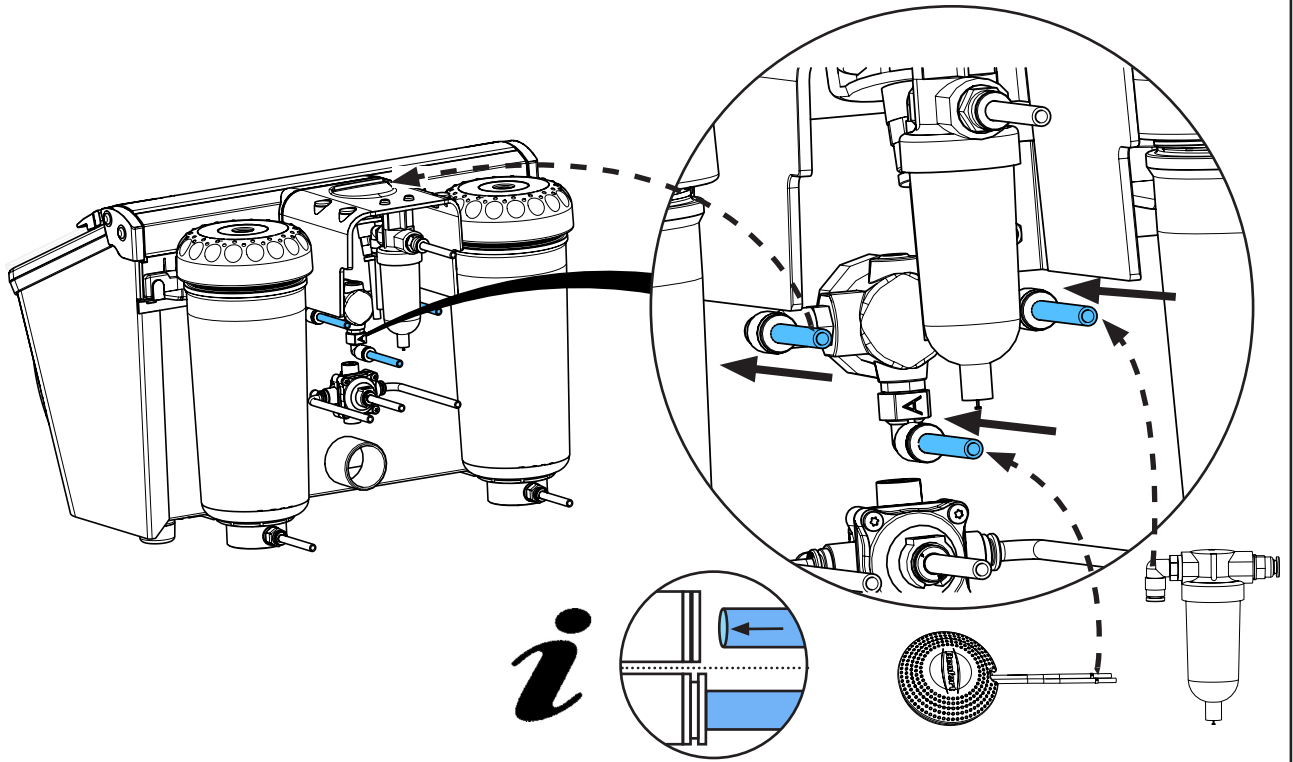

8.

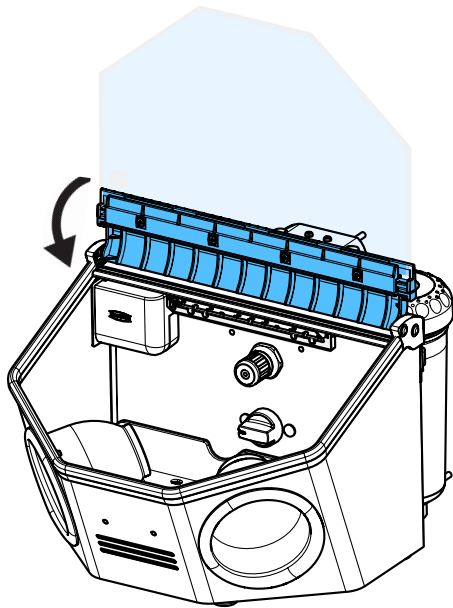

9.

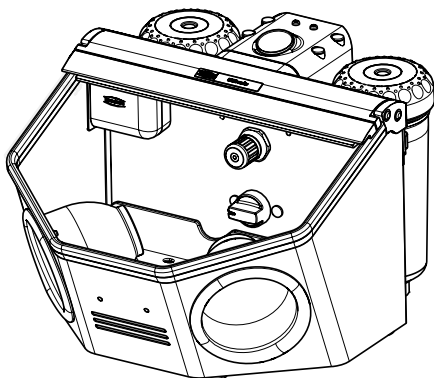





**Hochaktuell und ausführlich auf ...  
Up to date and in detail at ...  
Actualisé et détaillé sous ...  
Aggiornato e dettagliato su ...  
La máxima actualidad y detalle en ...  
Актуально и подробно на ...**

**[www.renfert.com](http://www.renfert.com)**

Renfert GmbH • Untere Gießwiesen 2 • 78247 Hilzingen/Germany  
Tel.: +49 7731 82 08-0 • Fax: +49 7731 82 08-70  
[www.renfert.com](http://www.renfert.com) • [info@renfert.com](mailto:info@renfert.com)

Renfert USA • 3718 Illinois Avenue • St. Charles IL 60174/USA  
Tel.: +1 6307 62 18 03 • Fax: +1 6307 62 97 87  
[www.renfert.com](http://www.renfert.com) • [info@renfertusa.com](mailto:info@renfertusa.com)  
USA: Free call 800 336 7422