



making work easy



EASY view+ 3D

Service Instructions

Made in Germany

20-5086 08112022

Inhalt

90003 7807

Haltearm • Boom arm • Bras de support • Braccio di supporto • Brazo de sujeción • Кронштейн

90003 7808

3D Monitor • 3D monitor • Écran 3D • Monitor 3D • Monitor 3D • Монитор 3D

90002 1891

Kopfteil • Head • Tête • Testata • Cabezal • Головка

90003 7818

Kleinteileset • Hardware kit • Petit matériel de montage • Kit di minuteria
Piezas pequeñas • Комплект мелких деталей

90003 7809

3D Brille • 3D goggles • Lunettes 3D • Occhiali 3D • Gafas 3D • Очки 3D

90003 7368

Maus • Mouse • Souris • Mouse • Ratón • Мышь

90003 3776

Tischauflage • Desk pad • Sous-main • Piano di appoggio • Apoyo para la mesa
Настольная подложка

90003 7810

Fußschalter • Foot switch • Pédale • Interruttore a pedale • Interruptor de pie
Педаль

90003 7900

HDMI-DVI Kabel • HDMI-DVI cable • Câble HDMI DVI • Cavo HDMI DVI • Cable HDMI DVI •
Кабель HDMI-DVI

90003 7883

Netzteil 3D Monitor • Power supply unit 3D monitor • Adaptateur secteur écran 3D
Alimentatore monitor 3D • Adaptador de corriente monitor 3D • Блок питания 3D монитора

90003 4769

Netzteil LED-Beleuchtung • Power supply LED lighting • Poste secteur LED lampe • Trasfor-
matore lampada LED • Bloque alimentación lampara LED • Блок питания со светодиодной
подсветкой

90003 5975

Netzkabel • Power cord • Câble alimentation • Cavo di rete • Cable de corriente
Сетевой кабель

90003 3758

Schutzglas • Protective glass • Verre de protection • Vetro protettivo • Cristal de protección •
Защитное стекло

90003 7903

HDMI-HDMI Kabel • HDMI-HDMI cable • Câble HDMI HDMI • Cavo HDMI HDMI
Cable HDMI HDMI • Кабель HDMI-HDMI

90003 7895

USB Stick • USB stick • Clé USB • Chiave USB • Lápiz de memoria USB
USB-накопитель

90003 7552

Universalnetzsteckerset • Universal power plug kit • Jeu universel de fiches secteur • Set spine
di rete universale • Kit de enchufes eléctricos universales • Сетевой штекер универсальный

90003 7916

Steckdosenleiste • Multiple socket • Bloc multiprise • Presa multipla • Regleta de enchufe •
3-кратный разъем

90003 7909

Netzwerkverlängerung • Network extension cable • Rallonge de câble réseau • Prolunga per
cavo di rete • cable alargador de red • Удлинитель сетевого интернет-кабеля

90003 3775

HDMI Video Capture • Карта видеозахвата HDMI Video Capture

90003 7778

Höheneinstellung inkl. Klemmschrauben • Height adjustment incl. clamping screw
Réglage de hauteur avec vis de serrage • Regolazione altezza inc. viti di bloccaggio
Ajuste en altura incl. tornillo de fijación • Регулятор высоты вкл. зажимные винты

90003 3759

LED-Modul • LED module • Module DEL • Modulo LED • Módulo LED • Модуль светодиодов

90003 3760




LED-Treiberleiterkarte • LED driver card • Circuit imprimé pilote DEL • Circuito stampato driver
LED • Circuito impreso de controlador LED • Плата драйвера светодиодов




making work easy

90003 7807


Haltearm • Boom arm • Bras de support • Braccio di supporto • Brazo de sujeci-
ón • Кронштейн

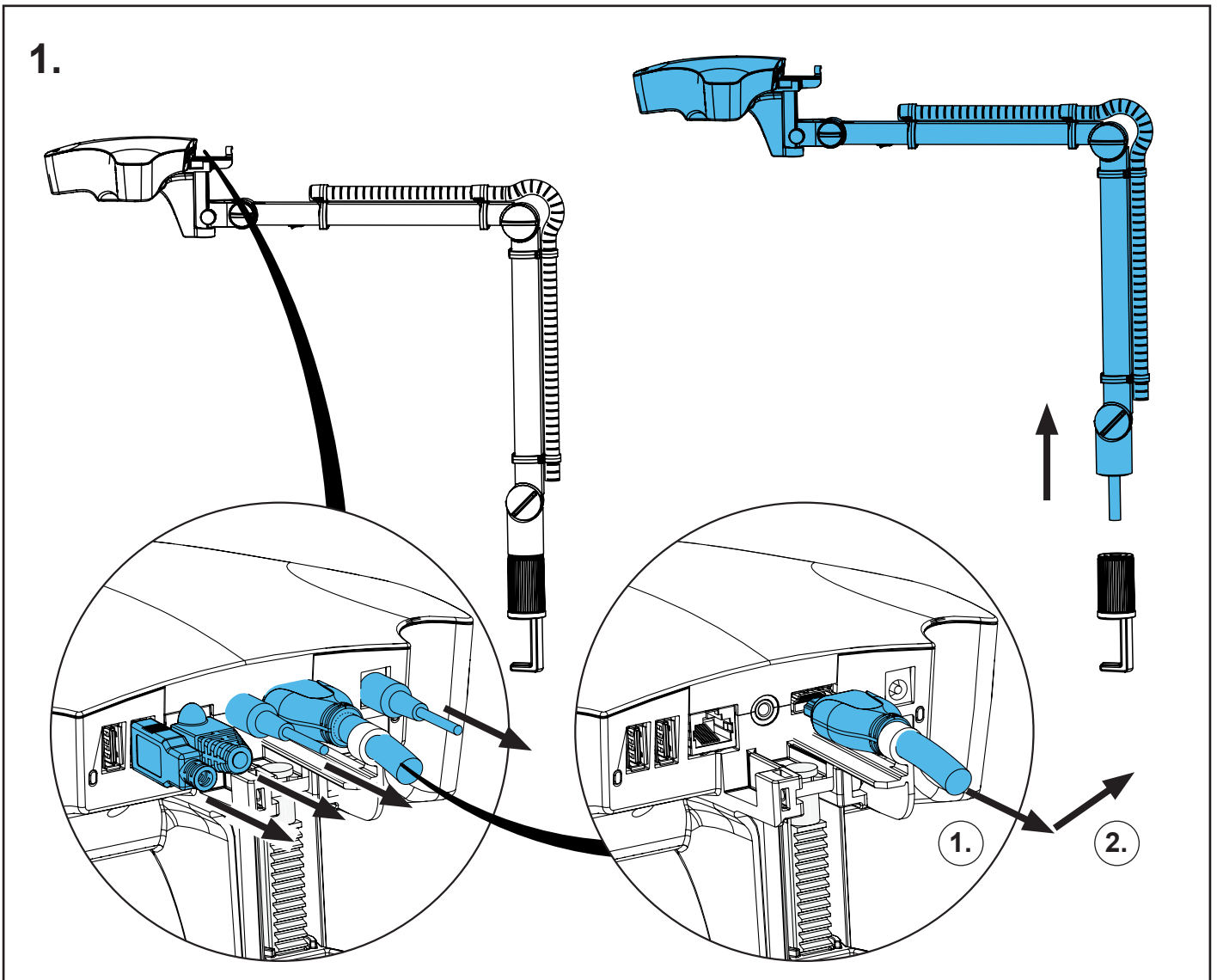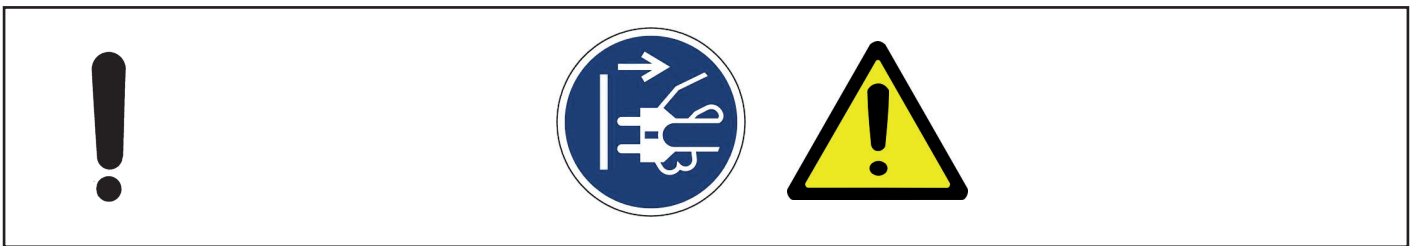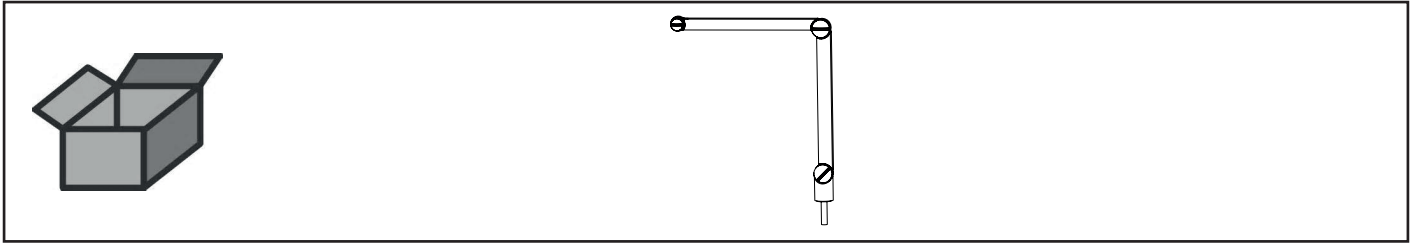


2.

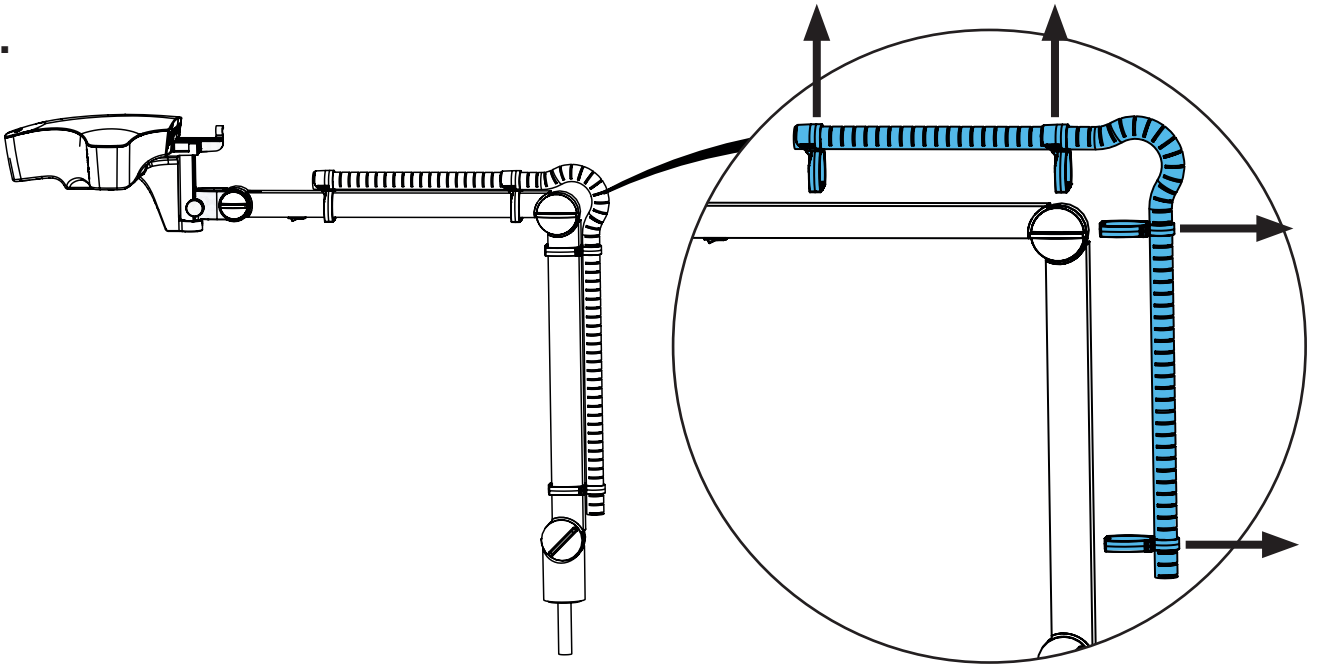


3.

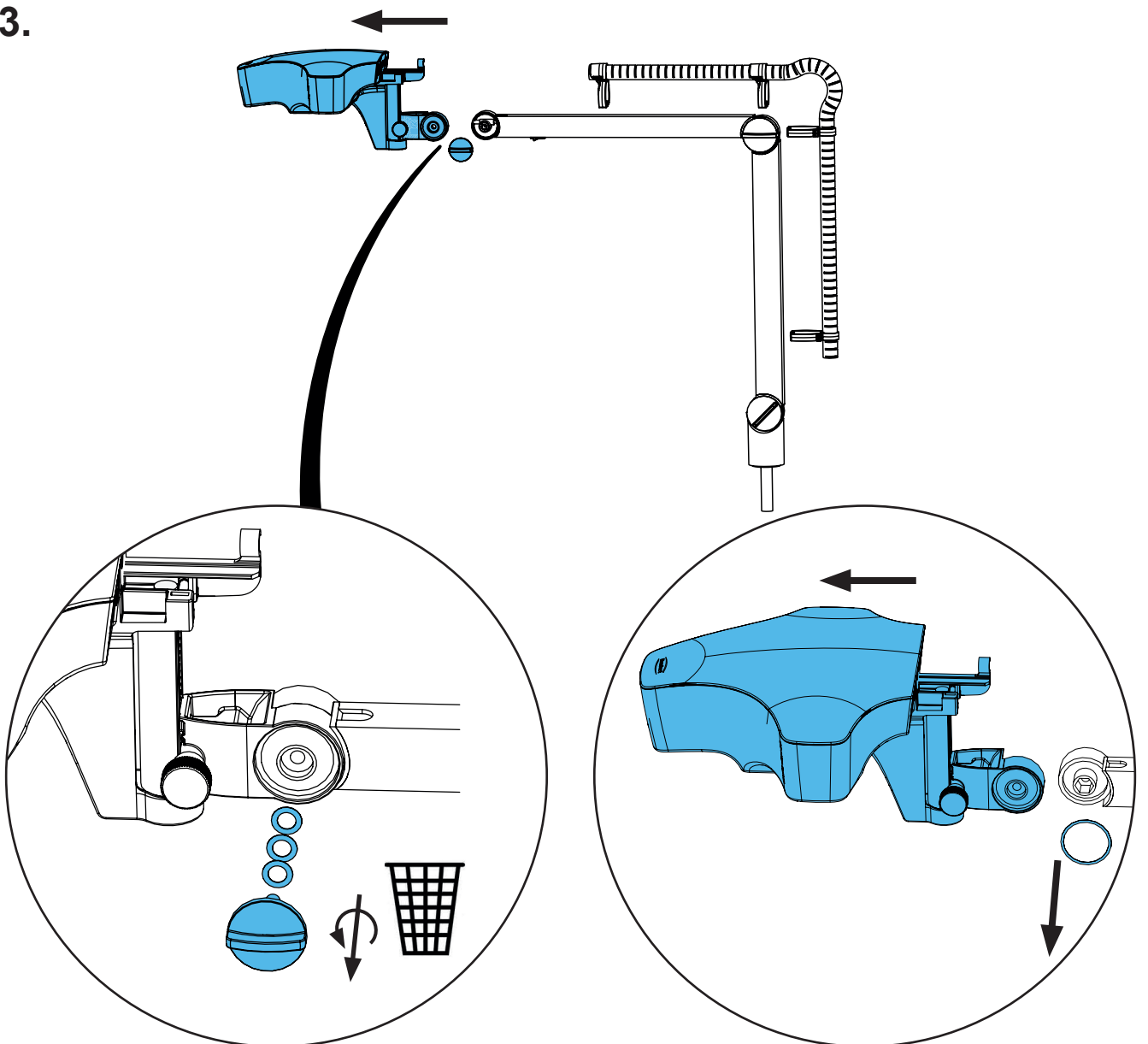


4.

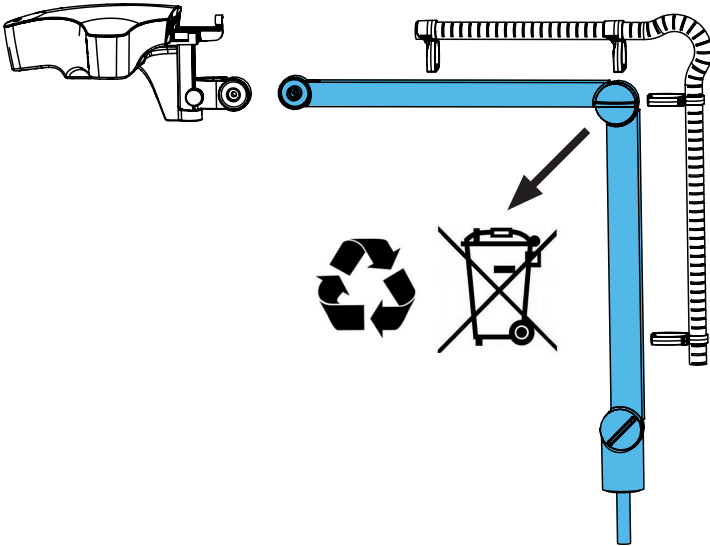


5.

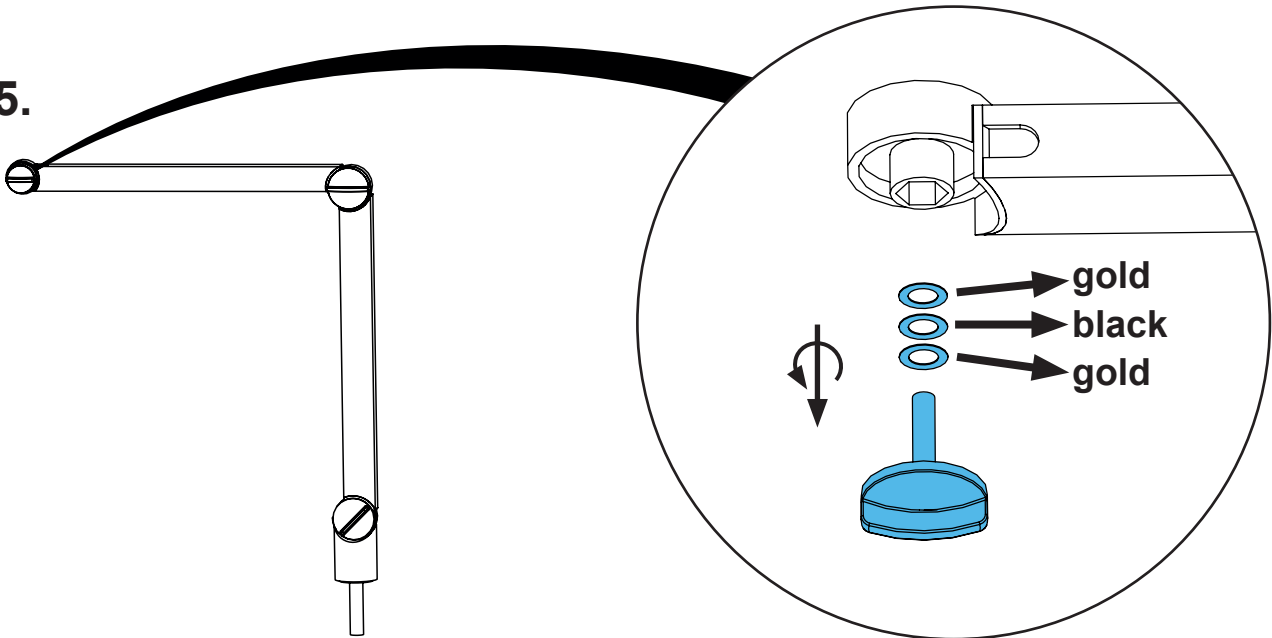


6.

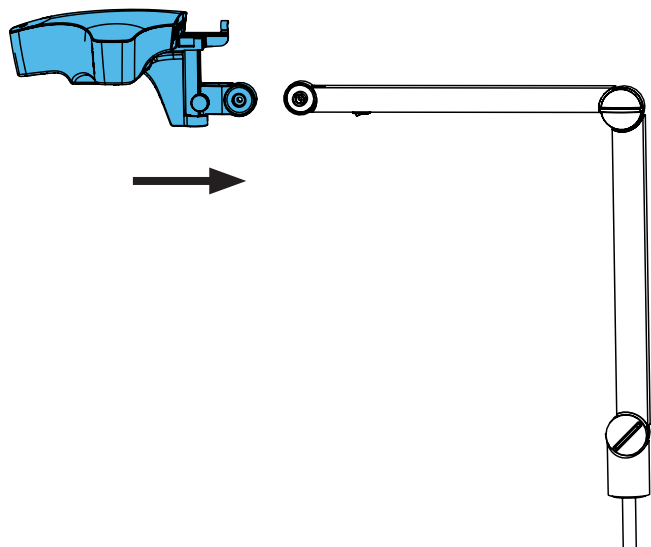


7.

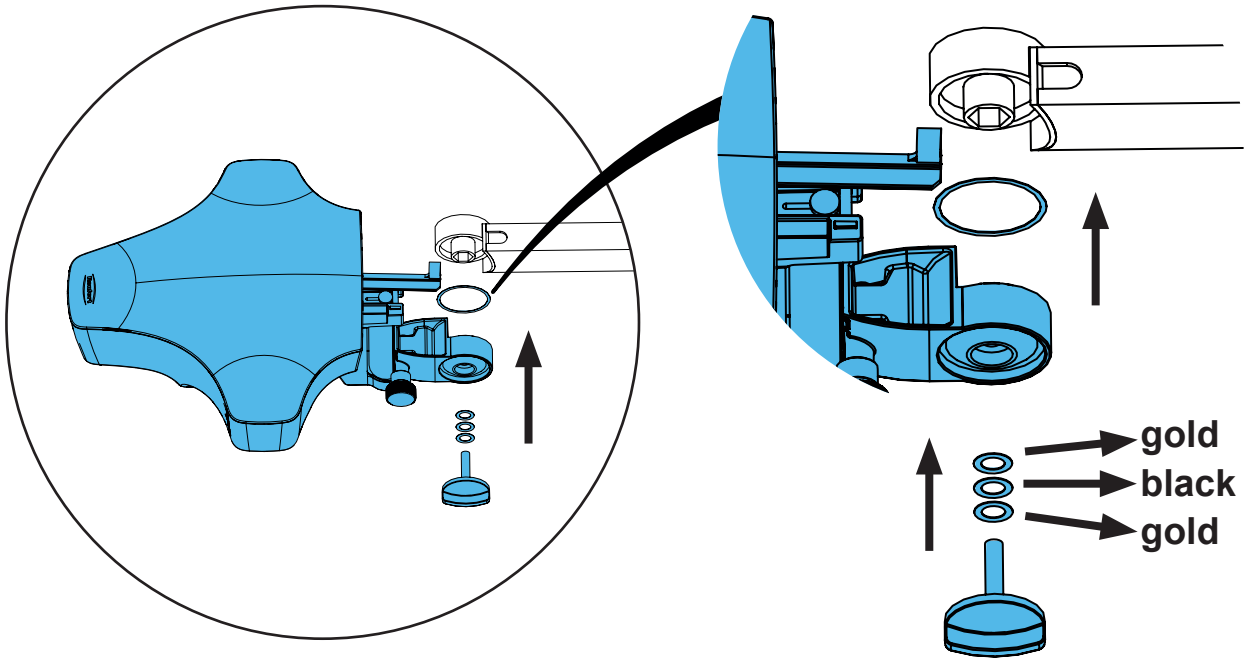

8.

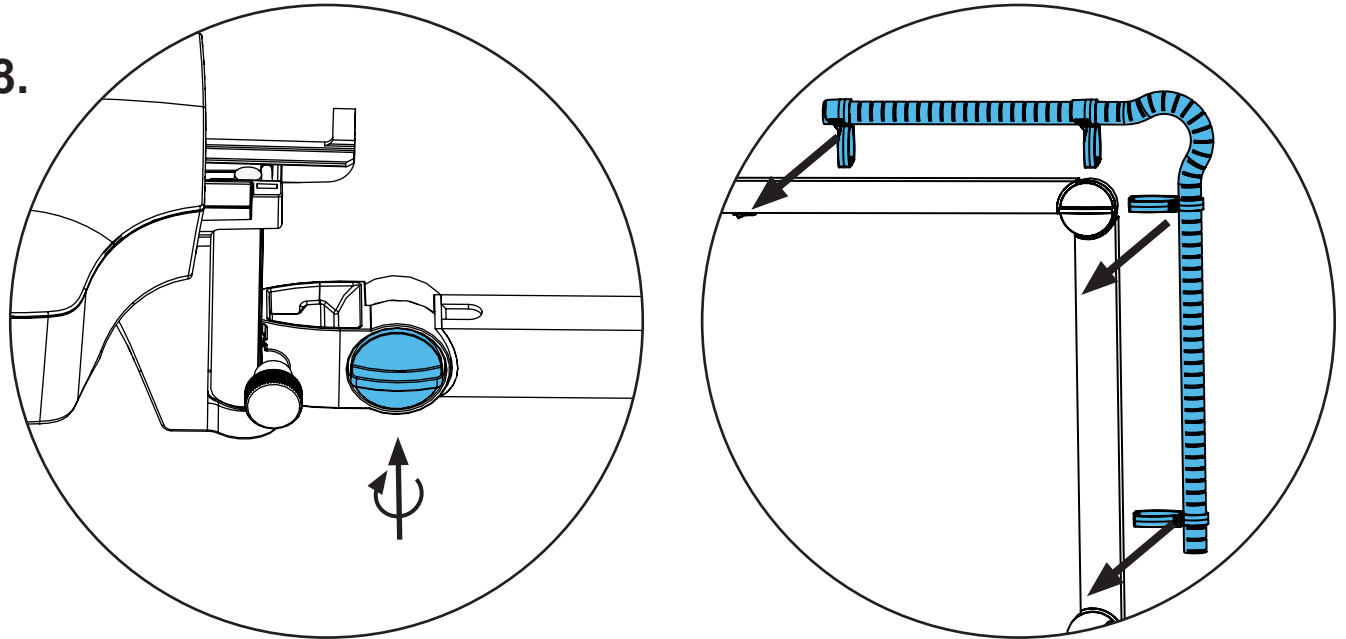

9.

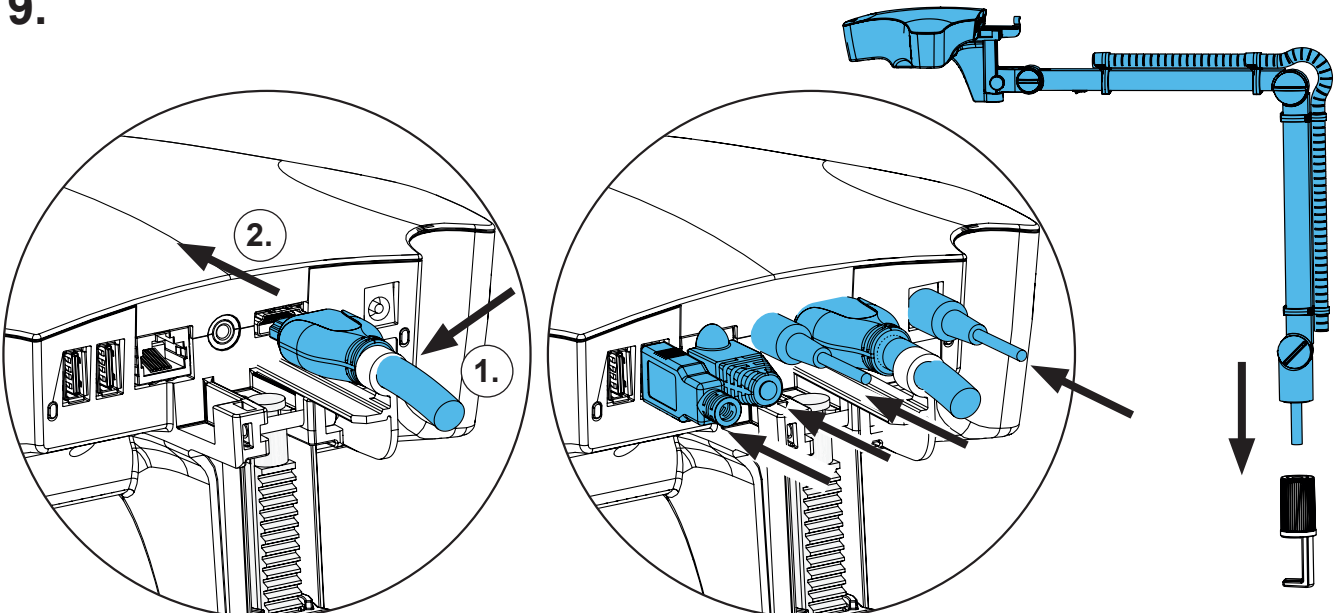






Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу


making work easy

90003 7808


3D Monitor • 3D monitor • Écran 3D • Monitor 3D • Monitor 3D • Монитор 3D

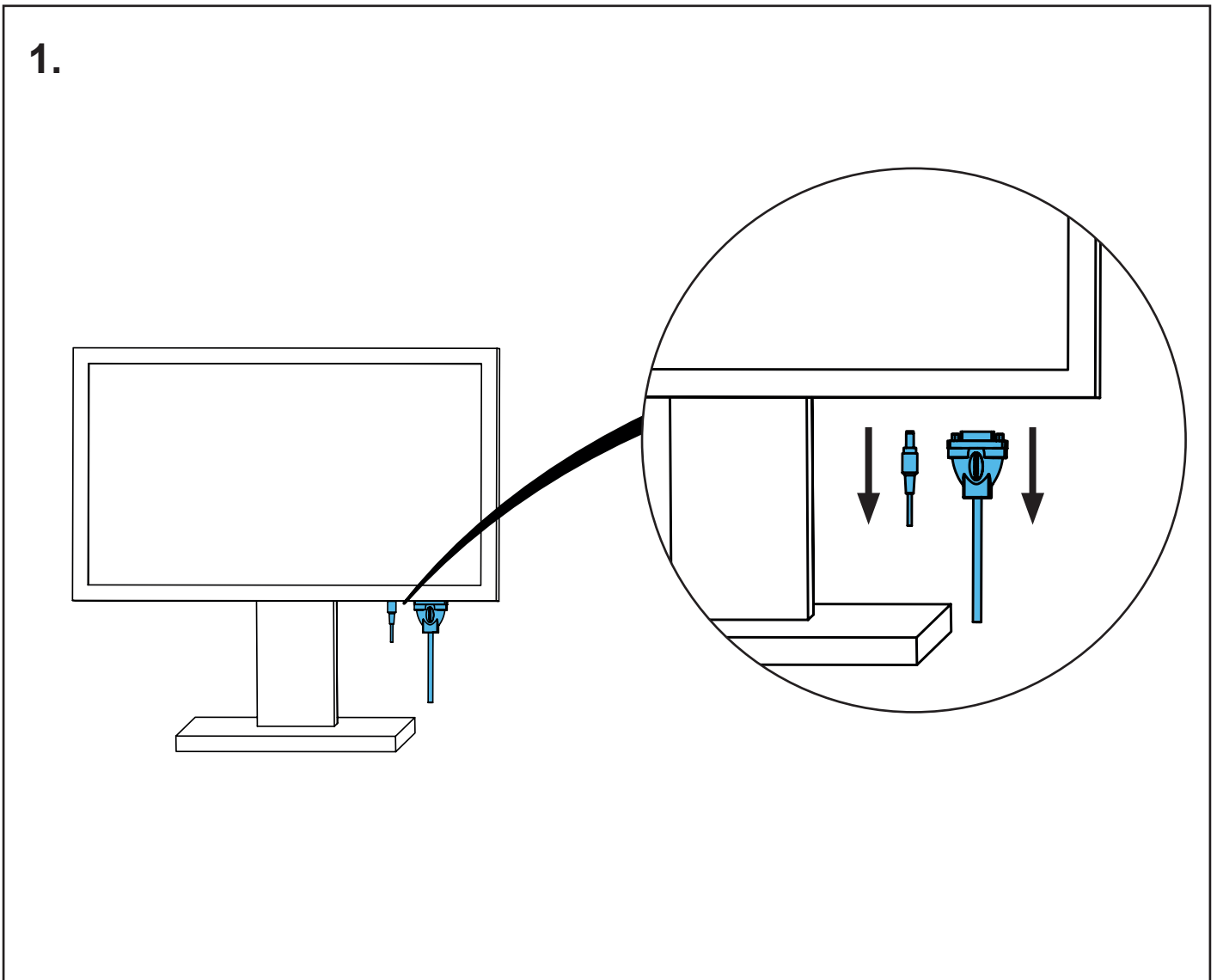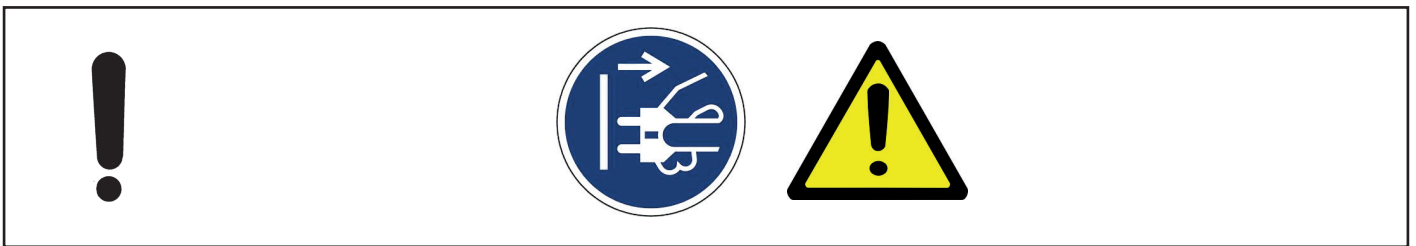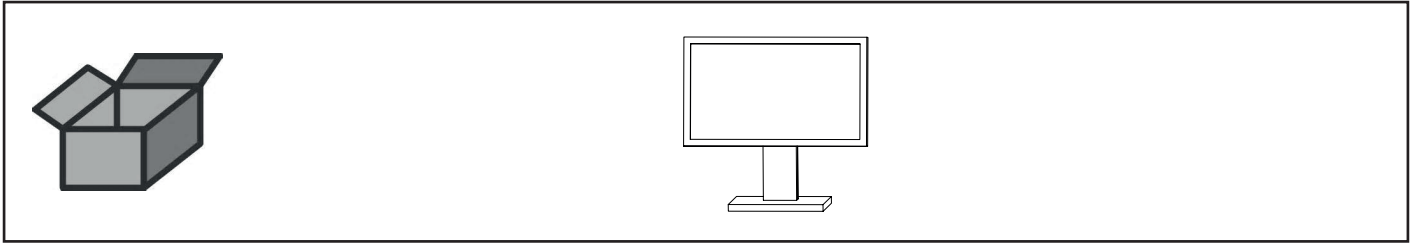

2.

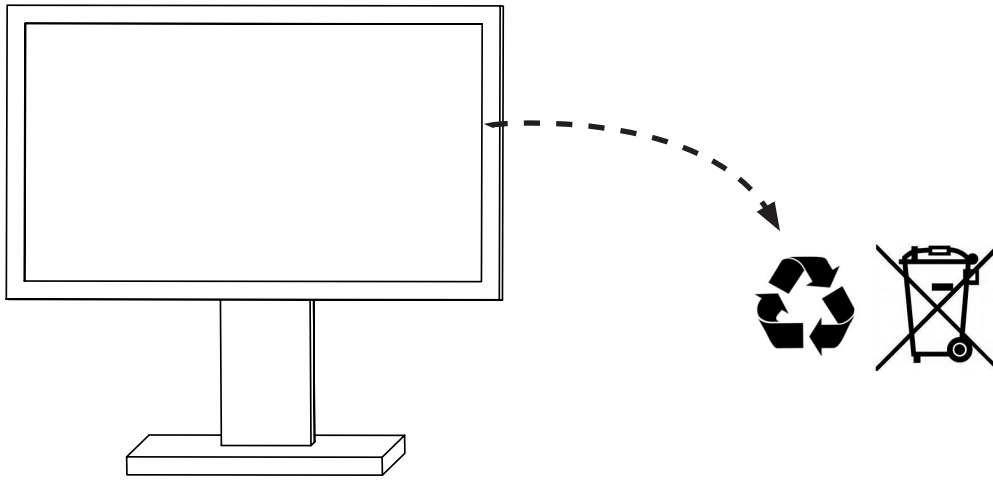

3.

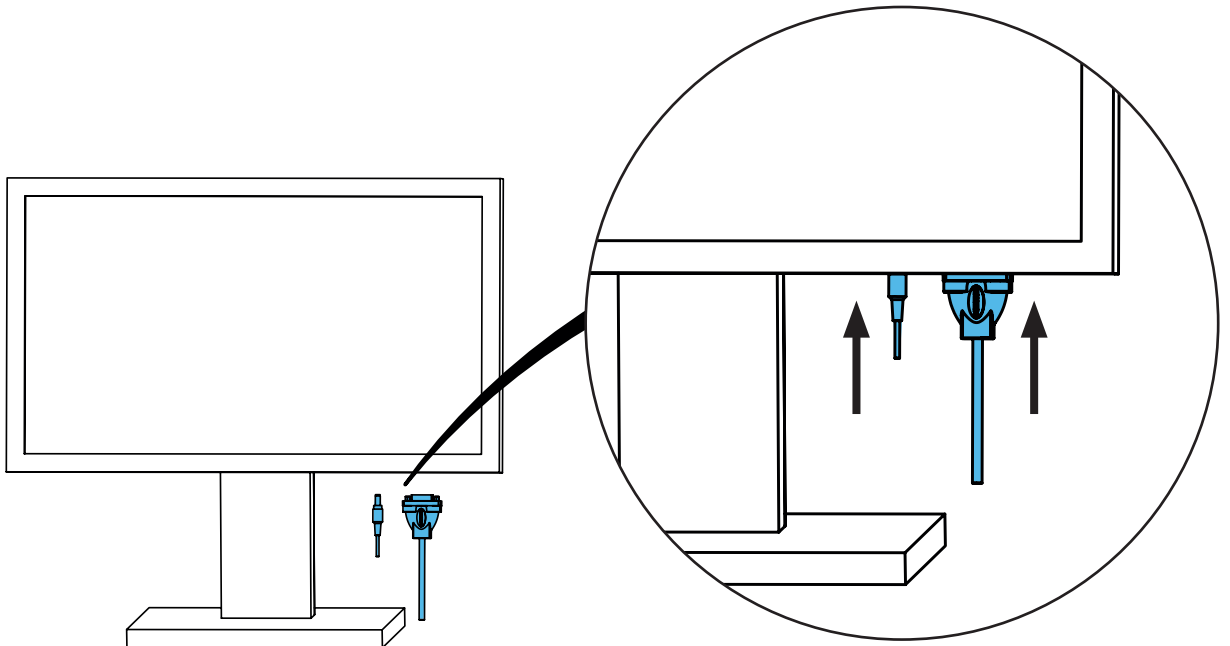






making work easy


Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу

90002 1891


Kopfteil • Head • Tête • Testata • Cabezal • Головка

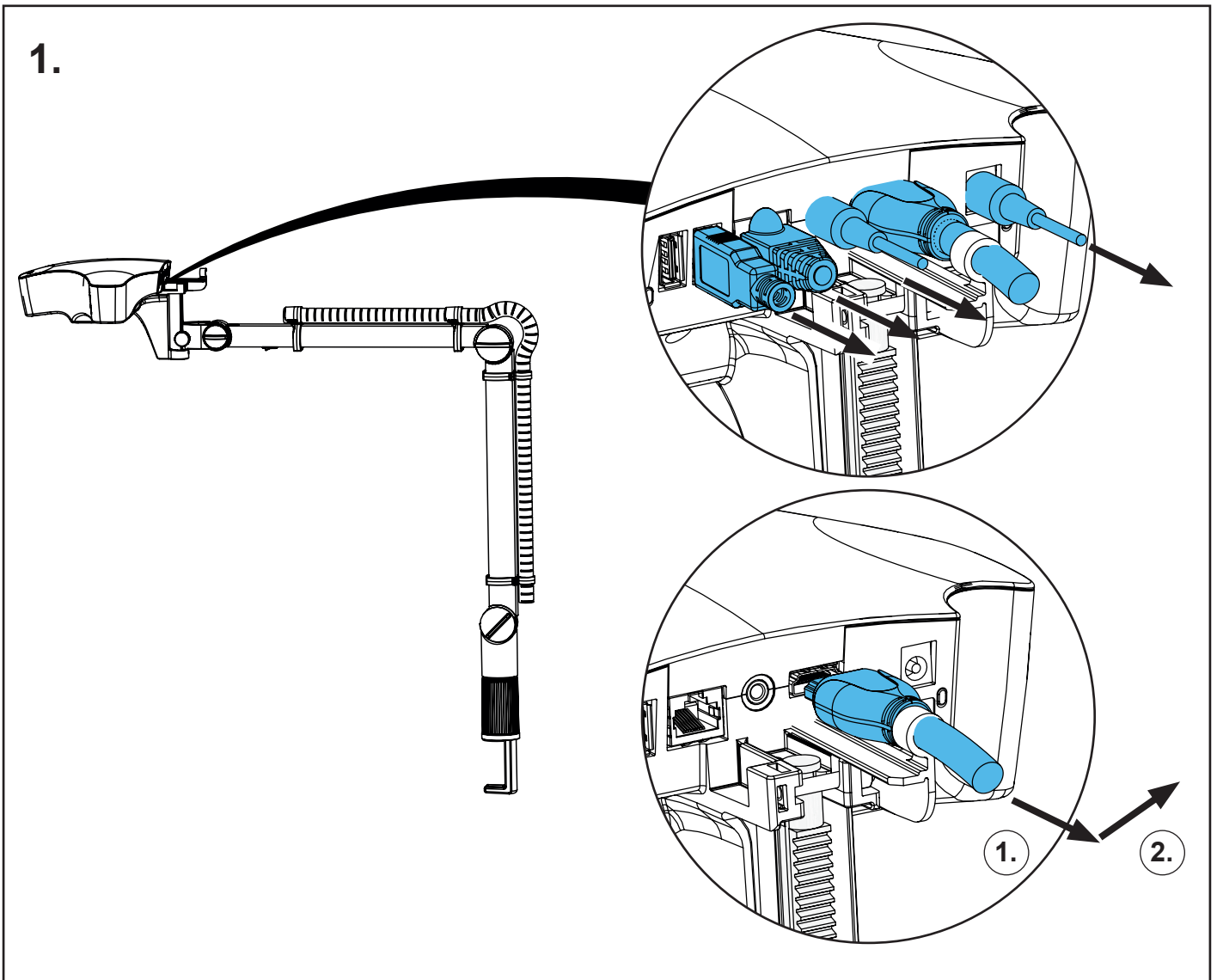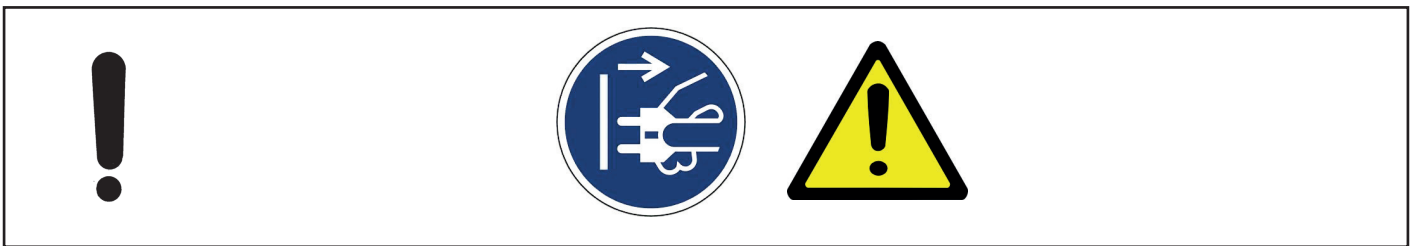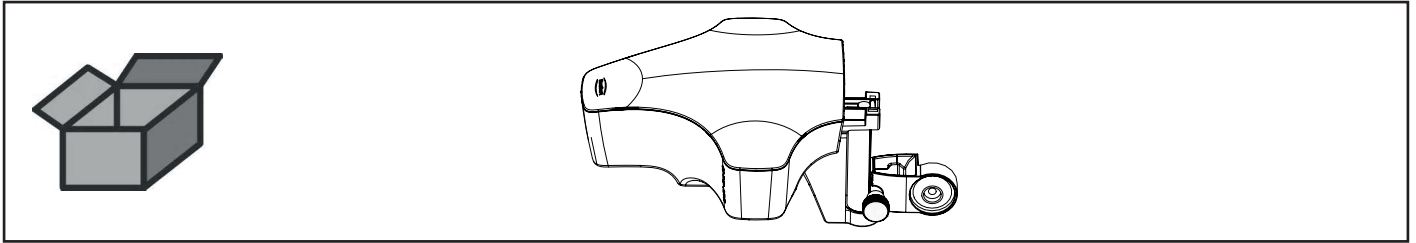


2

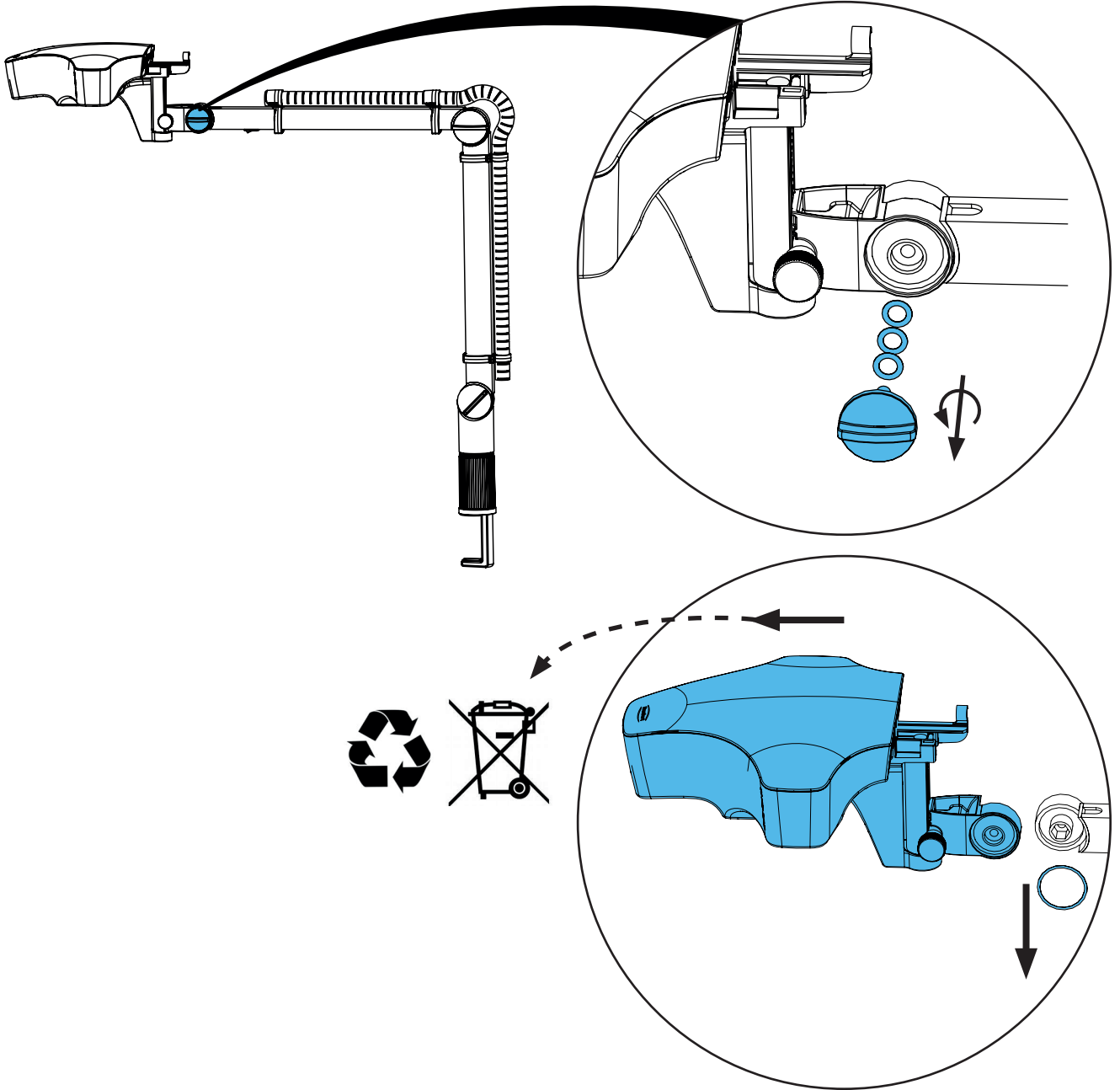


3.

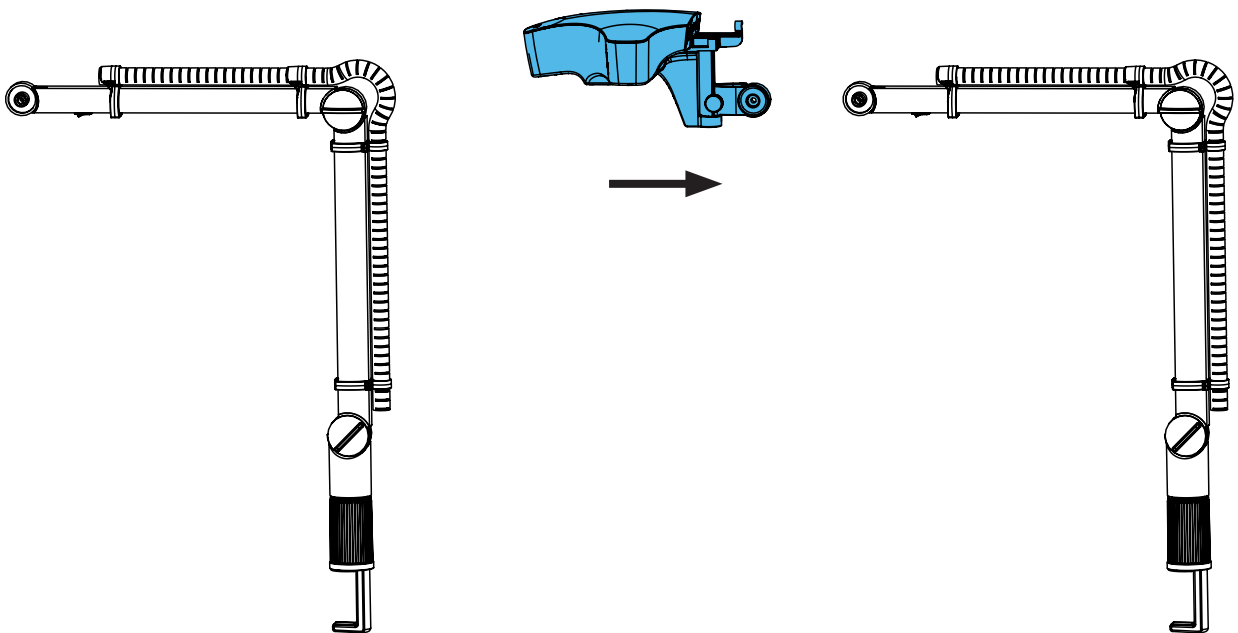

4.

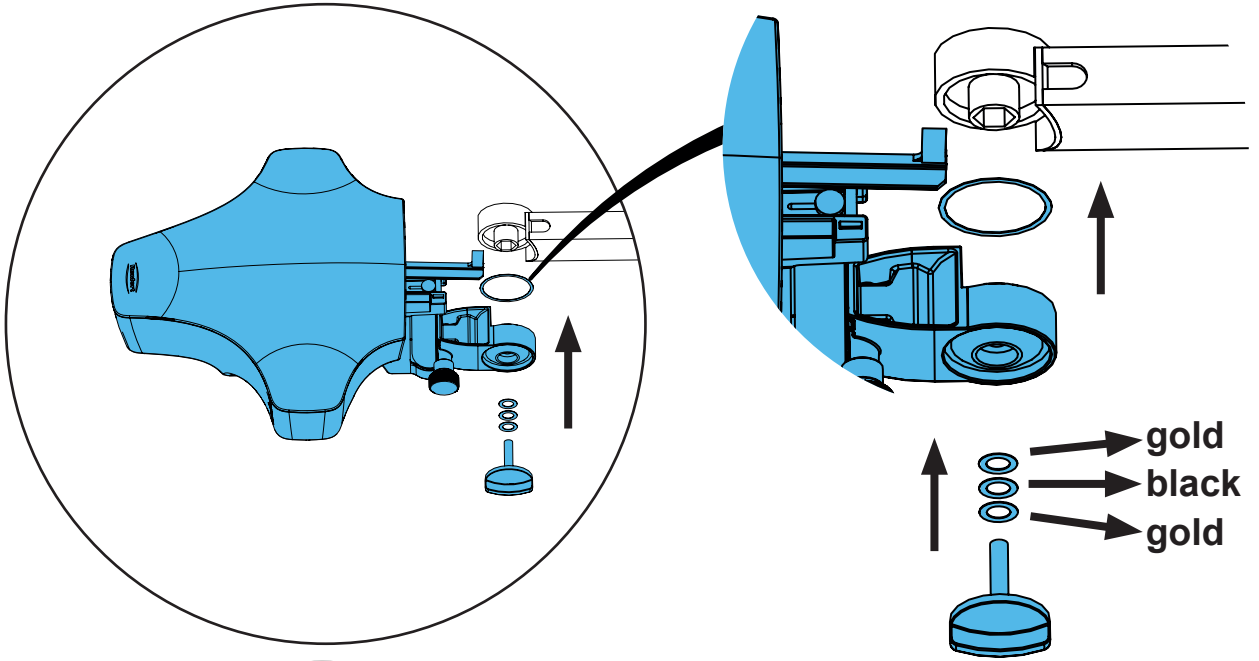

5.

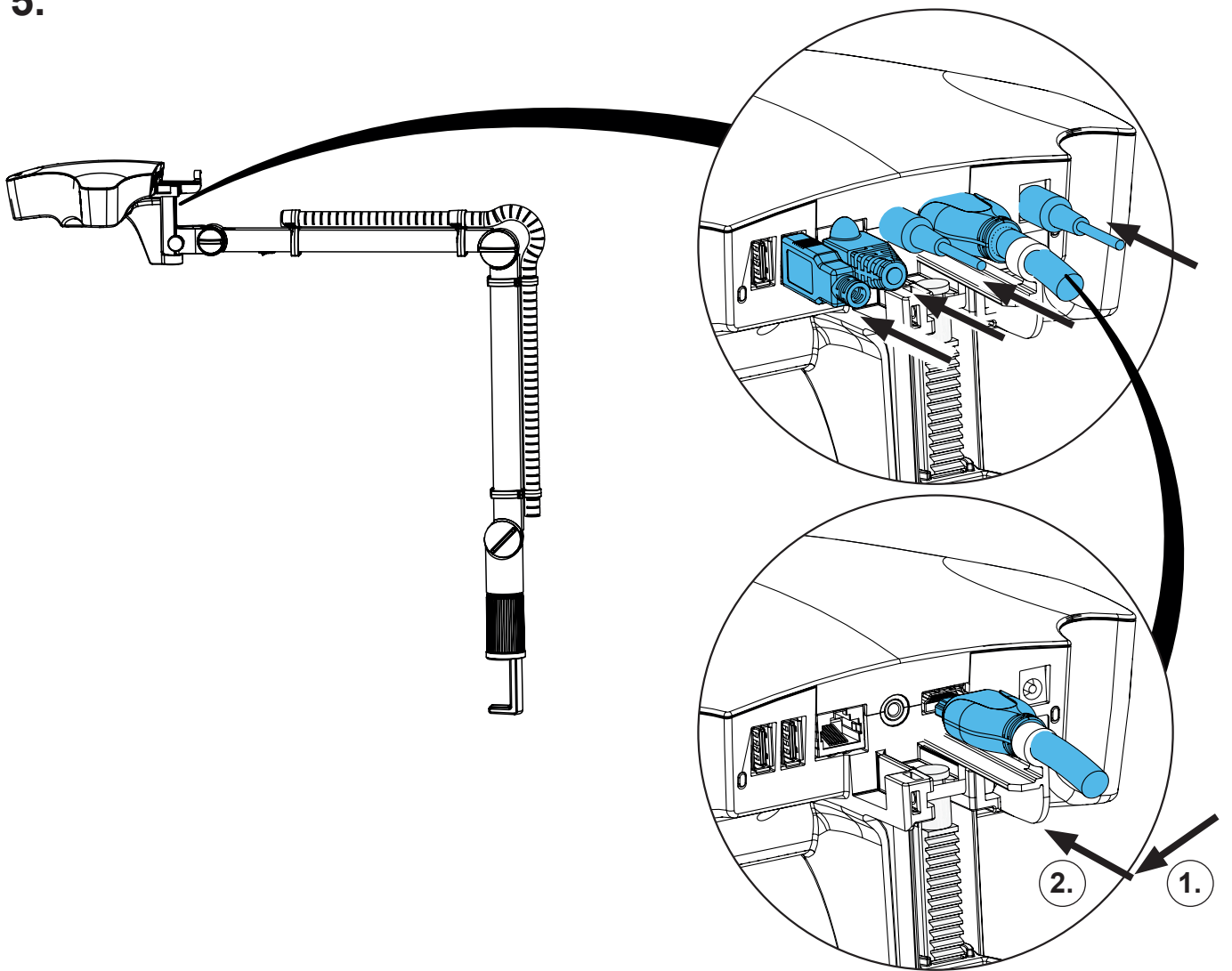










making work easy

Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по монтажу



90003 7818

Kleinteileset • Hardware kit • Petit matériel de montage • Kit di minuteria
Piezas pequeñas • Комплект мелких деталей

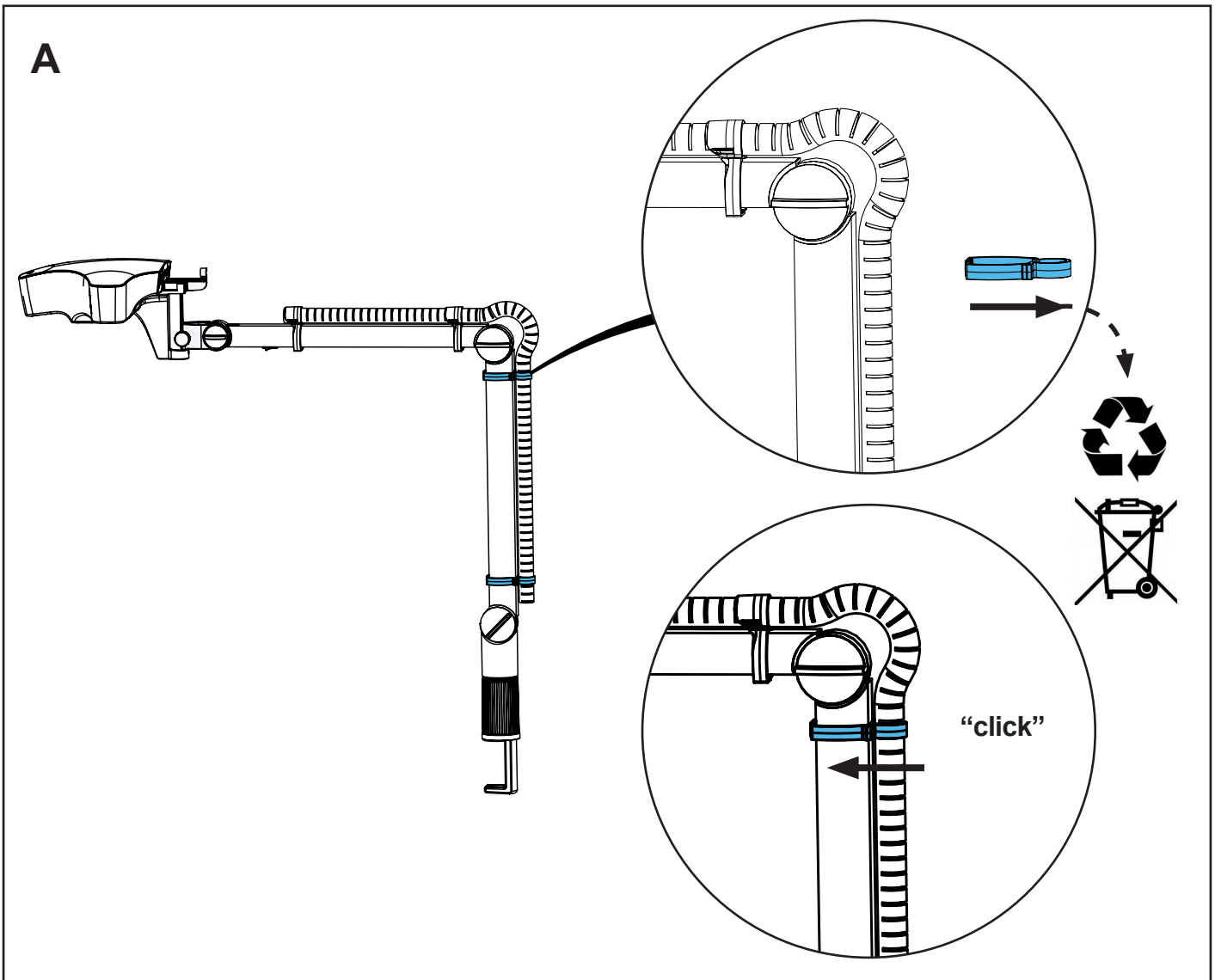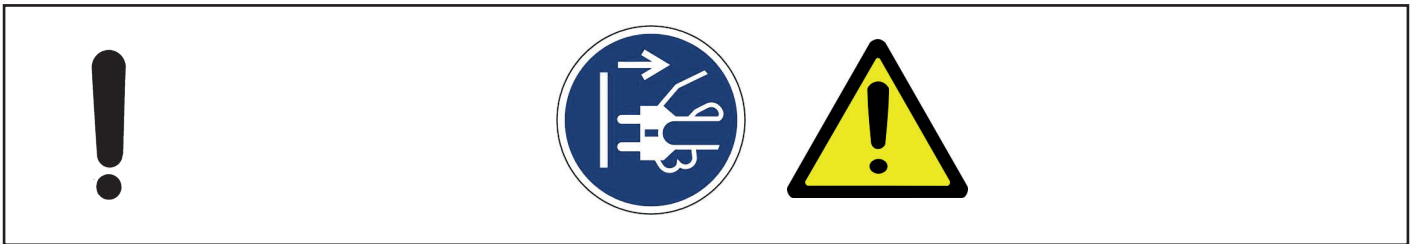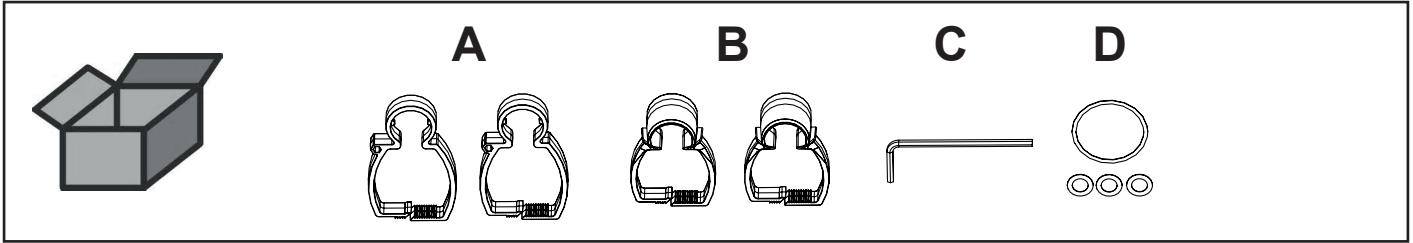


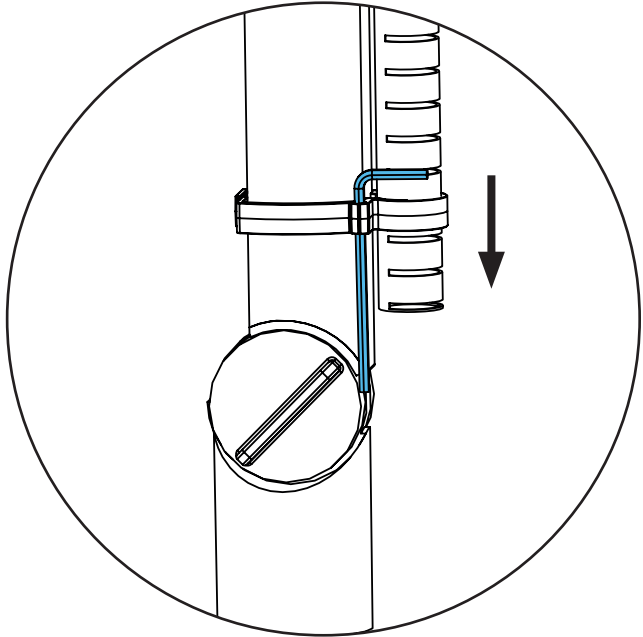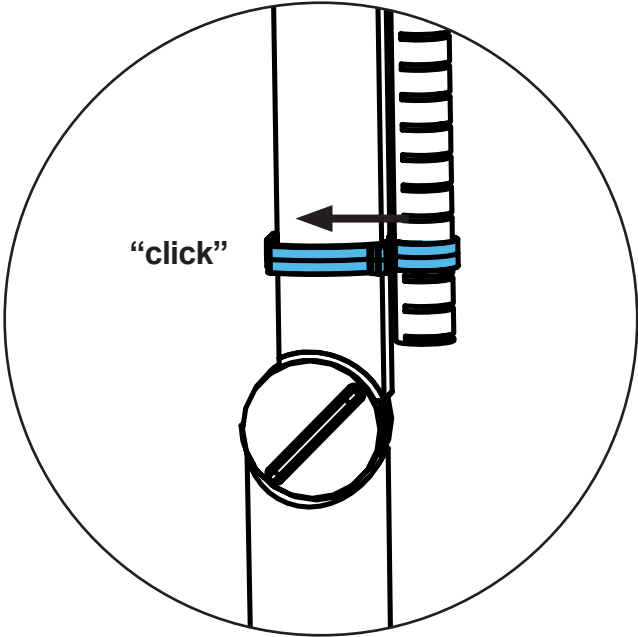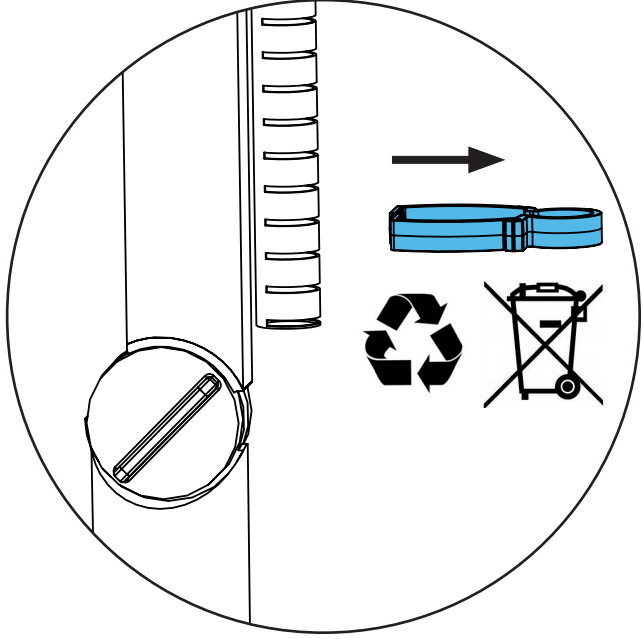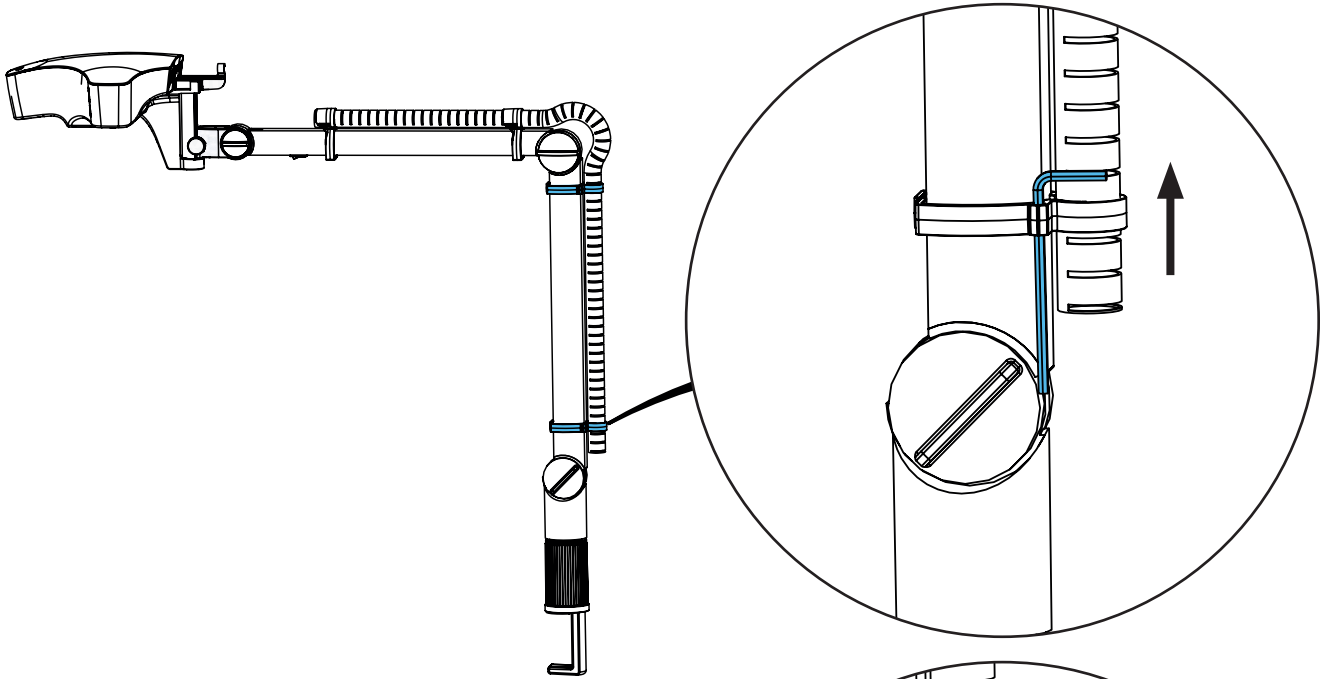



B

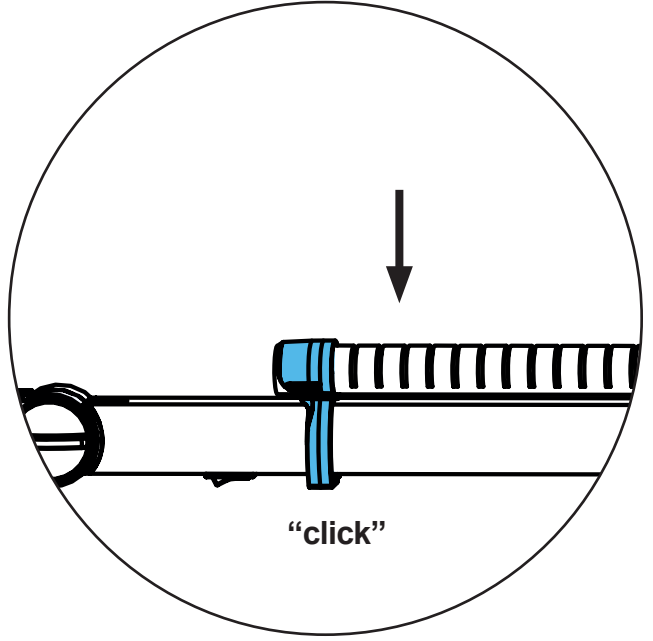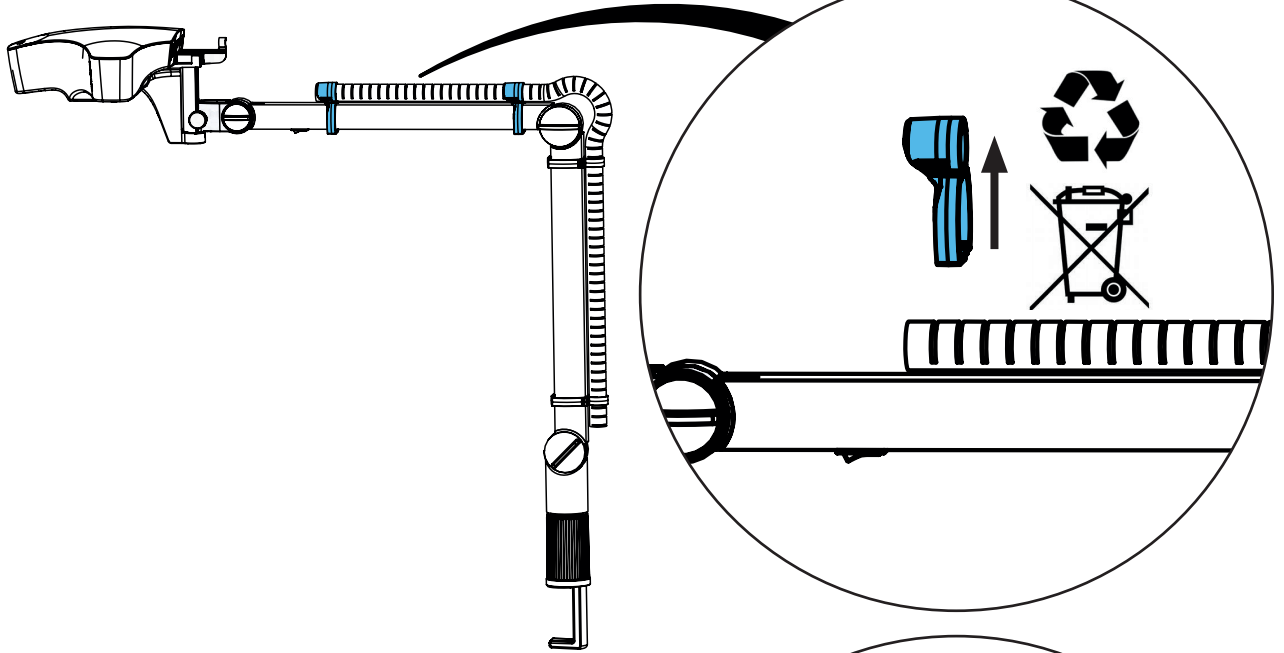

C

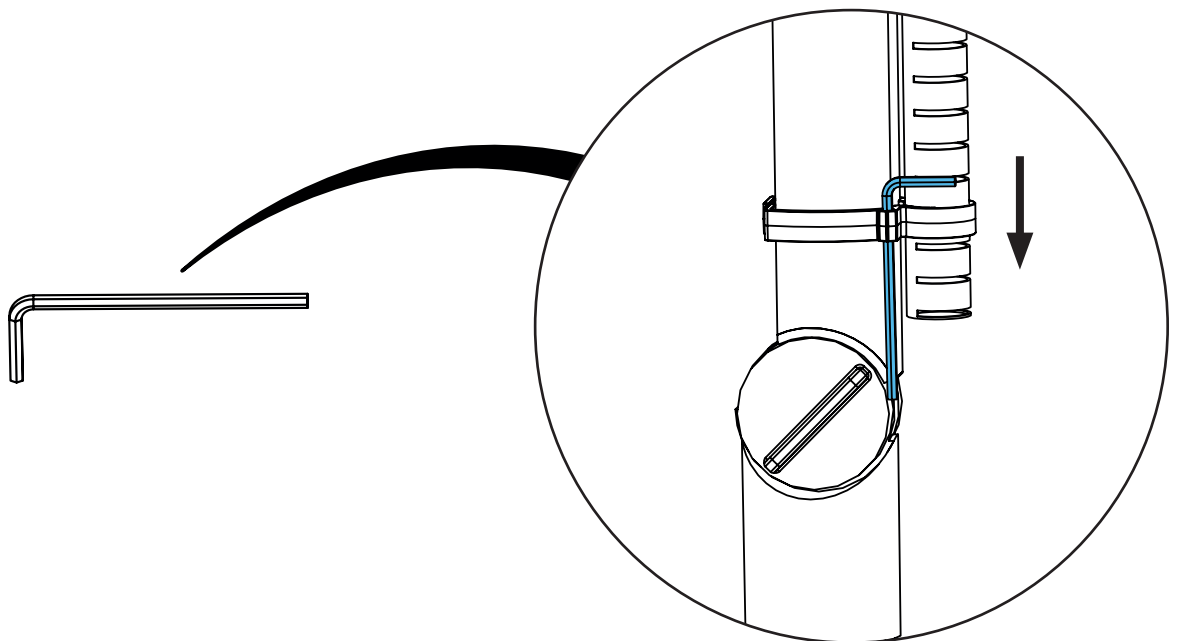

D

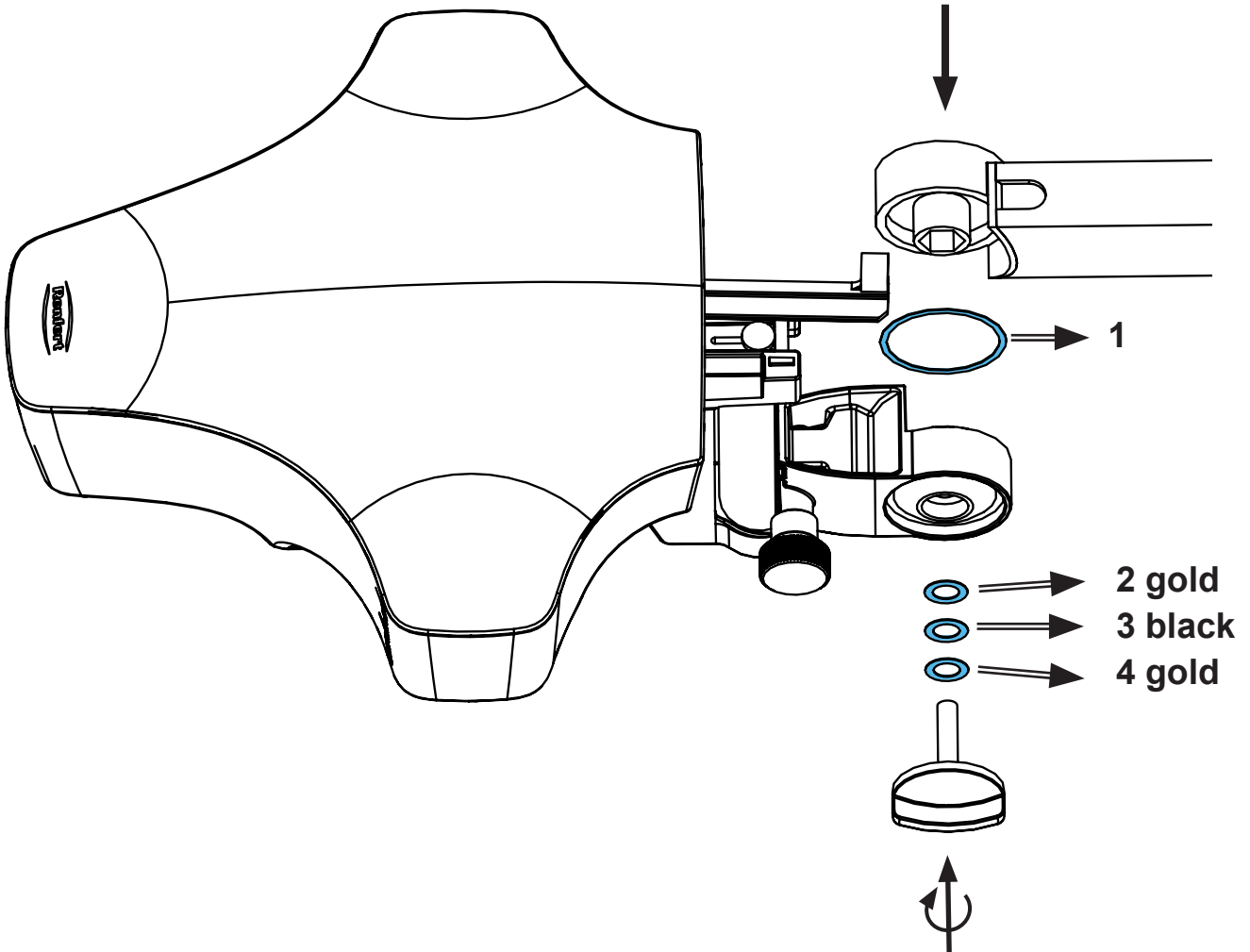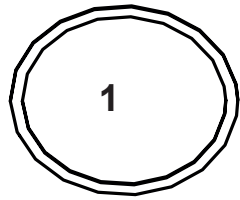





making work easy

Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу

90003 7809

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3D Brille • 3D goggles • Lunettes 3D • Occhiali 3D • Gafas 3D • Очки 3D

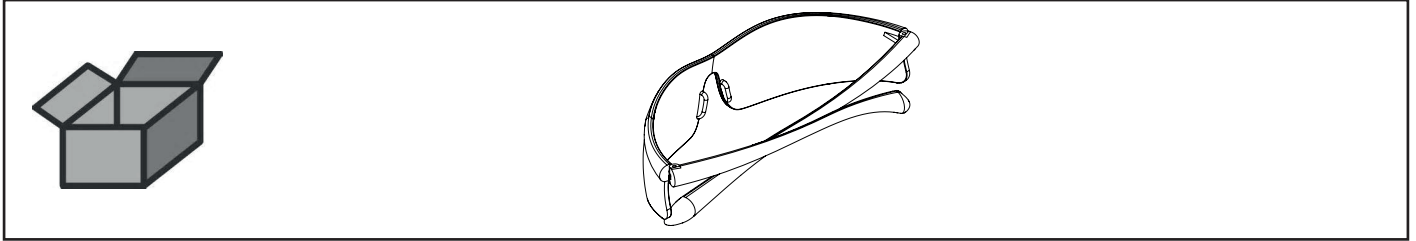







Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по монтажу






making work easy

90003 7368





Maus • Mouse • Souris • Mouse • Ratón • Мышь

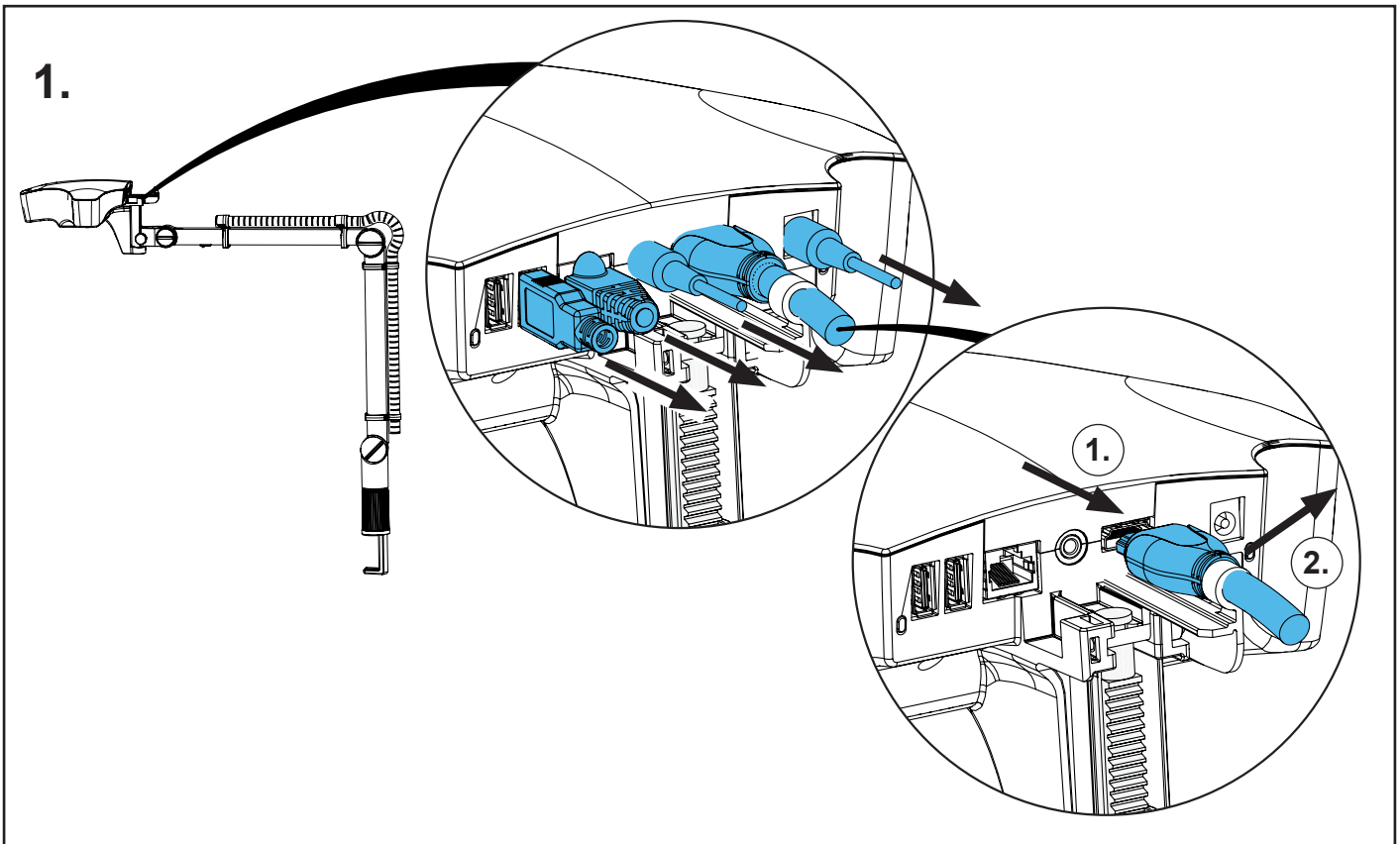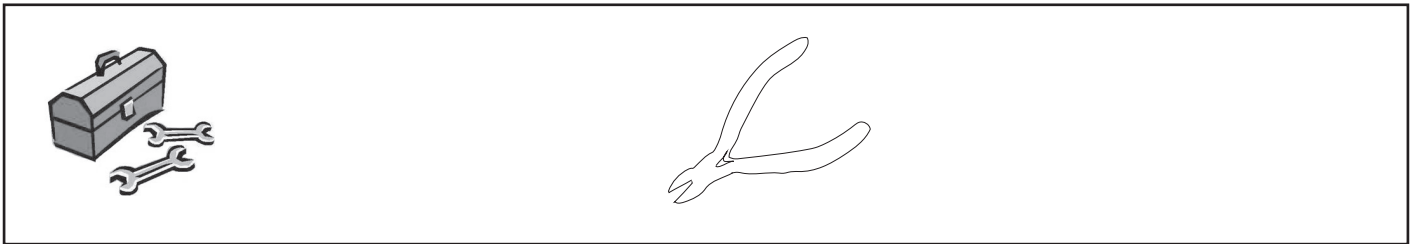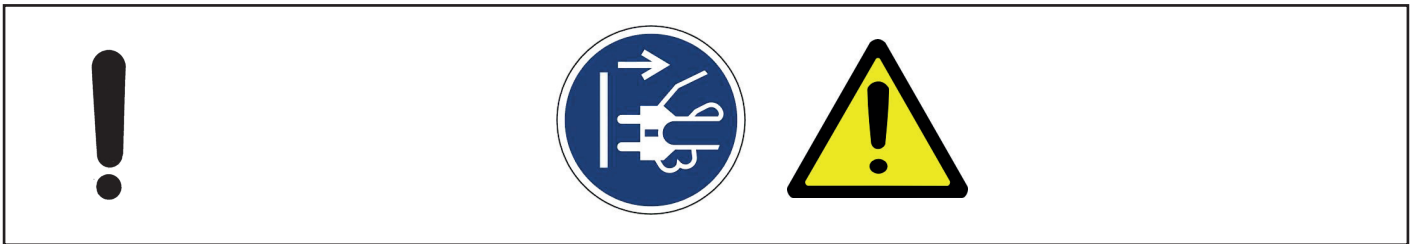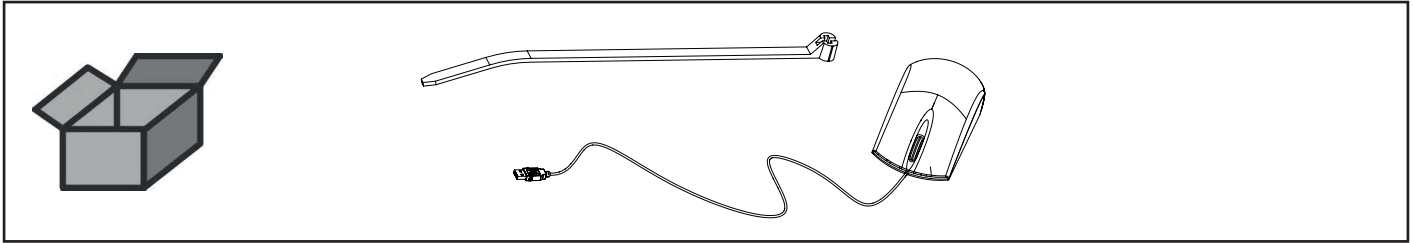




2.

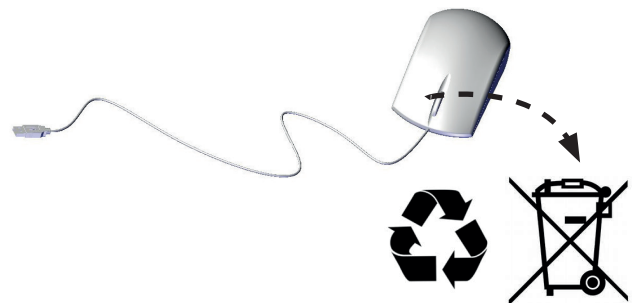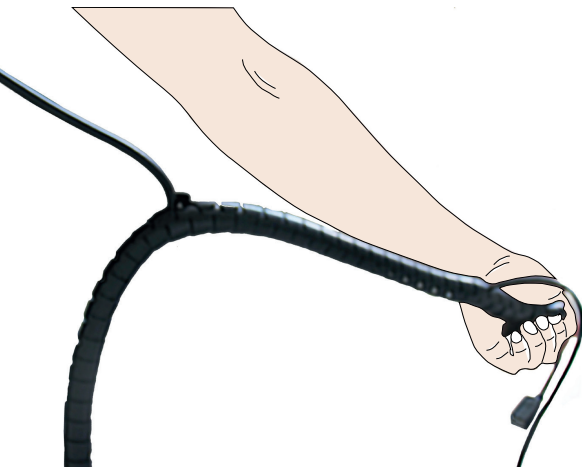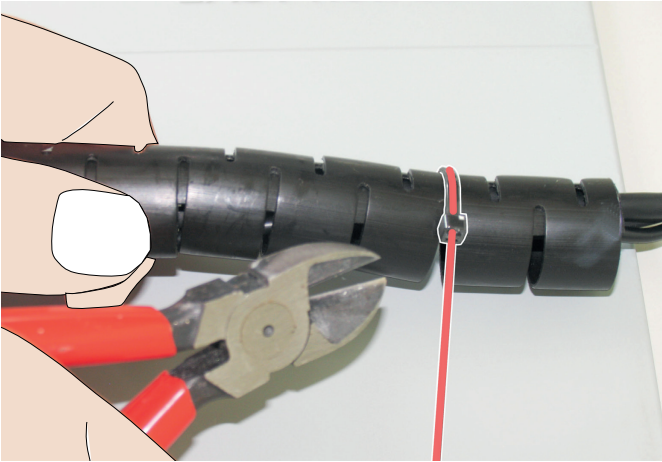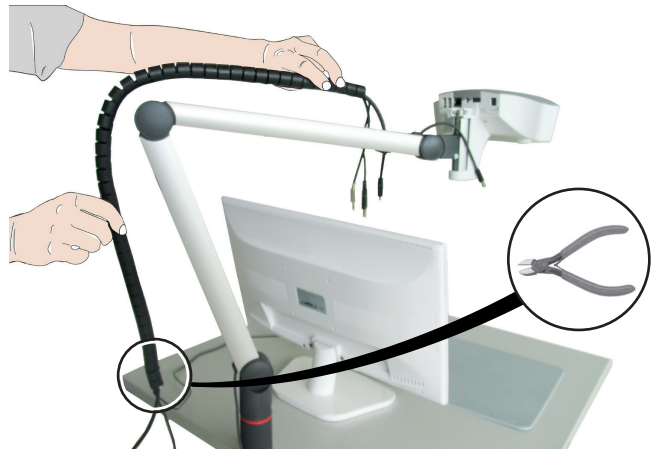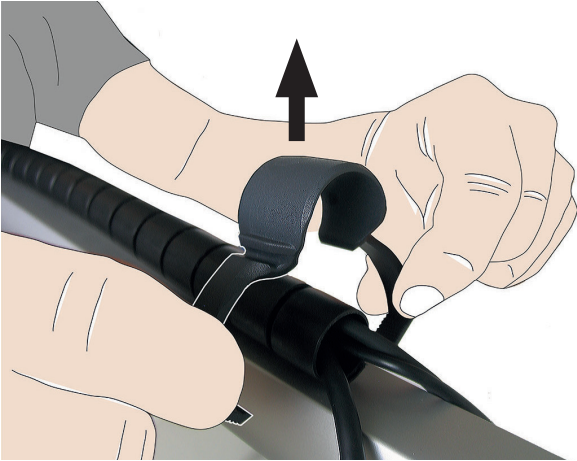

3.

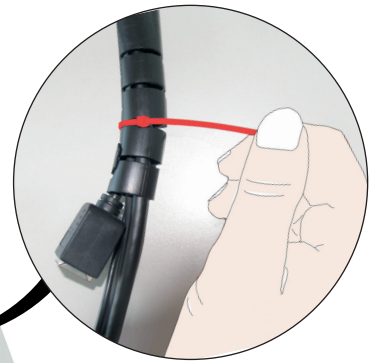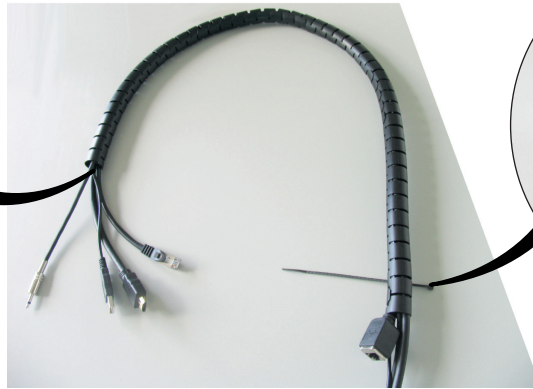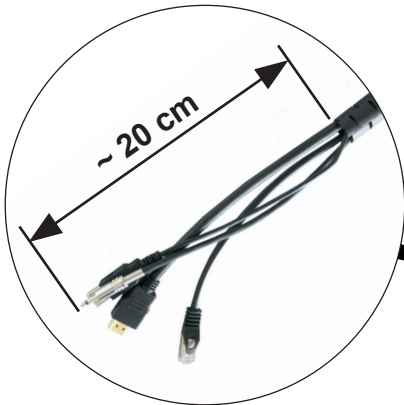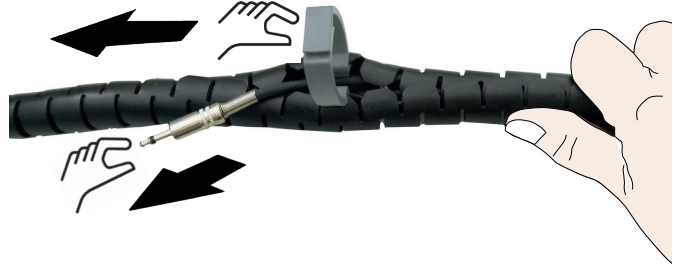

4.

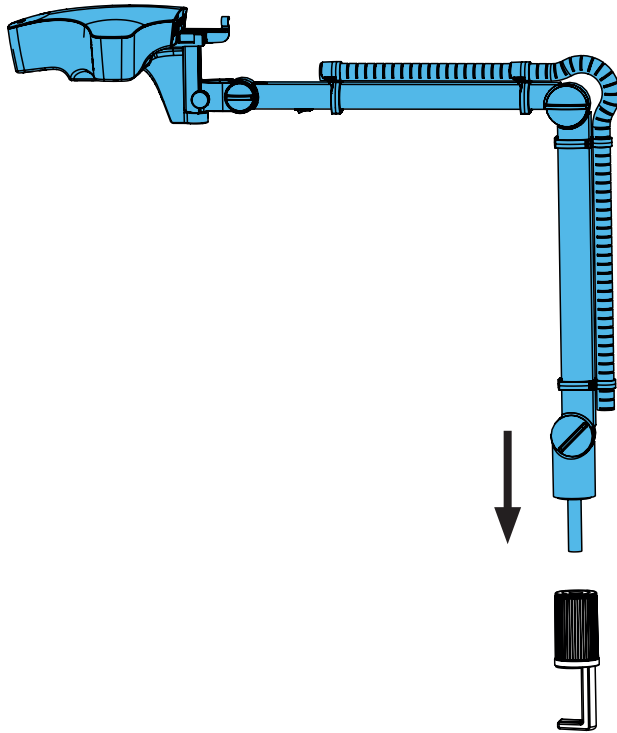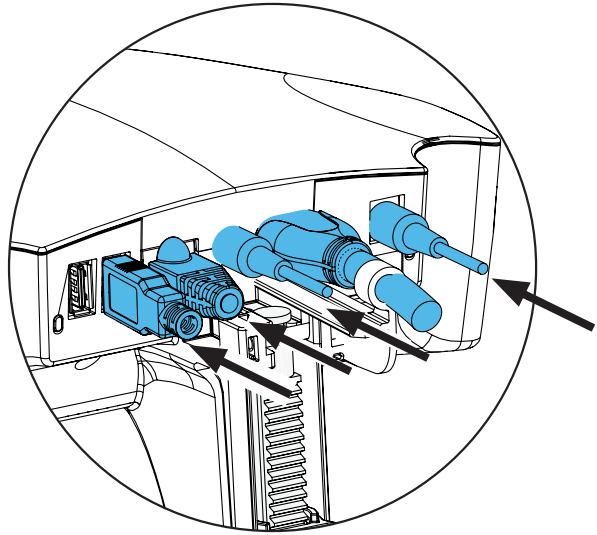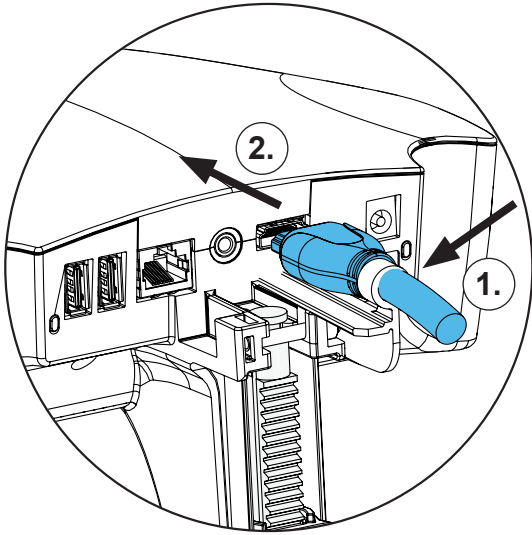






making work easy

Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу

90003 3776

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Tischauflage • Desk pad • Sous-main • Piano di appoggio • Apoyo para la mesa
Настольная подложка



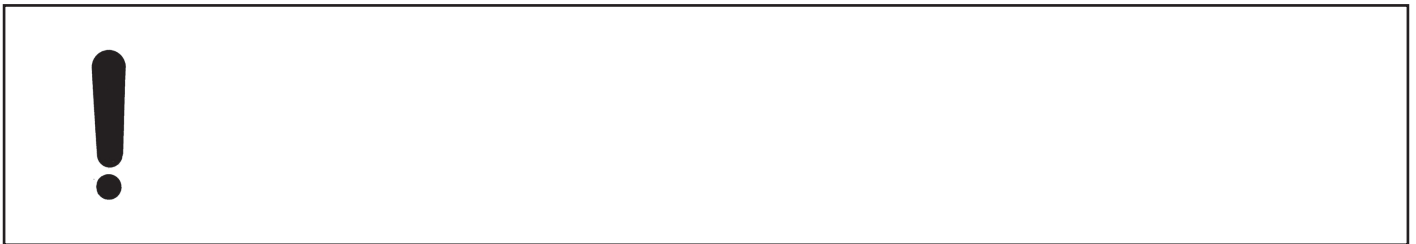
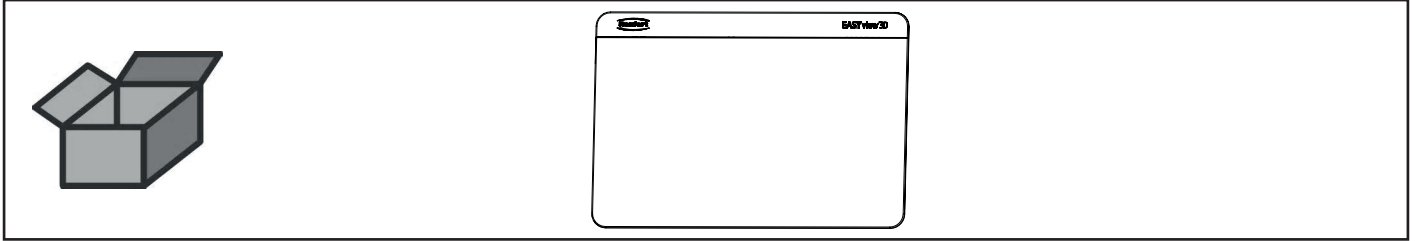






making work easy

90003 7810




Fußschalter • Foot switch • Pédale • Interruttore a pedale • Interruptor de pie
Педаль

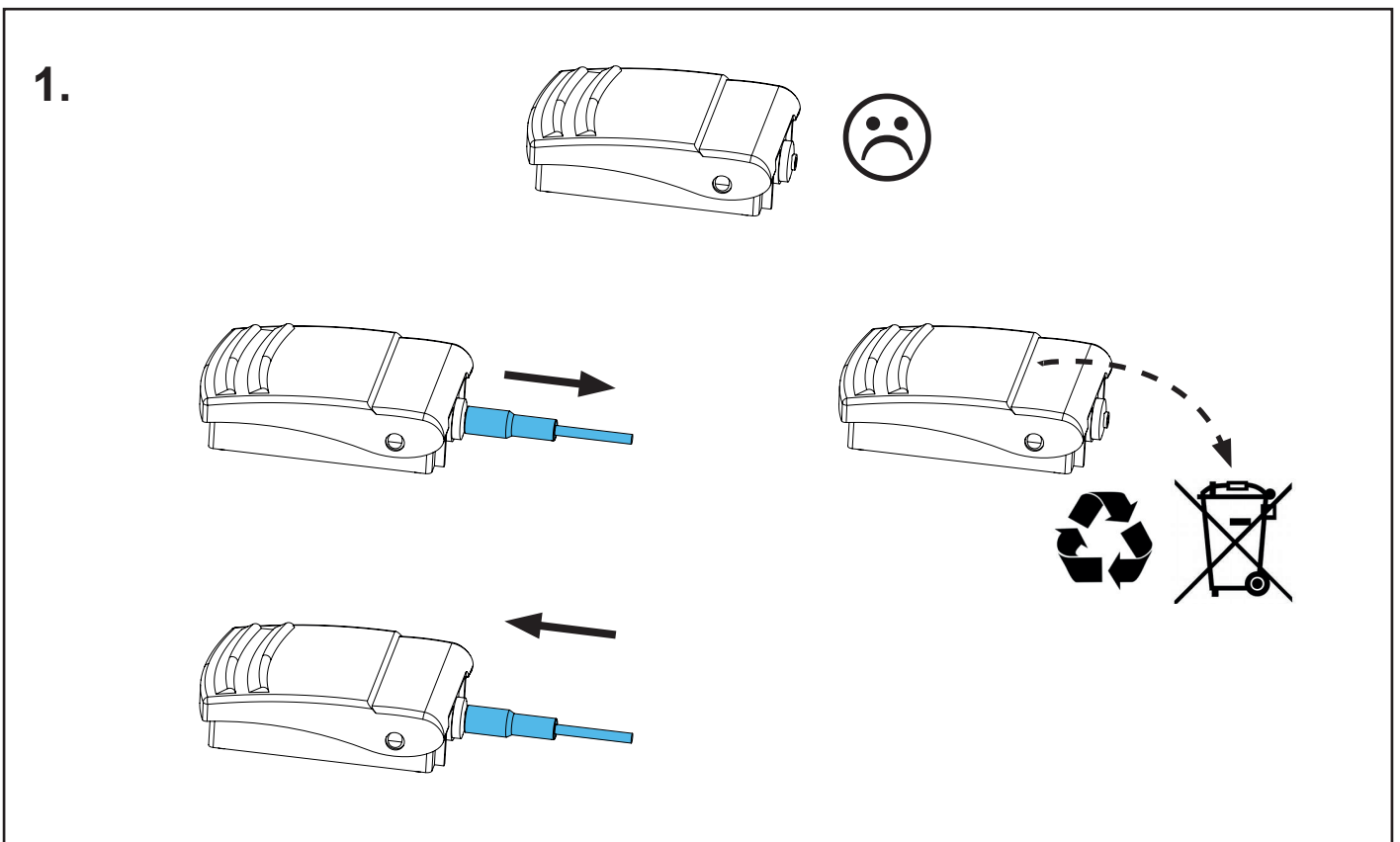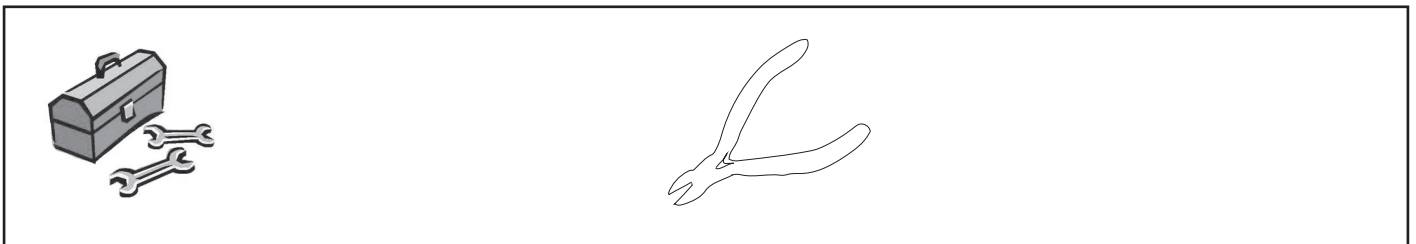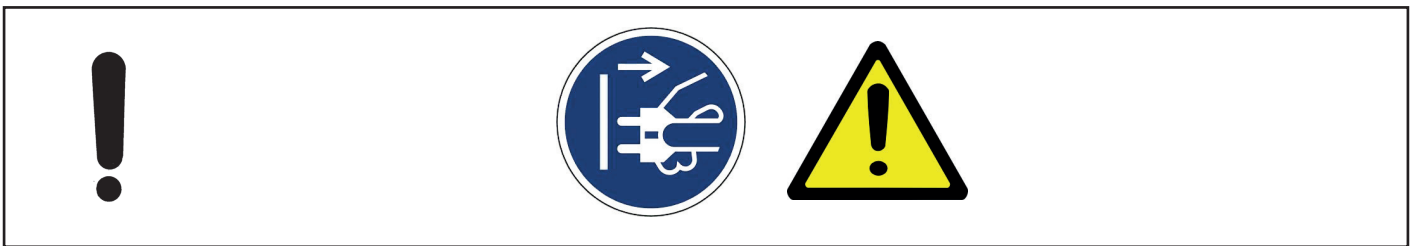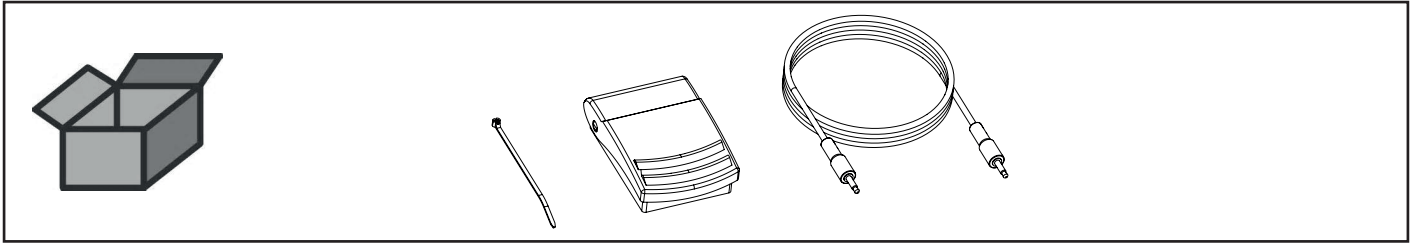


2.




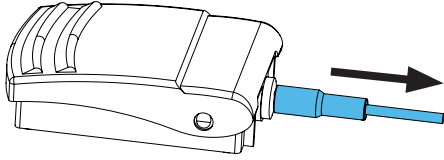
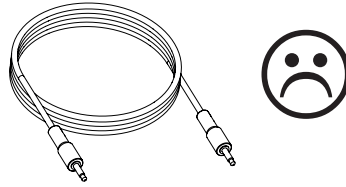





3.

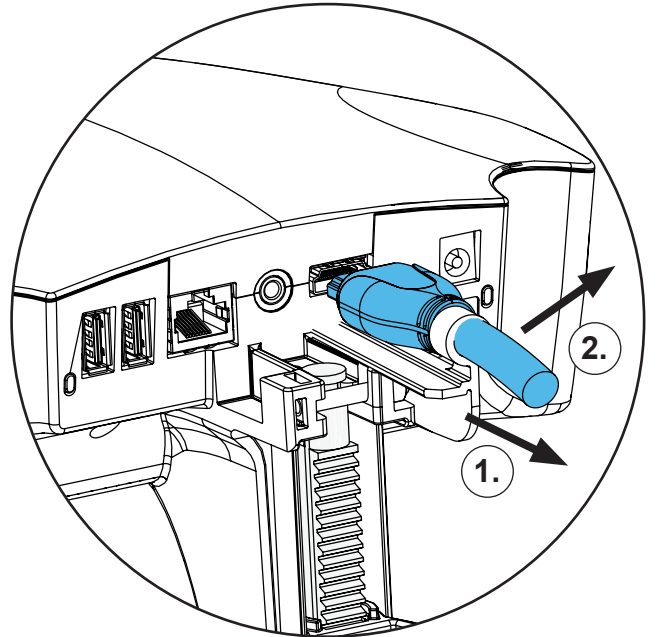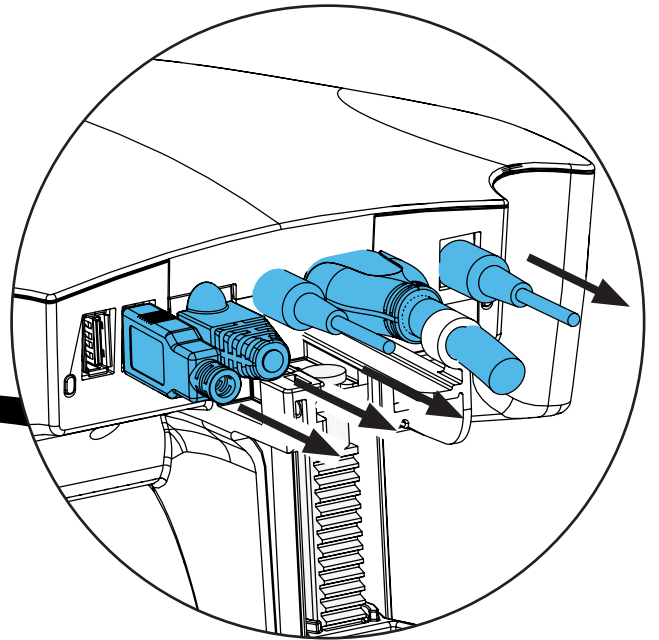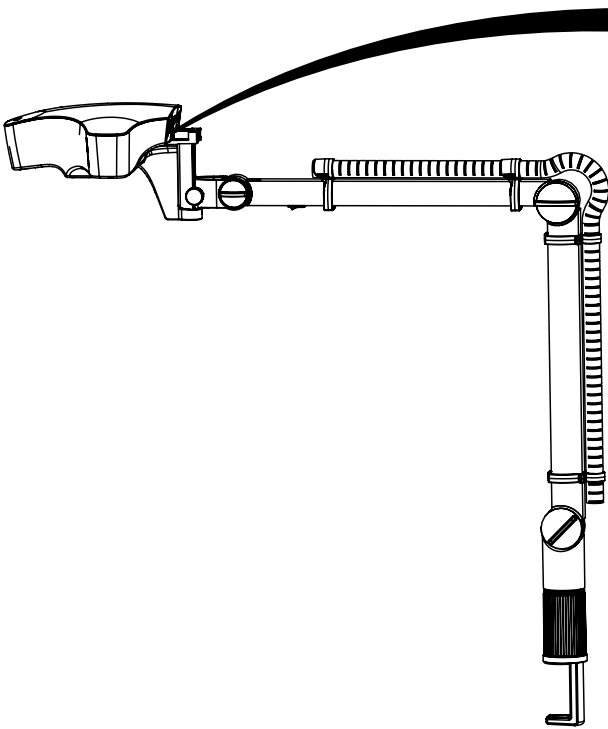



4.

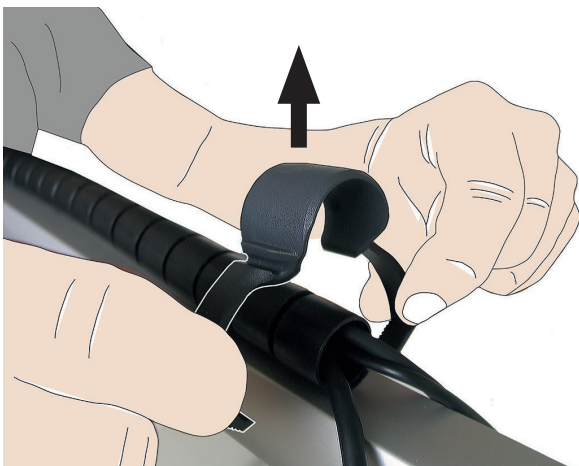

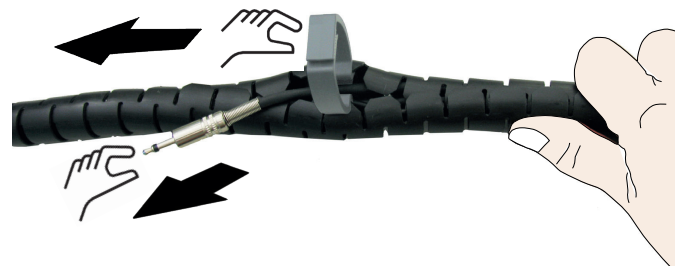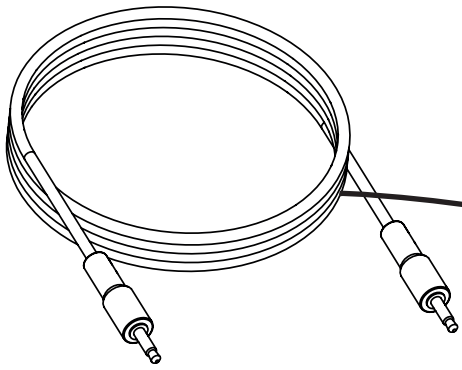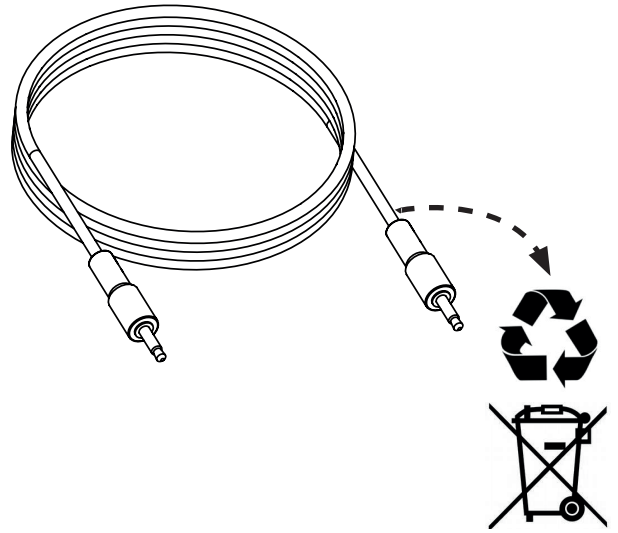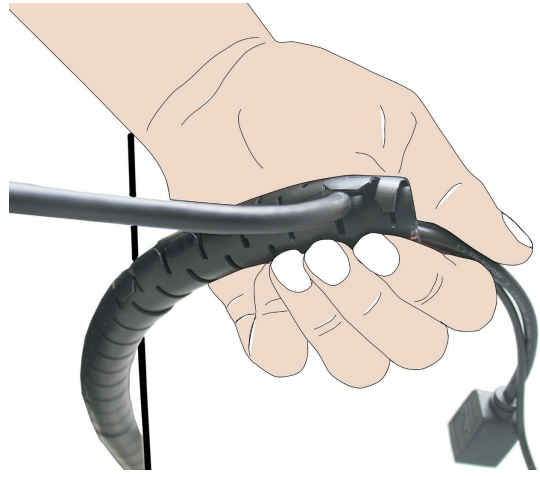
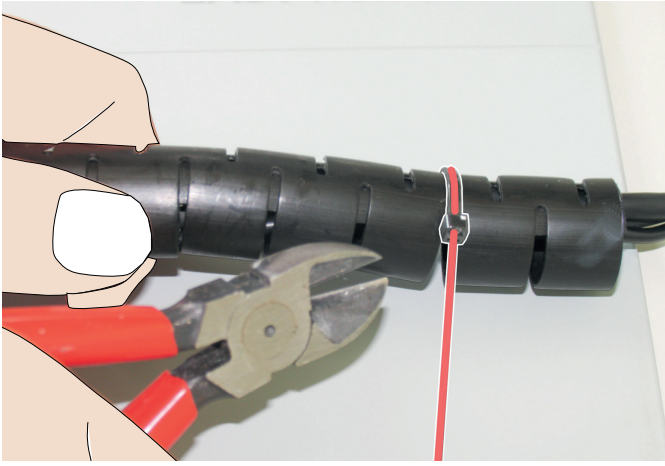


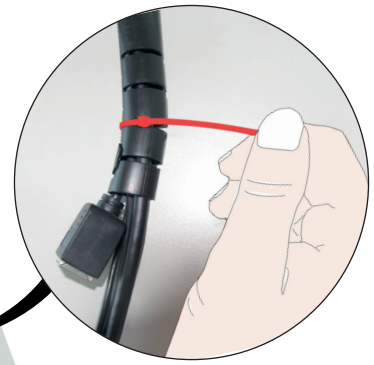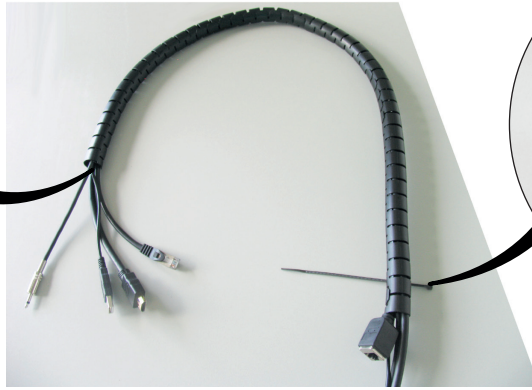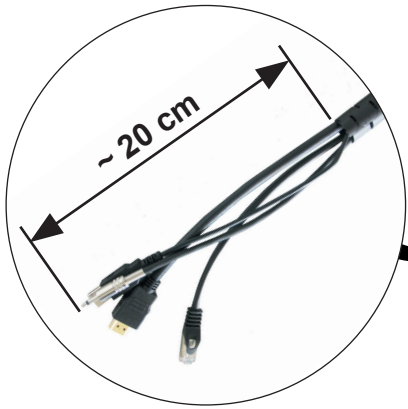

5.

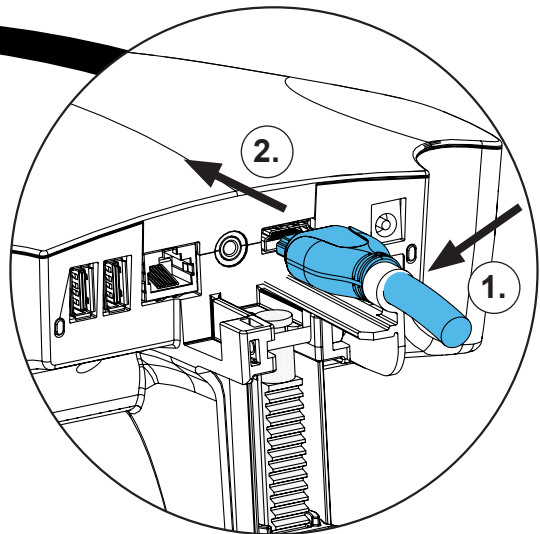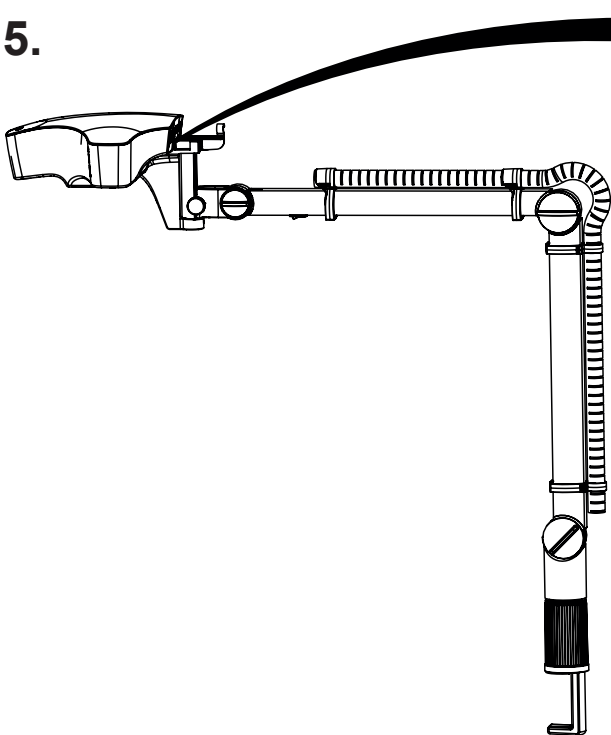

6.

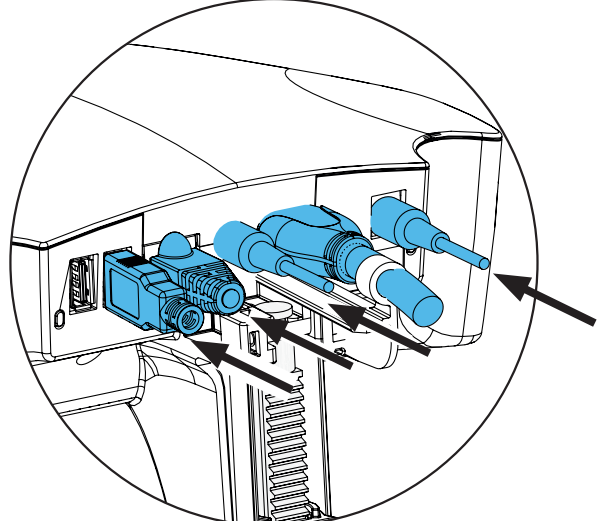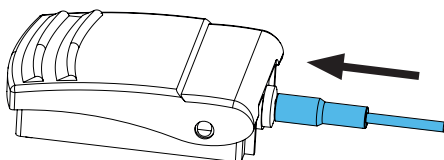







Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу


making work easy

90003 7900





HDMI-DVI Kabel • HDMI-DVI cable • Câble HDMI DVI • Cavo HDMI DVI • Cable
HDMI DVI • Кабель HDMI-DVI

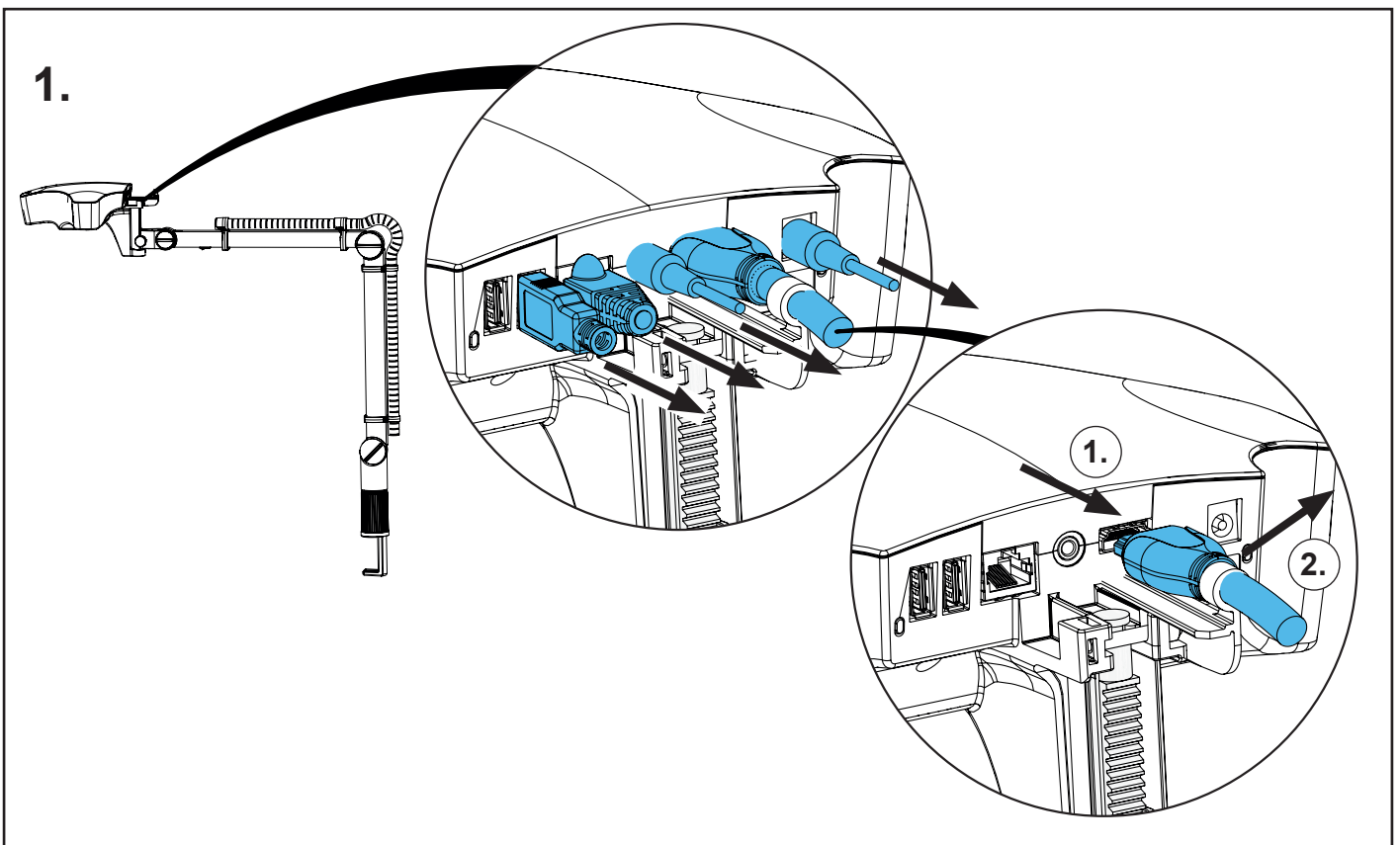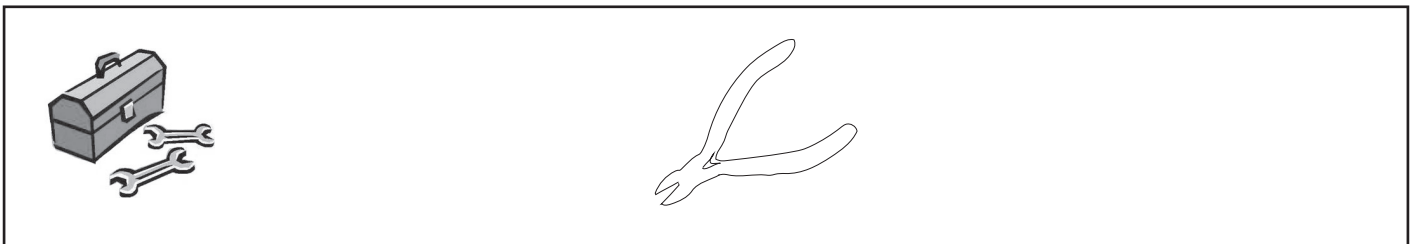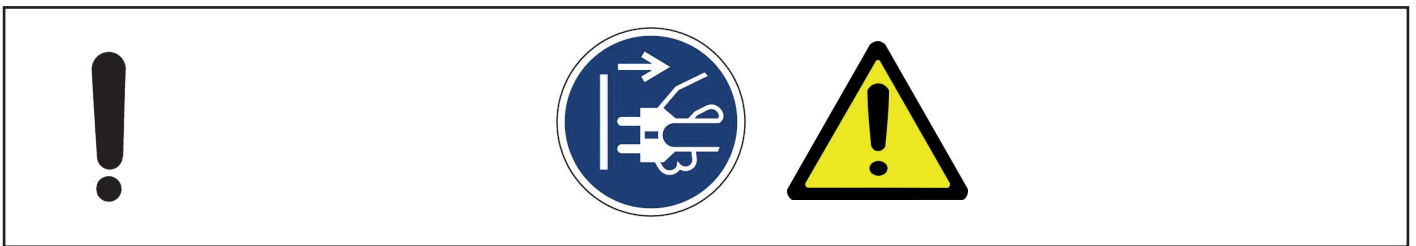
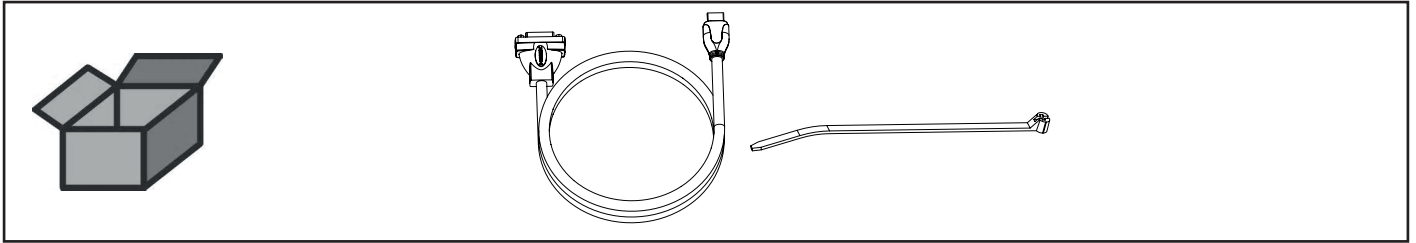


2.



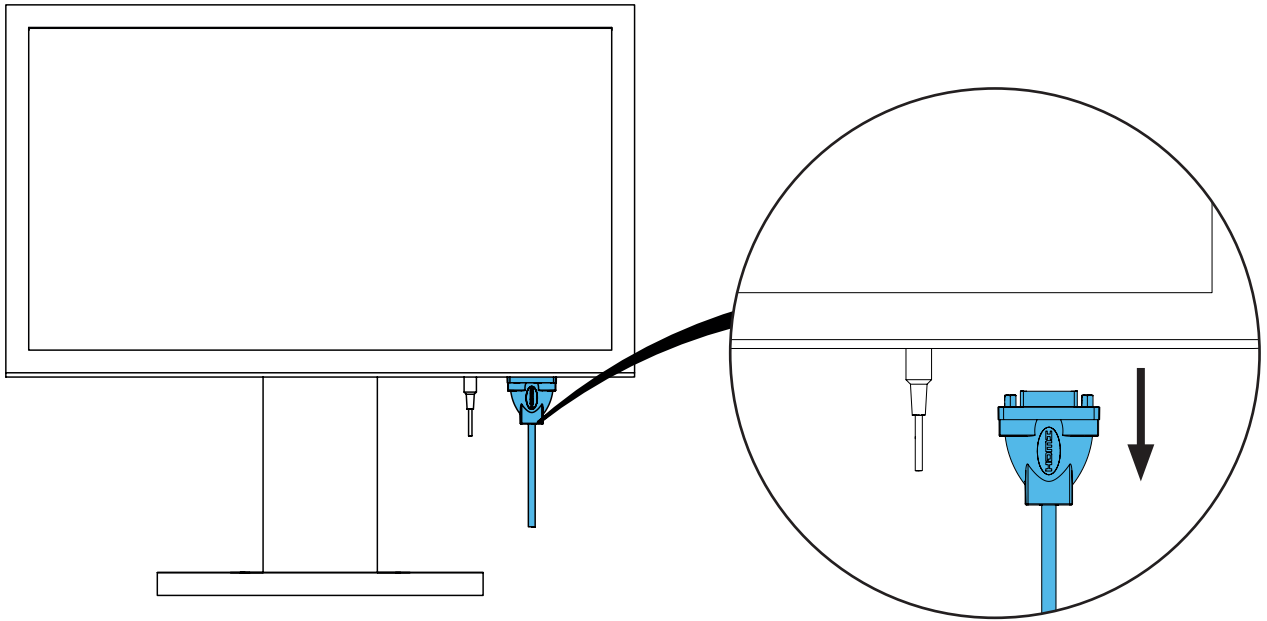


3.

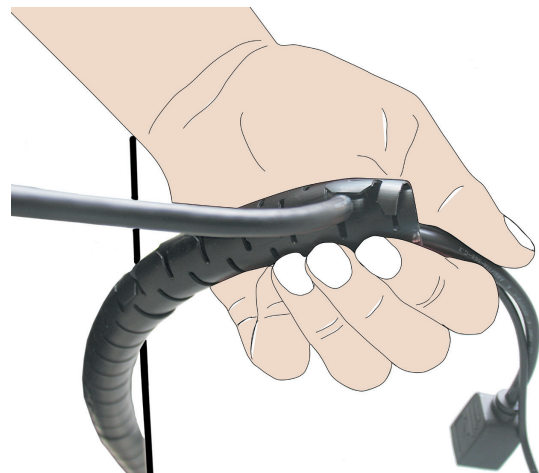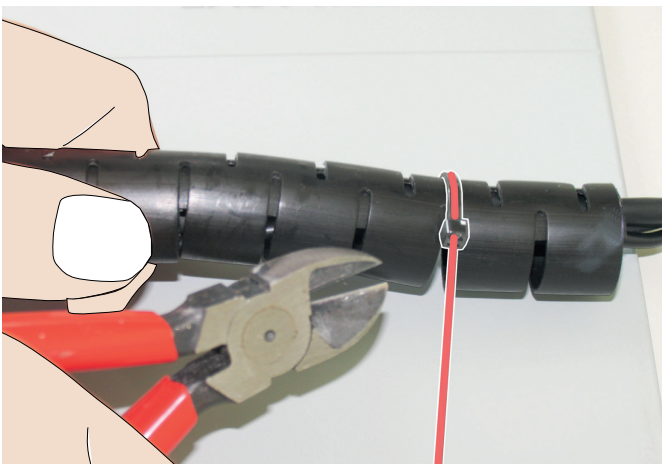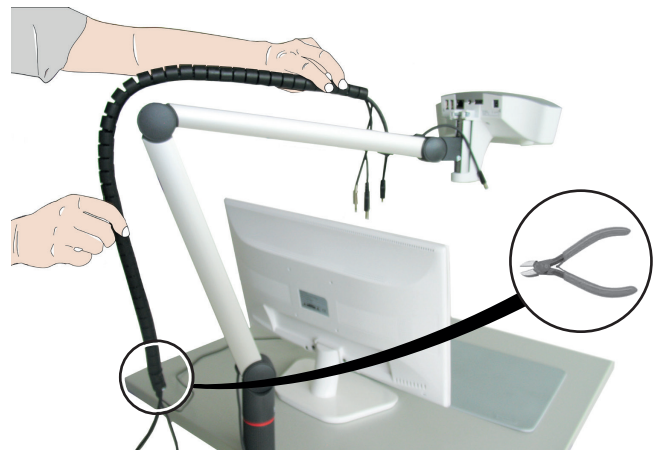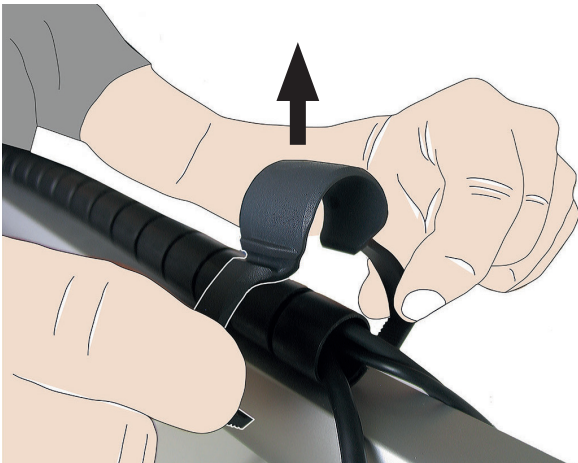

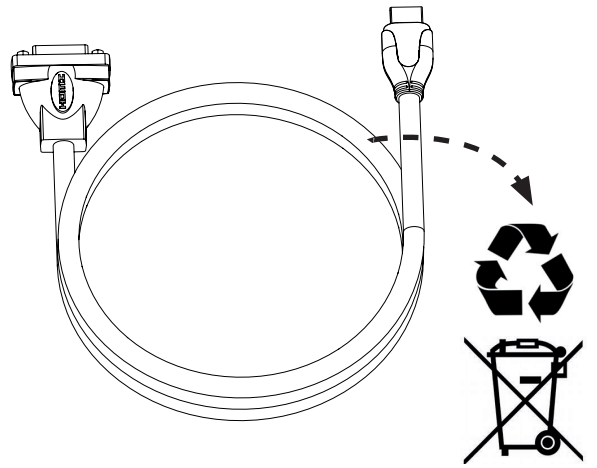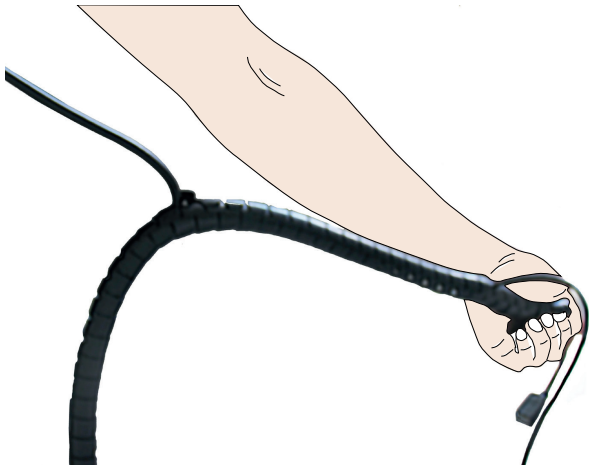


4.

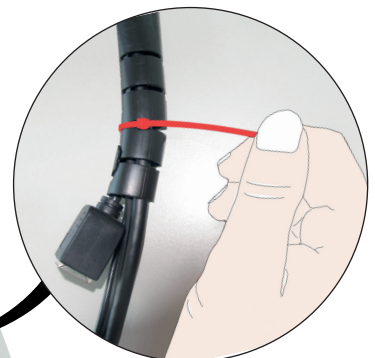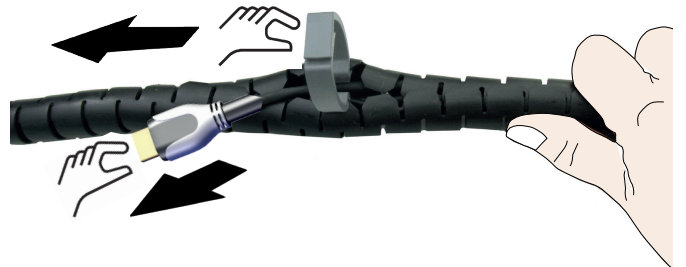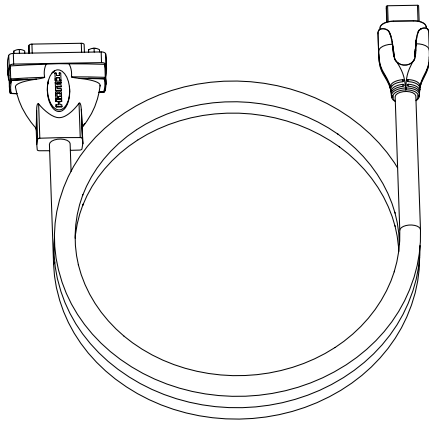


5.



6.

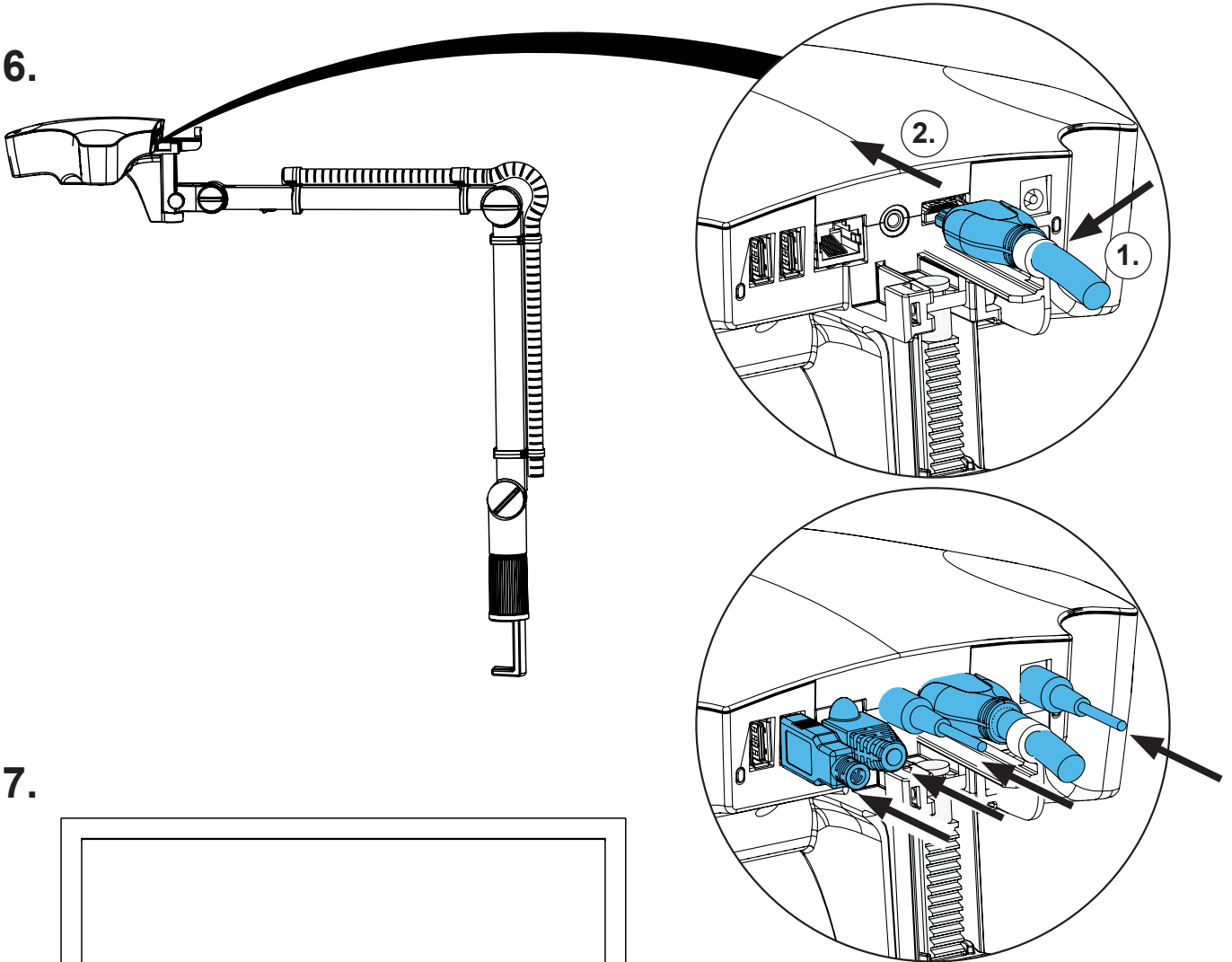

7.

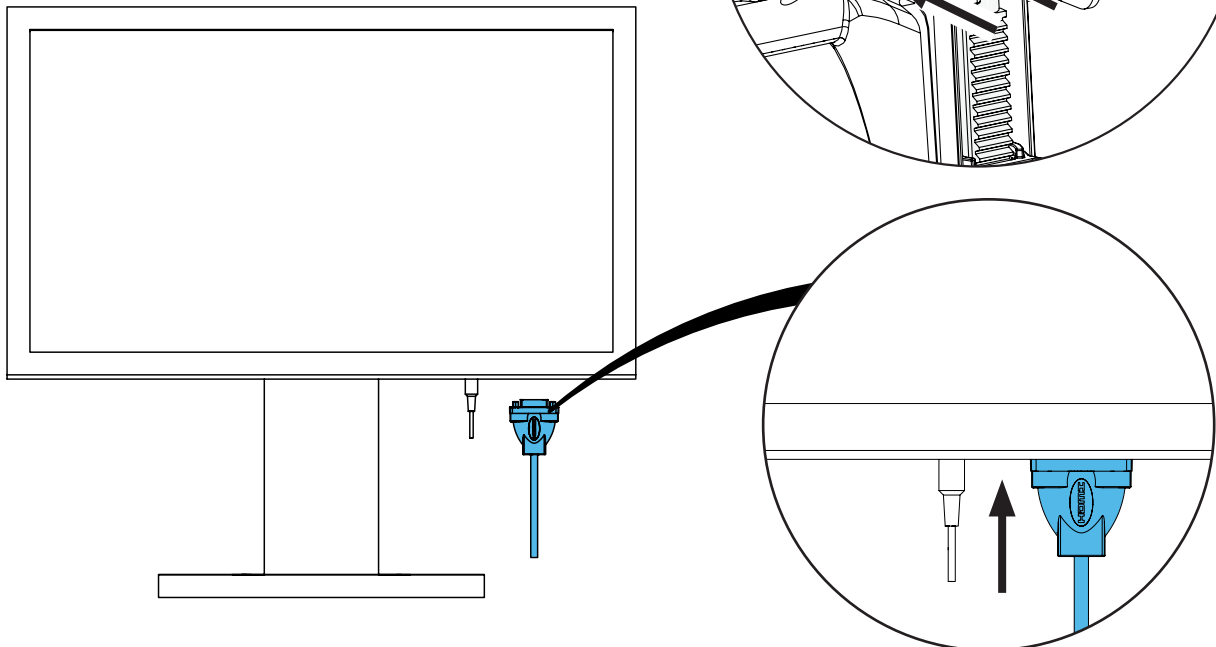






Montageanleitung • Assembly Instruction • Instruction d'installation
 Istruzioni l'istallazione • Instrucciones de montaje • Руководство по монтажу


making work easy

90003 7883

Netzteil 3D Monitor • Power supply unit 3D monitor • Adaptateur secteur écran 3D
 Alimentatore monitor 3D • Adaptador de corriente monitor 3D • Блок питания 3D монитора

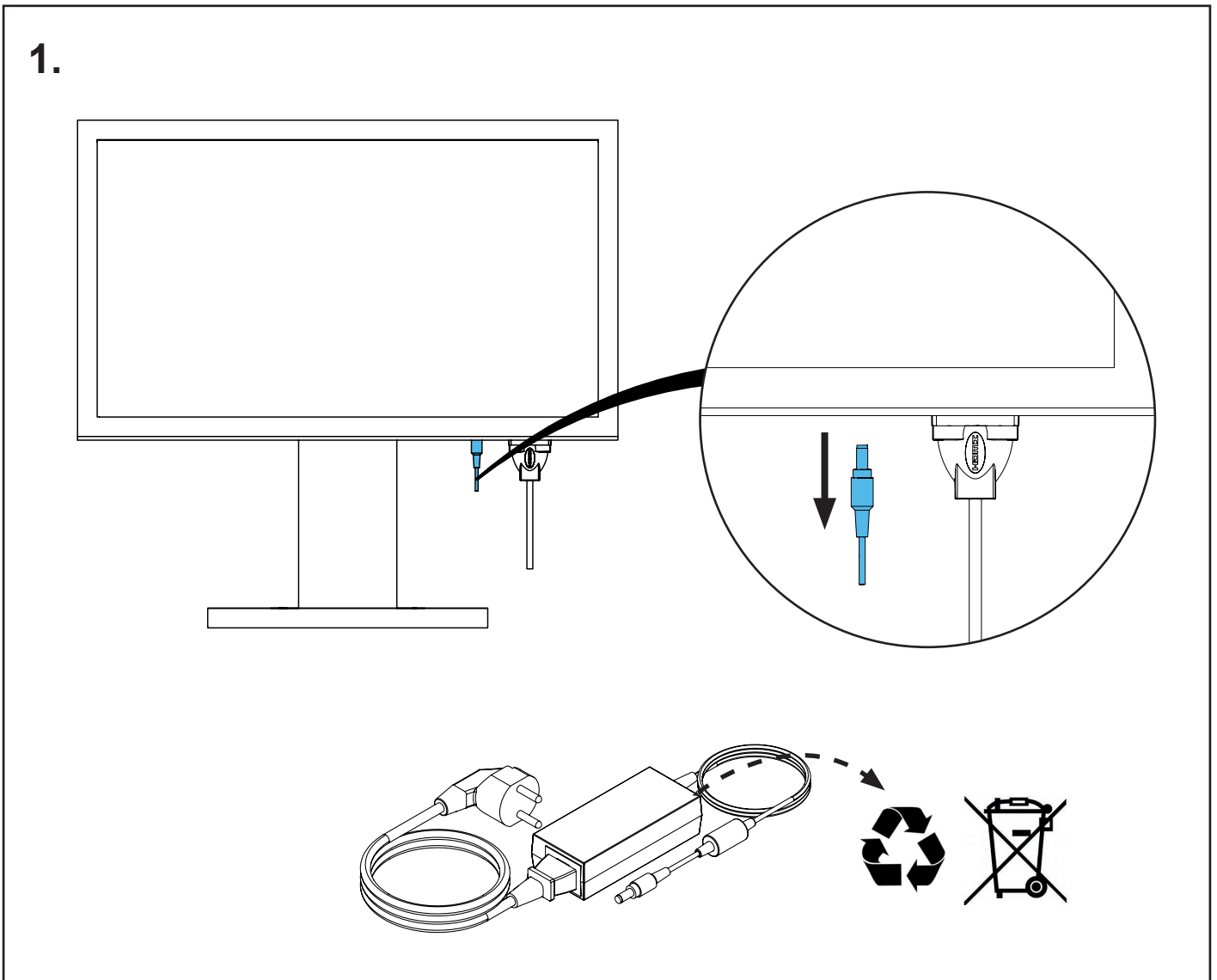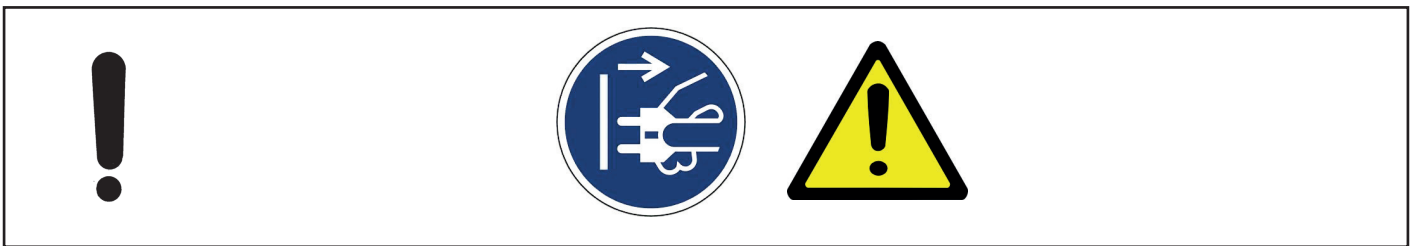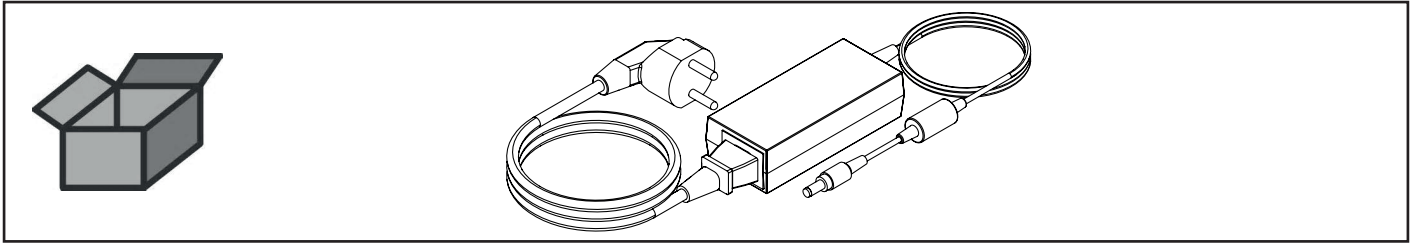

2.

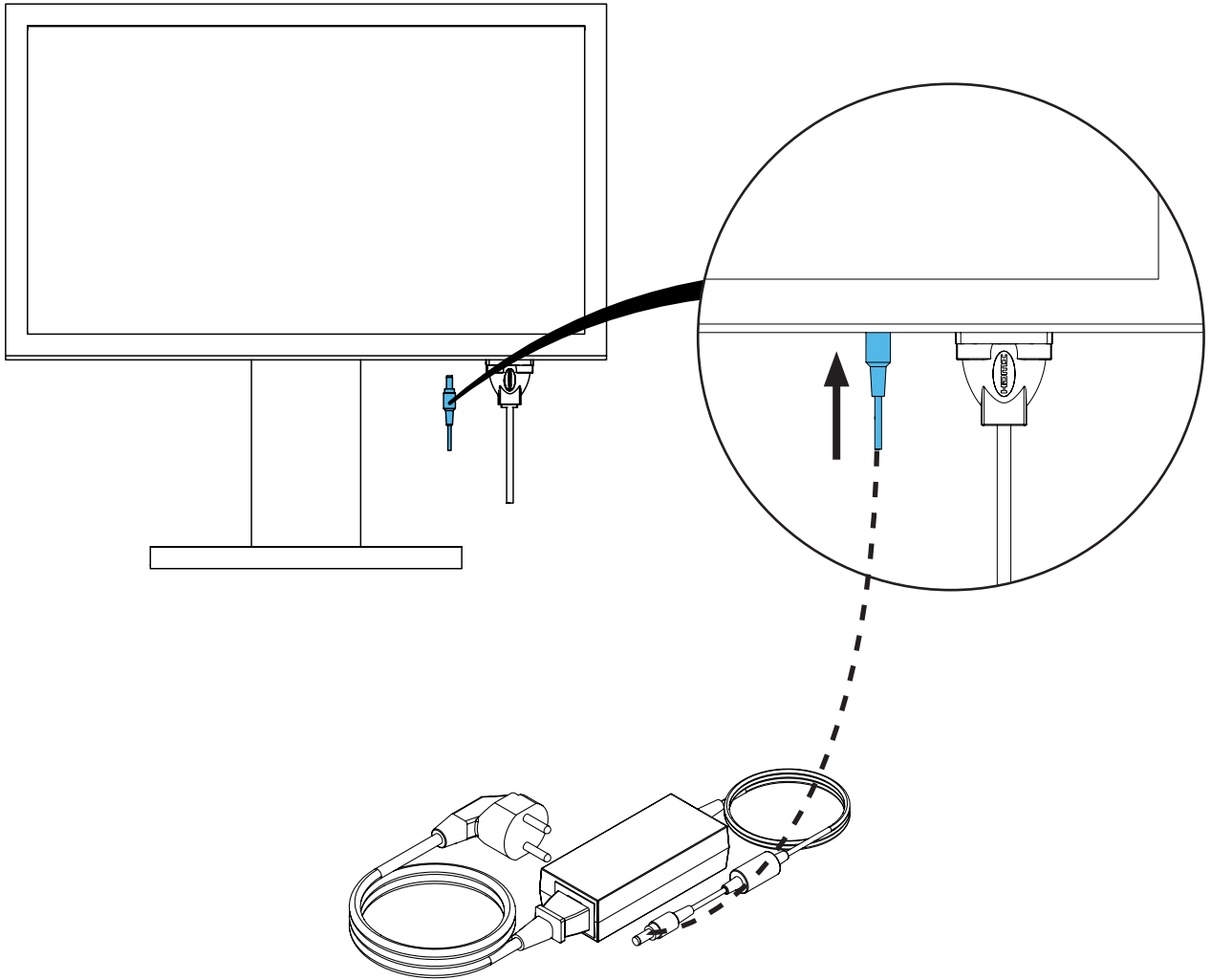






Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу



making work easy

90003 4769



Netzteil LED-Beleuchtung • Power supply LED lighting • Poste secteur LED lam-
pe • Trasformatore lampada LED • Bloque alimentación lampara LED • Блок
питания со светодиодной подсветкой

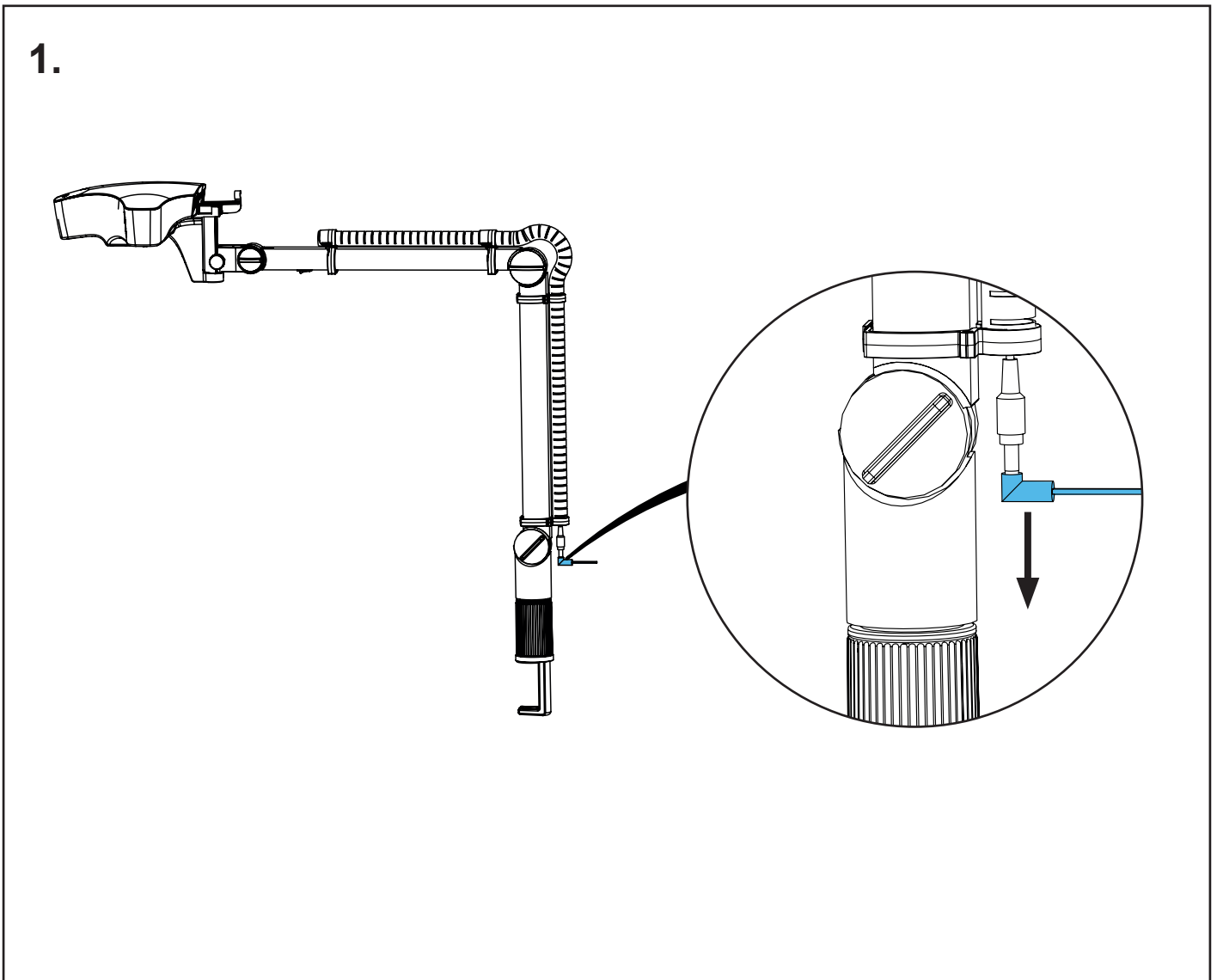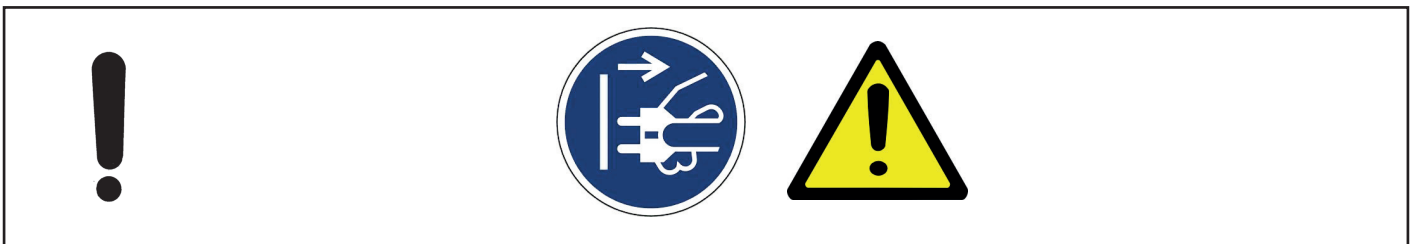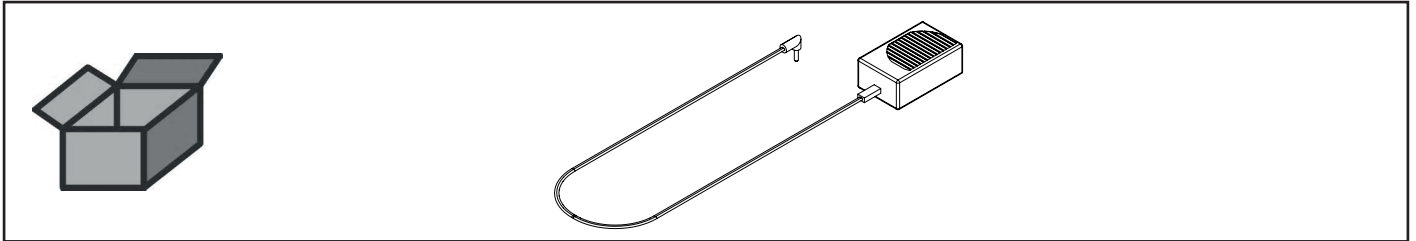


2.

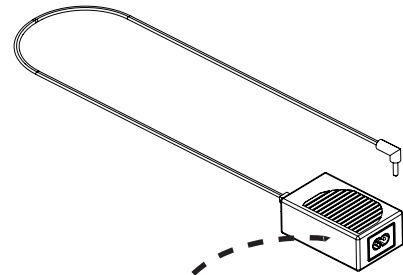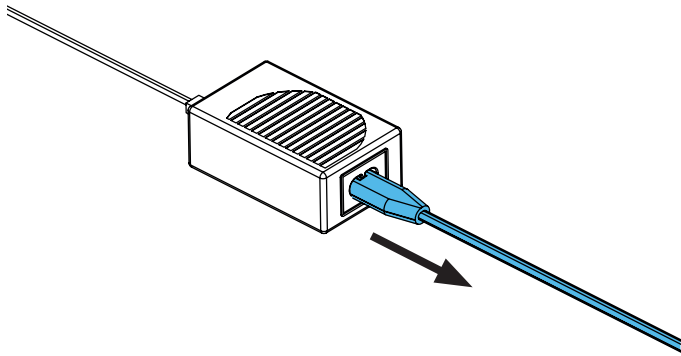

3.

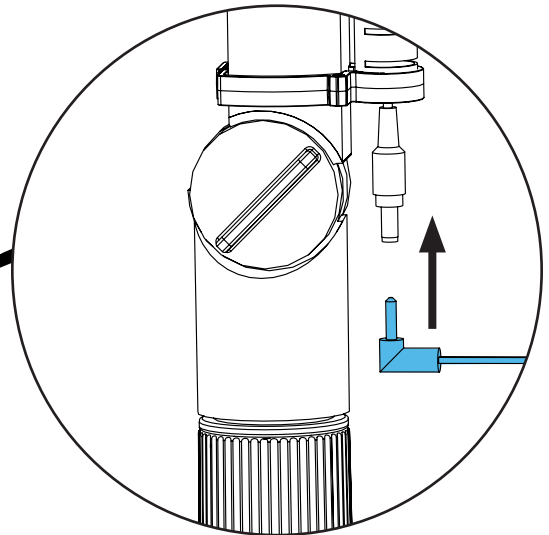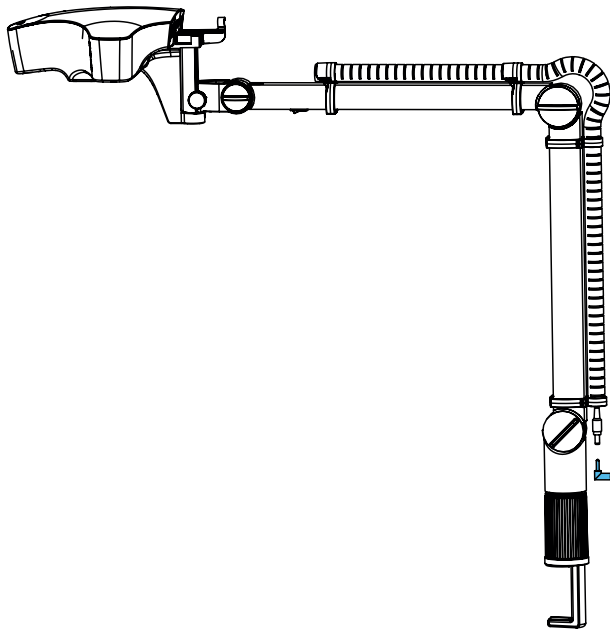

4.

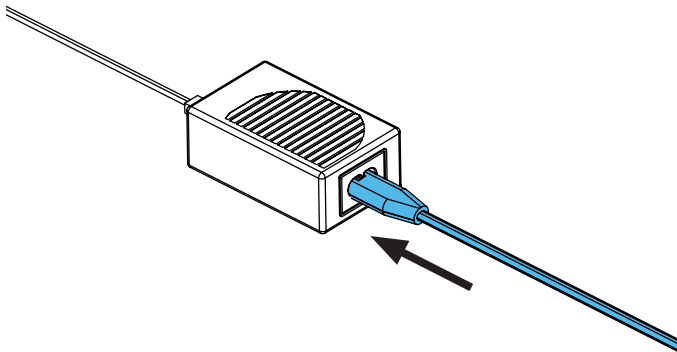






Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу


making work easy

90003 5975

Netzkabel • Power cord • Câble alimentation • Cavo di rete • Cable de corriente
Сетевой кабель

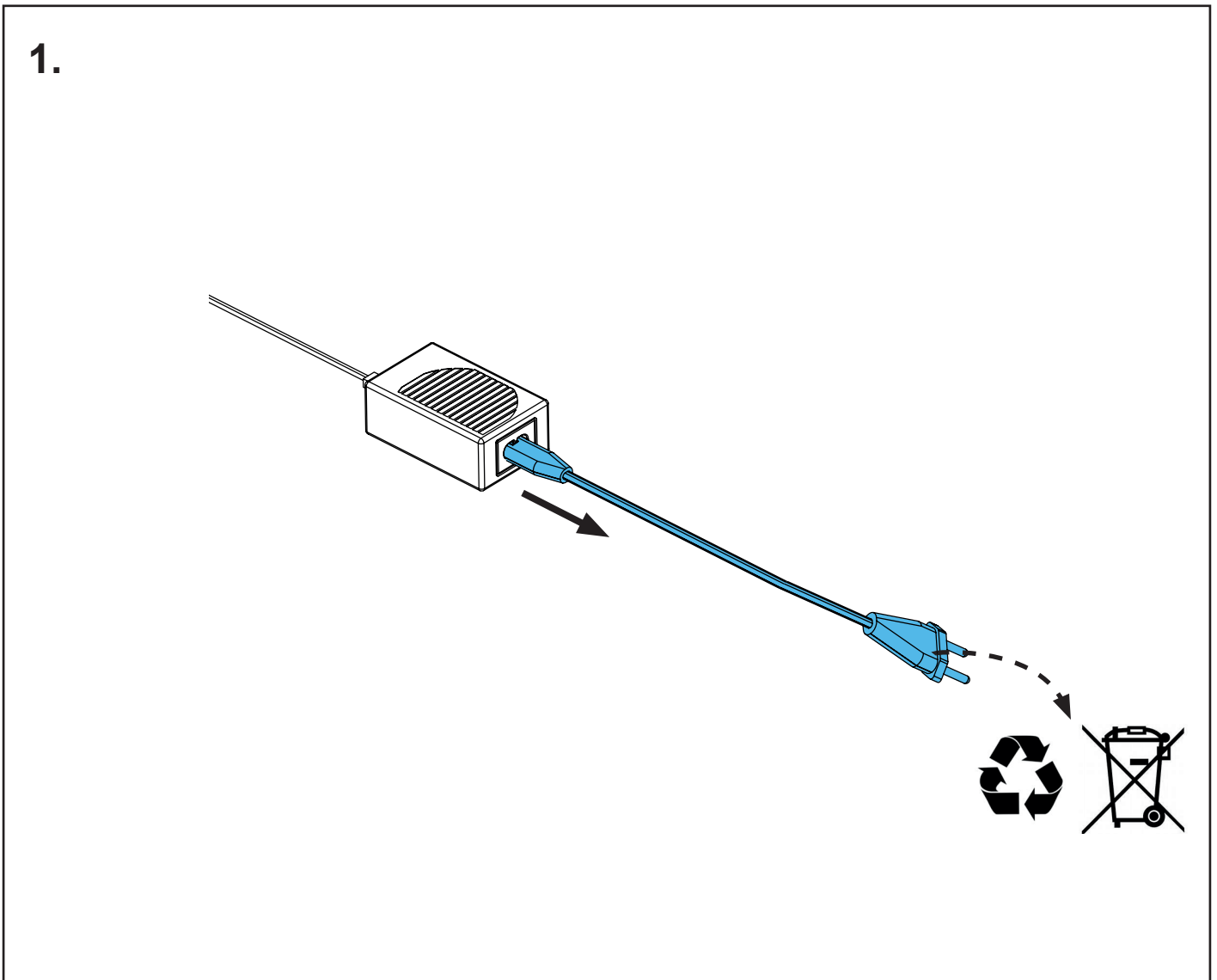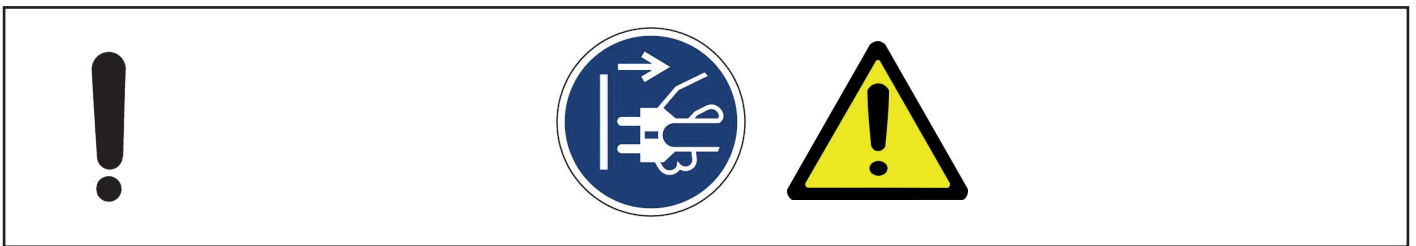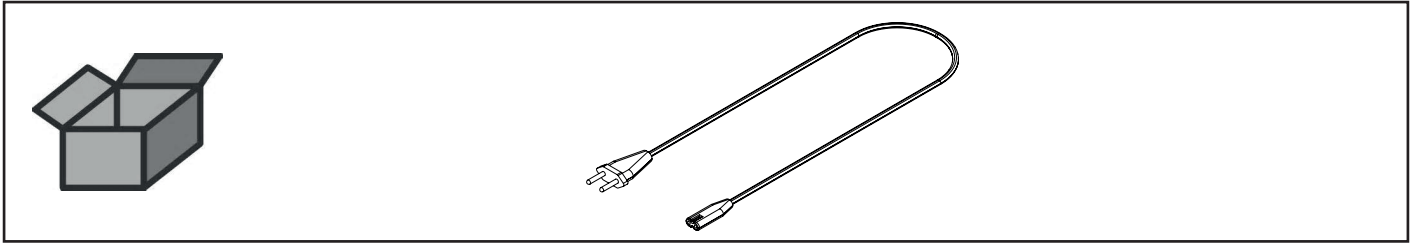

2.


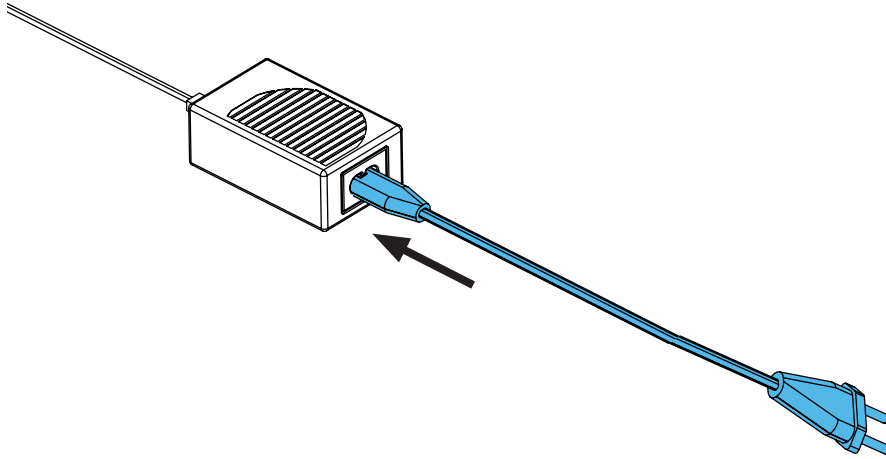





making work easy

90003 3758



Schutzglas • Protective glass • Verre de protection • Vetro protettivo • Cristal de
protección • Защитное стекло

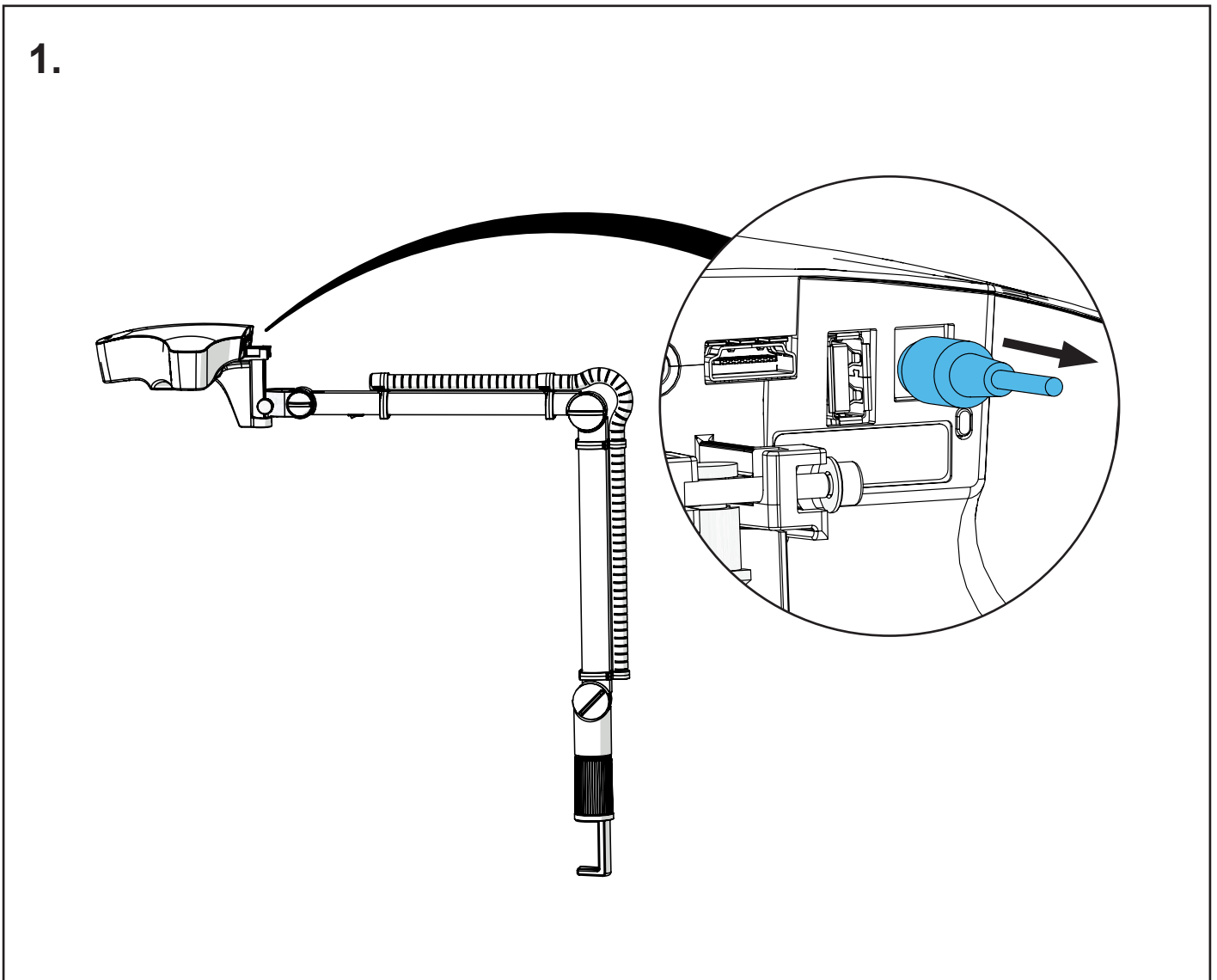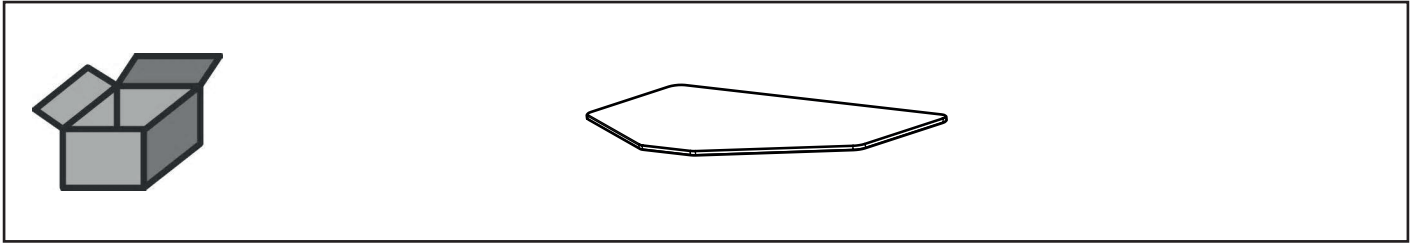




2.

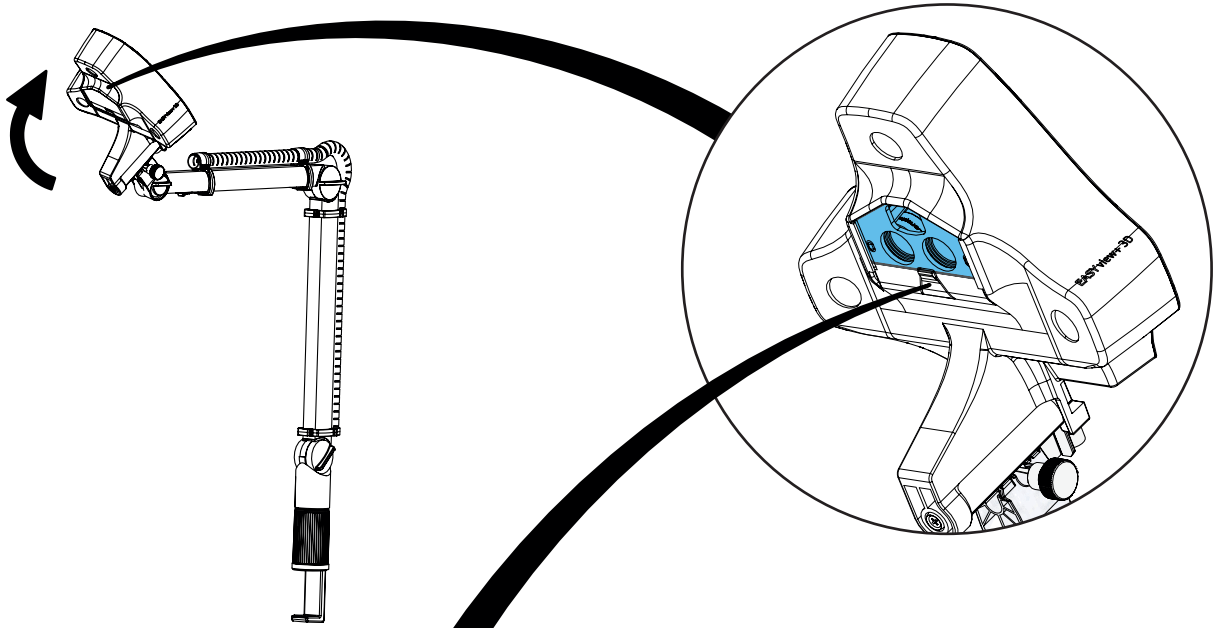


3.

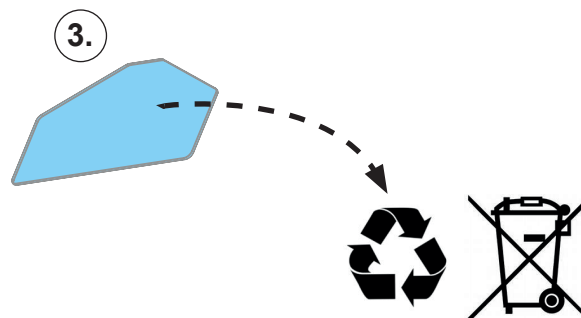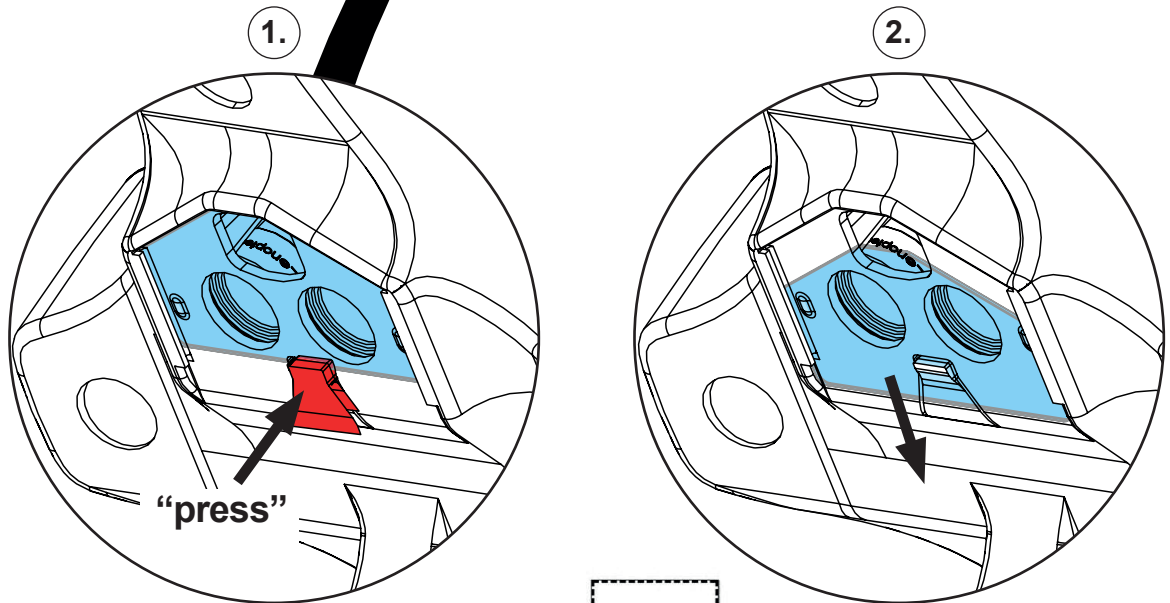


4.

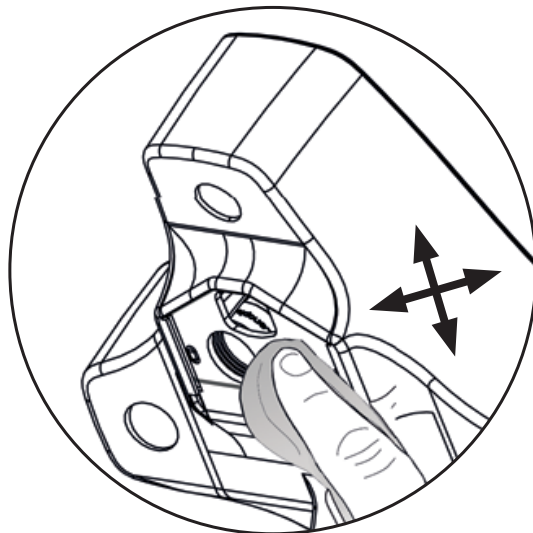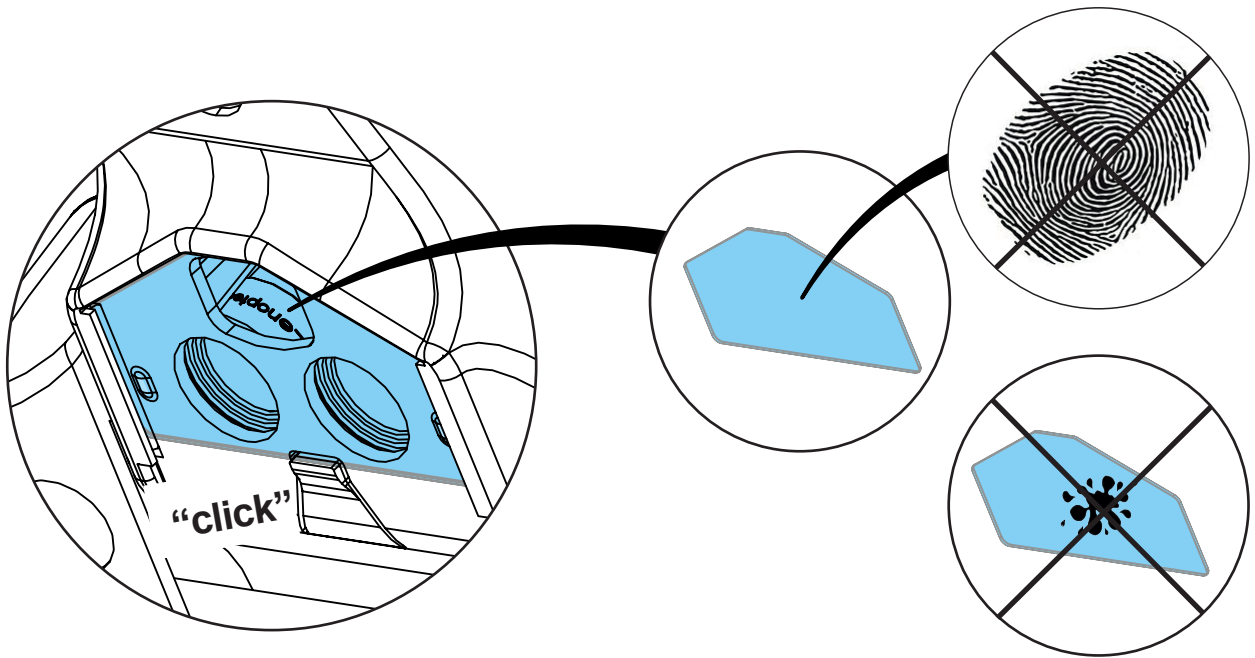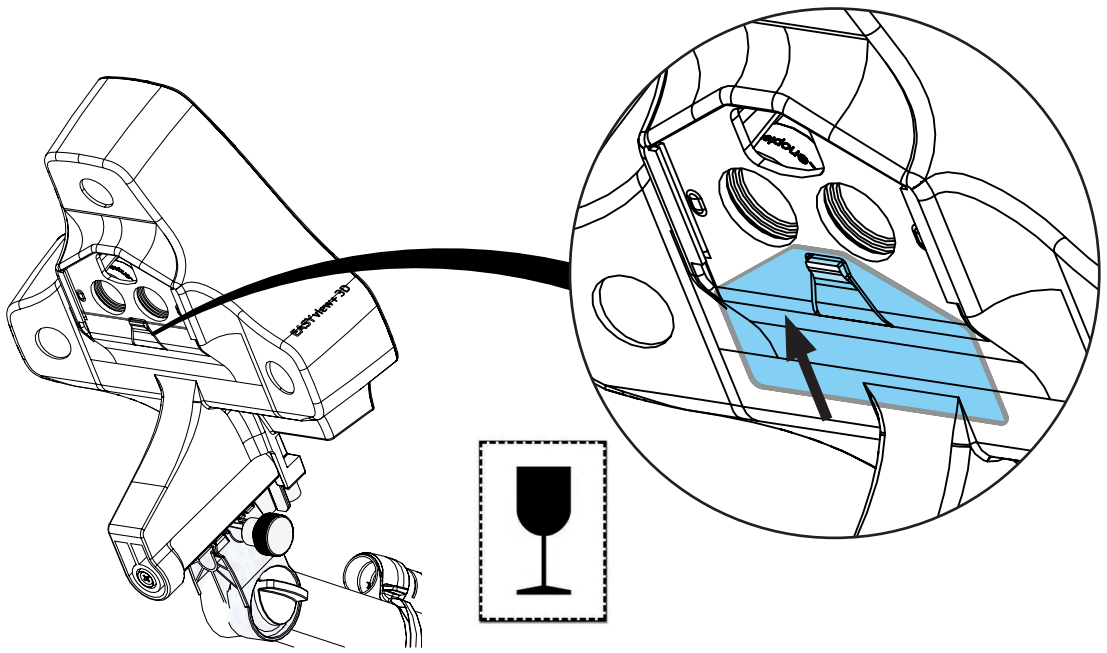

5.

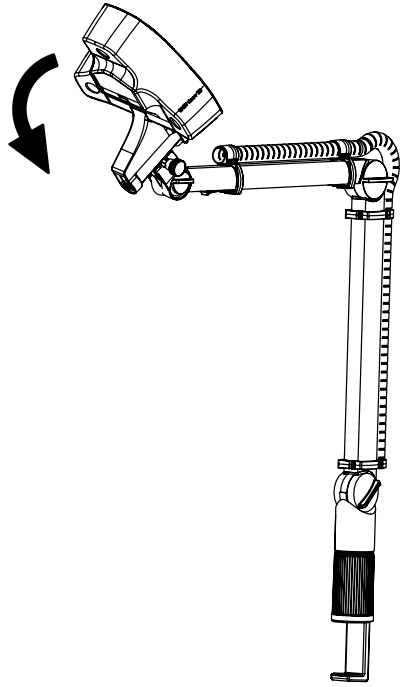

6.

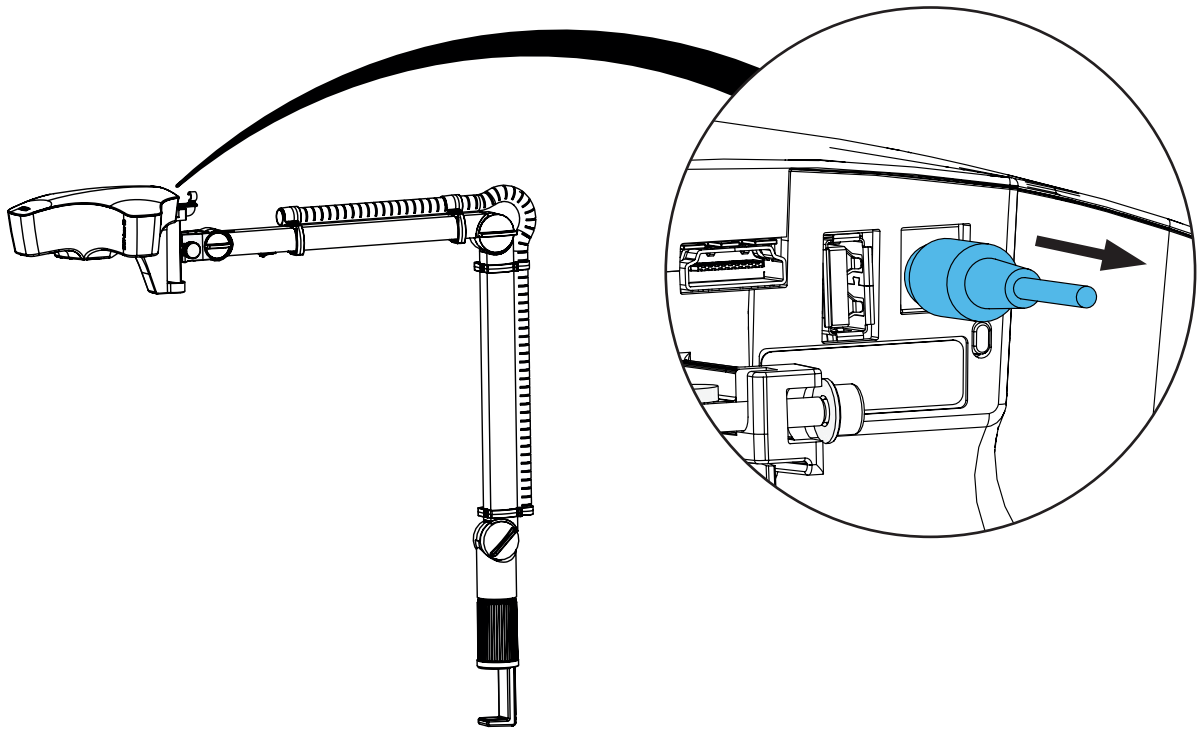







making work easy


Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по монтажу

90003 7903





HDMI-HDMI Kabel • HDMI-HDMI cable • Câble HDMI HDMI • Cavo HDMI HDMI
Cable HDMI HDMI • Кабель HDMI-HDMI

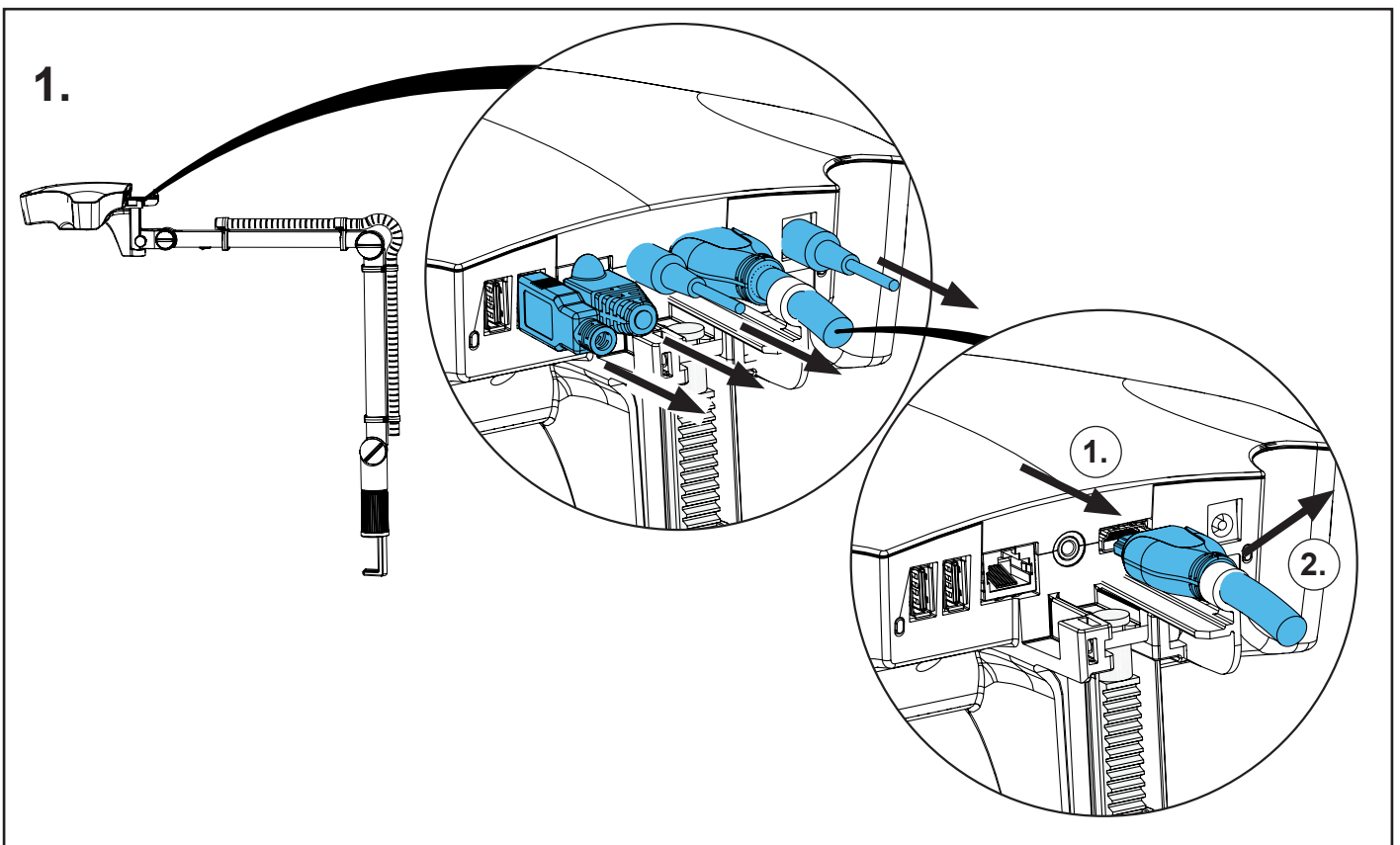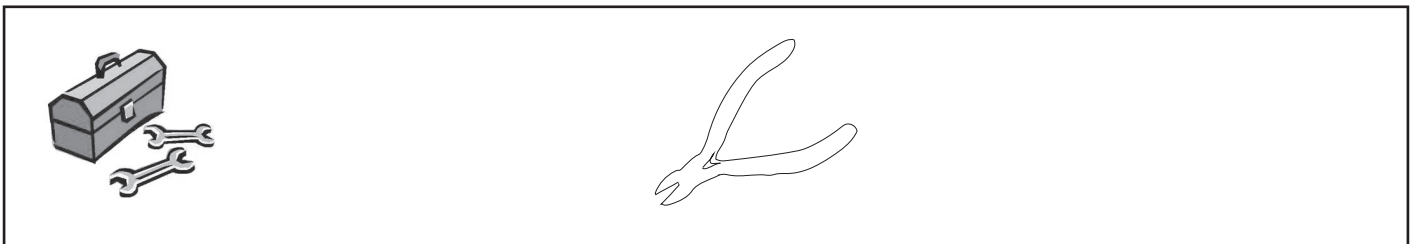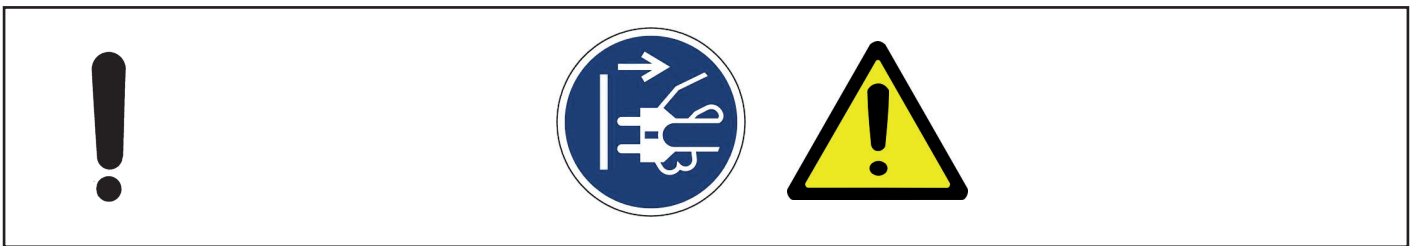
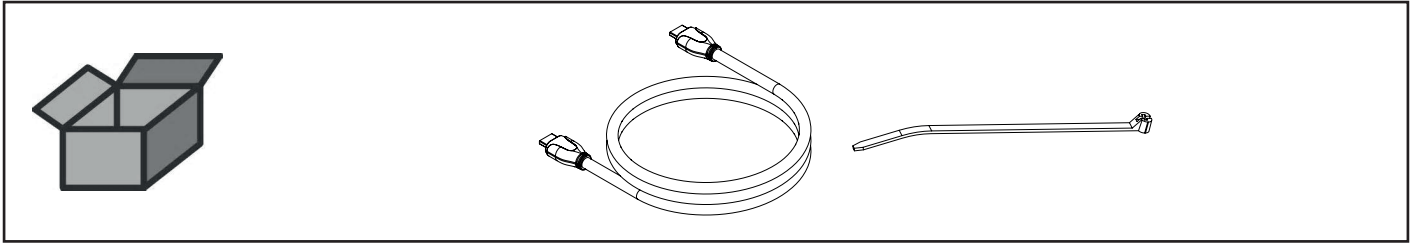


2.


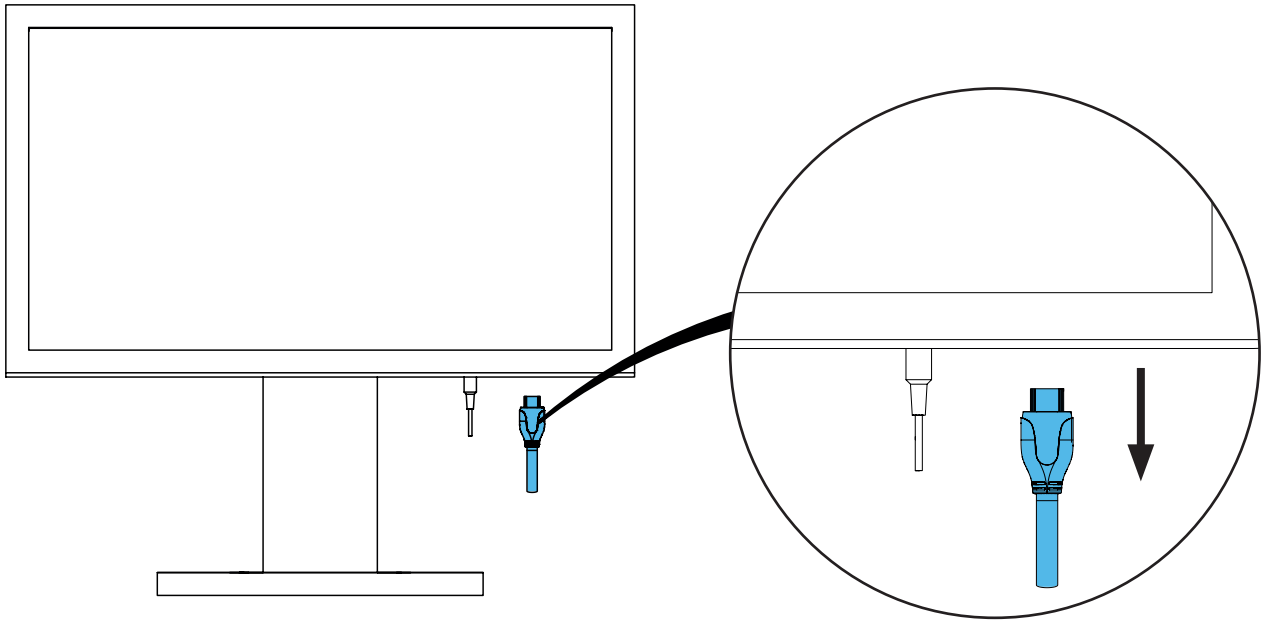


3.

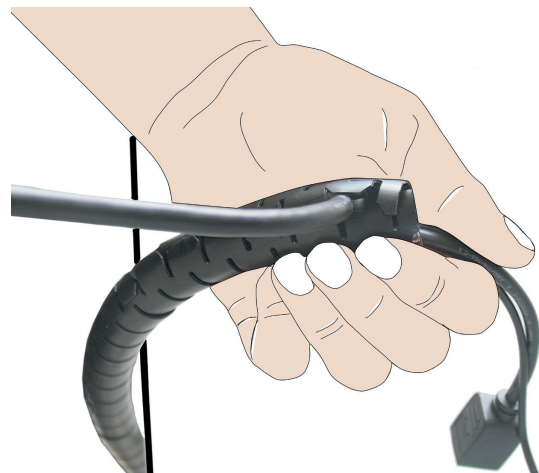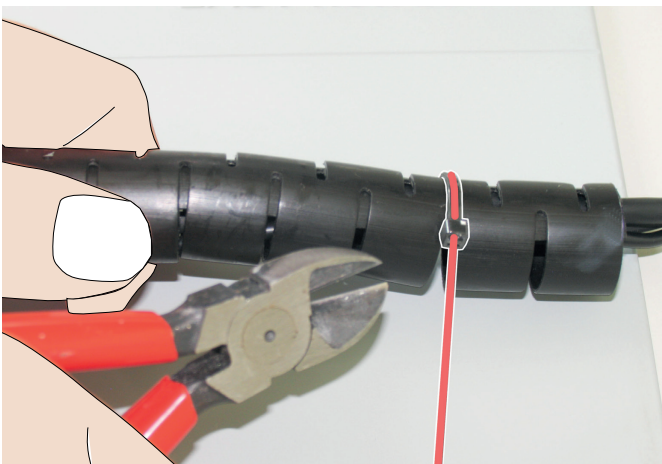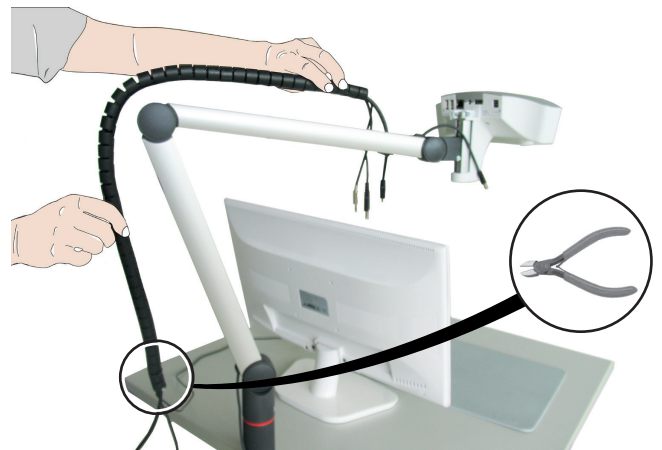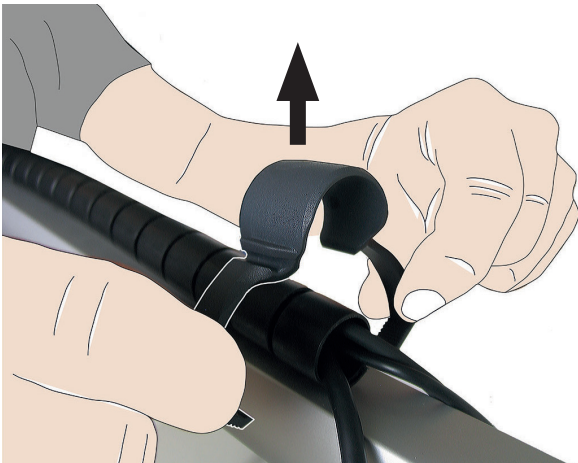

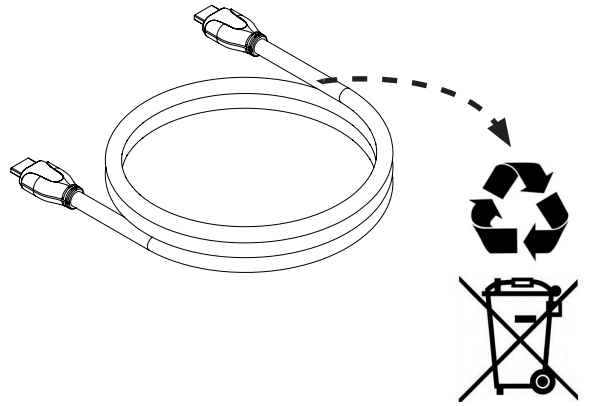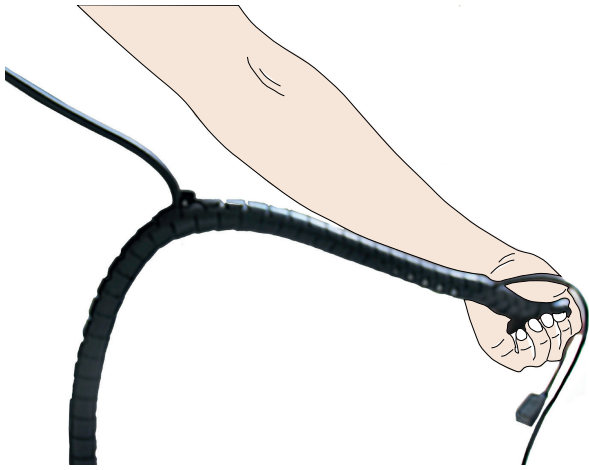


4.

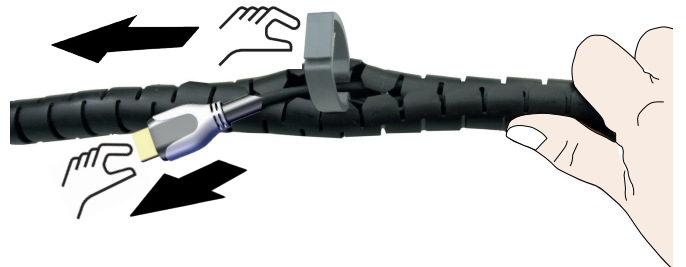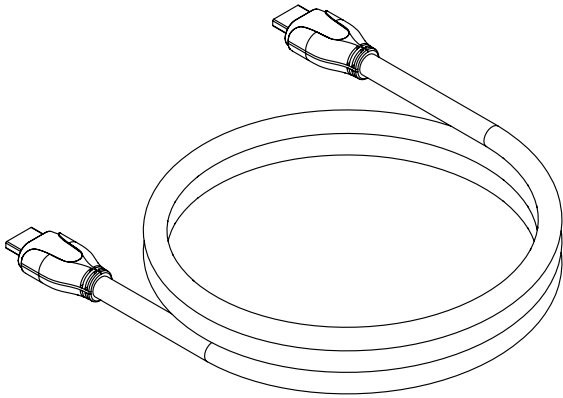


5.



6.

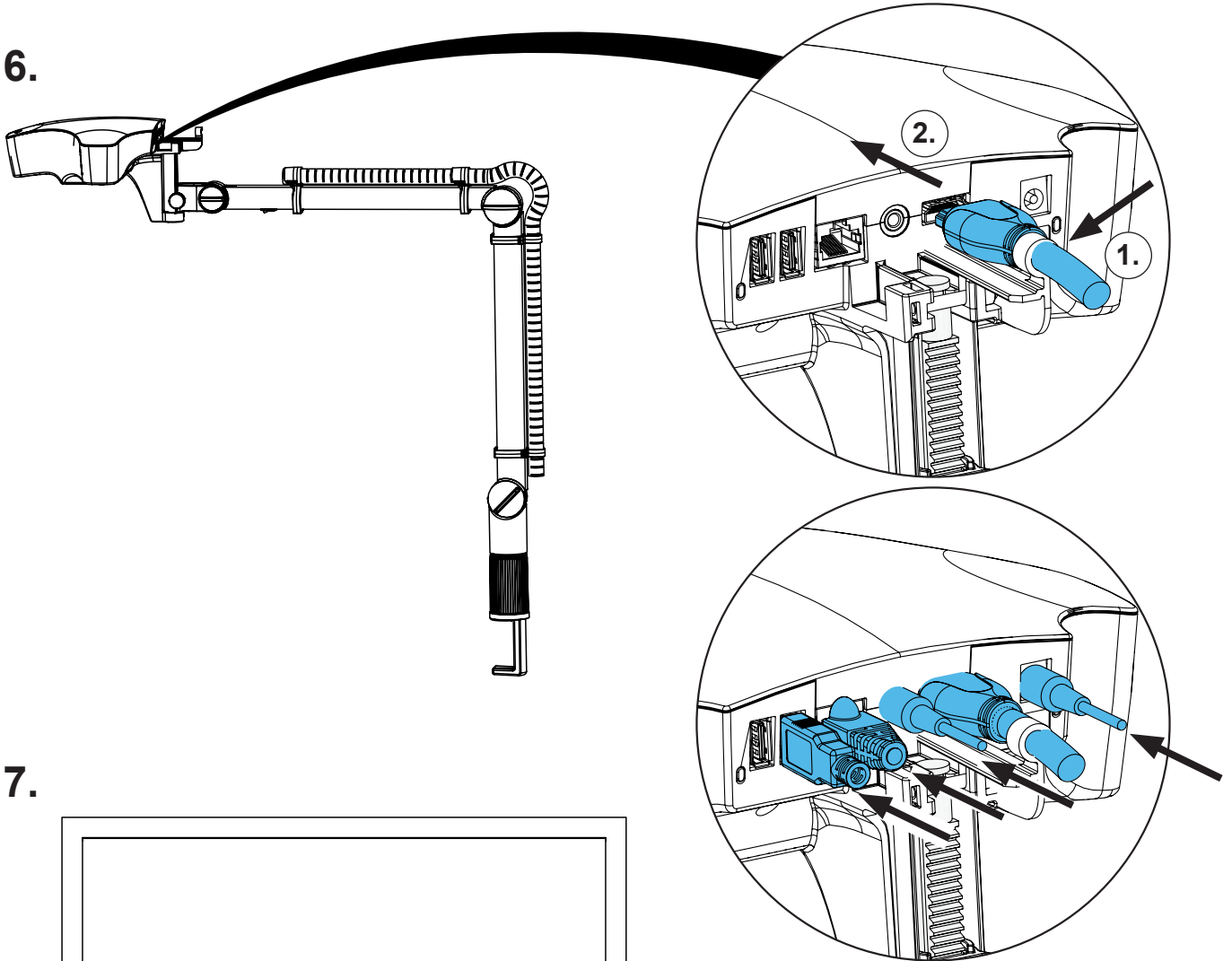

7.

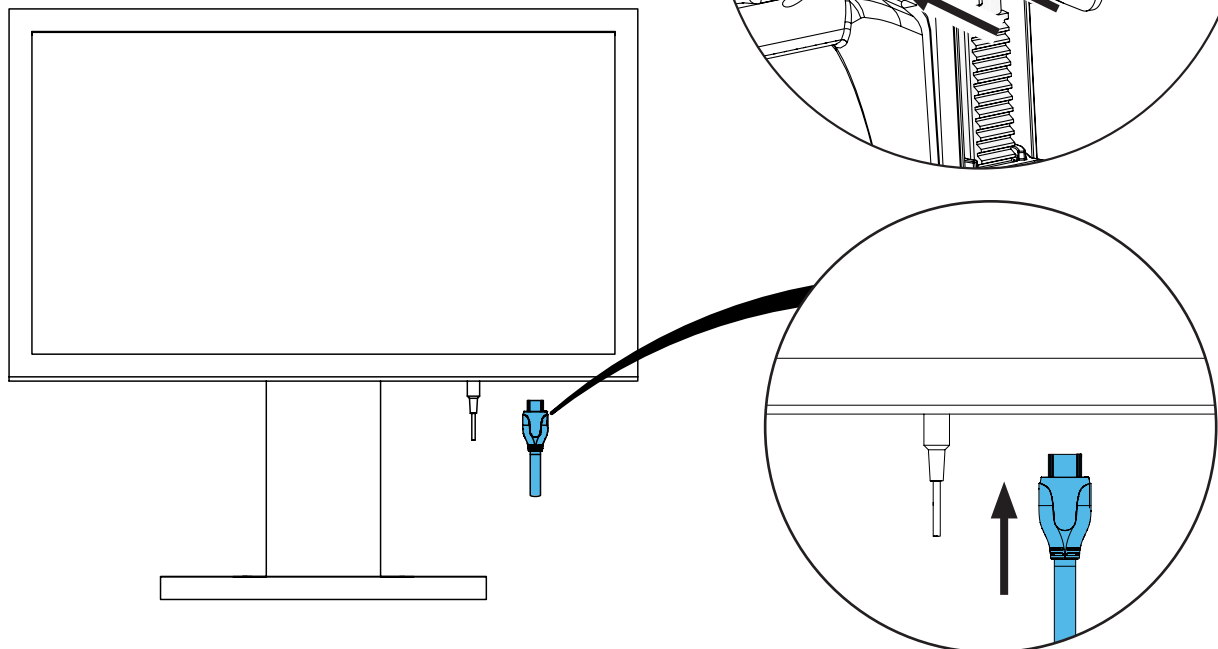






making work easy

Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу

90003 7895

USB Stick • USB stick • Clé USB • Chiave USB • Lápiz de memoria USB
USB-накопитель

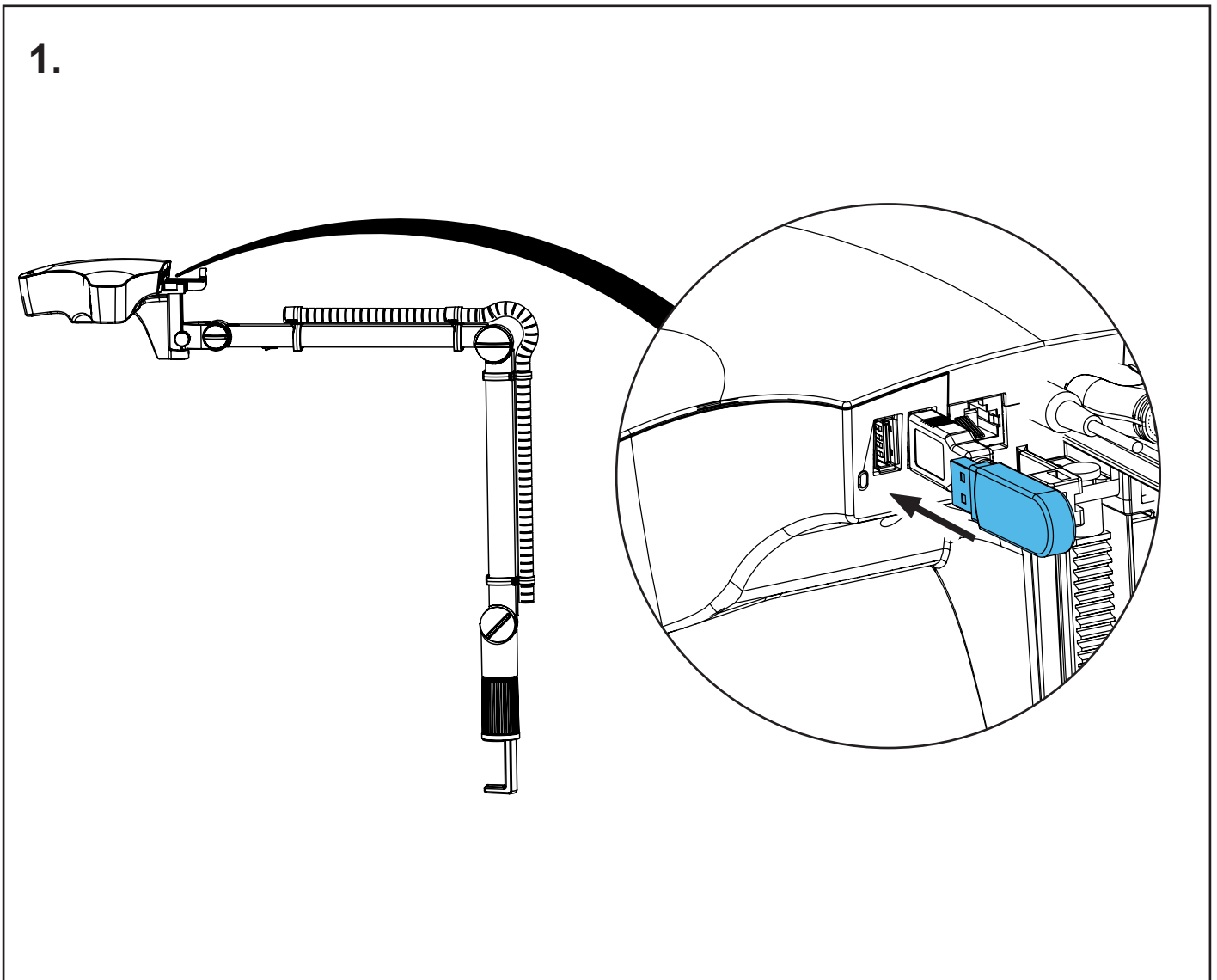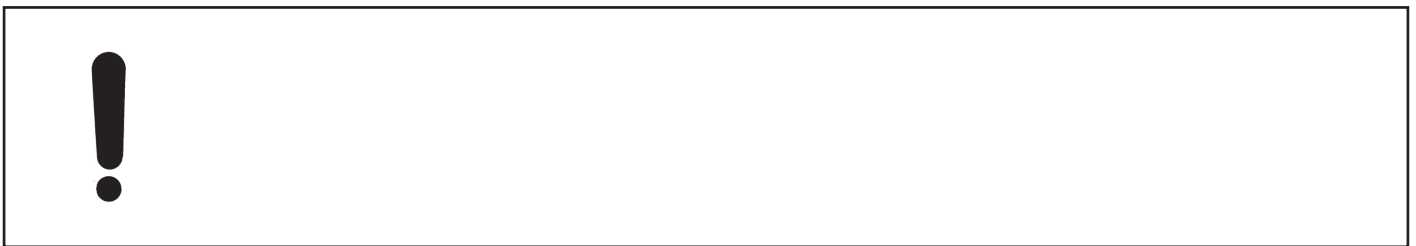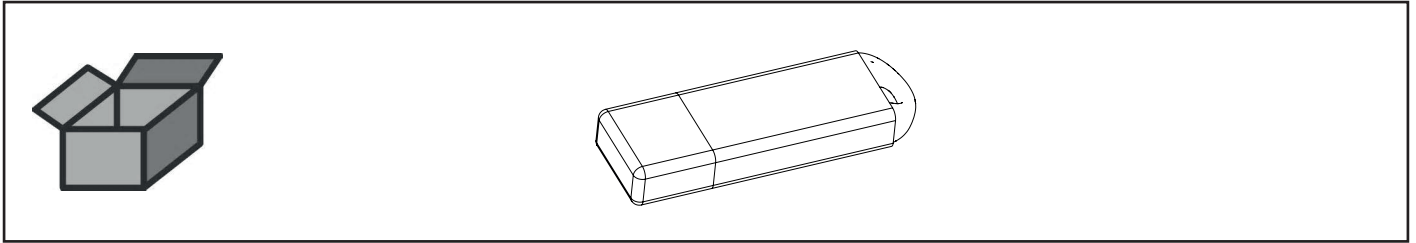





Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу

making work easy

90003 7552

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Universalnetzsteckerset • Universal power plug kit • Jeu universel de fiches
secteur • Set spine di rete universale • Kit de enchufes eléctricos universales •
Сетевой штекер универсальный

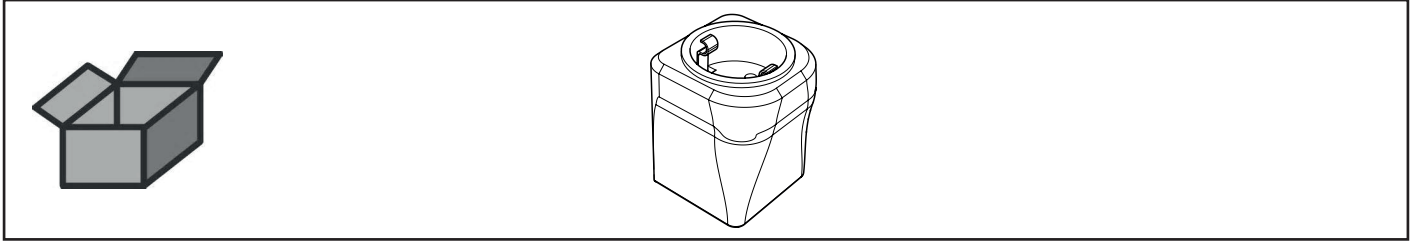





making work easy

Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу

90003 7916

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Steckdosenleiste • Multiple socket • Bloc multiprise • Presa multipla • Regleta
de enchufe • 3-кратный разъем

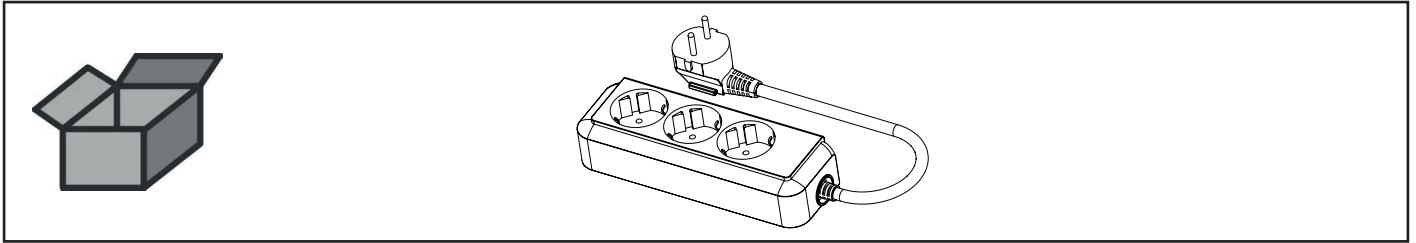







Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу


making work easy

90003 7909



Netzwerkverlängerung • Network extension cable • Rallonge de câble réseau
• Prolunga per cavo di rete • cable alargador de red • Удлинитель сетевого
интернет-кабеля

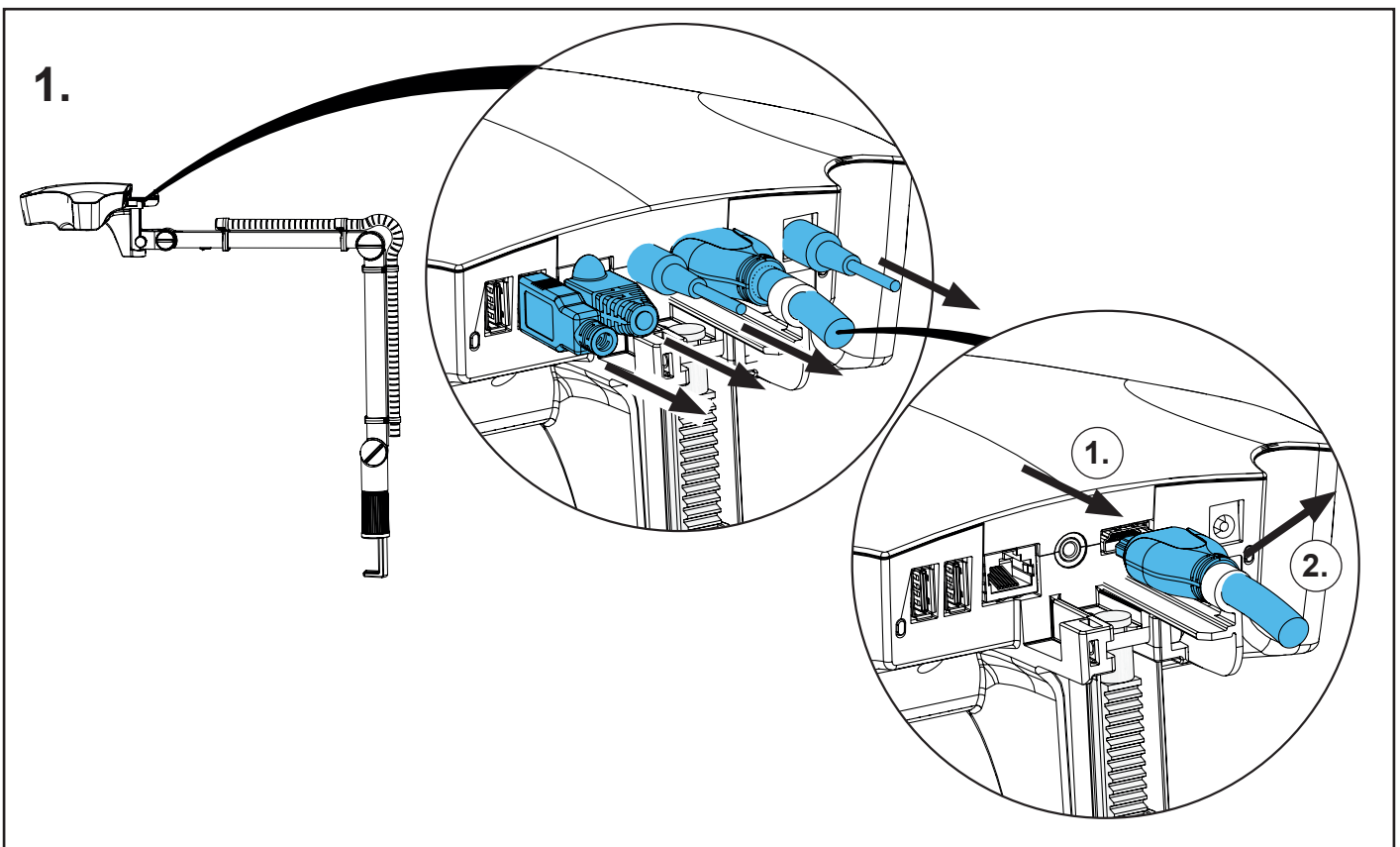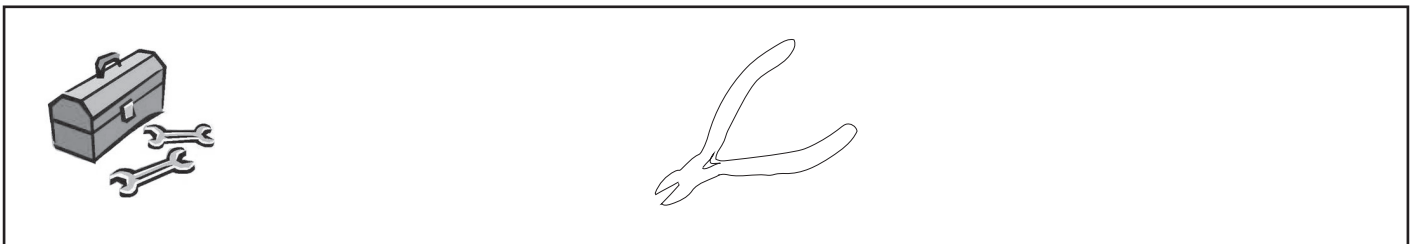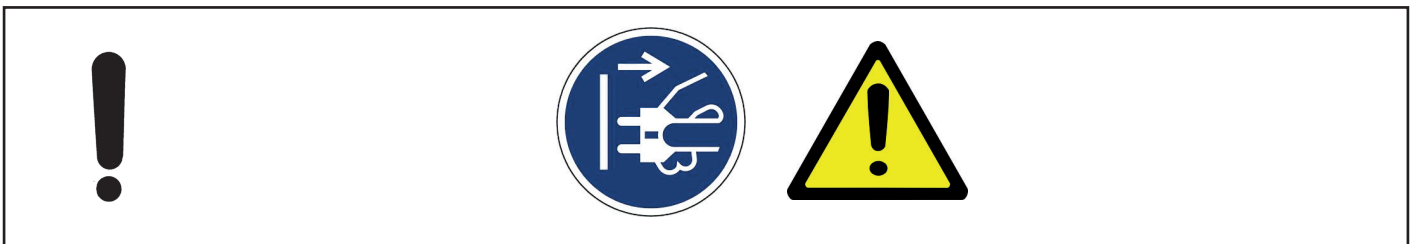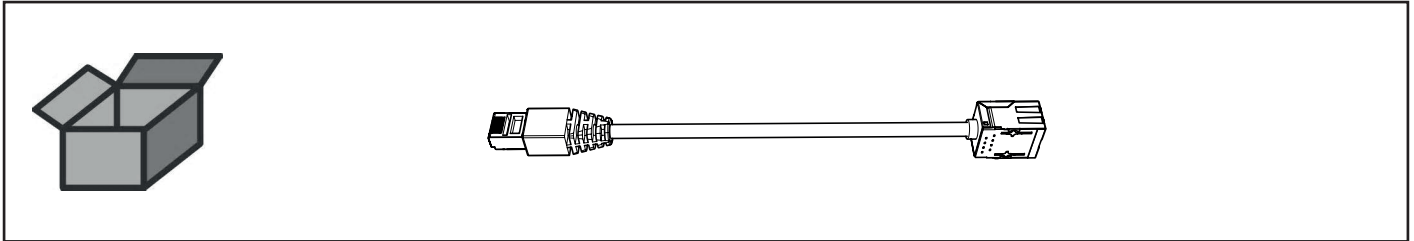



2.

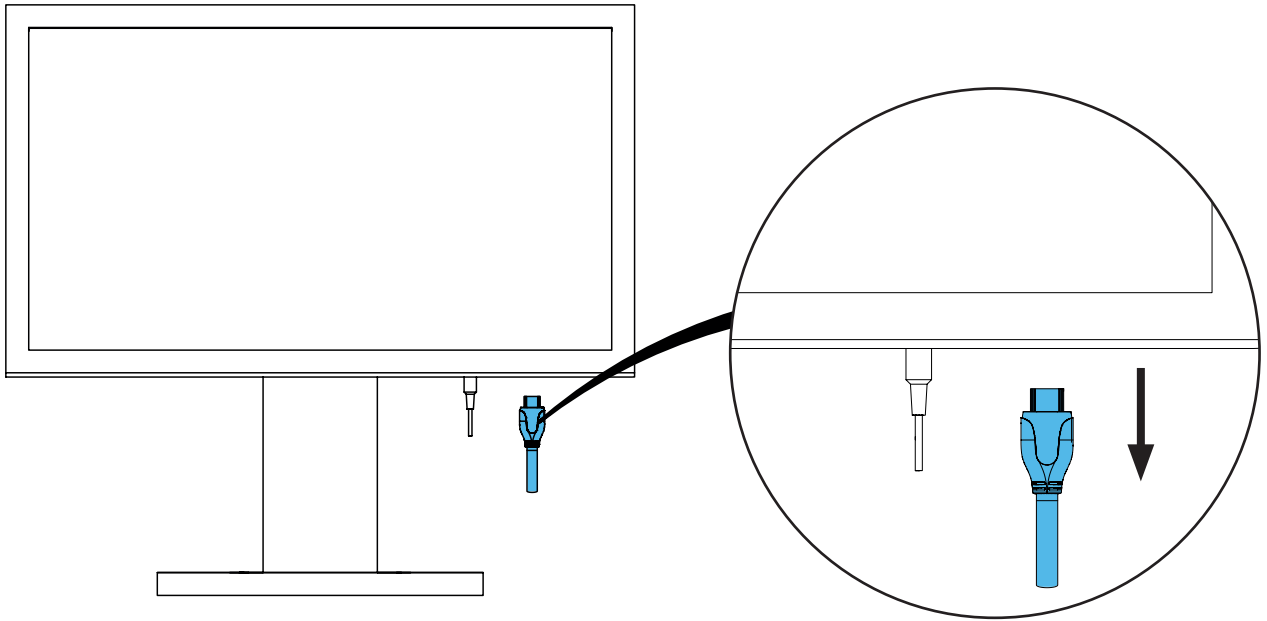




3.

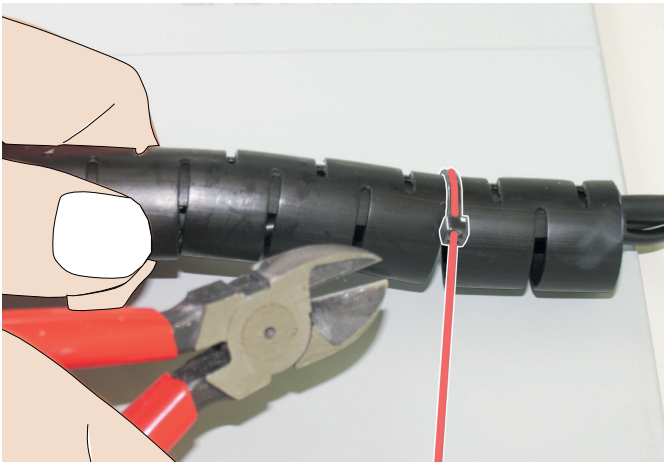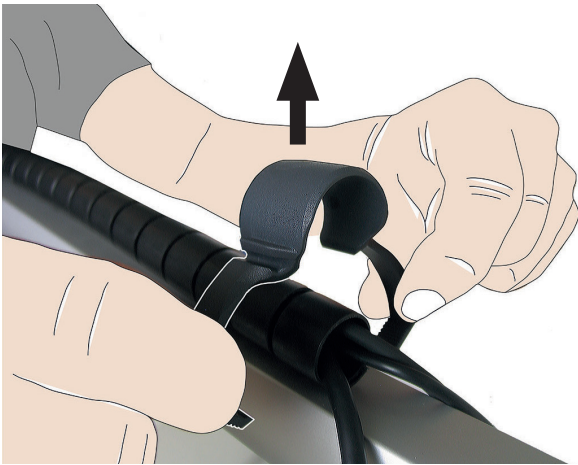


4.

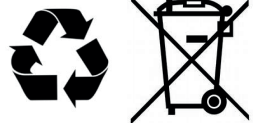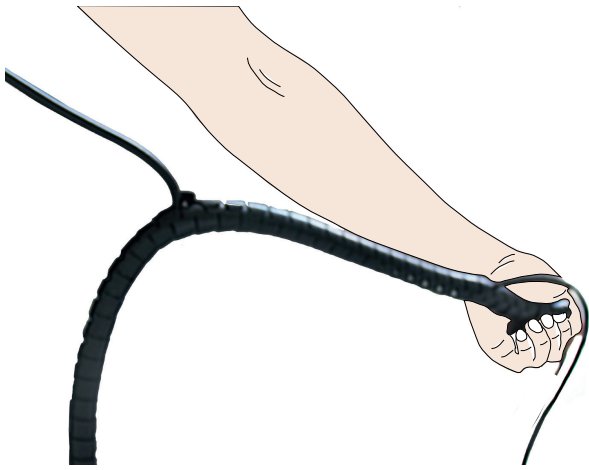

5.

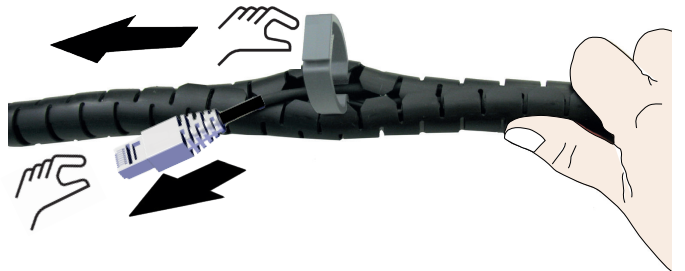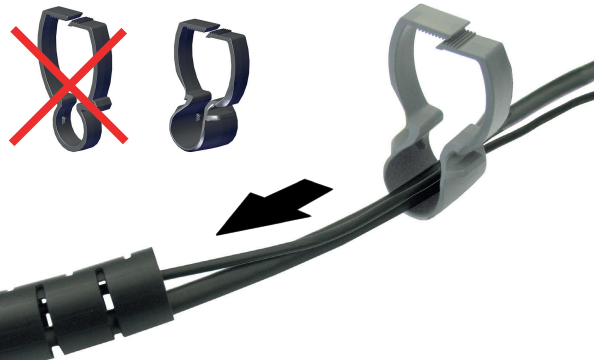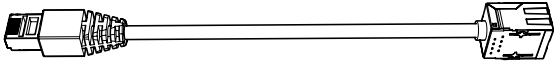


6.

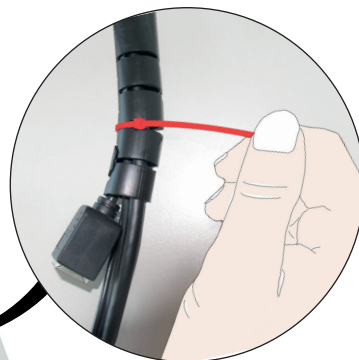


7.



8.

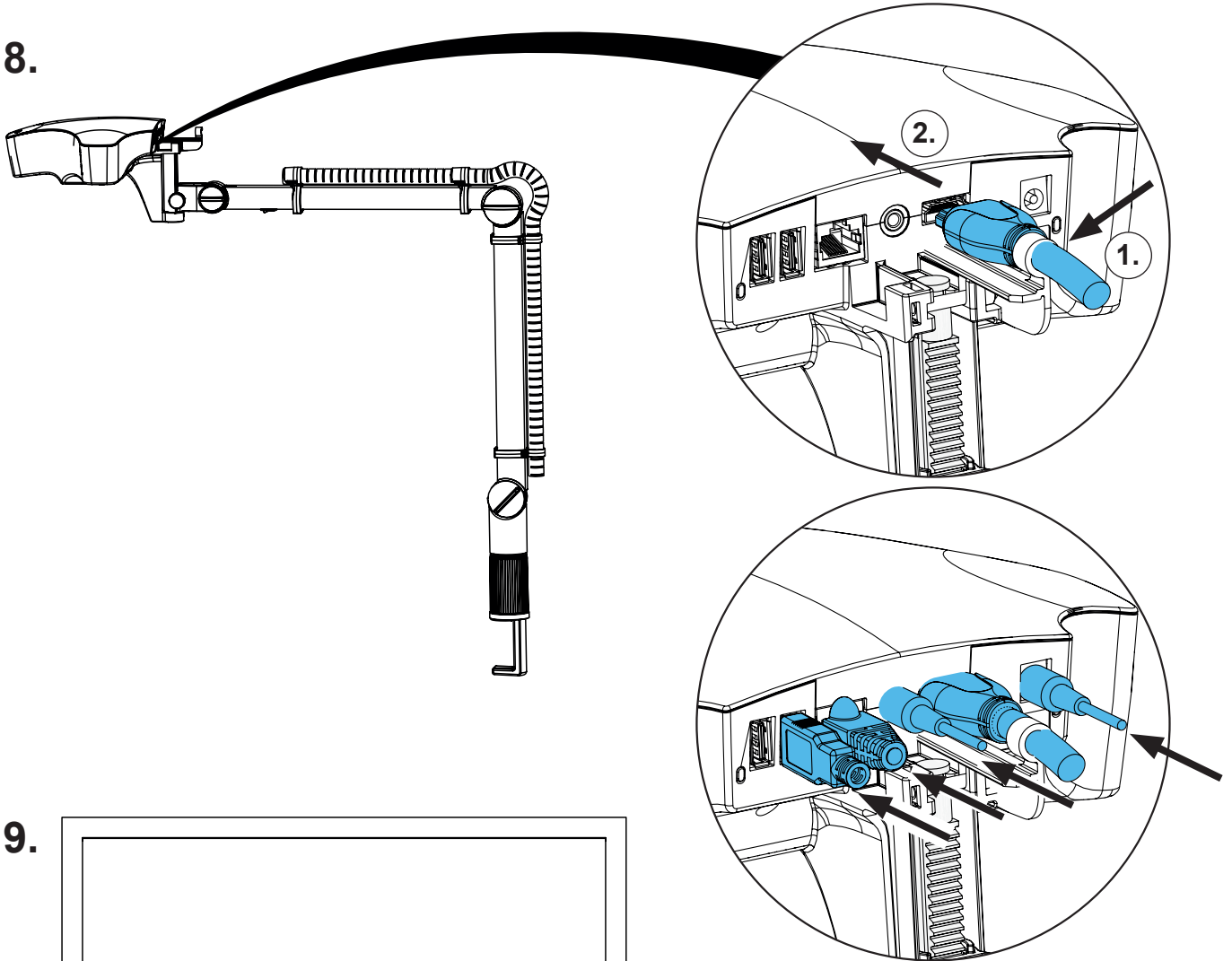

9.

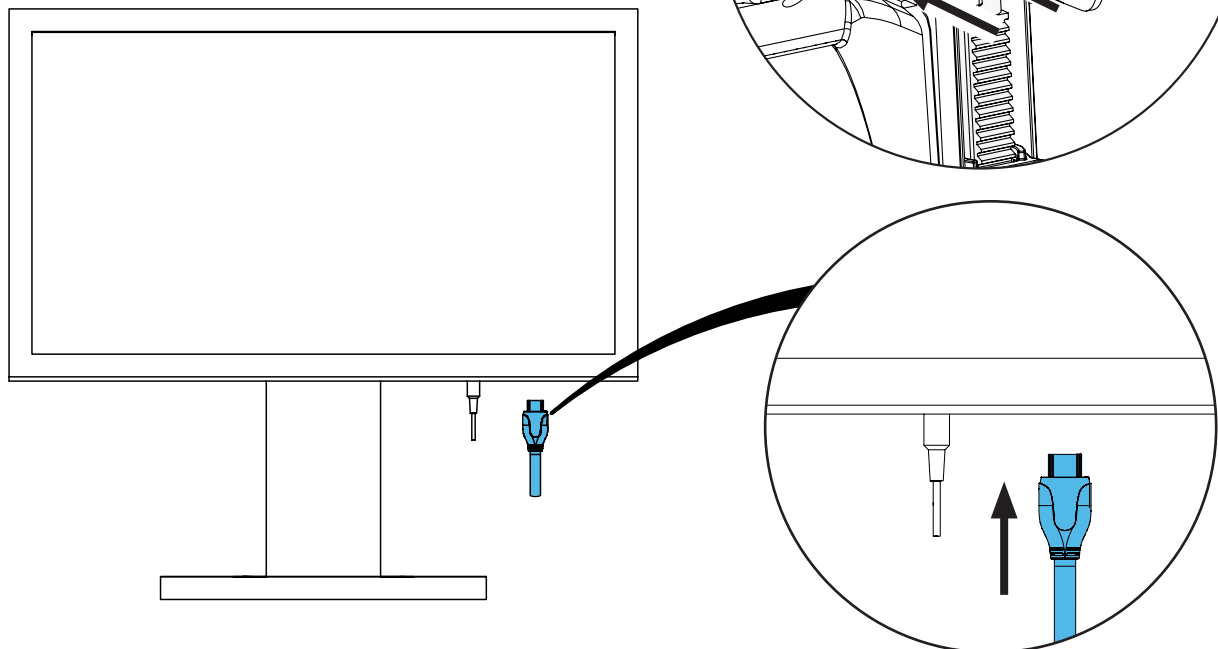





making work easy

Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по монтажу

90003 3775

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HDMI Video Capture • Карта видеозахвата HDMI Video Capture

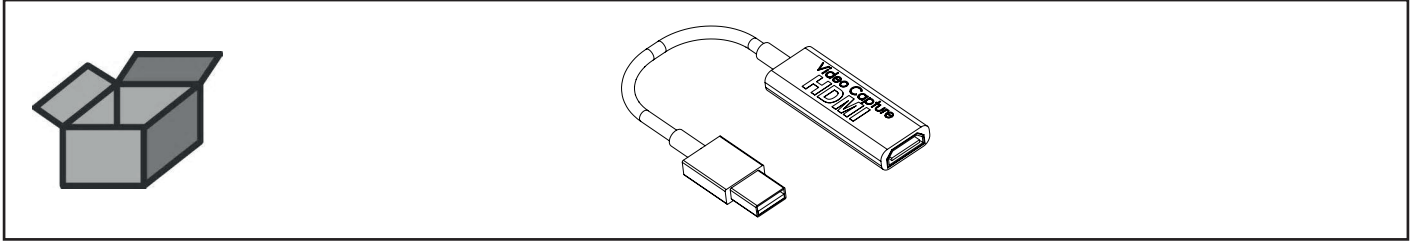







Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по монтажу


making work easy

90003 7778


Höheneinstellung inkl. Klemmschrauben • Height adjustment incl. clamping screw
Réglage de hauteur avec vis de serrage • Regolazione altezza inc. viti di bloccaggio
Ajuste en altura incl. tornillo de fijación • Регулятор высоты вкл. зажимные винты

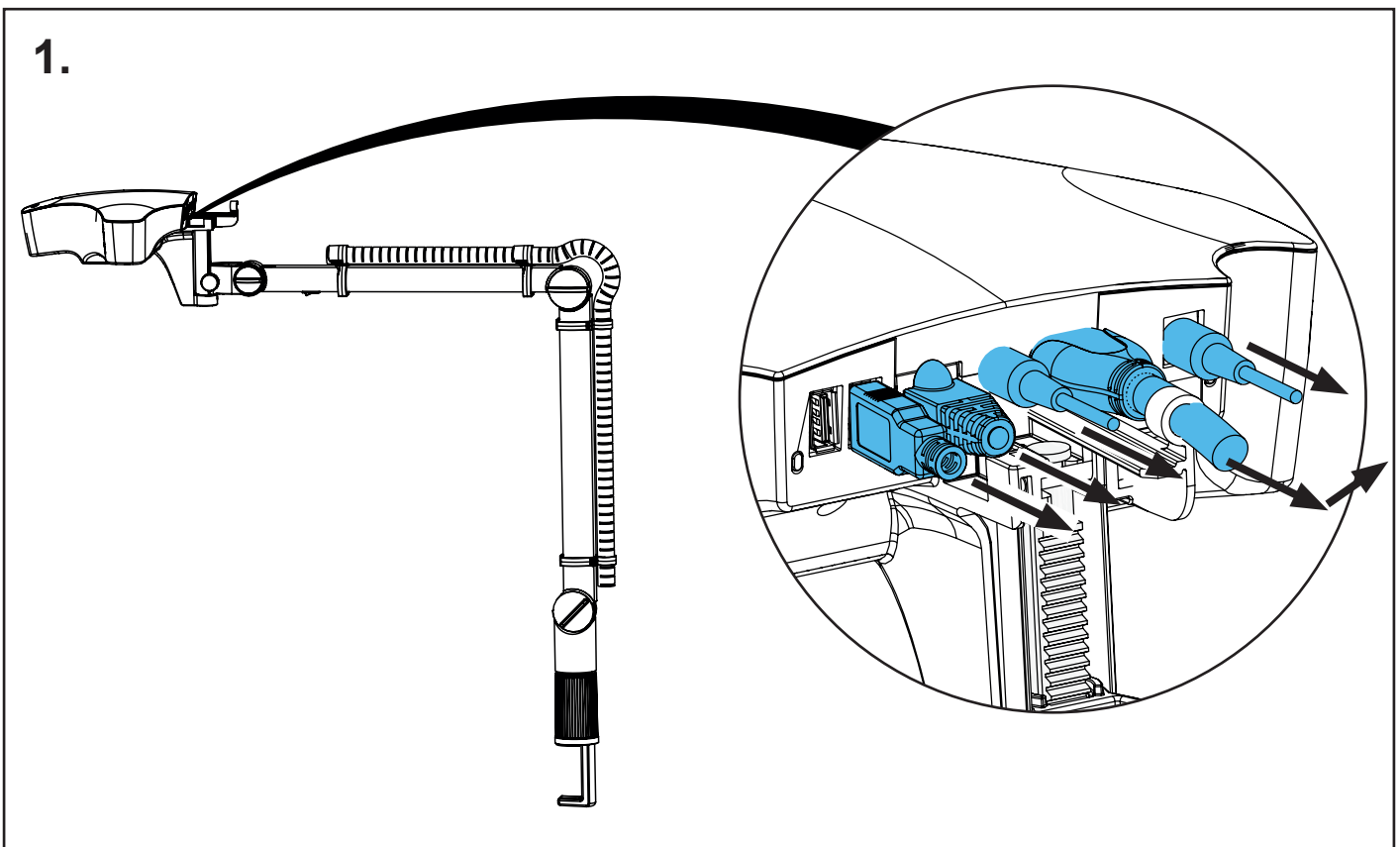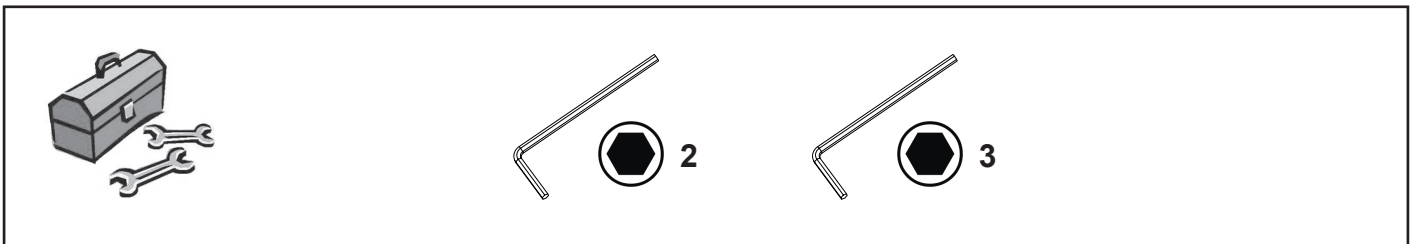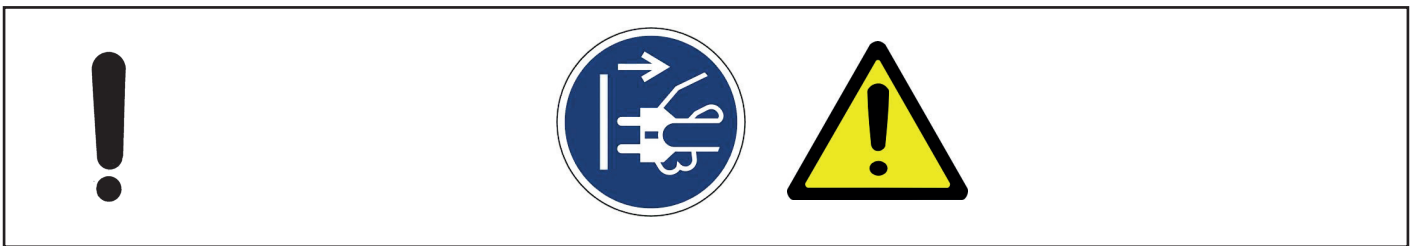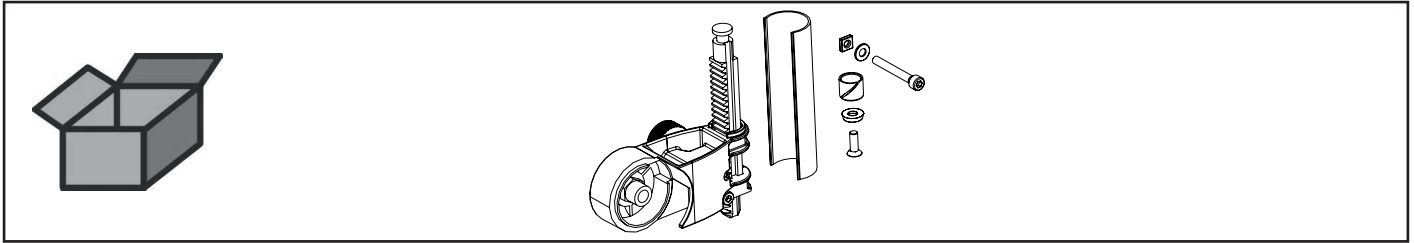


2.

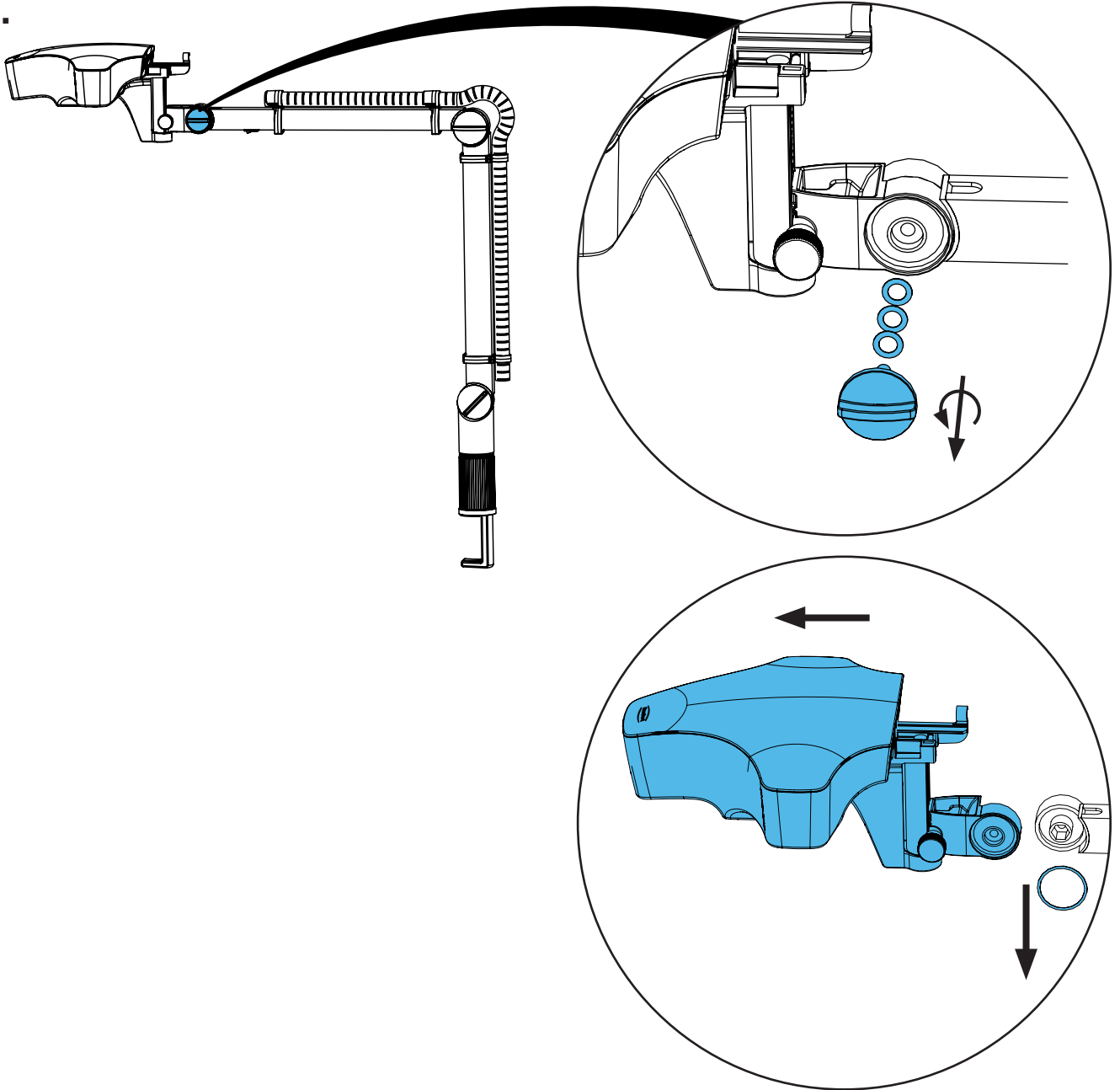


3.

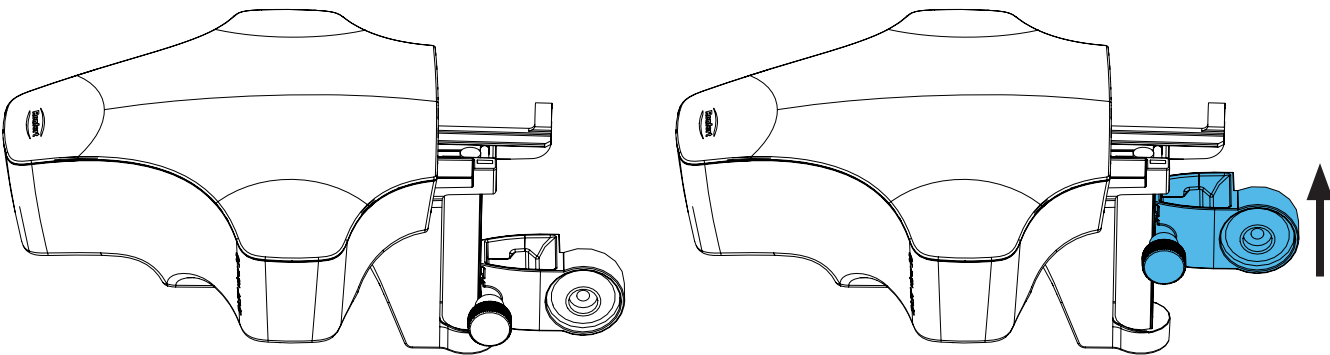


4.

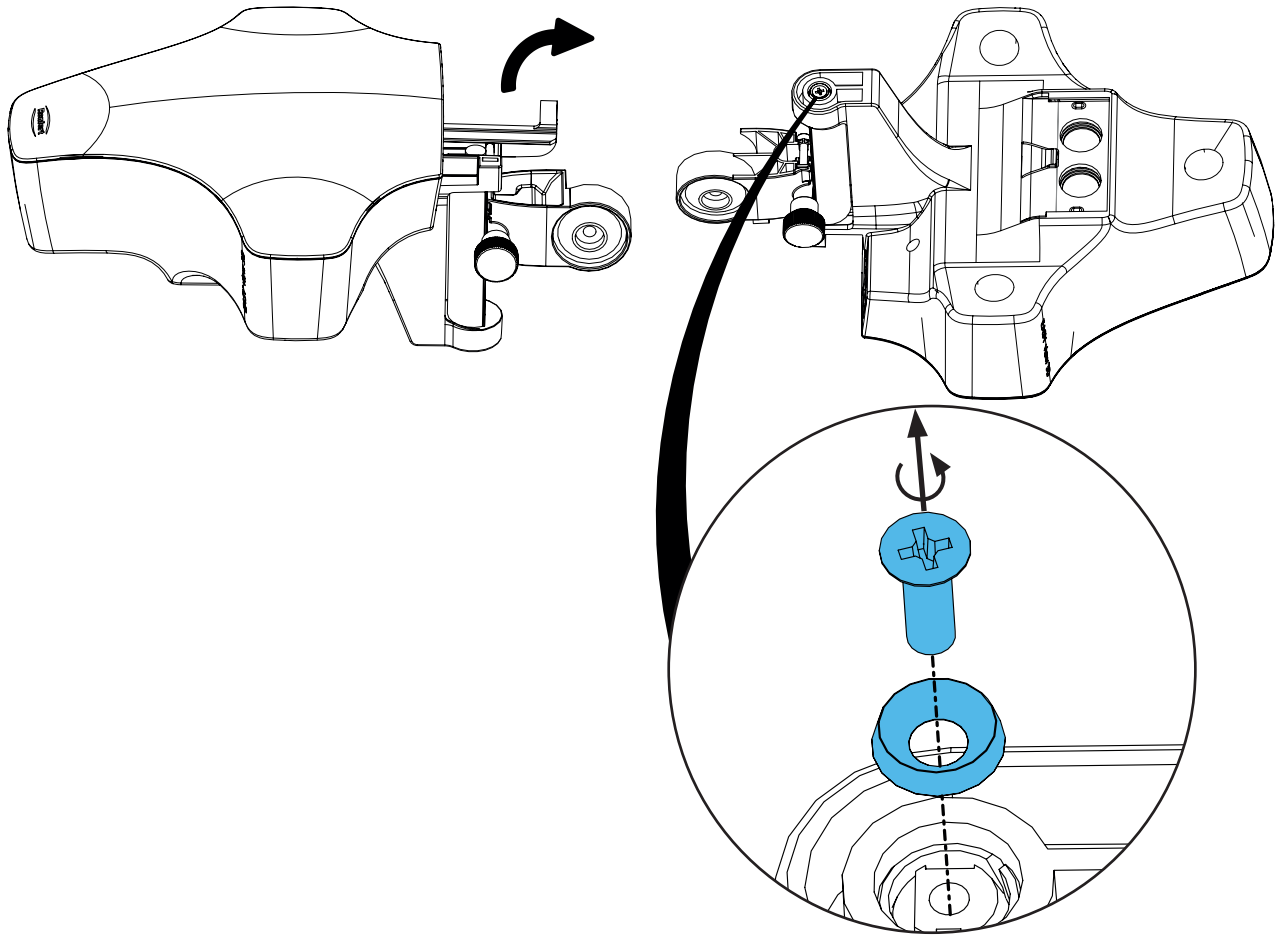


5.

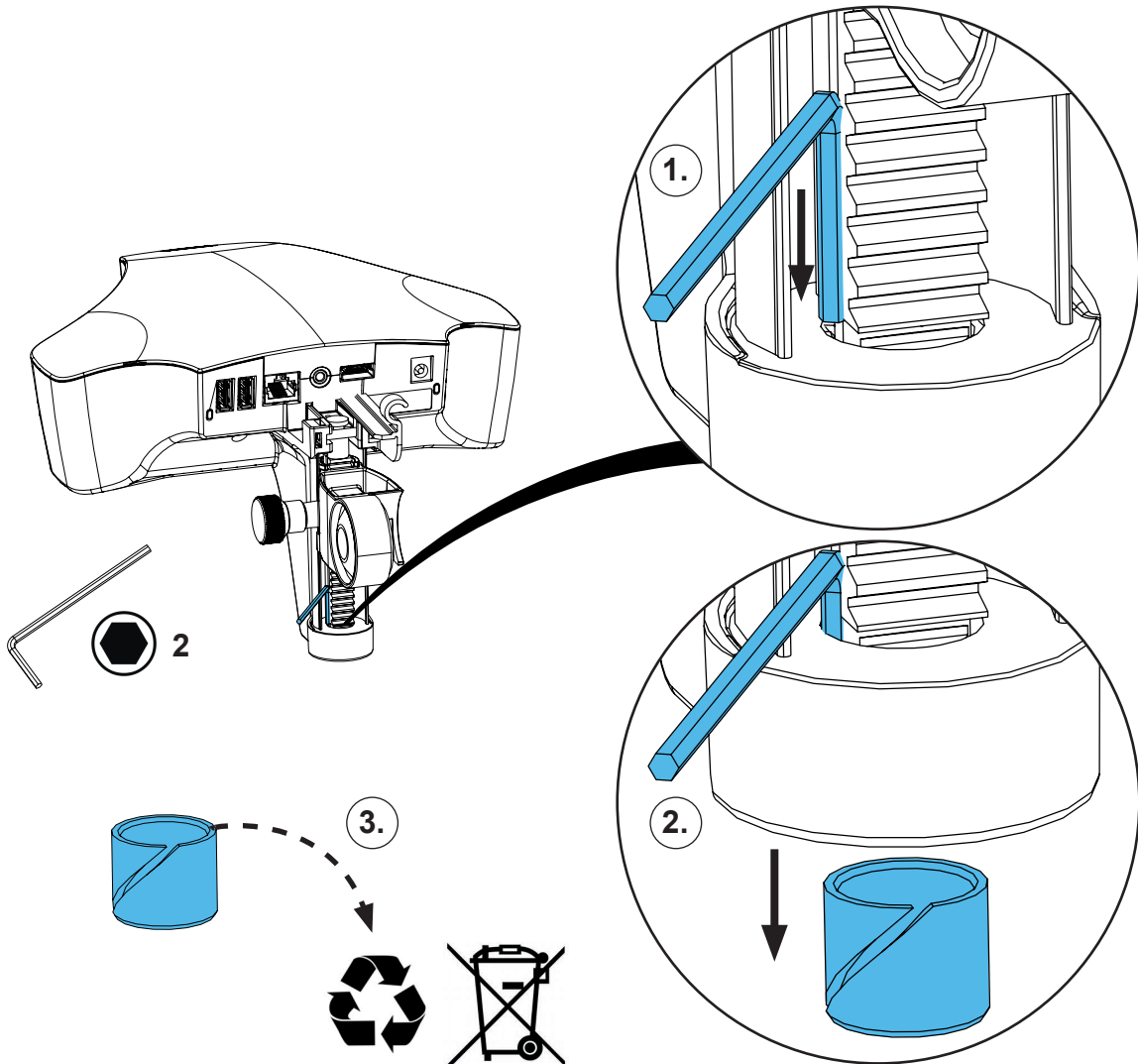


6.

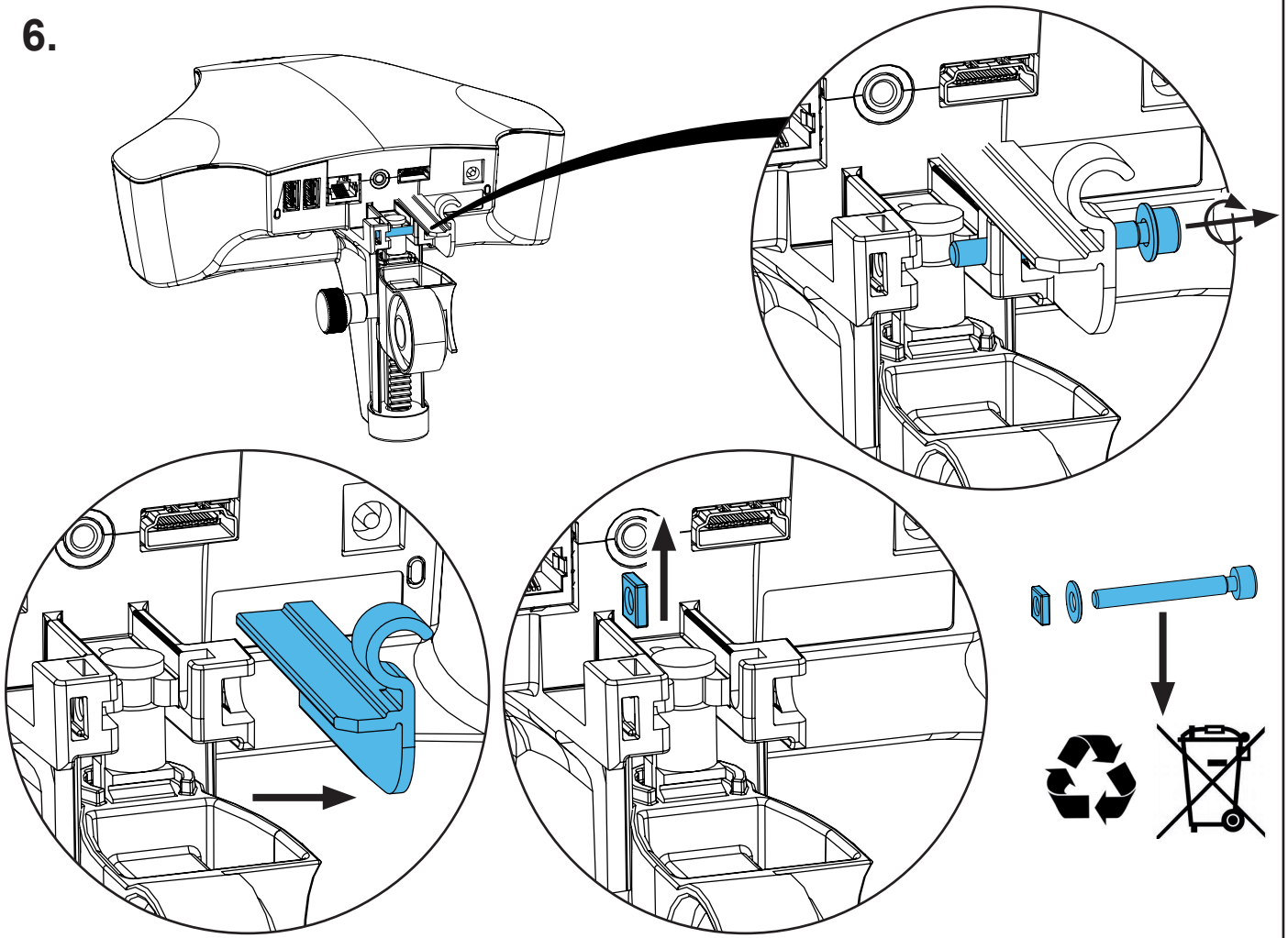


7.

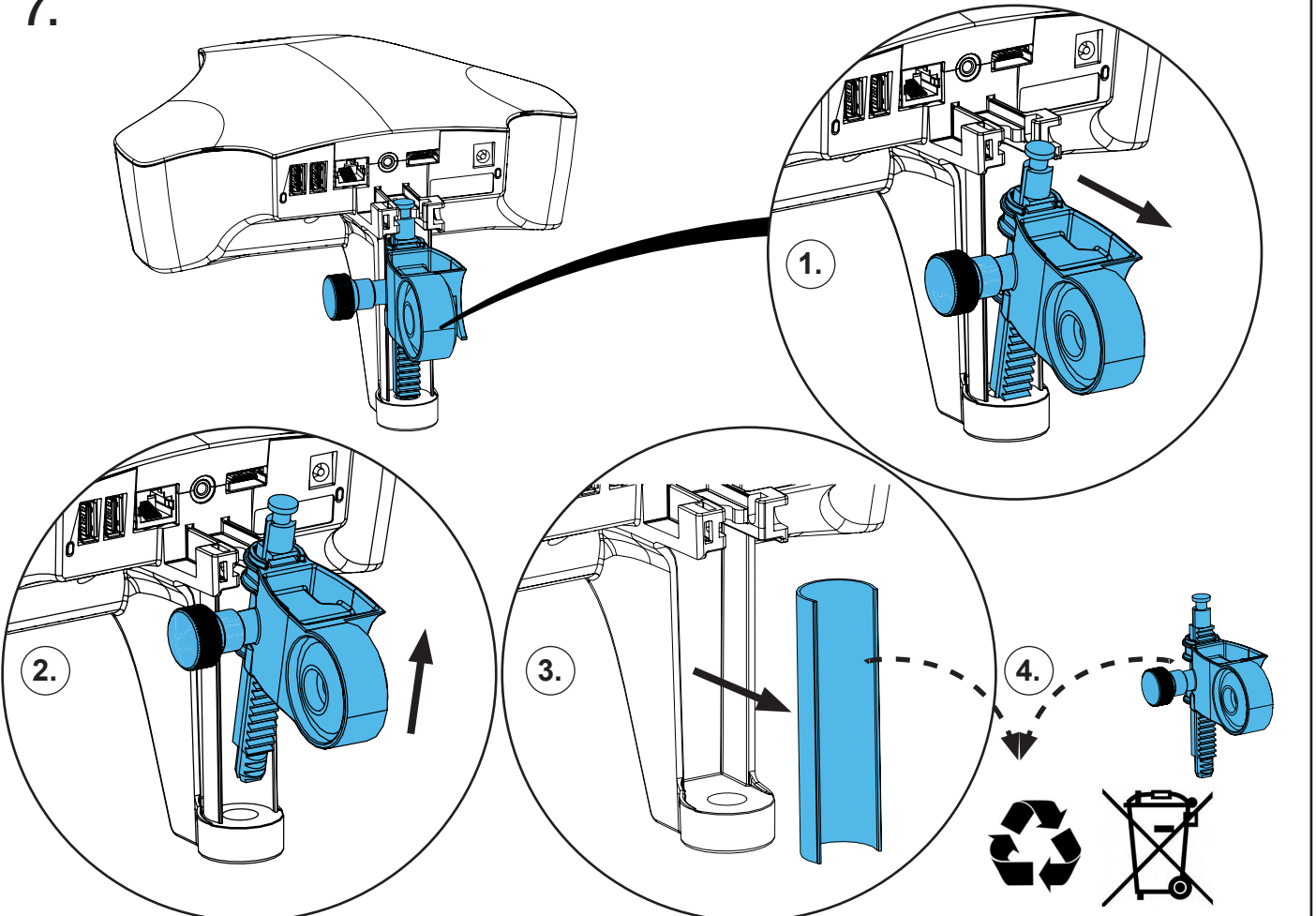


8.

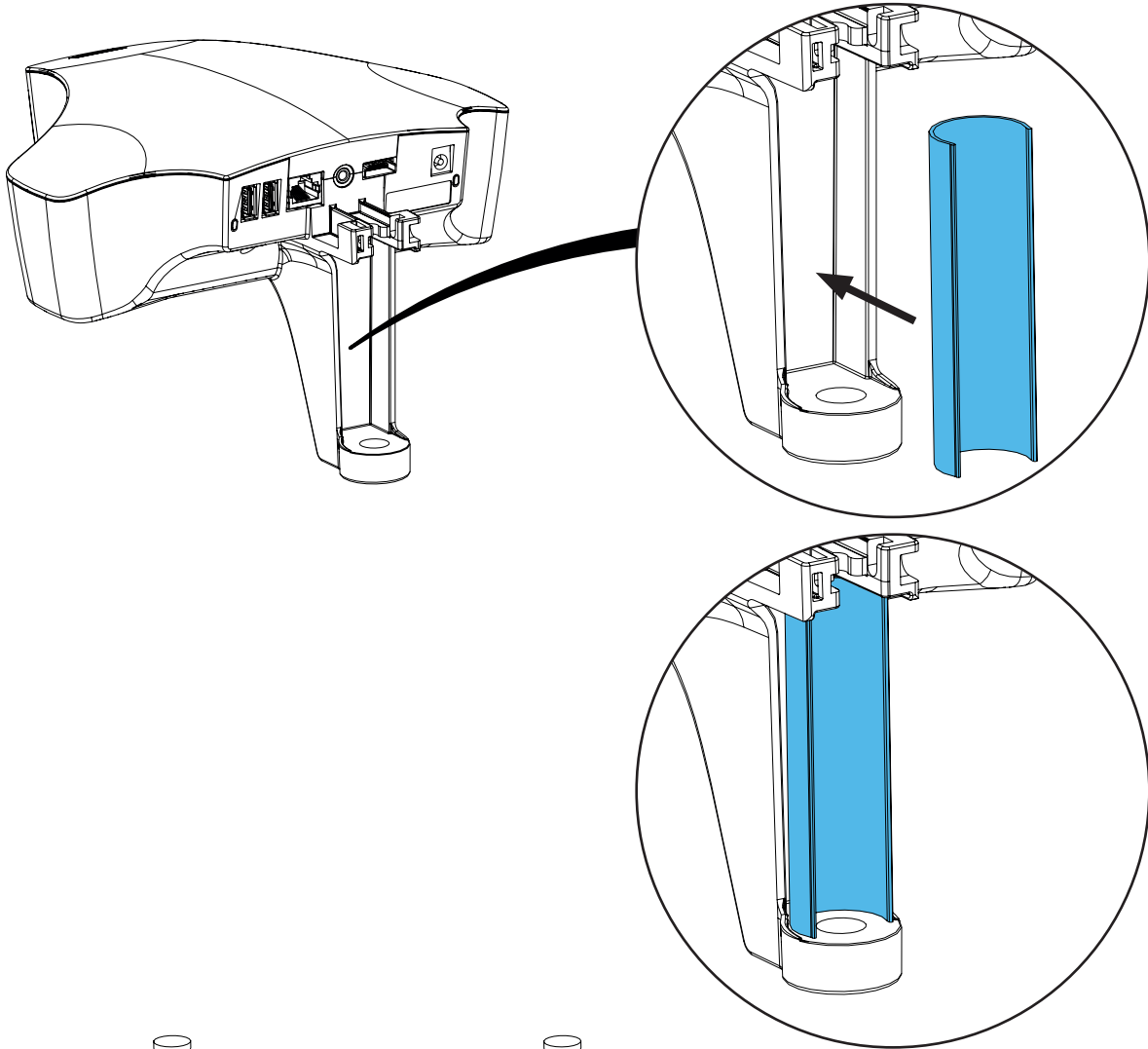



9.

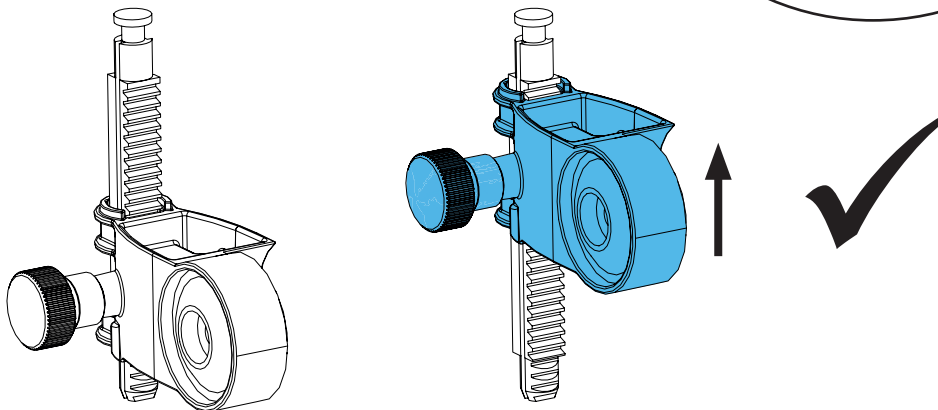



10.

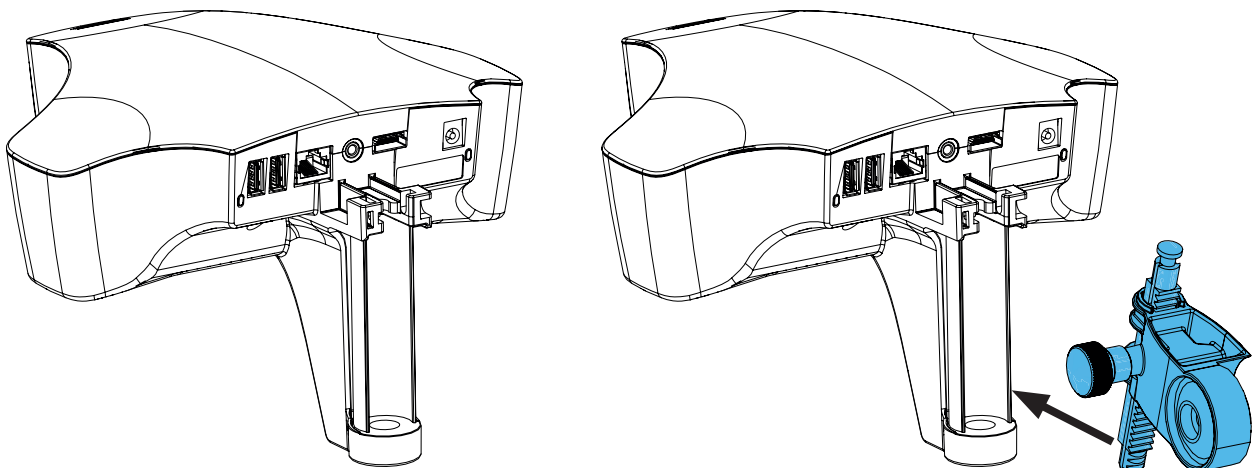





11.

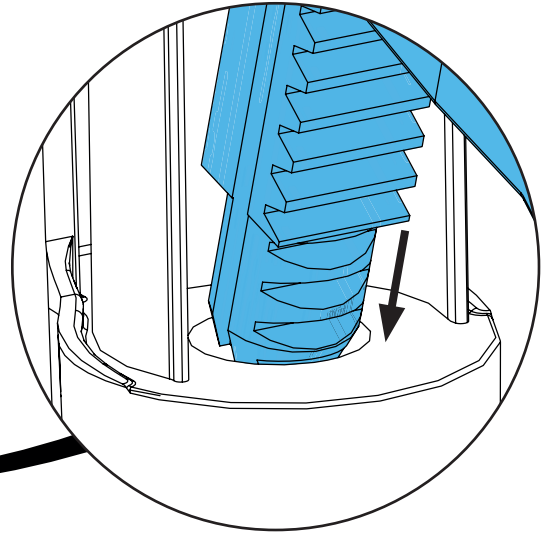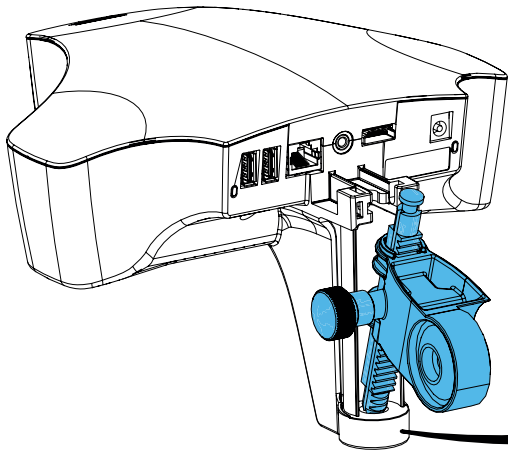





12.

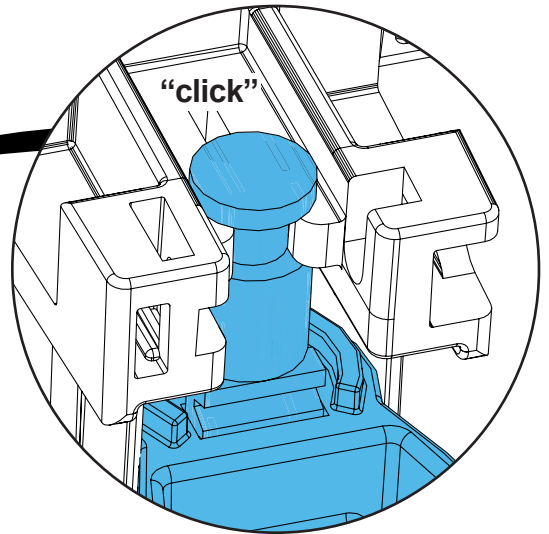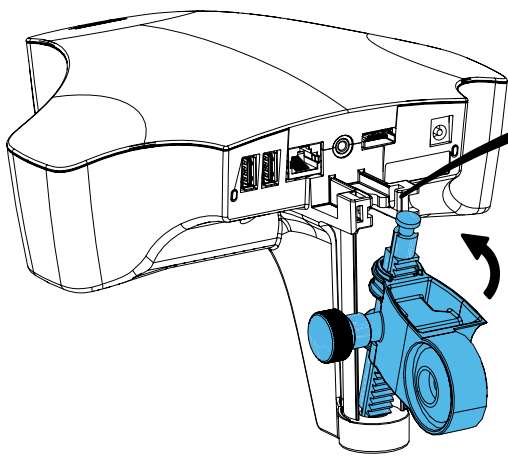

13.

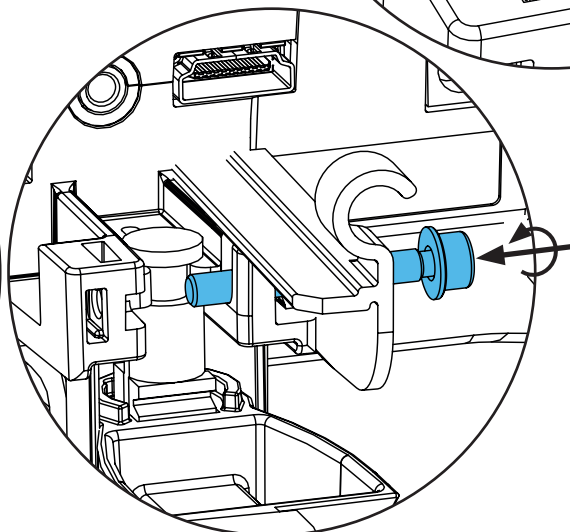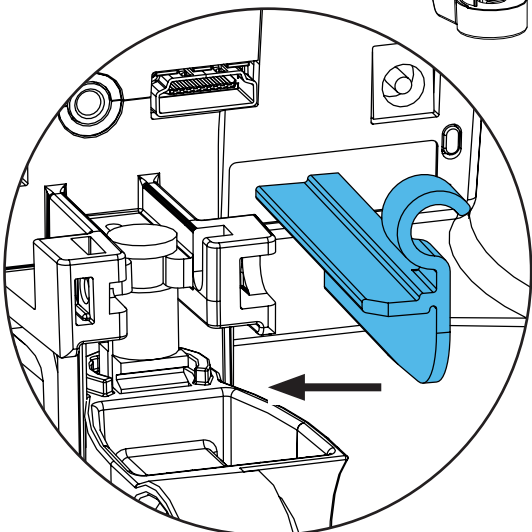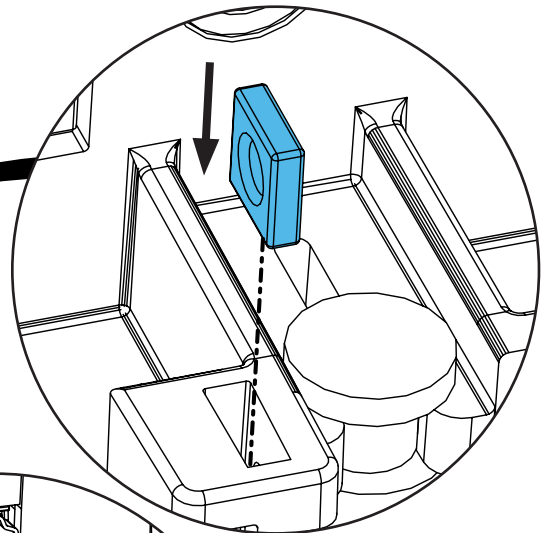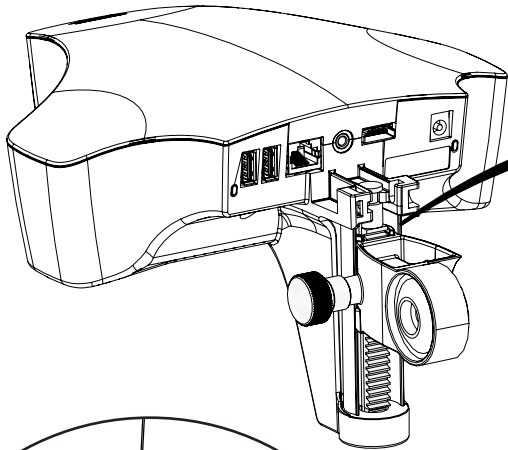


14.

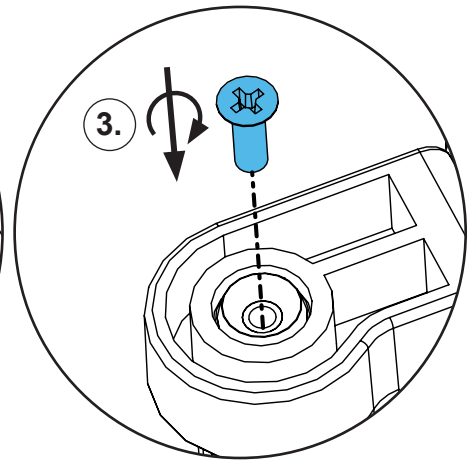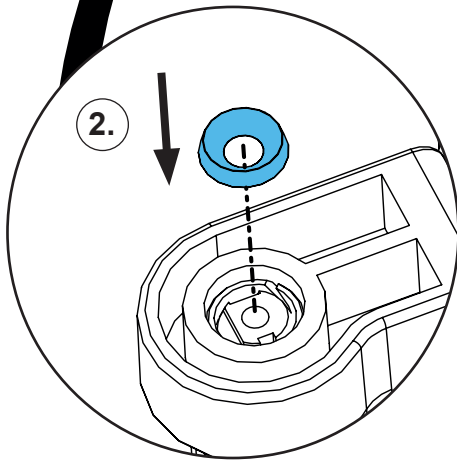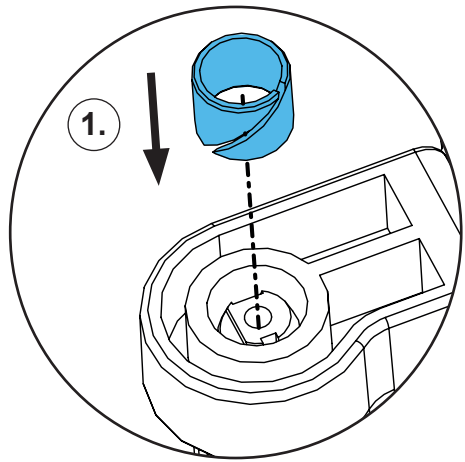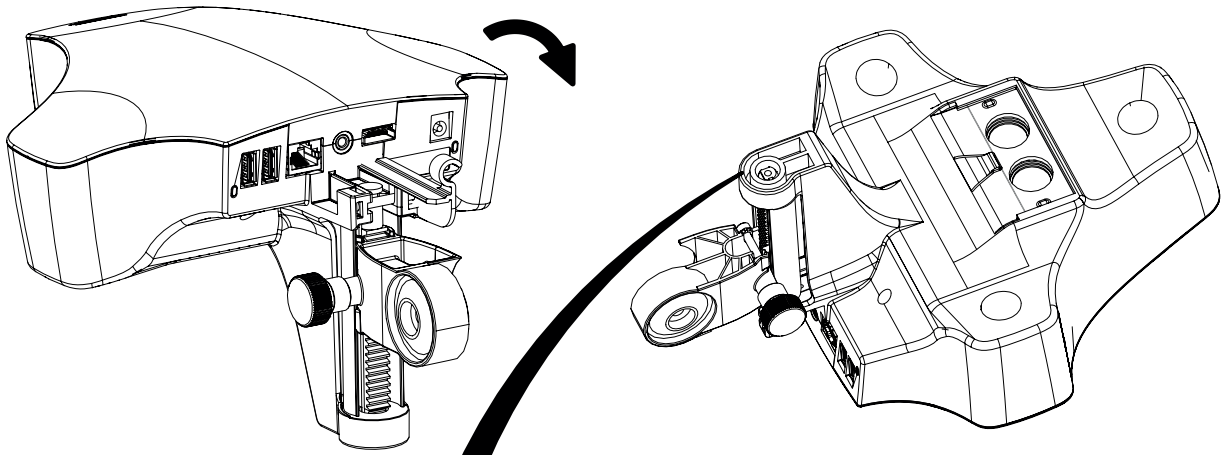


1,1 - 1,2 Nm


15.

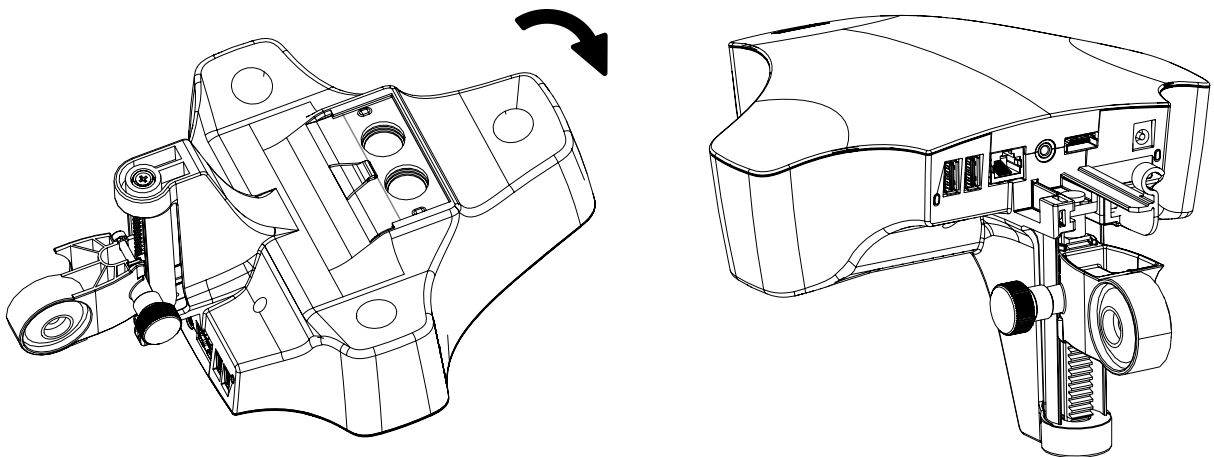

16.

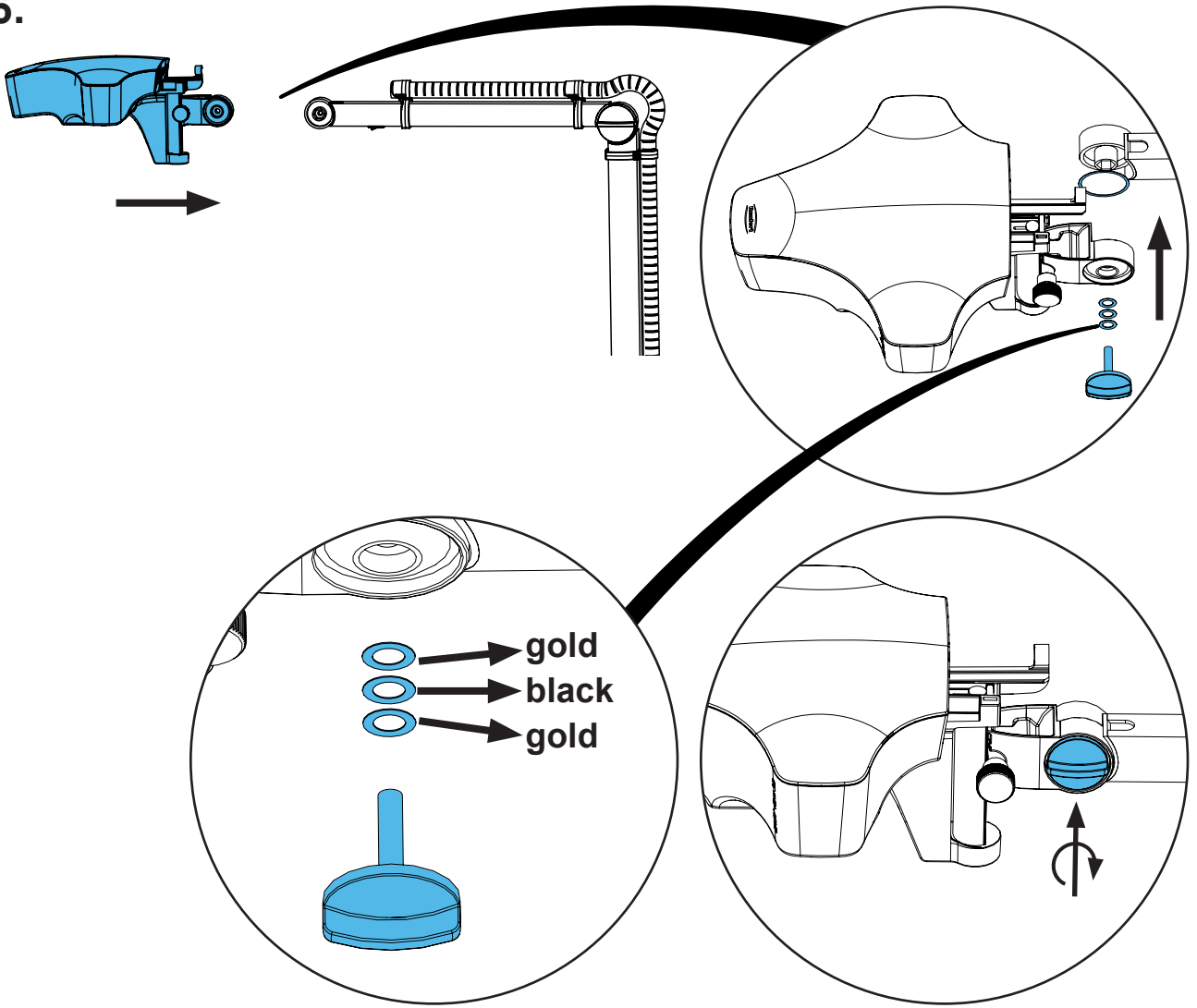

17.

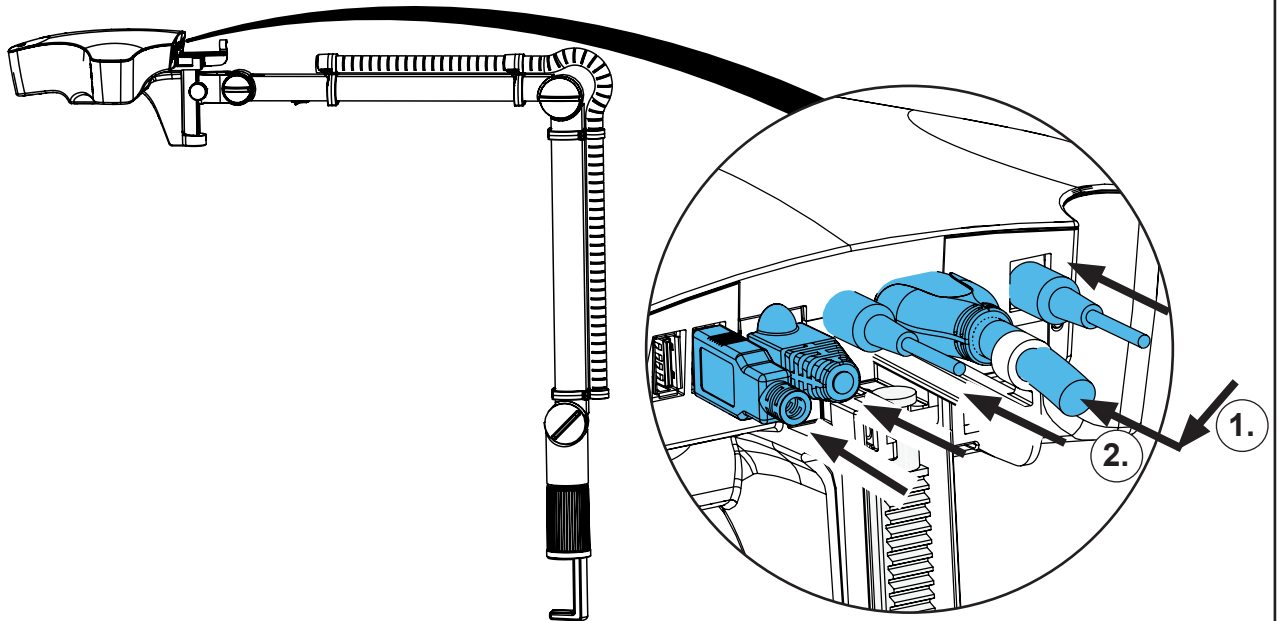






Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу


making work easy

90003 3759


LED-Modul • LED module • Module DEL • Modulo LED • Módulo LED •
Модуль светодиодов

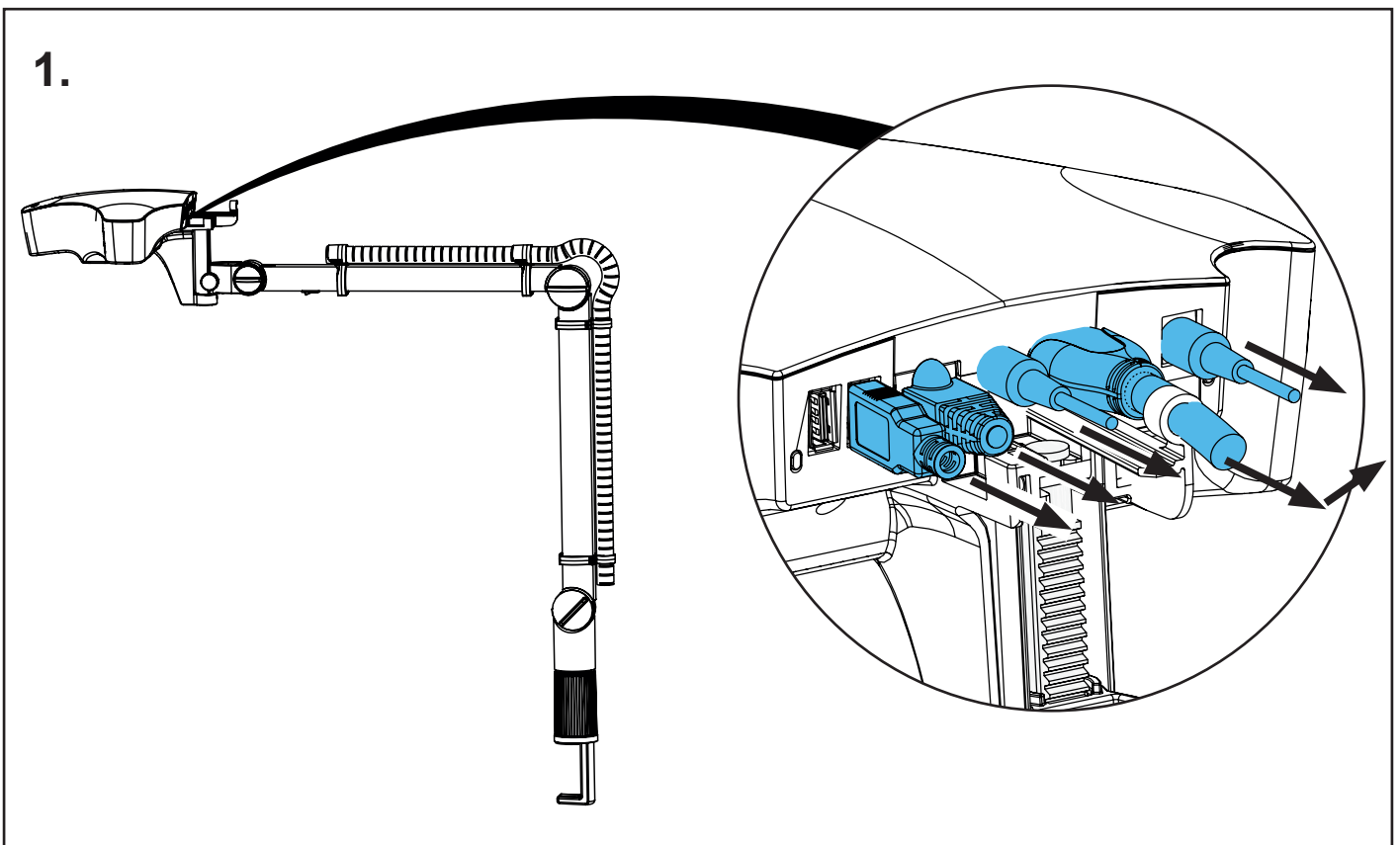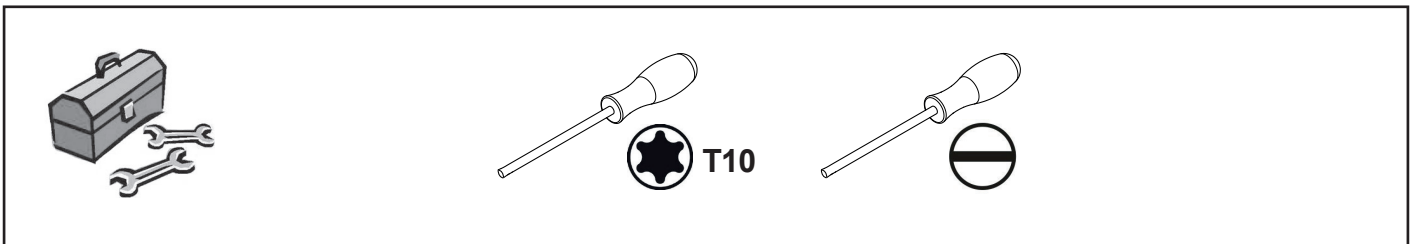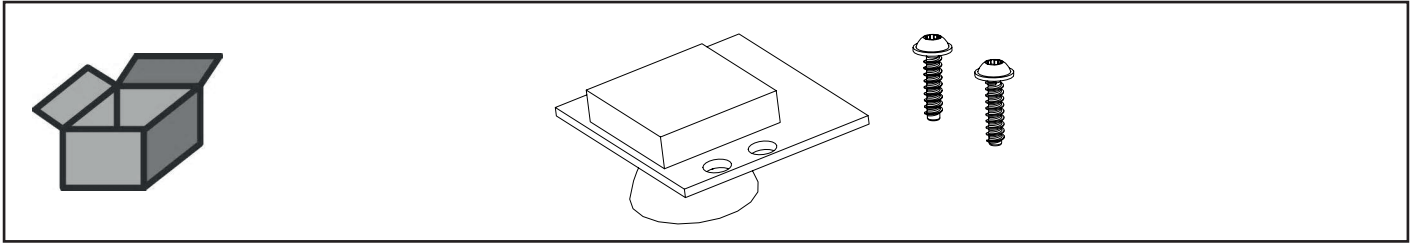


2.

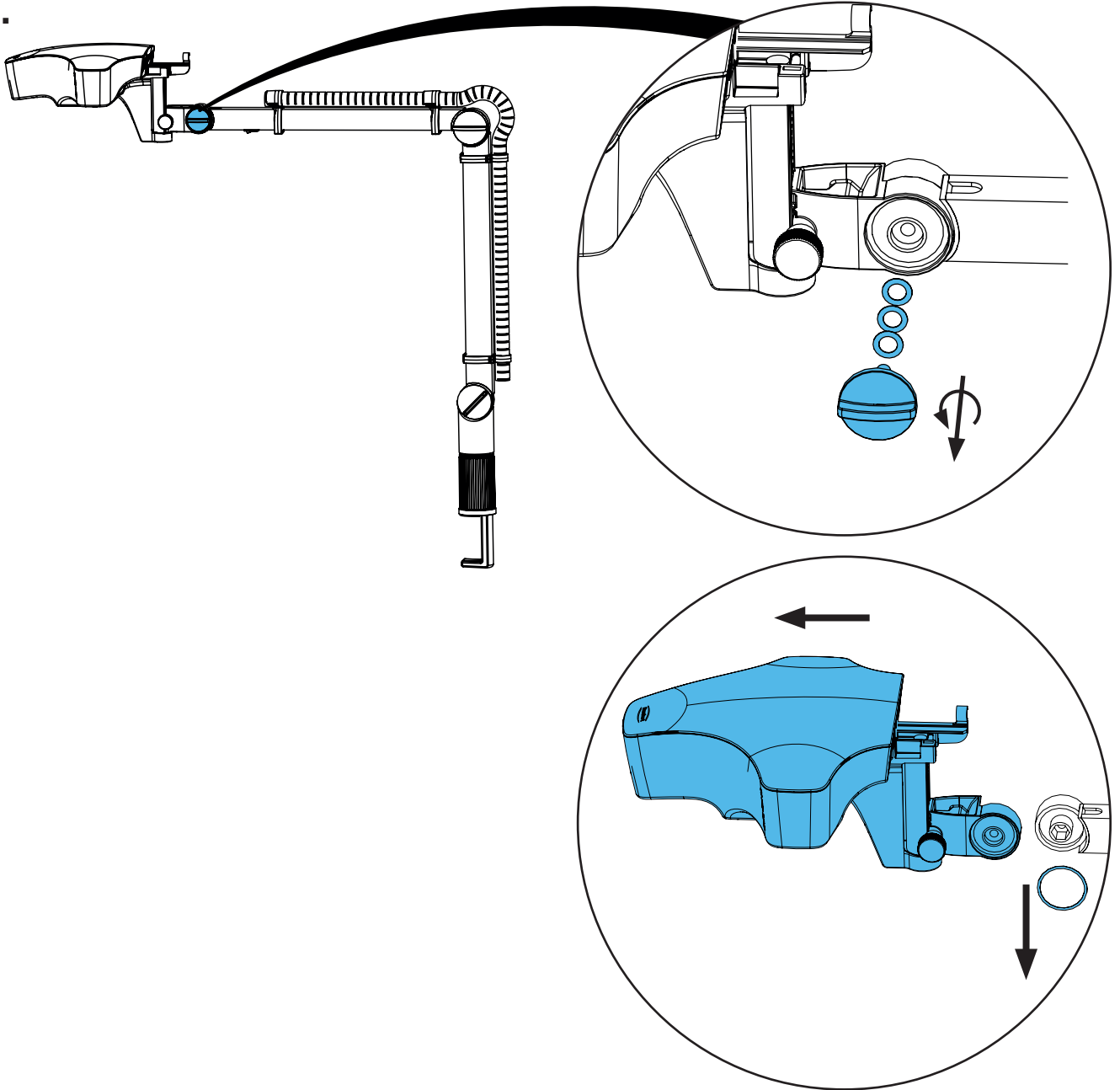


3.

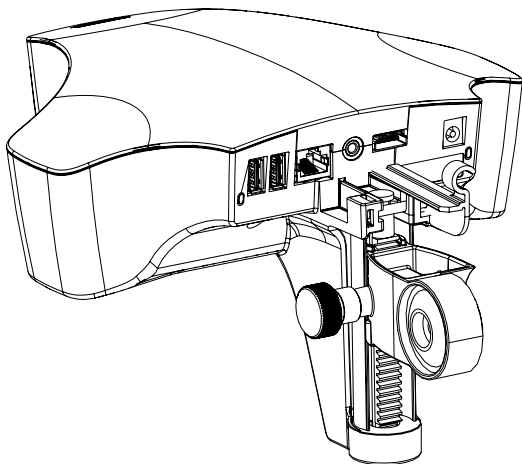





4.

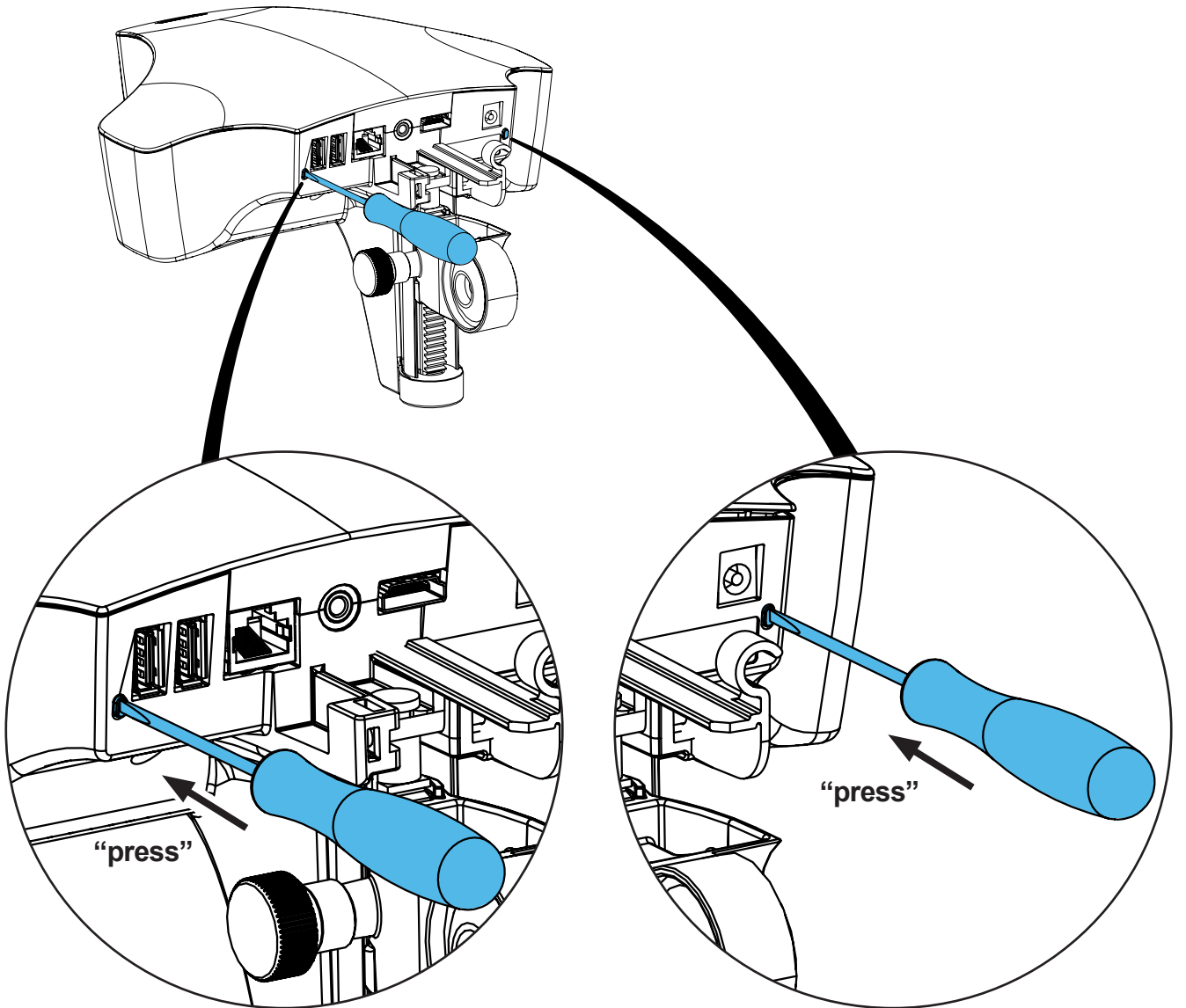


5.

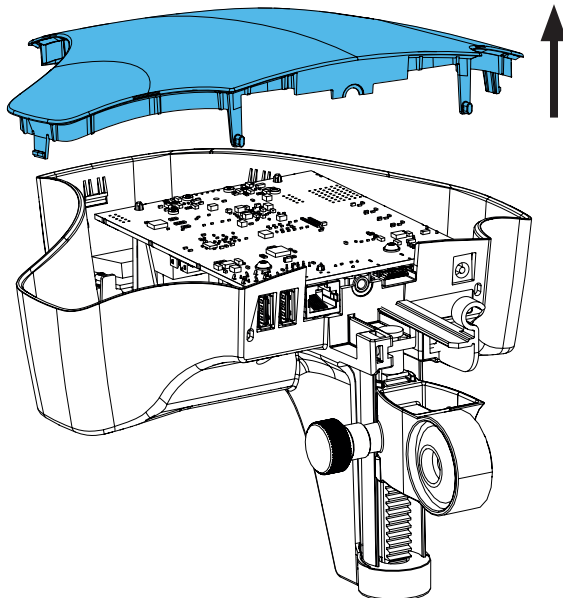




6.

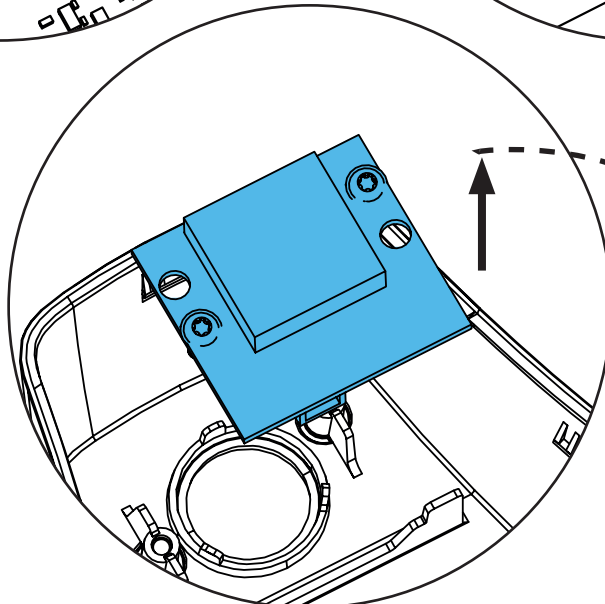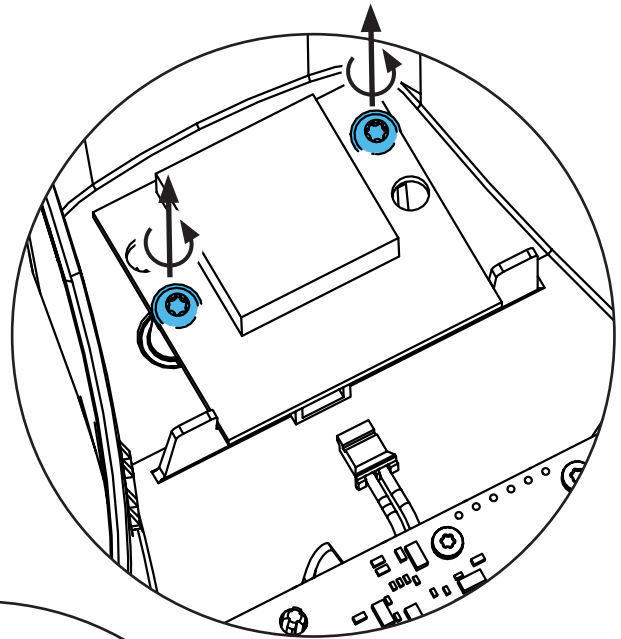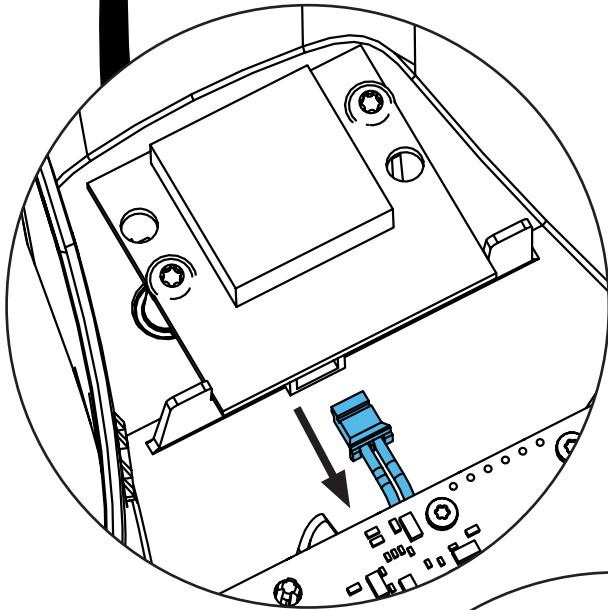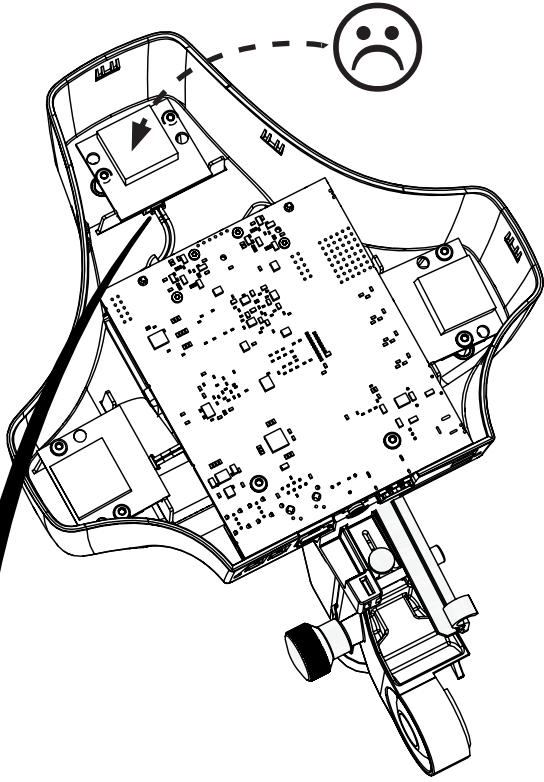

7.

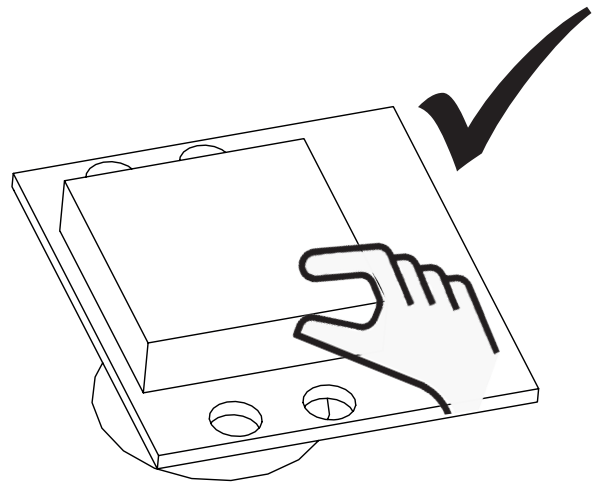



8.

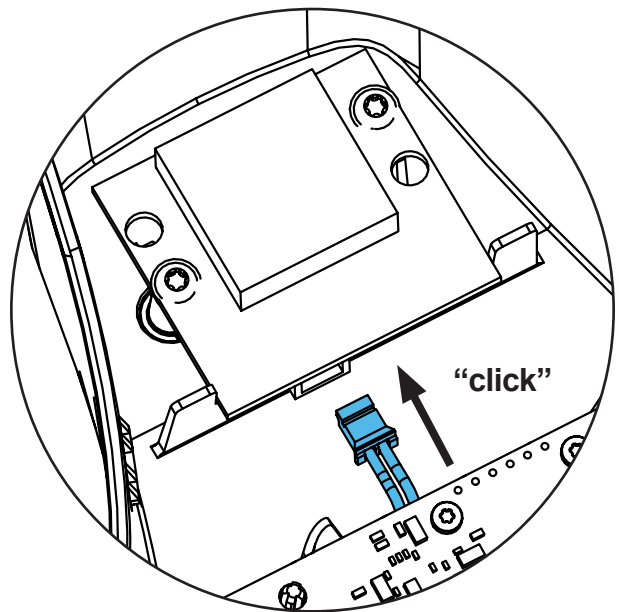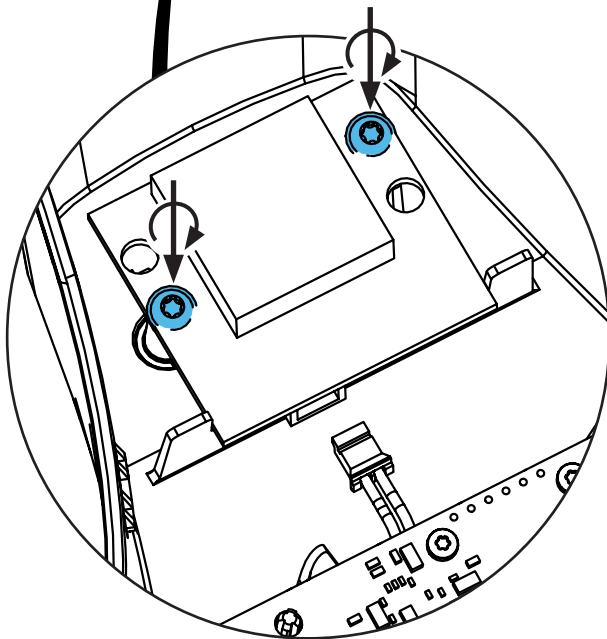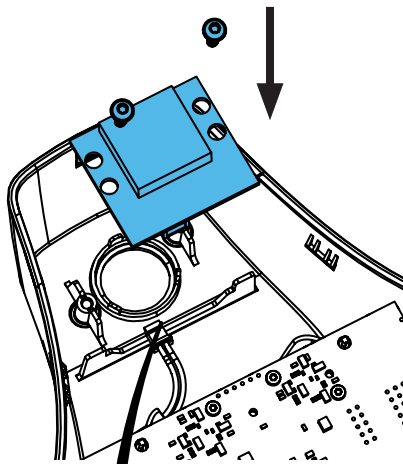


0,8 - 0,9 Nm




9.

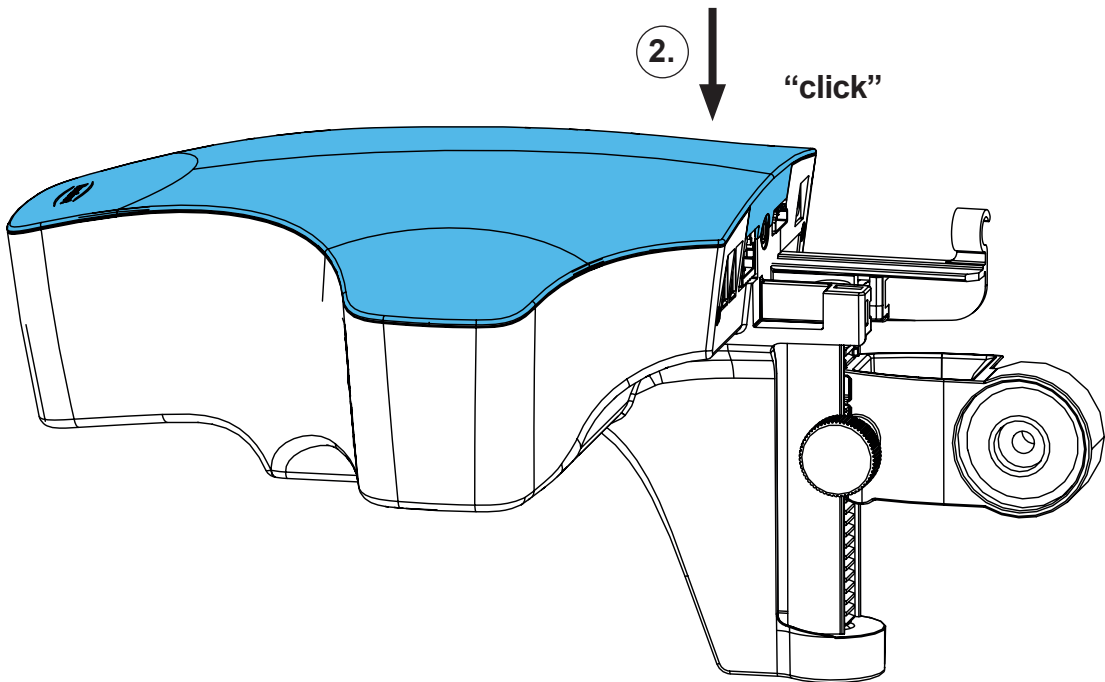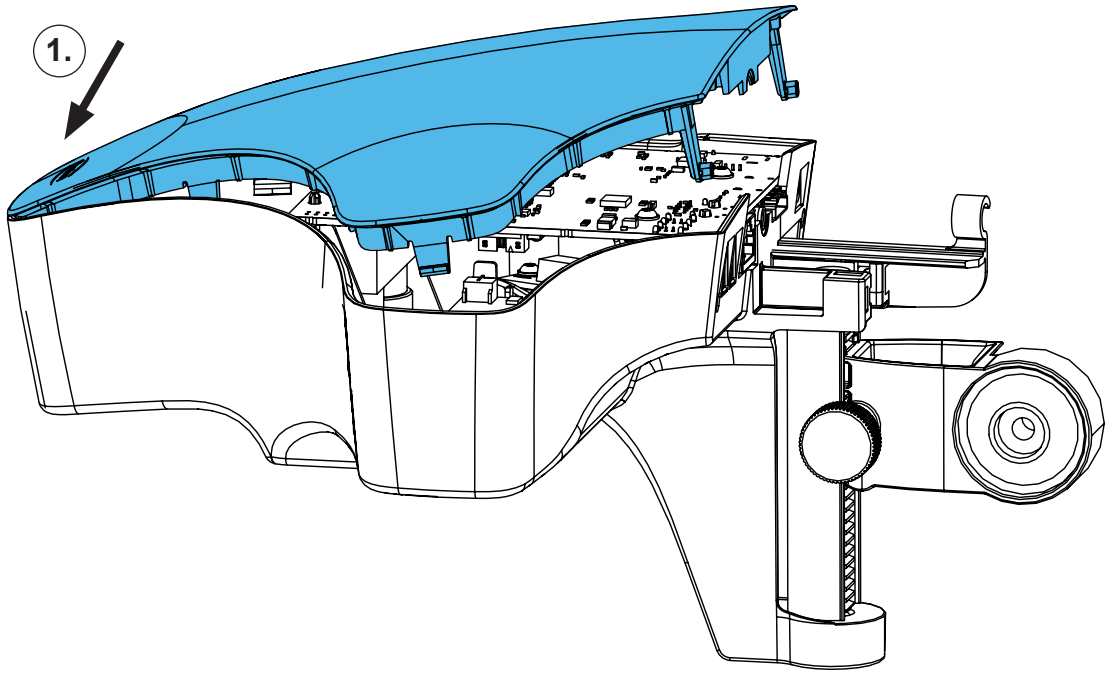

10.

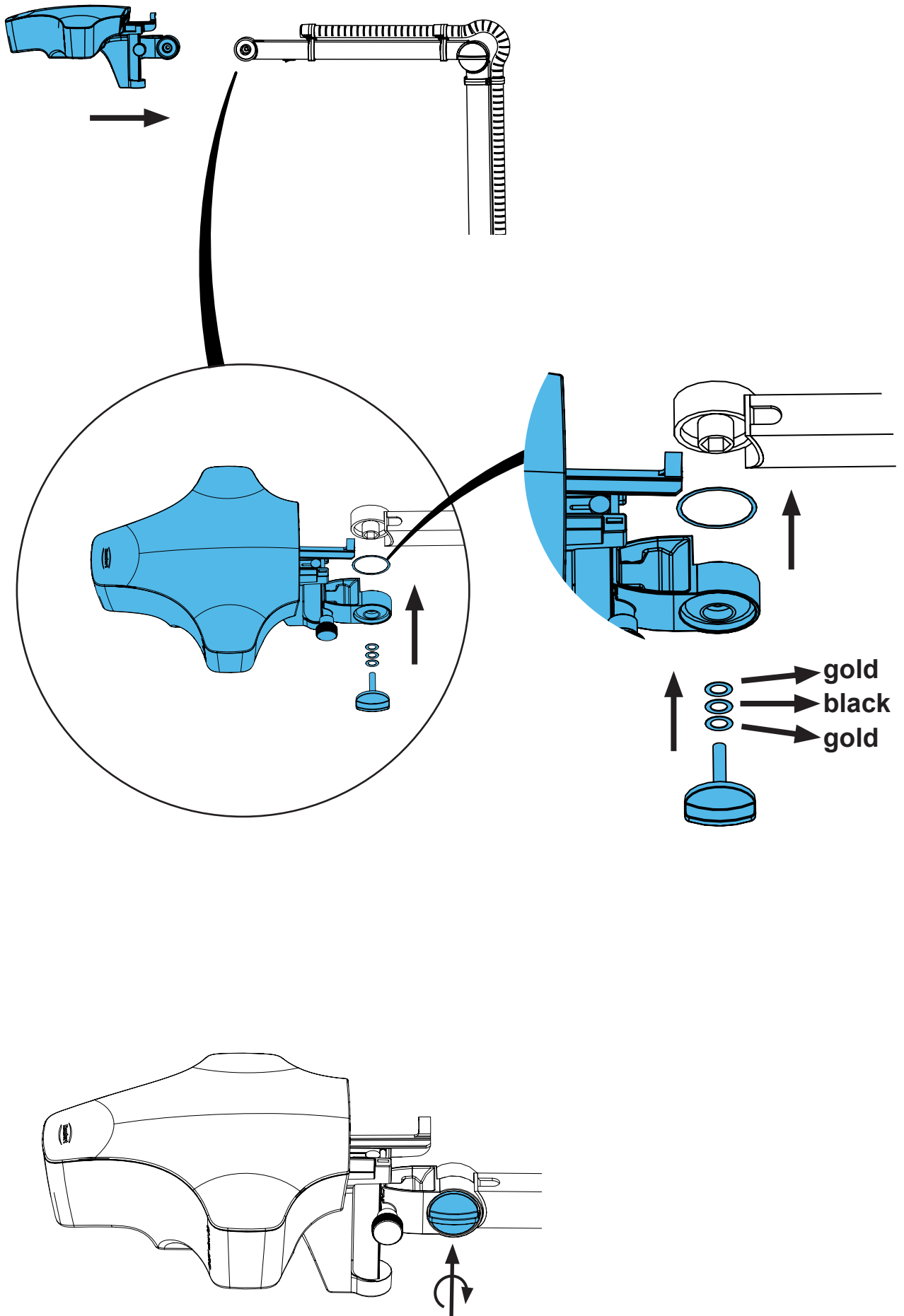

11.

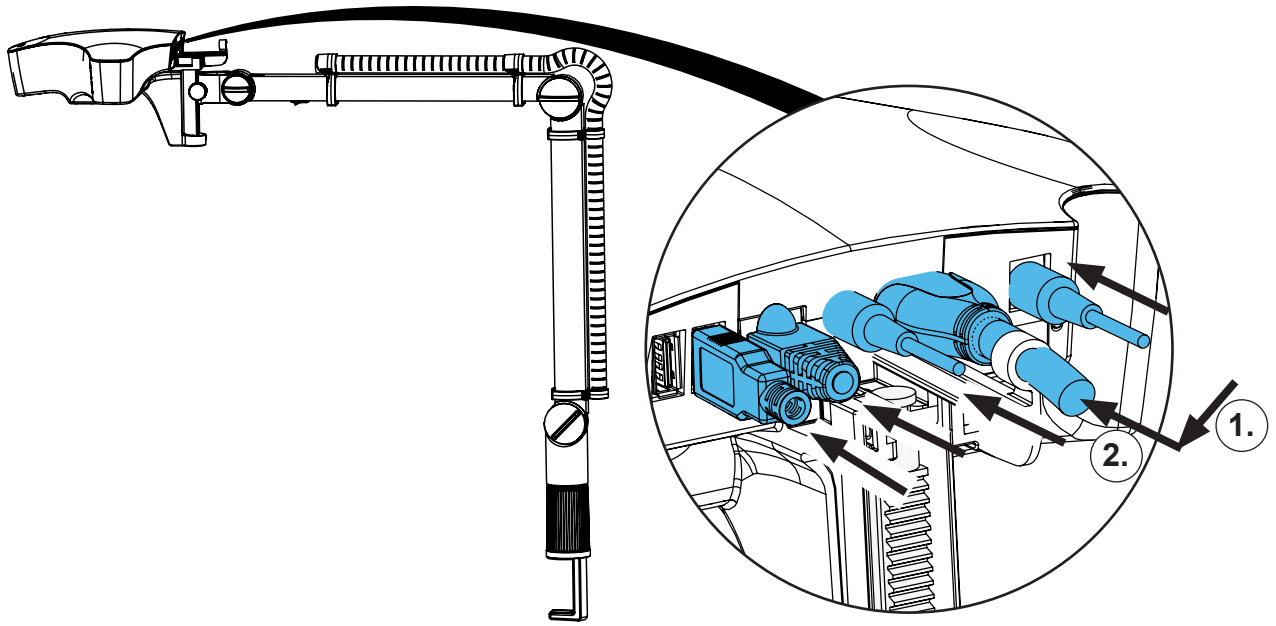






Montageanleitung • Assembly Instruction • Instruction d'installation
Istruzioni l'istallazione • Instrucciones de montaje • Руководство по
монтажу


making work easy

90003 3760


LED-Treiberleiterkarte • LED driver card • Circuit imprimé pilote DEL • Circuito
stampato driver LED • Circuito impreso de controlador LED • Плата драйвера
СВЕТОДИОДОВ

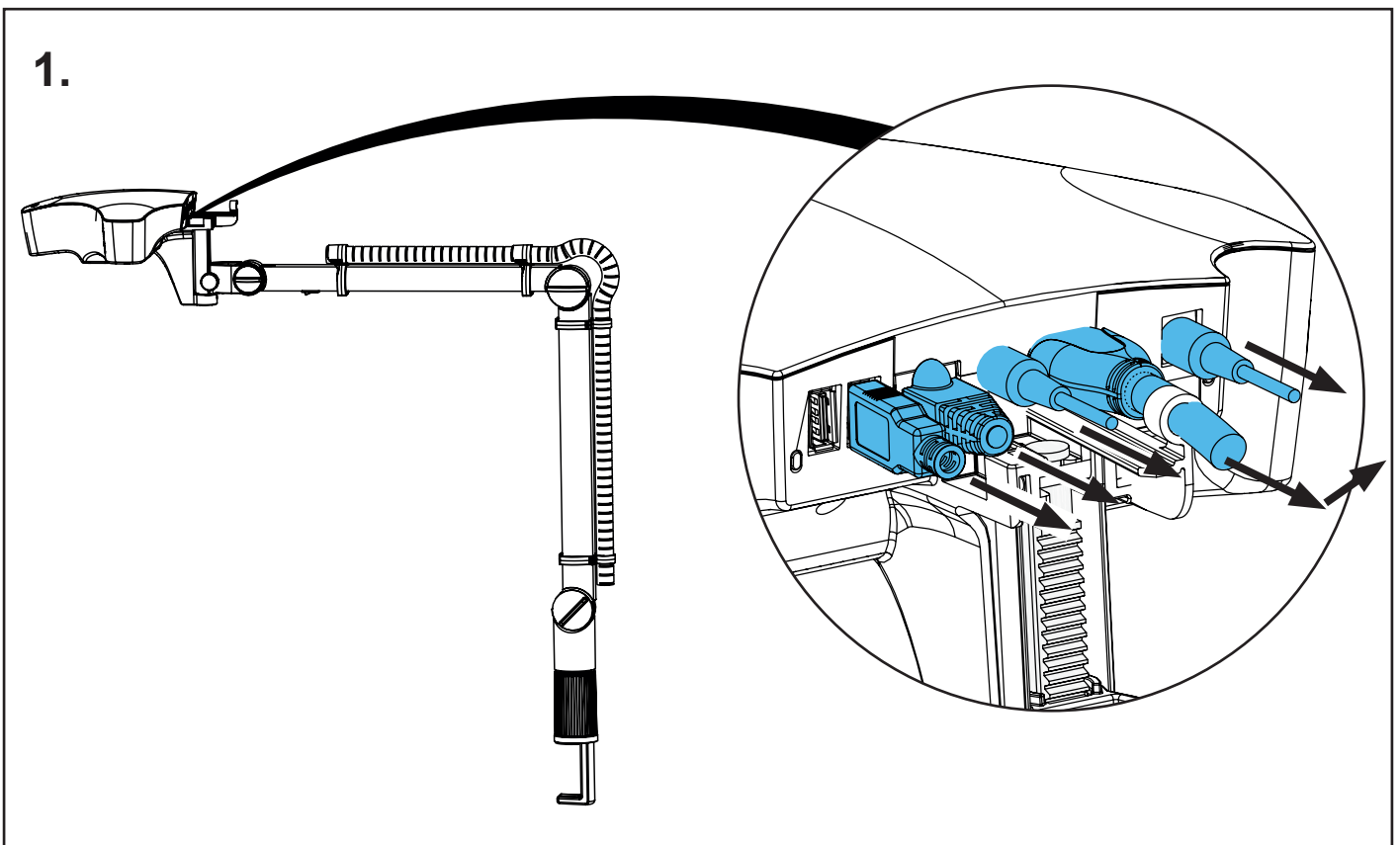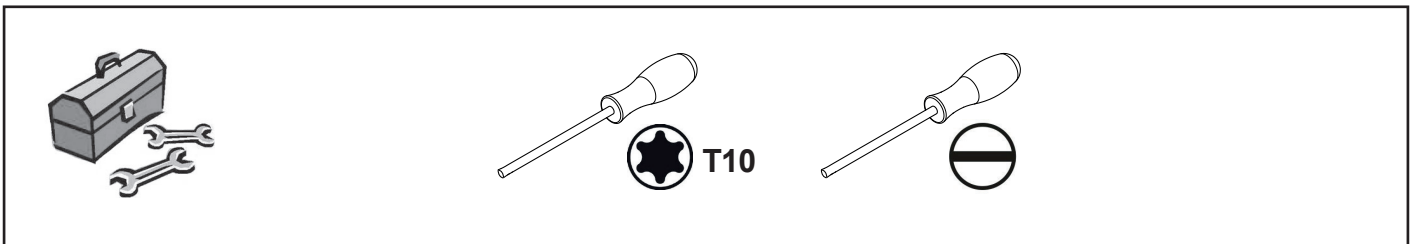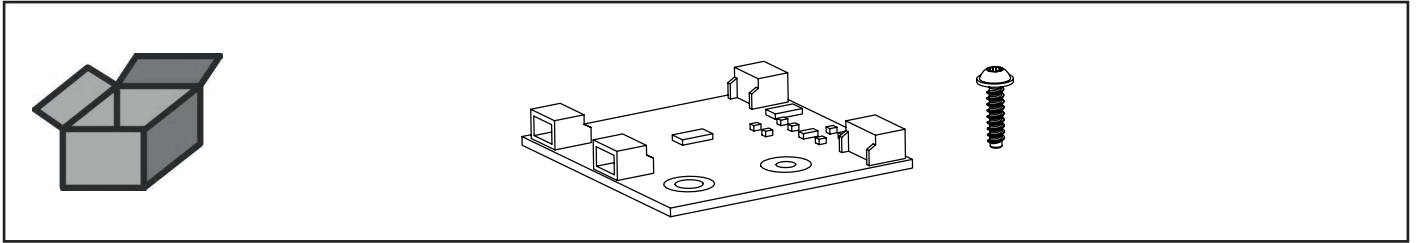


2.

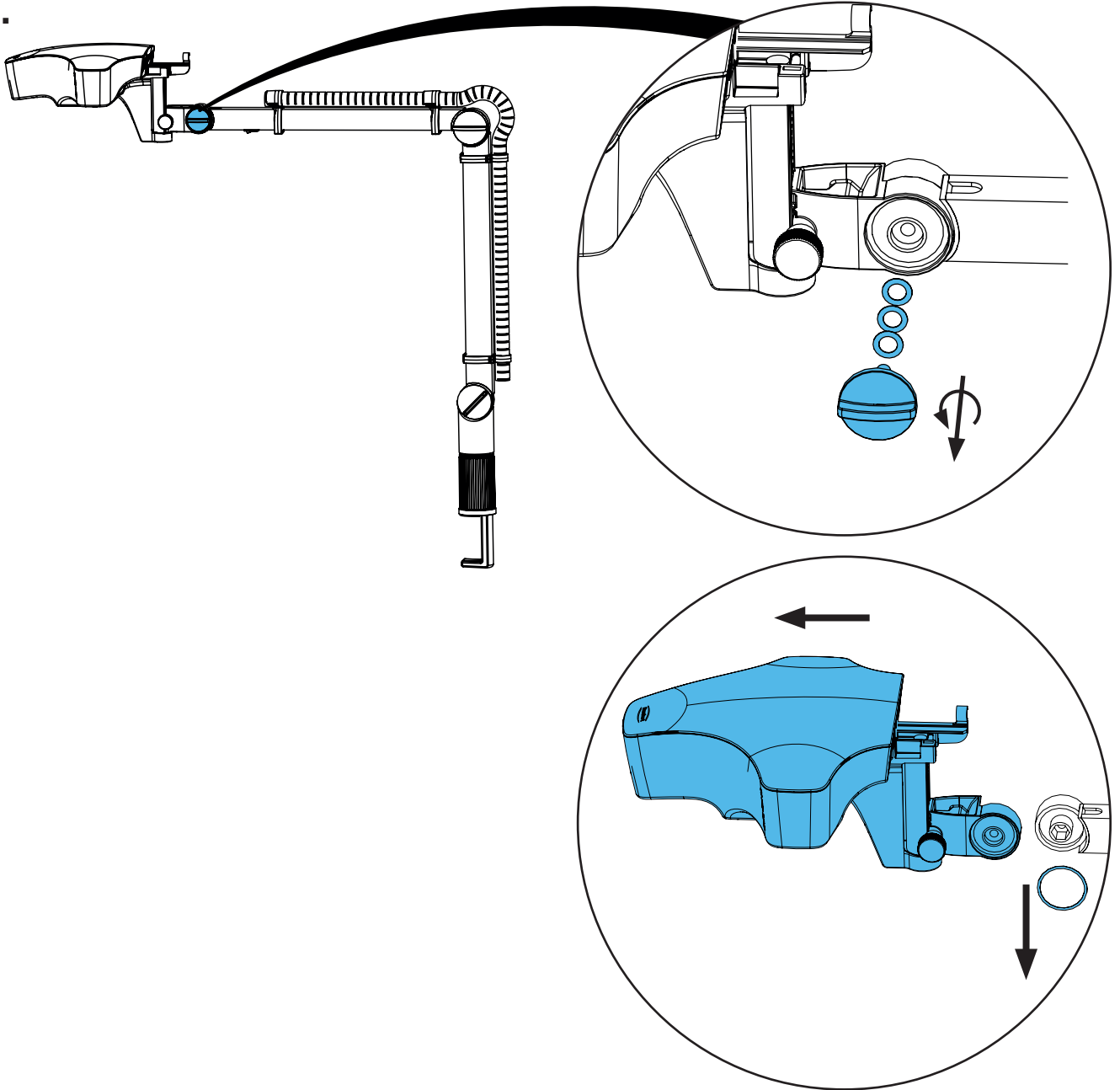


3.

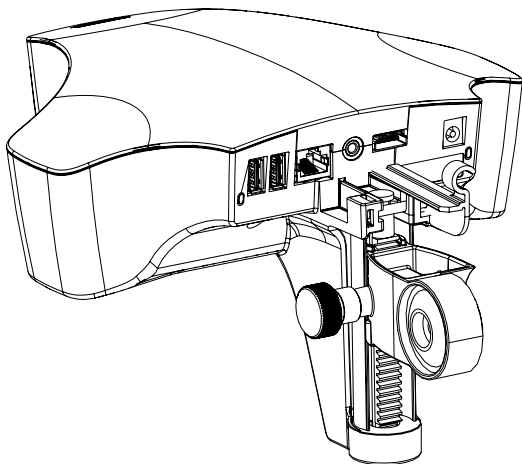


4.

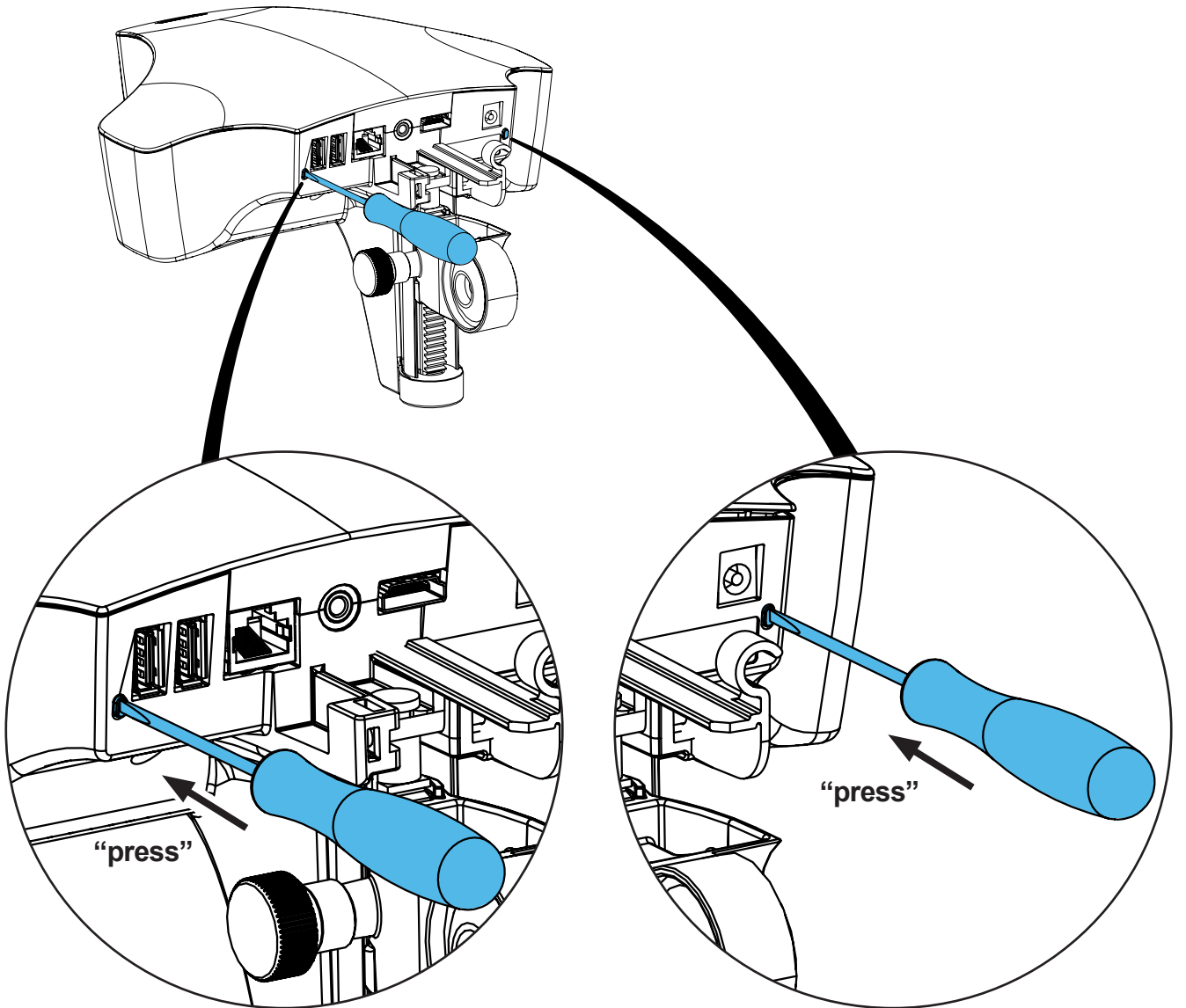


5.

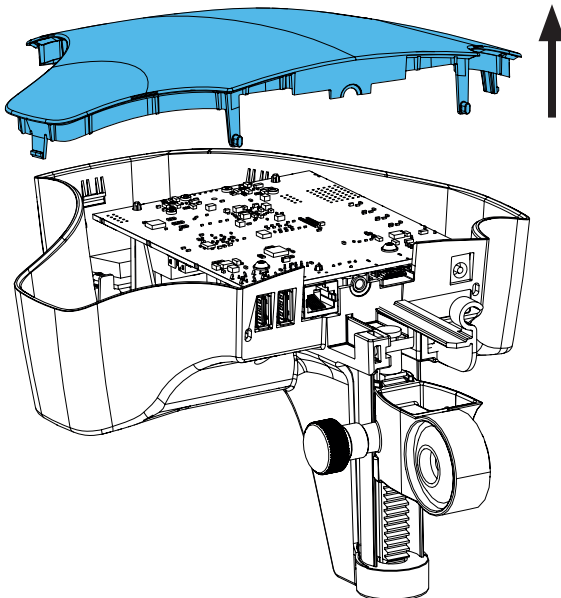


6.

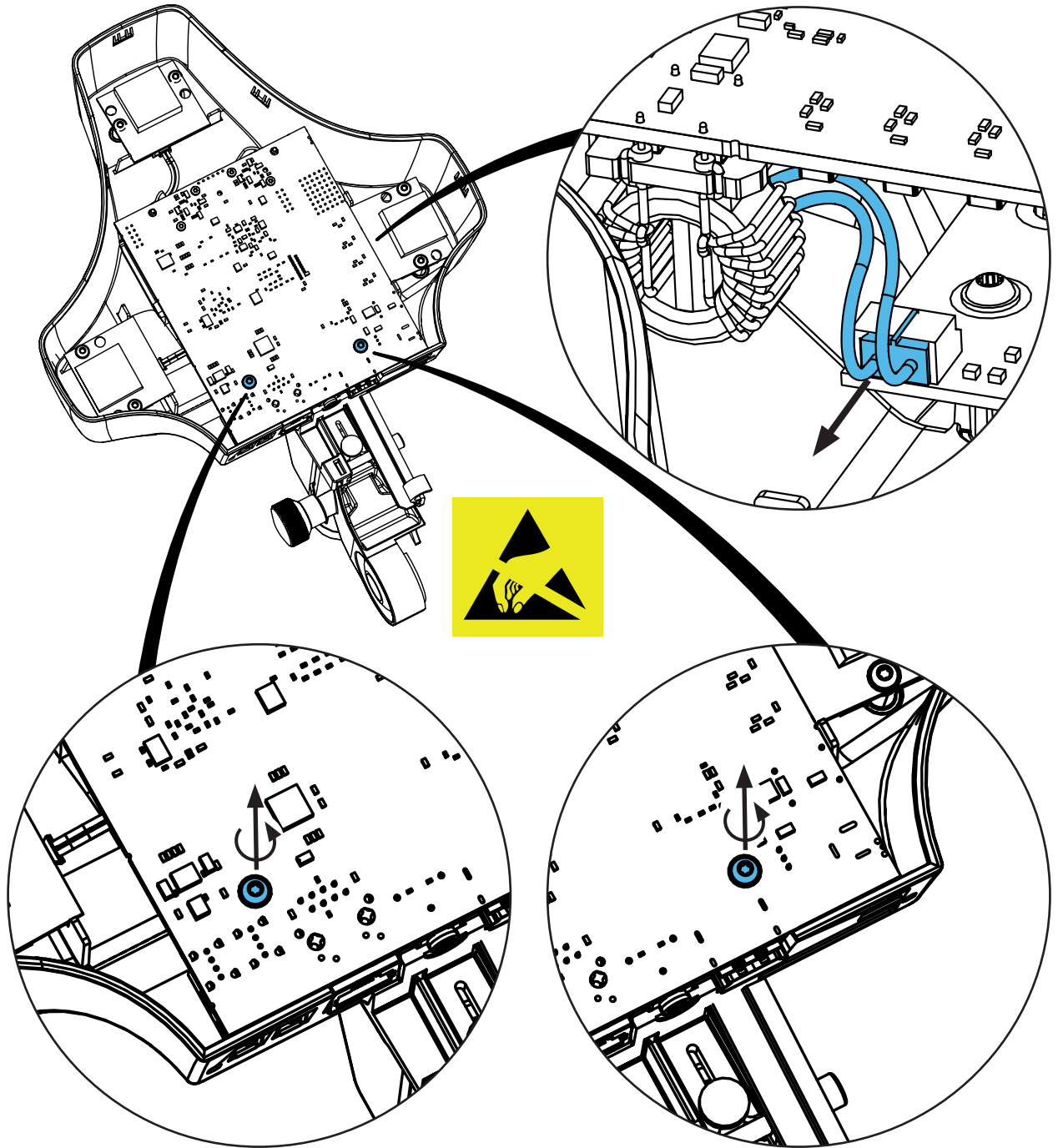


7.

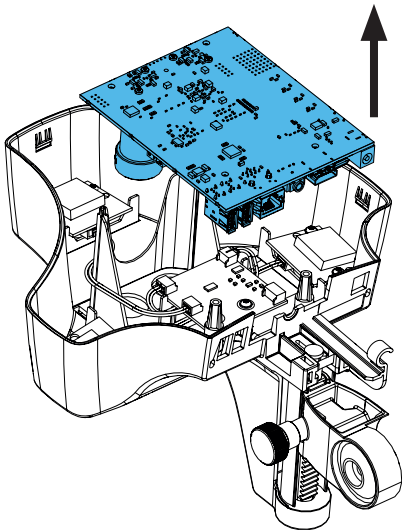


8.

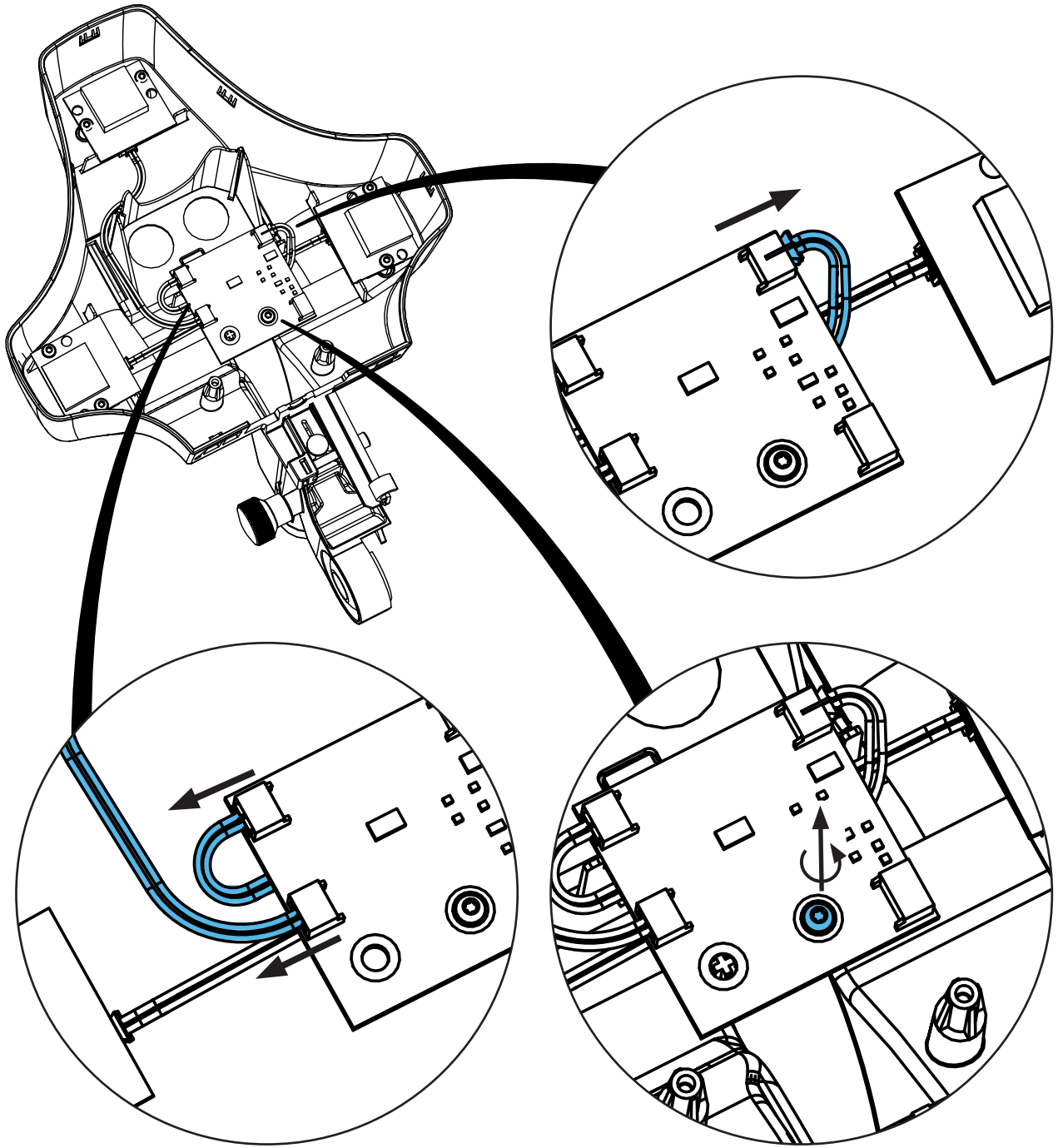


9.

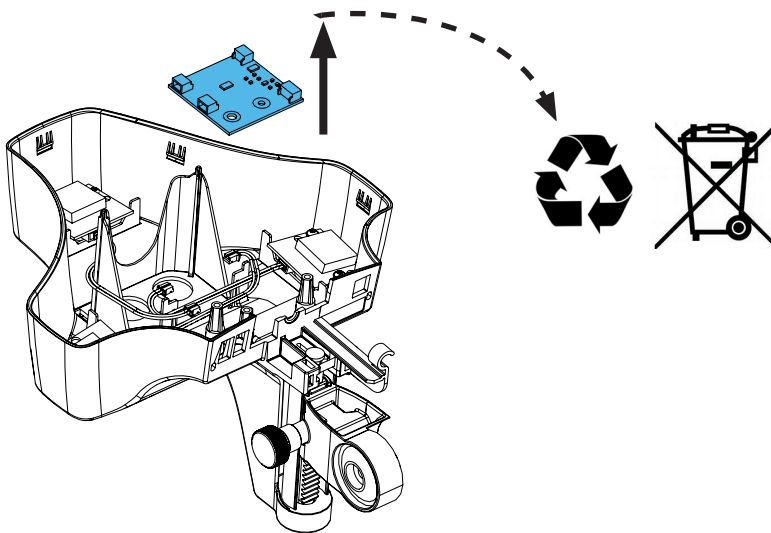



10.

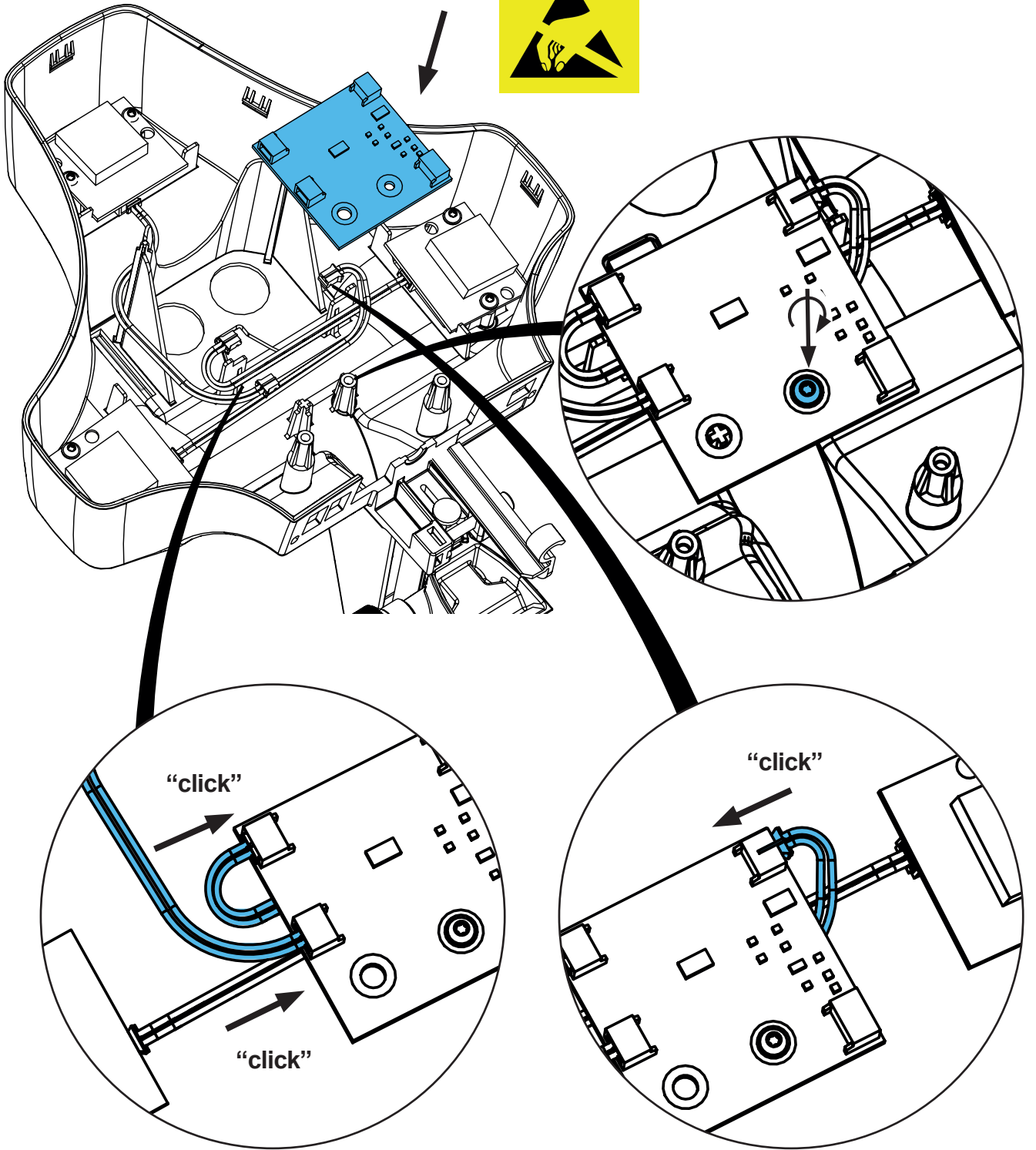


11.

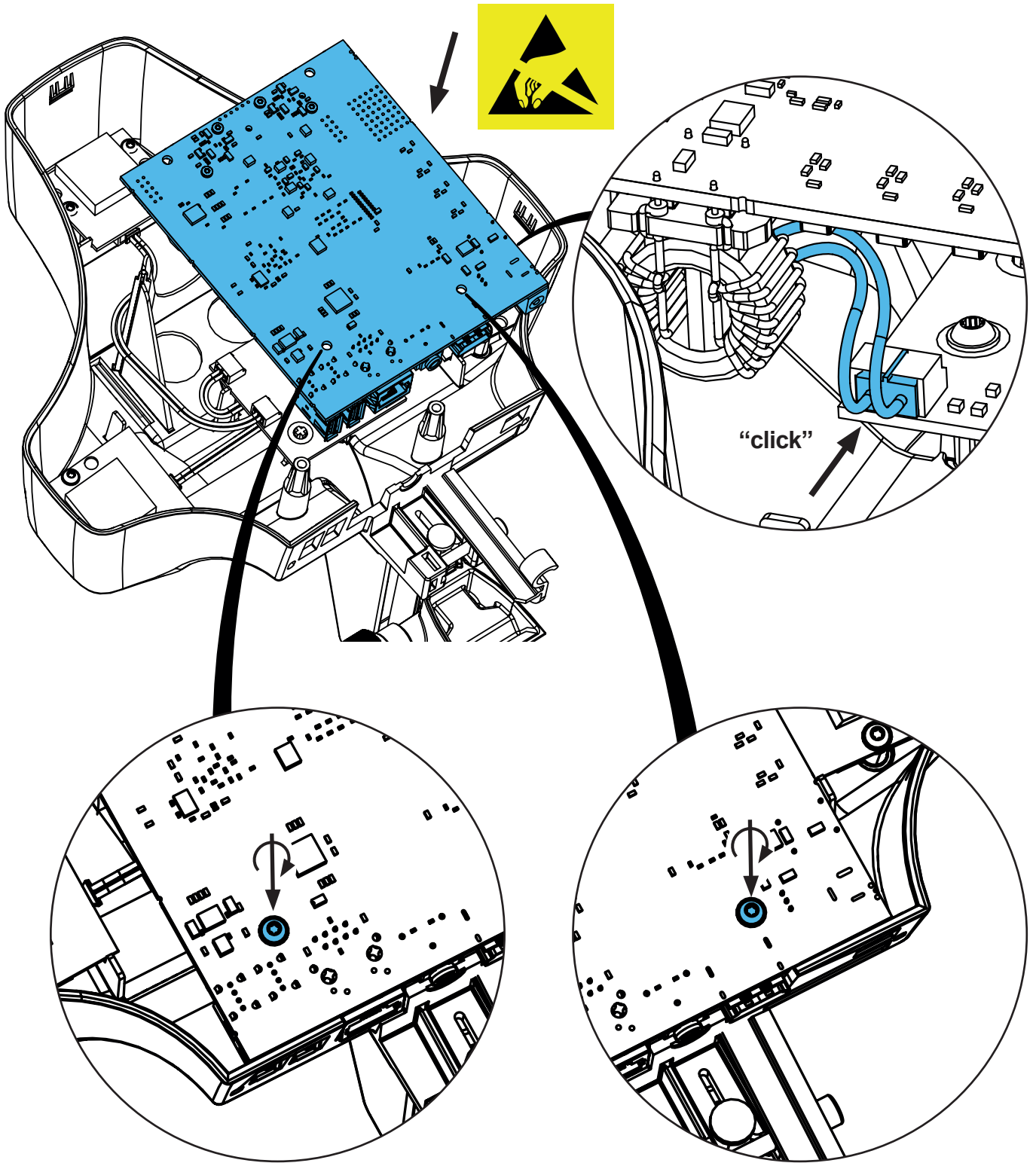

12.

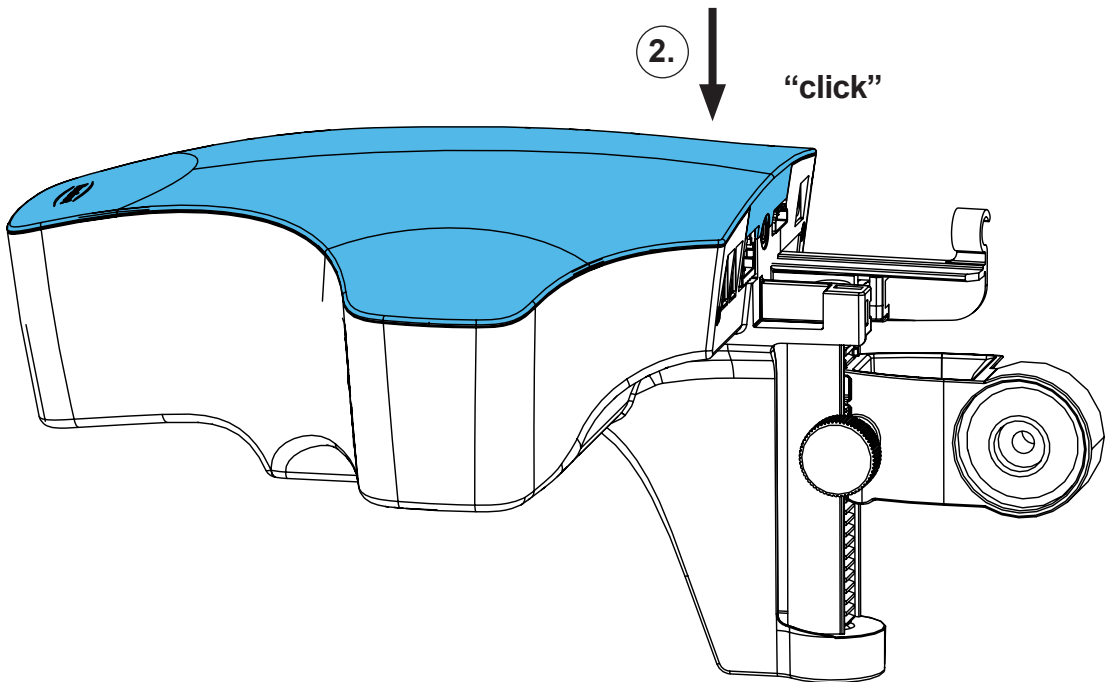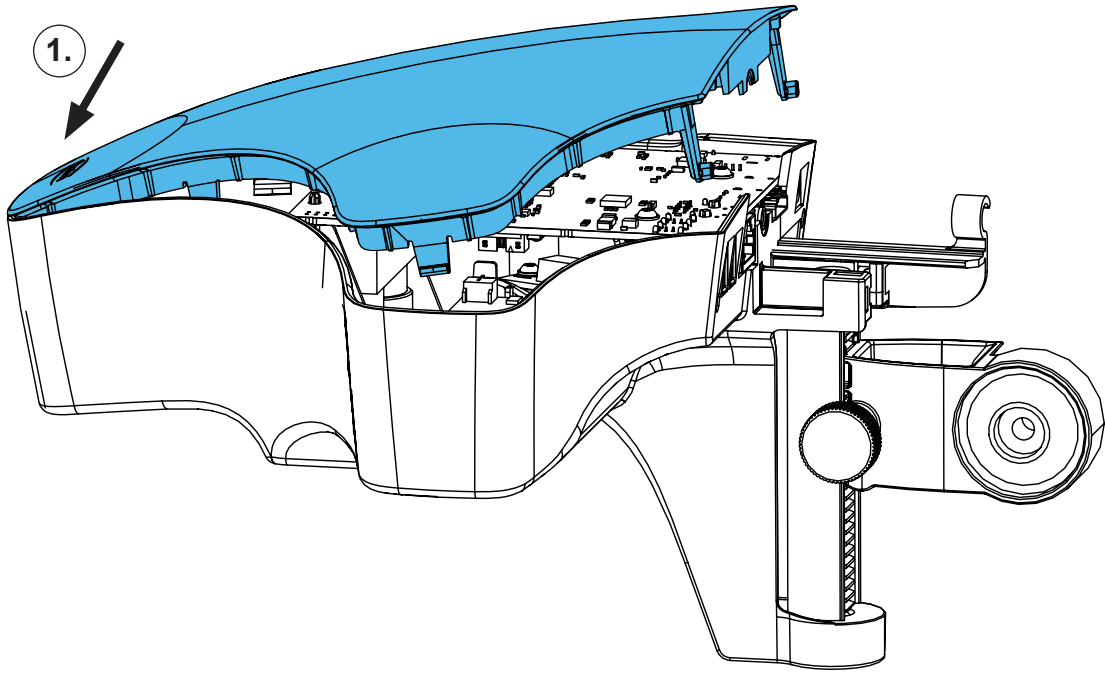

13.

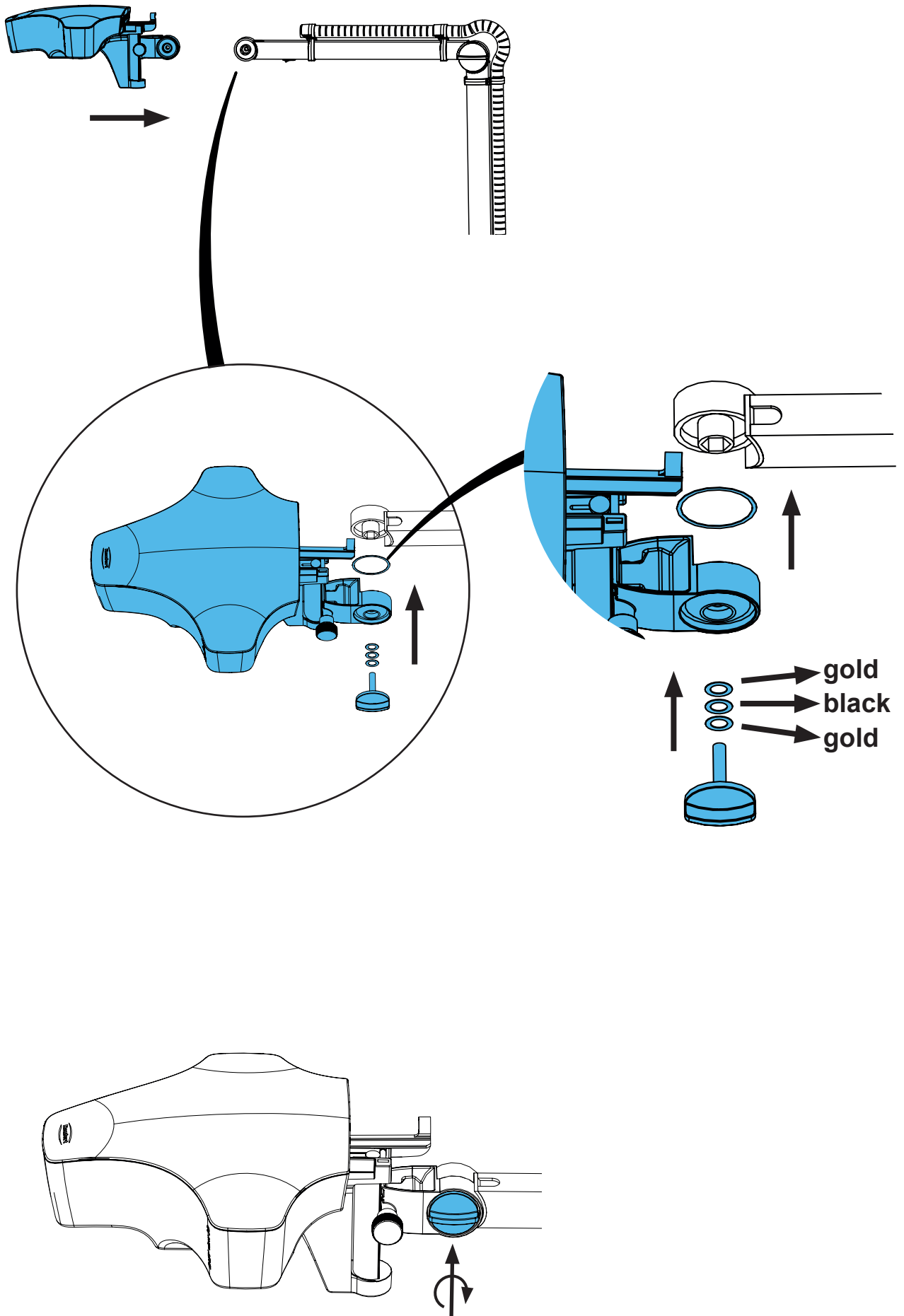

14.

